

Astell Autoclave Containers

Wire Baskets

Use these for vessels as they are the least expensive option. See picture in the brochure below.

Morrison Containers

Use these for discard runs where there is a possibility that the bags may burst or leak. Morrison containers are constructed of solid stainless steel with holes in the side to allow steam in (see picture below.) They have a solid base so that any spillages are contained. The solid base has a perforated false base above it so that if leaks occur, liquid will be taken away from the bag but contained below the false base. Bags can then be removed without exposure to the liquid.

New Swiftlock Autoclaves 90-330 Litres

Control System

The unique SECURE+ Micro-Computer controls the autoclave ensuring repeatable, safe and accurate sterilization time after time. A large LCD display gives an easy to read stage by stage report on cycle progress. Programs are factory set before delivery but re-programming by the user is extremely simple. Access to re-programming is PIN protected via the membrane touch pad and features simple question and answer stages.

Standard Selectable Parameters include:

- Delayed Start
- Timed Freesteaming Period (on non-vacuum units)
- Sterilization Time
- Sterilization Temperature
- Holdwarm On/Off
- Special Safety Valve Test Cycle
- Cooling Lock Release Temperature
- Profiled Overshoot Boost

Optional Selectable Parameters include:

- Assisted Air Cooling
- Load Sensed Process Timing
- Advanced Water Cooling
- Pulsar Free Steaming (on non-vacuum units)
- Pulsed Pre-Vacuum/Vacuum Drying (Front Loading Units Only)
- Air Ballast

New

Vacuum Swiftlocks are now available with the new

'LogiColor' Colour Touch Screen Controller.

Please see full details on page 8, and also ordering details covered in OPTIONS, pages 11-13.



Accessories

A wide selection of baskets, containers, load support plates and shelving systems are available for all machines giving optimum capacity for numerous applications.

ACCESSORIES		
90-330 Litre Autoclaves (dimensions in mm)		
AAN036	S/Steel Wire Basket (400 dia x 400h)	Top Loading Models
AAN042	S/Steel Wire Basket (400 dia x 230h)	Top Loading Models
AAN044	S/Steel Wire Basket (400 dia x 210h)	Top Loading Models
AAN048	Container with Perforated Base (380 dia x 400h)	Top Loading Models
AAN056	Morrison Container (390 dia x 500h)	Top Loading Models
AAN058	Morrison Container (390 dia x 355h)	Top Loading Models
AAN300	Rectangular Morrison (290 x 330 x 280)	Front Loading Models
AAN002	S/Steel Load Support Plate	Top Loading Models
AAN080	Container Tray (279 x 279 x 127)	Front Loading Models
AAN316	Additional Shelf Kit (2 Pos) 120 Litre	Front Loading Models
AAN317	Additional Shelf Kit (2 pos) 150 Litre	Front Loading Models
AAN318	Additional Shelf Kit (3 Pos) 230 Litre	Front Loading Models
AAN319	Additional Shelf Kit (3 Pos) 275 Litre	Front Loading Models
AAN320	Additional Shelf Kit (3 Pos) 330 Litre	Front Loading Models

FOR ALL OPTIONS SEE PAGES 11-13



MORRISON DISCARD CONTAINER



WIRE BASKET