


EggPunch

Pneumatic Egg-Opener

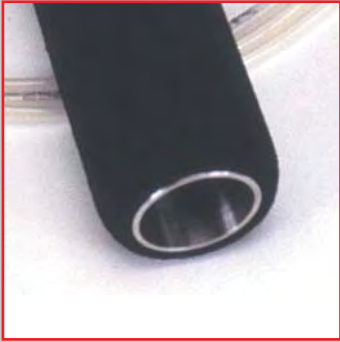

Quick & Clean

EggPunch **Pneumatic egg-opener**

- **opening of 60 embryonated chicken eggs per minute**
- **clean and circular fracture**
- **sterilizable**
- **handheld**

The EggPunch pneumatic egg-opener was designed for rapid and clean opening of embryonated chicken eggs by using air pressure.

For operation, simply hold the EggPunch in your hand and put your thumb on the trigger button. Pushing the button triggers the tool and produces a clean, circular fracture (Ø 25 mm) in the shell of embryonated chicken eggs.

Note:

In the moment when the circular punching blade meets the egg's shell, some particles of lime are released.

Convenient

The cap of the egg may be lifted off easily with a sharp knife.

The lower air sac membrane is exposed, intact and readily accessible.

With brief practice, you will be able to quickly and efficiently open 60 or more chicken eggs per minute.



Easy

The pneumatic egg-opener may be sterilized by suitable disinfectants or by heating in dry air.

It should not be totally immersed in liquids.

The EggPunch comes with approx. 1.8 metres of tubing to connect to a compressed air source.

The air pressure may range from 3 to 6 bar (50 to 100 psi).

The circular punching blade is subject to wear and has to be replaced by the user.

The safety nut securing the blade is a disposable suited for one-time use only.

Technical Data

Blade diameter	25 mm
Outer diameter	28 mm
Compressed air range	3 to 6 bar (50 to 100 psi)
Weight	approx. 1 kg

Ordering Information

EggPunch	Cat-No.
Pneumatic egg-opener	3.684 002

Spare parts

Punching blade including disposable safety nut	3.984 020
--	------------------

schuett **biotec**.de

Rudolf-Wissell-Straße 13, D-37079 Göttingen, Germany
Fon +49 (0) 551/5 04 10-0, Fax +49 (0) 551/5 04 10-99
info@schuett-biotec.de
www.schuett-biotec.de