


## MINI BUT MIGHTY LAB INSTRUMENTS

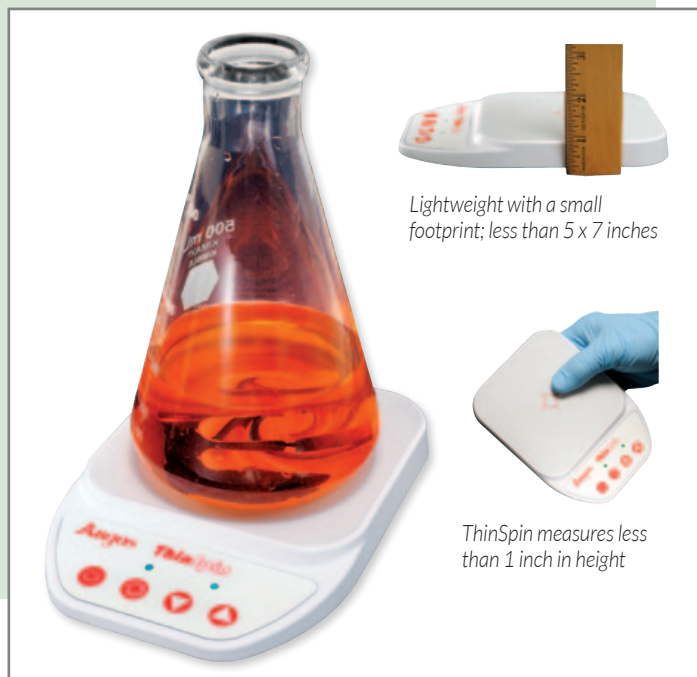

Lightweight with a small footprint; less than 5 x 7 inches

ThinSpin measures less than 1 inch in height

### ThinSpin™ Magnetic Stirrer

The Argos ThinSpin™ Magnetic Stirrer is an ultra-thin, light-weight piece of equipment that can be programmed to mix and stir between 15 and 1500 rpm. The preset speed is 15 rpm, but speed can be easily adjusted in increments of 15 rpm. Each time the speed is changed the handy LED light will flash. For added versatility the spin direction can be alternated clockwise or counter clockwise by pressing the reverse key.

The low-profile and sleek design of the ThinSpin enables it to be used and stored in tight spaces such as in fume hoods or in cabinets. The unit weighs only 11 oz., and has a tiny footprint of 175 x 125 x 20.5 mm. The working platform has dimensions of 110 x 110 mm. The maximum load is 2.0 kg.

This small but mighty instrument can mix and stir up to 500 mL of laboratory liquids, solutions and media. The top plate of the ThinSpin is water-proof and resistant to most laboratory chemicals.

The ThinSpin is CE marked, and has an IP rating of DIN EN 60529. Unit includes a 2 year warranty.

### Ordering Information

Cat. No.	Description	Qty.
S2001	ThinSpin magnetic stirrer (220/240V)	1/each

### Vortamix™ Mini Vortexer

The Argos Vortamix Mini Vortexer is an unbelievably compact, yet powerful instrument engineered to mix tubes quickly and easily. The vortexing motion is activated by placing a tube on the gasket cap and then pressing down lightly. The innovative design of the gasket cap allows centrifuge and test tubes up to 30 mm in diameter to be mixed with very little effort – including 50mL conical tubes! The standard unit comes with a red rubber gasket cap – blue, green and purple colored caps are offered separately.

Despite its small size and weight (dimensions: 95 x 85 mm/weight: .55 kg), the Vortamix offers an excellent mixing speed of 3,000 rpm. The orbital diameter is 4.5mm. The ambient temperature range of the unit is 5 - 40°C.

The Vortamix is CE marked and has an IP rating of DIN EN 60529. Unit includes a 2 year warranty.

### Ordering Information

Cat. No.	Description	Qty.
V0001	Gasket cap, red	1/each
V0002	Gasket cap, blue	1/each
V0003	Gasket cap, green	1/each
V0004	Gasket cap, purple	1/each



Use with a variety of tube sizes

Colored gasket caps available

For tubes from 0.5 - 50mL



**Wolflabs**

# Wolf Laboratories Limited

[www.wolflabs.co.uk](http://www.wolflabs.co.uk)

Tel: 01759 301142

Fax: 01759 301143

[sales@wolflabs.co.uk](mailto:sales@wolflabs.co.uk)



**Use the above details to contact us if this literature doesn't answer all your questions.**

**Pricing on any accessories shown can be found by keying the part number into the search box on our website.**

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

