ThinSpin Magnetic Stirrer

Operator's Manual
Manuel de l'opérateur
Bedienungshandbuch
Manual del operador
Manuale d'uso
オペレーターマニュアル

ARGOS
TECHNOLOGIES *manufacturing solutions for science*

Argos Technologies, Inc.
customer-service@argos-tech.com
Toll Free: 800-886-8675
International: 847-622-0456
www.argos-tech.com

CE **RoHS**
2011/65/EU

Introduction

Thank you for your purchase of the Argos Technologies ThinSpin Magnetic Stirrer. This unit was designed for mixing and stirring laboratory liquids, solutions and media up to 500mL. The small footprint requires little laboratory space, while the low profile allows for easy storage between use.

Package Contents

- ThinSpin Magnetic Stirrer
- Pin Jack Power Supply (US or International)
- Instruction Manual
- Warranty Card

Materials & Specifications

Model	S2000 (US), S2001 (International)
Speed	15 – 1500 rpm
Capacity	500mL Water
Working Platform	110 x 110 mm
Load	Max. 2.0 kg
Ambient	5-40°C, ≤ 80% RH
Weight	0.3 kg
Power Supply	AC100-240V, 50/60Hz
Power	6W
Dimensions (DxWxH)	175 x 125 x 20.5 mm
IP Rating (DIN EN 60529)	IP 45
Certifications	CE Marked

This device complies with RoHS 2011/65/EU

Maintenance, Care and Safety Precautions

CAUTION


The ThinSpin Magnetic Stirrer contains magnets and may generate a magnetic field. Use caution around data storage material (including credit cards), cardiac pacemakers or any other device that is sensitive to magnetic fields.

Safety Precautions

- When mixing or stirring liquids or solutions, wear suitable personal protection to prevent injury from splashing of liquids, hazardous, toxic fumes or gases, or the ejection of spin or stir bars.
- Use extra caution when mixing:
 - Combustible liquids
 - Flammable materials
- Only use the power supply adapter supplied with the unit.
- Do not overfill mixing vessel to avoid splashing while stirring.
- Place the mixing container on the center of the top plate.
- With prolonged use, the unit may heat up.
- Discontinue use or reduce speed when:
 - The mixing container moves or vibrates on the top plate.
 - The unit emits a loud, audible noise during operation.
 - Liquid splashes out of the mixing container.
- Clean the top plate and stir bar after each use.

Cleaning & Care

- Prior to cleaning, disconnect the power supply from the outlet.
- Clean the unit with a soft cloth and a non-corrosive liquid. Isopropyl Alcohol can be used.
- Do not allow any liquid to get into the ThinSpin Magnetic Stirrer during cleaning.
- Do not submerge unit into water!
- Ensure the unit is completely dry prior to use.
- There are no serviceable parts inside the ThinSpin Magnetic Stirrer.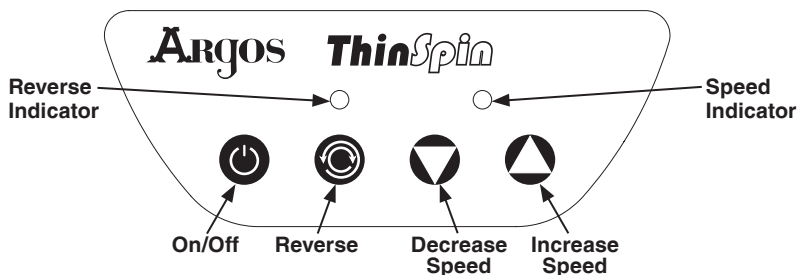

Operating Instructions

- Place the ThinSpin Magnetic Stirrer on a flat, stable surface.
- Connect the pin jack adapter to the ThinSpin base unit prior to inserting the power cord into an electrical socket.
- Place the stir-bar into the mixing container first.
- Fill the mixing container with the liquid to be mixed.
- Place the mixing container in the center of the ThinSpin Magnetic Stirrer top plate.
- Turn the unit “ON” with the On/Off key.
- The initial preset speed is 15 rpm. Each time the up or down key is pressed, the speed is increased or decreased by 15 rpm. When changing speeds, the LED light will flash.
- The speed setting will remain programmed into the unit when the unit is turned off or unplugged.
- Spin directions can be alternated clockwise and counter clockwise by depressing the reverse key. When selected, the reverse LED light will turn on and the spin-bar will reverse direction every 30 seconds. The spin-bar will pause 5 seconds between each change in direction.

Warranty Information

ThinSpin Magnetic Stirrer 2 Year Warranty

Argos ThinSpin Magnetic Stirrer is warranted to be free of defects in materials and workmanship for a period of 24 months from the date of purchase, providing that:

- The unit has not been opened or attempted to be opened by the end user
- The user ensures that it has been carefully cleaned and decontaminated
- The unit has been operated according to the instructions and not in any improper way.

In order to qualify for warranty repair; the customer must return the warranty card supplied with the unit that includes all of the information required to validate proof of purchase. Should warranty service be required, repair or replacement will be provided by Argos without charge. Freight shipping the unit to Argos will be the responsibility of the purchaser. There are no other warranties or guarantees, implied or otherwise expressed including any warranties of application or merchantability.

Repair Service


Should your ThinSpin Magnetic Stirrer require service or repair, please contact Argos Technologies customer service. They will issue a Return Authorization Number and advise on the proper packaging and method of shipment.

Customer Service: in US: 800-886-8675 outside US: 847-622-0456

**Argos Technologies, Inc.
1551 South Scottsdale Court
Suite 200
Elgin, IL 60123
Attn: Technical Service**

Be sure to list your RA number on the shipping label.

Please use the same numbers listed above, should you require any technical assistance or support.

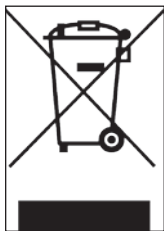This equipment is labeled with the crossed out wheeled bin symbol to indicate that the equipment must not be disposed of with unsorted waste. It is your responsibility to correctly dispose of the equipment by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical or radioactive contamination. So as to protect from health hazards the persons involved in the disposal or recycling of products.

For more information about where you can drop off your waste equipment, please contact your local dealer from whom you originally purchased the product.

By doing so, you will help to conserve our natural resources and you will ensure that your waste products are recycled in a manner that protects human health and protects the environment.

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