Bibbyjet Pro

(Pipettierhelfer/Pipette Controller)

Bibby Scientific Ltd

Beacon Road
Stone
Staffordshire
ST15 0SA
United Kingdom
E-Mail: info@bibby-scientific.com
Internet: www.bibby-scientific.com

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Operating Manual

Please read carefully before use,
and follow all operating and safety
instructions!

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Please read the following carefully!

Safety Instructions

This instrument may sometimes be used with hazardous materials, operations, and equipment. It is beyond the scope of this manual to address all of the potential safety risks associated with its use in such applications. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Every user must read and understand this operating manual before operation.
2. Follow general instructions for hazard prevention and safety instructions; e.g., wear protective clothing, eye protection and gloves.
3. Observe all specifications provided by reagent manufacturers.
4. Never use the instrument in an atmosphere with a danger of explosion. Highly flammable liquids must not be pipetted.
5. Use the instrument only for pipetting liquids, with strict regard to the defined limitations of use. Observe operating exclusions (s. page 7)!

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Warning!

Improper use of the instrument or the battery (short-circuit, mechanical damage, overheating, etc.) may cause the explosion of the battery.

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+ 10 °C bis + 40 °C

Application

The instrument is designed to assist the filling and dispensing of graduated and volumetric pipettes of glass or plastic in the volume range of 0.1 ml to 200 ml with a suction tube outer diameter < 9,2 mm for measuring liquids. If the instrument is used correctly, the pipetted liquid will only contact the pipette.

Working- and recharging temperature:

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Limitations of use

The instrument is designed for pipetting liquids, observing the following physical limits:

- +10 °C to +40 °C (50 °F to 104 °F)
(of instrument and reagent)
- Vapor pressure up to max. 500 mbar.
Aspirate slowly above 300 mbar, in order to prevent the liquid from boiling.
- Density up to 9.0 g/cm³

Operating exclusions

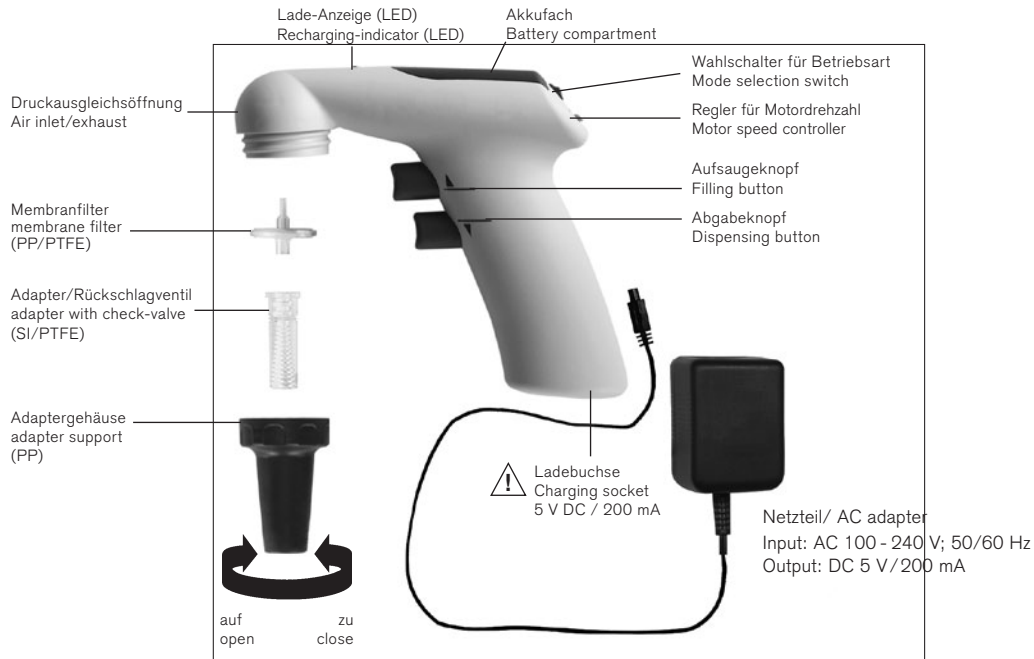
Never use the instrument with liquids whose vapors attack the materials silicone or EPDM.

The instrument is not designed for Pasteur pipettes.

Warning!


Never use or recharge the instrument in an atmosphere with danger of explosion. Highly flammable liquids (e. g., ether, acetone and other liquids with a flash point below 0 °C) must not be pipetted.

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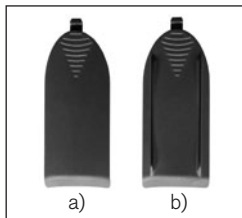Items supplied

- Pipette Controller
- Nickel-metal hydride battery
- 2 battery compartment covers
 - a) smoothly rounded
 - b) with ridges to place the instrument inverted on the bench
- AC adapter
- 2 spare membrane filters, 0.2 µm
- wall support
- Operating Manual

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Battery Pack, NiMH 2.4 V/730 mAh

Storage conditions

Store instrument and accessories in a cool, dry place.

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Getting started

1. Grip battery plug by the wire, confirm alignment of the plug and insert it in the socket. Push it firmly into the socket until you can hear it snap into place. Place battery in compartment.

2. Select battery compartment cover:
 - a) smoothly rounded
 - b) with ridges for resting the instrument inverted on the bench (> page 14)

3. Close battery compartment.
4. Recharge the battery (charging time is 4 hours).

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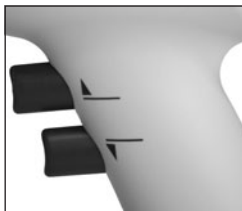
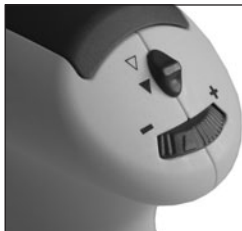
▼ **Blow-out**
(with motor power)

Adjust Motor speed

Maximum motor speed is continuously adjustable from maximum (+) to minimum (-) by turning the thumbwheel. We recommend using the maximum speed setting for larger volume pipettes and a slower speed setting for lower volume pipettes.

Pipetting speed control

The amount of pressure on the pipetting buttons provides continuously variable speeds up to the rate set by the speed controller. Press upper pipetting button for filling and lower button for dispensing.

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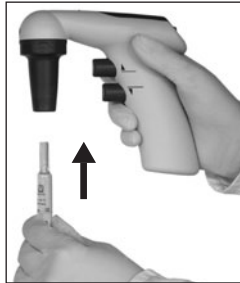
Pipetting

Attach the pipette

Hold the pipette as near to its upper end as possible, and carefully insert it into the adapter until it fits tightly.

Warning!

Be sure that the pipette fits tightly into the adapter. Never use force. Thin pipets are particularly liable to break. Avoid the risk of injury!

Once the pipette has been securely attached, always hold the instrument in a vertical position, tip down.

Filling

Immerse the pipette tip into the liquid. Slowly depress the upper pipetting button and fill the pipette so that the meniscus is slightly above the mark desired.

Attention!

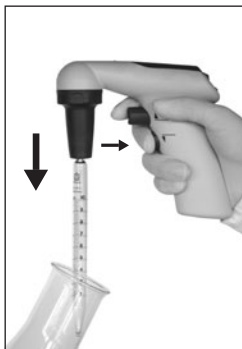
Take care to avoid overfilling of the pipette.

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Adjust the volume

Use suitable lint-free tissue to wipe the pipette tip. Slowly press the lower button and dispense liquid until the meniscus is adjusted exactly to the desired volume.

Dispensing

Slowly press the lower pipetting button. The dispensing speed is dependent on the amount of pressure on the pipetting button, as well as the setting of the speed controller, when the mode switch is in the motor powered delivery position.

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Wall support/ Inverted bench placement

Place the instrument in the wall support for storage.

Mounting the wall support

The wall support is attached with snap tapes to permit easy removal for cleaning purposes. To use the snap tapes, the mounting surface must be smooth, clean and grease-free. Prepare the surface for mounting by cleaning with a 50% solution of isopropanol/water. Alternatively, the wall support can be attached to the wall with two screws.

If the cover with ridges is mounted, instrument can be placed inverted on the bench.

Attention!

Never place the instrument inverted on the bench with a pipette inserted.

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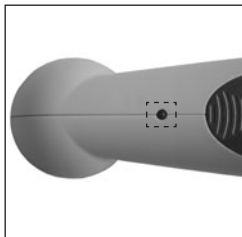
Recharging the battery

Warning!

Use original manufacturer's AC adapter only!
Using a different AC adapter can damage both the instrument and the AC adapter, and will void the warranty.

Confirm the voltage of the charger matches the voltage of your electrical outlet.

The power supplied to the AC adapter should differ by no more than $\pm 10\%$.

One full charge of the nickel-metal hydride battery allows 8 hours of non-stop pipetting. A flashing LED indicates the need for recharging.

Recharging indicator

LED off:

- battery charge is sufficient

LED flashes slowly:

- remaining battery charge is sufficient for a maximum of two hours of pipetting; charging of battery is recommended

LED flashes quickly:

- finish pipetting activity and charge battery

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Recharging procedure

1. Temperature range for recharging:
+ 10 °C to + 35 °C
2. When the AC adapter is initially connected, the LED on the instrument lights continuously.
3. After 4 hours of charging, the AC adapter switches to a pulsed, long-term charging mode (indicated by the LED flashing every 15 seconds).

Note:

- The charging system is designed to prevent the battery from overcharging and minimizes the lazy-battery-effect. To maintain maximum battery capacity, it is best to charge the battery only when the slowly-flashing LED indicates charging is needed.
- The instrument can be used during recharging.
- If the LED does not light up during charging see chapter 'Troubleshooting' (➤ page 25).

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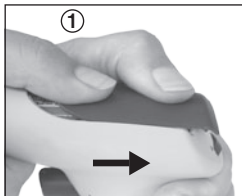
Replacing the battery

Warning!

Only use original manufacturer's accessories (➤ page 29). The battery must not be replaced by non-rechargeable or rechargeable batteries of other manufacturers.

Use of the wrong batteries or improper use of the instrument (e.g., short-circuit, mechanical damage, overheating, etc.) may cause the batteries to explode.

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- To open the battery compartment press at the embossed arrow, position ①, and simultaneously slide cover back.

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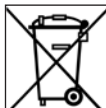
- Remove the battery.
- Grip battery plug by the wire and pull it gently out of the socket.
- Grip the plug of the new battery by the wire and push it firmly into the socket until you can hear it is snapping in.
- Insert new battery and close battery compartment.

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Disposal

The adjoining symbol means that storage batteries and electronic devices must be disposed of separately from household trash (mixed municipal waste) at the end of their service life.

- According to the Directive 2002/96/EC of the European Parliament and of the Council on Waste Electrical and Electronic Equipment (WEEE) of 27 January 2003, electronic equipment requires disposal according to the relevant national disposal regulations.
- Batteries contain substances that can have harmful effects on the environment and human health. Therefore according to the Directive 2006/66/EC of the European Parliament and the Council on Waste Batteries of 6 September 2006 batteries require disposal according to the relevant national disposal regulations. Dispose of batteries only when completely discharged.

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Warning!

Do not short-circuit the battery to discharge it!

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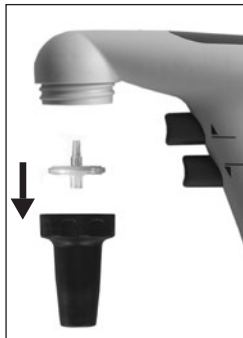
Cleaning/Autoclaving

Occasionally wipe the housing with a damp cloth. When properly used, the instrument is maintenance-free.

Membrane filter, adapter and adapter support are autoclavable at 121 °C (250 °F), 2 bar absolute (30 psi) with a holding time of at least 15 minutes according to DIN EN 285.

Replacing the filter and cleaning the adapter

Only necessary if a pipette has been overfilled or the suction capacity is impaired.

Warning!

Wear protective gloves and eye protection. Avoid the risk of contamination!

1. Pull out the pipette.
2. Unscrew the adapter support.
3. Pull out the membrane filter.

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4. Remove the adapter from the adapter support by pushing it upwards. The check valve remains in the adapter!

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5. To clean adapter with built-in check valve, carefully rinse it by using an appropriate compatible solvent (e.g., water) in a wash bottle. Blow out the liquid and allow to dry completely.

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6. Place the new membrane filter with the thicker end facing downwards into the adapter (with the smaller conical side up).

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7. The entire pipette adapter incl. membrane filter is autoclavable at 121°C.

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Note:

Not every membrane filter is suitable.

Only use genuine manufacturer's recommended accessories.

Each instrument will be supplied with a 0.2 µm membrane filter (red color code). The membrane filter is autoclavable up to 5 times at 121 °C. It is the user's responsibility to ensure effective autoclaving.

The filter with 0.2 µm pore size (for cell culture) is sufficient to meet class B tolerances.

For class A accuracy, it is necessary to use a 3 µm membrane filter. This filter has a larger pore size and thus permits better airflow for gravity-delivery.

(Spare parts ➤ page 29).

8. Reassemble the instrument in reverse order and carry out leak test.

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(> Seite 24)

Leak Test

When the meniscus has been set, liquid should not drip out of the pipette until the valves are activated.

Should the pipette drip, see chapter 'Troubleshooting'

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Troubleshooting

Trouble	Possible Cause	Action to be taken
Suction capacity impaired	<ul style="list-style-type: none"> - Motor speed is reduced - Check valve closed - Filter or check valve dirty 	<ul style="list-style-type: none"> - Increase motor speed - Set mode selection switch for "power delivery" and press lower button for blow out (➤ p. 13) - Clean and/or replace check valve (➤ p. 20-22)
Pipette drips	<ul style="list-style-type: none"> - Filter not properly positioned - Adapter or pipette damaged 	<ul style="list-style-type: none"> - Insert filter properly (➤ p. 21) - Replace adapter or pipette (➤ p. 17-18)
Pipette not held properly	<ul style="list-style-type: none"> - Adapter dirty or damaged 	<ul style="list-style-type: none"> - Clean the adapter, allow it to dry and/or replace it (➤ p. 20-23)
After connecting the AC adapter, the LED-display of the instrument flashes or does <u>not</u> light up	<ul style="list-style-type: none"> - Plug of the battery not firmly pushed in - Wrong AC adapter - Battery/ AC adapter defective - Instrument defective 	<ul style="list-style-type: none"> - Firmly push in the battery plug (➤ p. 10) - Only use original AC adapter - Replace battery/AC adapter - Send the instrument in for repair (➤ p. 26)
Low battery charge	<ul style="list-style-type: none"> - Battery problem 	<ul style="list-style-type: none"> - Replace battery (➤ p. 17-18)
Instrument doesn't work	<ul style="list-style-type: none"> - Internal error (carry out reset) 	<ul style="list-style-type: none"> - Pull out battery plug and press a pipetting button for 10 seconds. Then firmly push in the battery plug again. (➤ p. 10)

If recommendations in the Troubleshooting section do not solve current problems, contact your authorized dealer or the manufacturer directly.

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Return for Repair

Attention!

For safety reasons, only clean/decontaminated instruments can be checked/repared.

- Therefore: Clean and decontaminate the instrument carefully.
- Complete the “Declaration on Absence of Health Hazards” (ask your supplier or manufacturer for forms) and send the instrument to the manufacturer or supplier and describe:
 - the nature of the problem
 - the liquid used.

Shipment is at the risk and the cost of the sender.

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Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operation or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual. We are not liable for damage resulting from any actions not described in the operating manual or if non-original parts have been used.

Subject to technical modification without notice. Errors excepted.

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Ordering Data

Bibbyjet Pro - pipette controller with AC adapter,

incl. battery, 2 battery compartment covers, 2 spare
membrane filters (0.2 µm), wall support and
operating manual

with AC adapter for	Cat.No.
UK/Ireland 230 V/50 Hz	PC 2000
Europa (Continental) 230 V/50 Hz	PC 2001
Without AC adapter	PC 2002

Adapter support, battery compartment cover, pipetting keys, mode selection switch and controls are royal blue, instrument body is grey.

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Akku-Pack NIMH
2,4 V / 700 mAh **Best.-Nr.
PC 2000/B**

Accessories and spare parts

**Silicone adapter
with non-return valve** **Cat. No. PC 1011**

Spare membrane filter 0.2 µm, sterile
(red color code)
1 pc. in blister pack **Cat. No. PC 1012**

Spare membrane filter 0.2 µm, non-sterile
10 pc. in PE-bag **Cat. No. PC 1013**

Spare membrane filter 3 µm, non-sterile
1 pc. in blister pack **Cat. No. PC 1014**

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This sign certifies that the product meets the requirements of the EC directive and has been tested according the specified test methods.