

Instruction Manual

CertoClav Vacuum Pro 45

Steam Steriliser Model STE-45-T



Dear customer,

Thank you for the trust you have placed in us by purchasing this autoclave.

CertoClav, a medium-sized family company, has specialised in the manufacture of table-top autoclaves for more than 60 years.

This autoclave has been manufactured and tested according to strict requirements. However, please read through this instruction manual carefully.

The long-lasting functionality and value retention of your autoclave depend primarily on the maintenance of the device.

Instruction Manual CertoClav Vacuum Pro 45

CertoClav

Valid for model STE-45-T

Issue 1 June 2019

Responsible for content:

CERTOCLAV Sterilizer GmbH

Peintner Straße 10

A-4060 Leonding

Austria

Email: support@certoclav.com

www.certoclav.com

© 2019 CertoClav

Last change: 11.06.2019

Document: BA_VACPRO45_D.doc / Revision: 8

Subject to technical changes

Symbolic images: Images may differ from the original

CERTOCLAV

Since 60 years.


Table of Contents

1	General Information	6
1.1	Intended use	6
1.2	Environmental protection	6
1.3	Customer Service	6
1.4	Ordering spare parts and special accessories	7
1.5	Guarantee	7
2	Meaning of the Symbols on the Autoclave	8
3	Safety Information	9
4	Non-autoclavable Substances	10
5	Initial Start-up	11
5.1	Device description	11
5.2	Device dimensions	11
5.3	Scope of supply	12
5.4	Before initial start-up	12
6	Installation	13
6.1	Setting up the autoclave	13
6.2	Unlocking the autoclave door	14
6.3	Touchscreen assembly	14
6.4	Installing accessories	16
6.4.1	Connecting the protocol printer	16
6.4.2	Connecting the barcode reader	16
6.4.3	Connecting the label printer	16
6.4.4	Connecting the CertoServer	17
6.5	Pressure sensor calibration	17
6.6	Validation and calibration	18
7	Sterilisation Programs	19
7.1	Default settings	19
7.2	Editing program parameters	20
8	Basic Operation	21
8.1	Switching on the touchscreen	21
8.2	Switching off the touchscreen	21
8.3	Standby Mode for the touchscreen	21
9	Instructions for Daily Use	22
9.1	Preparations before use	22
9.2	Replenishing with distilled water	22
9.3	Opening the autoclave door	23

9.4	Steps to start a sterilisation process	24
9.4.1	Selecting the sterilisation program on the screen	24
9.4.2	CertoTrace – Documentation and load management	25
9.4.3	Step-by-step video assistant	26
9.4.4	Loading the autoclave	26
9.4.5	Closing the autoclave door	26
9.4.6	Material tests and automatic repetitions	27
9.4.7	Status screen	27
9.5	Unloading the autoclave	28
9.6	Emptying the used water tank	29
9.7	PC software	29
9.7.1	Functions of the CertoCloud	30
9.7.2	Display of sterilisation protocols via web browser	31
10	Control Panel of the Autoclave	32
10.1	Login Menu	32
10.2	Support Menu	34
10.2.1	Tutorial videos	34
10.2.2	Send email to CertoSupport	35
10.2.3	TeamViewer, remote support and share screen	36
10.3	Protocols Menu	38
10.3.1	Viewing protocols	38
10.3.2	Printing protocols	38
10.3.3	Example of a protocol printout	39
10.3.4	Print barcode	39
10.3.5	Tracing marked products	40
11	Settings	41
11.1	User administration	41
11.2	Network settings	41
11.2.1	Connecting to a WLAN network	41
11.2.2	Connecting to a LAN network	42
11.2.3	Network communication	42
11.2.4	Connecting to the CertoServer	43
11.2.5	Connecting to the CertoCloud server	44
11.3	Sterilisation settings	45
11.3.1	Customising sterilisation programs	45
11.3.2	Material tests and automatic repetitions	45

11.3.3	CertoTrace — Your support during material tests	47
11.3.4	Label designer	47
11.3.5	Real-time protocolling to USB	48
11.3.6	Exporting all protocols to a USB	48
11.3.7	Exporting all protocols to an SD card	49
11.3.8	Exporting all protocols via a network	50
11.4	System settings	51
11.4.1	Changing the date and time	51
11.4.2	Installing updates via the Internet	51
11.4.3	Installing updates via USB	52
11.4.4	Installing updates via SD card	53
11.4.5	Restoring the default settings	53
11.4.6	Changing the language	54
11.5	Configuration of SMS and email notifications	54
11.6	Sensor calibration	55
11.7	GLP and GMP data	55
11.7.1	Header	55
11.7.2	Autoclave designation	55
11.7.3	Project designation	55
11.7.4	Advanced GLP and GMP settings	56
11.8	Locking settings	56
12	Error Messages	57
13	Maintenance	59
14	Transport and Storage	60
14.1	Preparations	60
14.2	Emptying the water containers	60
14.3	Transport conditions	60
14.4	Storage conditions	60
15	Technical Data	61
16	Appendix	62
16.1	Appendix 1: Structural diagram	62
16.2	Appendix 2: Circuit diagram	63
16.3	Appendix 3: CE Declaration of Conformity for Autoclave	63
16.4	Appendix 4: CE Declaration of Conformity for Safety Valve	65

1 General Information

	<p>Dear customer,</p> <p>Before putting your autoclave into operation for the first time, please read through this instruction manual and follow its information and instructions. Do not use another instruction manual.</p> <p>This instruction manual applies only to the CertoClav Vacuum Pro Series autoclaves with model designation STE-45-T.</p> <p>Keep this instruction manual within easy reach for later use or for subsequent owners.</p>
---	--


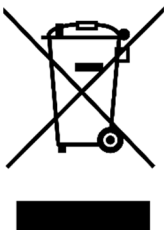
1.1 Intended use

This autoclave is intended for laboratory use. For example for the sterilisation of liquids, solids, waste and textiles as well as swabs, glasses, petri dishes, porous materials, powder, hollow bodies and instruments. Containers sealed with compressed air must not be sterilised. This autoclave has not been developed for private use.

Operation and maintenance of the autoclave may only be performed by trained personnel.

Comply consistently with all safety regulations in this instruction manual.

1.2 Environmental protection

	<p>The packaging materials are recyclable. Please do not dispose of the packaging in domestic waste but instead hand it over for recycling.</p>
	<p>Old devices contain valuable recyclable materials, which should undergo recycling. Please therefore dispose of old devices via suitable collection systems.</p>

1.3 Customer Service

If you have any questions or problems regarding the autoclave, please contact your dealer or the manufacturer by email to support@certoclav.com.

1.4 Ordering spare parts and special accessories



Use original accessories only and original spare parts for maintenance, which are approved by the manufacturer, so as not to impair the safety of the device.

You can obtain spare parts and accessories from your dealer or from the manufacturer.

1.5 Guarantee

The guarantee is valid for 2 years. The claim to free service shall be nullified if one of the following conditions is not met:

- The damage has been caused by incorrect installation.
- The damage has been caused by negligent conduct.
- The damage has been caused by improper operation.
- The damage has been caused by force majeure such as fire.

2 Meaning of the Symbols on the Autoclave



ATTENTION!



PROTECTIVE EARTH CONNECTION



HOT SURFACE, KEEP AWAY



MANUFACTURER



COMMUNAUTÉ EUROPÉENNE



SERIAL NUMBER



ENVIRONMENTAL PROTECTION



3 Safety Information

Besides the information in this instruction manual, the general safety and accident prevention regulations of the legislation must be observed.



Please read carefully! To avoid personal injury, fire or damage to the autoclave, please pay attention to the following safety rules:

- ⚠ Do not install the autoclave anywhere where the power supply cannot be disconnected quickly.
- ⚠ Operate the autoclave with the supply voltage indicated on the type plate of the autoclave only.
- ⚠ **DANGER** Use a socket with protective conductor and make sure that the protective conductor is well earthed.
- ⚠ Do not touch the mains plug or socket with wet hands.
- ⚠ Do not strain, bend or twist the cables. Do not place any heavy objects on the cables.
- ⚠ Do not block the door of the autoclave.
- ⚠ Do not cover the ventilation openings on the rear of the autoclave.
- ⚠ If you notice any unusual odours or noises during operation, disconnect the autoclave immediately from the power supply and contact your dealer or the manufacturer by email to support@certoclav.com.
- ⚠ **WARNING** If you do not operate the autoclave for a longer period, please disconnect it from the power supply and empty the internal water tank.
- ⚠ Do not modify the device.
- ⚠ The autoclave must not be installed anywhere where there is a risk of explosion.
- ⚠ Do not autoclave containers sealed with compressed air, as these containers could implode or explode.
- ⚠ Fill the water tank with distilled or demineralised water only.
- ⚠ Please fill the bottles and other containers with liquid to maximum half only.
- ⚠ Only remove containers with liquids from the chamber as soon as a liquid temperature of less than 80 °C is measured. There is a risk of boiling retardation of the liquid.
- ⚠ If the pressure is relieved rapidly during autoclaving of liquids owing to a program termination, there is a risk boiling retardation of the liquid. Therefore please always wait until the temperature of the medium has reached ≥ 80 °C before opening the autoclave.
- ⚠ The autoclave must not be transported full.
- ⚠ The autoclave must not be opened with force.

4 Non-autoclavable Substances

Never use the autoclave to sterilise the following materials or alkaline solutions. The sterilisation of such substances can cause an explosion, corrosion of the sterilisation chamber, pipes or wear to the seals.

Explosive substances

- Nitroglycol, nitroglycerin, nitrocellulose and other explosive nitric acid esters
- Trinitrobenzene, trinitrotoluene, picric acid and other explosive nitro compounds
- Peracetic acid, methyl ethyl ketone peroxide, benzoyl peroxide and other organic peroxides

Combustible substances

- Metallic lithium, potassium, sodium, yellow phosphorous, phosphorous sulphide, red phosphorous
- Celluloids, calcium carbide (carbide), calcium phosphide, magnesium powder
- Aluminium powder, magnesium powder, metal powders apart from aluminium powder
- Sodium dithionite (or sodium hydrosulphite)

Oxidising agents

- Potassium chlorate, sodium chlorate, ammonium chlorate and other chlorates
- Potassium perchlorate, sodium perchlorate, ammonium perchlorate and other perchlorates
- Potassium peroxide, sodium peroxide, barium peroxide and other inorganic peroxides
- Potassium nitrate, sodium nitrate, ammonium nitrate and other nitrates
- Sodium chlorite and other chlorites
- Calcium hypochlorite and other hypochlorites

Flammable substances

- Ethyl ether, benzine, acetaldehyde, propylene oxide, carbon disulphide and other substances whose flashpoint is in the range from -30 to 0 °C
- Methanol, ethanol, xylene, benzyl acetate (or amyl acetate) and other substances whose flashpoint is between 0 and 30 °C
- Kerosene, gas oil, turpentine oil, isopentyl alcohol (or isoamyl alcohol), acetic acid and other substances whose flashpoint is between 30 and 65 °C
- Flammable gas (hydrogen, acetylene, ethylene, methane, ethane, propane, butane and other substances, which work at a temperature of 60 °C at an absolute pressure of 0.2 bar)



If a salty solution is spilled inside the chamber, remove all liquid from the chamber and wipe up the liquid around the door seal. Clean the door seal with distilled or demineralised water.



5 Initial Start-up

Check whether all parts are provided when unpacking. If any parts are missing or you discern transport damage during unpacking, please contact your dealer or the manufacturer immediately by email to support@certoclav.com.

5.1 Device description

This autoclave is an automatic steam steriliser with vacuum pump and steam generator. Depending on the program and setting, the sterilisation temperature is between 105 °C and 134 °C. Solid materials and liquids can be sterilised. The autoclave has a flexible media sensor in order to measure and control the temperature directly in the medium. The program parameters and sterilisation programs are displayed on the front LCD screen as well as the mobile touchscreen. If a problem is discerned during the sterilisation process, the autoclave automatically issues an error message. If the boiler pressure or boiler temperature are too high, the autoclave will automatically disconnect the power supply and thus guarantee the safety of the operator. The autoclave is equipped with an internal water tank for fresh and used water.

5.2 Device dimensions

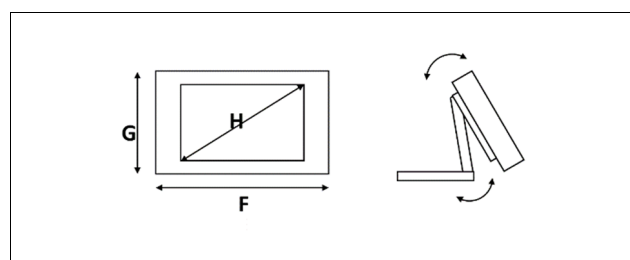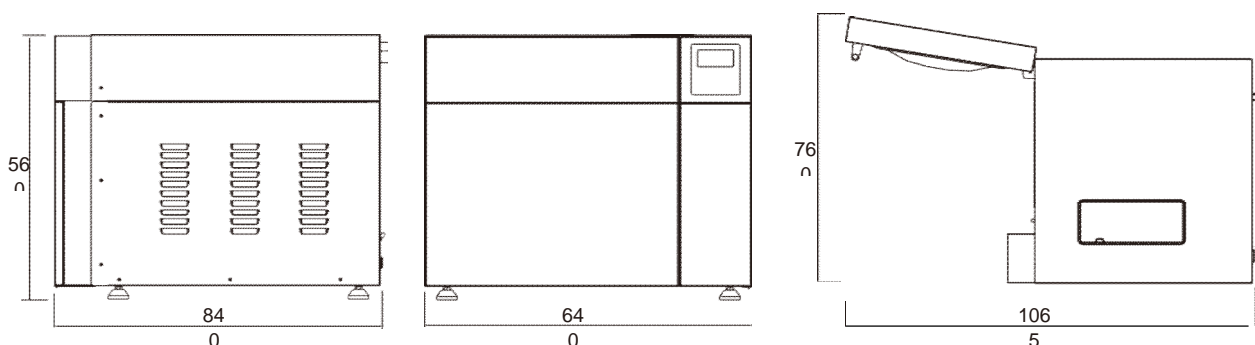

	F	G	H
Touchscreen	200 mm	128 mm	178 mm

5.3 Scope of supply

Standard scope of supply

- CertoClav Vacuum Pro 45
- Declaration of Conformity
- Instruction Manual
- 1 power cable for the autoclave
- 1 waste water hose
- 1 mobile touchscreen
- 1 power cable for the touchscreen
- 1 base for the touchscreen with four fastening screws
- 1 interface box for the touchscreen
- 1 data cable (2 m length)
- 1 tray set with 3 screws

Recommended accessories

Product	Order number
CertoServer for automatic backups with browser-based interface	8500508
CertoCloud Premium licence for the online autoclave management platform www.certocloud.com	8500212
Baseplate for bottles	8500404A
Wire basket including drip pan	8500406
Distillation device for water	8500482
Log printer	1300187
Thermal paper	8500496
Label printer	8500502
Barcode reader	8500503
Direct thermal labels	8500507

5.4 Before initial start-up


Unpacking the autoclave	This autoclave is delivered in a wooden crate. The metal tabs on the bottom of the crate can be opened with a large slotted screwdriver. After this the wooden crate can be raised and removed.
Location	Install the autoclave in a dry location free of dust. The air humidity should be between 30 – 60 % and the ambient temperature 5 – 26 °C. Make sure that the table is suitable for the weight of the autoclave.
Scope of supply	The touchscreen and remaining accessories are in the autoclave chamber when delivered.

6 Installation

6.1 Setting up the autoclave

If you wish to place the autoclave on a laboratory table, check whether the table is suitable for the weight of the autoclave.

For problem-free operation, the autoclave must be aligned by adjusting the device feet using spirit level in direction inclined backwards so that condensation can drain off in the drainage port at the rear end of the autoclave. The autoclave should be installed horizontally in a vertical direction.

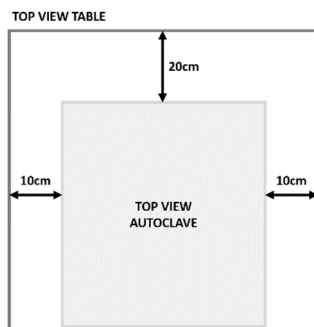


Please leave at least 10 cm free space around the autoclave. Leave at least 20 cm free space at the ventilation slits on the right of the autoclave. Make sure that the room is well ventilated



Keep the specified distance to the surrounding areas. Failure to comply can lead to heat accumulation. This could impair the function of the autoclave and shorten the service life of the pressure pump and result in prolonged program times.

Connect the power cable to the device socket on the rear of the autoclave.


Note: Check whether the data on the type plate of the autoclave correspond to the data for the power source. Check whether the safety switch of the autoclave is in the upper position. Switch on the autoclave at the rocker switch.

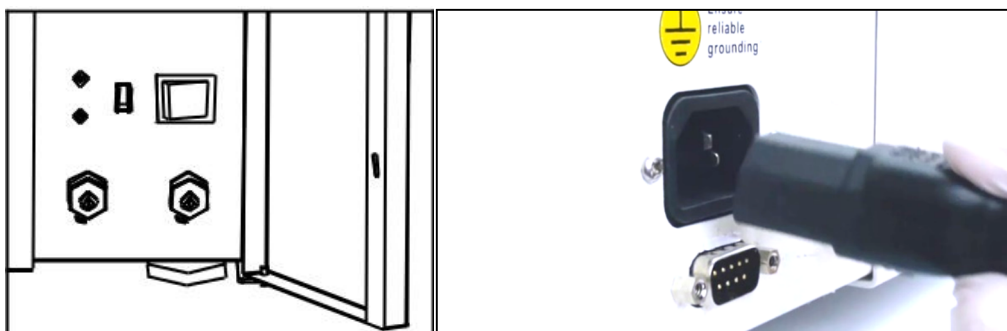

6.2 Unlocking the autoclave door

The touchscreen and remaining accessories are in the autoclave chamber when delivered.

To open the autoclave door, please hold the UP arrow for about 3 seconds until you hear the electronic interlock humming. The door now opens the you ca remove the content and continue installation.

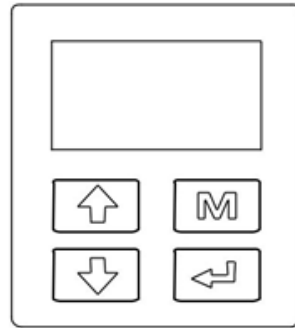

6.3 Touchscreen assembly



Mount the base with the four screws supplied on the touchscreen.



After assembling the base, the touchscreen should appear as shown in the figure opposite.



Connecting the touchscreen to the interface box.




Connecting the power cable of the touchscreen to the interface box.



Place the touchscreen on an even stable surface only. You can also place the watertight touchscreen on the autoclave or next to it.

Connect the autoclave to the RS232 cable at the COM4-port of the extension box to the RS232-port on the rear of the autoclave.

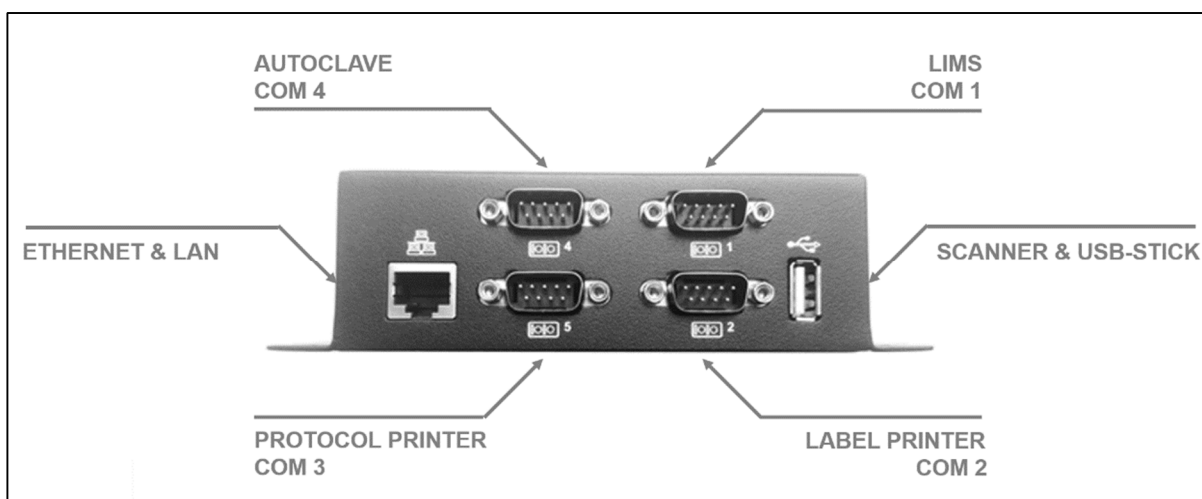

Switch on the touchscreen.

If a login window appears on the screen after several seconds, the touchscreen is correctly connected to the autoclave.

If a notification indicating a connection failure with the autoclave appears on the screen, the touchscreen and autoclave are not connected correctly. Check whether the autoclave is switched on and connected to the COM4-port of the interface box.

6.4 Installing accessories

Optionally available accessories.

6.4.1 Connecting the protocol printer



Connect the protocol printer (thermal or dot-matrix printer) to the COM1-port of the interface box. You can find the connecting cable in the box for the protocol printer. Follow the instructions in the protocol printer manual.

6.4.2 Connecting the barcode reader



Connect the barcode reader to the USB port of the interface box. You can find the connecting cable in the box for the barcode reader. Follow the instructions in the barcode reader manual.

6.4.3 Connecting the label printer



Connect the label printer to the COM2-port of the interface box. You can find the connecting cable in the box for the label printer. Follow the instructions in the label printer manual.

6.4.4 Connecting the CertoServer




If you do not have an Internet connection but wish to use the CertoCloud, we recommend connecting the local CertoServer to the autoclave.

6.5 Pressure sensor calibration

As the atmospheric pressure depends on the installation site, you should recalibrate the pressure sensor after changing the installation site.

Calibration method:

The autoclave is calibrated via the Service screen on the front of the autoclave. The Calibration Menu should only be operated by trained personnel. To access the Calibration Menu, hold down the UP and DOWN arrow keys simultaneously for 10 seconds while switching on the autoclave. To calibrate the pressure sensor to atmospheric pressure, open the door of the autoclave and confirm P-ZERO with the M button.

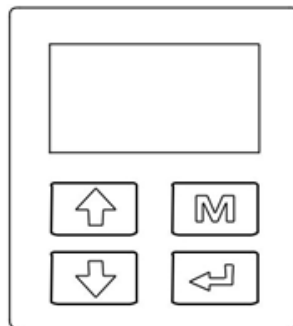

If the pressure sensor has not been calibrated properly after installing the autoclave, a situation may occur where the autoclave fails to start a program.

6.6 Validation and calibration

All sensors of the autoclave have been validated and calibrated at the factory. The autoclave has a G ¼ access on the side panel of the chamber for calibration purposes. We recommend using battery powered sensors such as those made by EBRO. If a calibration on site is necessary after installation, ask your dealer or contact the manufacturer directly by email. To obtain instructions for calibrating the autoclave, please ask your dealer for IQ/OQ documentation or send an email to support@certoclav.com.




The temperature sensors should be calibrated by trained technical personnel only.

To access the calibration menu, hold down the UP and DOWN arrow keys simultaneously for 10 seconds while switching on the autoclave. This menu lets you display and edit the current offset values of the chamber temperature probe, heating temperature probe, steam generator temperature probe or media temperature probe.

7 Sterilisation Programs

7.1 Default settings

Program	Vacuum phases	Sterilisation time [min]	Sterilisation temperature [°C]	Sterilisation pressure [kPa]	Vacuum retention time [min]	Drying time [min]
134°C SOLID	1	4	134	210	-	5
VACUUM TEST	1	-	-	-75	10	15
121°C WRAPPED	3	20	121	110	-	5
134°C WRAPPED	3	4	134	210	-	5
B&D TEST	3	3.5	134	210	-	-
HELIX TEST	3	3.5	134	210	-	-
134°C PRION	3	-	134	-75	15	-
121°C LIQUID	-	30	121	110	-	-

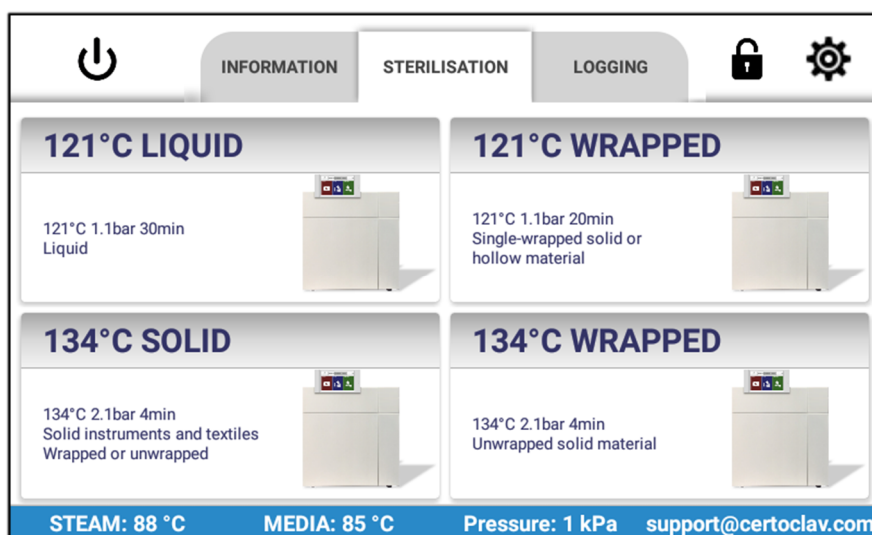

The SOLID program should only be used for unwrapped solids. Hollow bodies of type A and B or porous materials cannot be sterilised with this program.

The VACUUM TEST program performs a pressure leak test.

The WRAPPED programs are suitable for the sterilisation of porous material, hollow bodies of type A and B, packaged waste and instruments. Both programs do not differ in end result. If the load is temperature sensitive and cannot therefore be sterilised at 134°C, please select the 121°C program.

The B&D TEST or HELIX TEST program tests hollow bodies of type A. The programs are almost identical. The only difference involves the fact that the B&D TEST program is a B&D test package, while the HELIX TEST is used as a test device. (PCD: Process Challenge Device).

The 134°C PRION program is specially designed for the sterilisation of prions on all kinds of surface.

The 121°C LIQUID program is used for the sterilisation of liquids in bottles and other open vessels. We recommend using the integrated flexible media sensor for this program. The media sensor is immersed in the liquid medium or in a reference glass with distilled water during the sterilisation. The autoclave now adjusts the retention time during sterilisation so that the medium retains the required sterilisation temperature for at least 30 minutes. If you do not use the media sensor the media temperature cannot be recorded during sterilisation.



Never place pressure-sealed bottles, bags or containers without pressure equalisation mechanism into the autoclave. These could explode due to overpressure during cooling.

Only fill the glasses to maximum half their capacity.

Please perform a vacuum test each month. If the program is not completed successfully, please contact your dealer or the manufacturer by email (support@certoclav.com).

All programs can be customised by the user.

7.2 Editing program parameters

To edit the program parameters, press and hold the M key for 5 seconds while switching on the autoclave at the rocker switch.


Use the arrow keys to toggle between the menu entries.

Use the M key to confirm and save an entry.

Use the Enter key to cancel the process. A menu for editing the program parameters opens. Select the desired program to change the retention time and drying time after the sterilisation phase.


Retention time: 1-59 minutes.

Drying time: 1-19 minutes. Setting parameter N to 0 deactivates the drying.

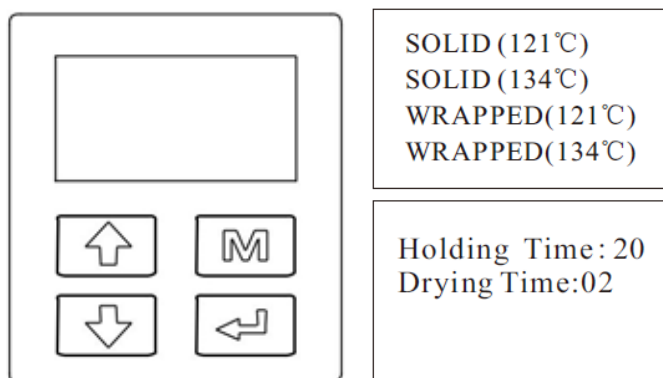


8 Basic Operation


Click and swipe

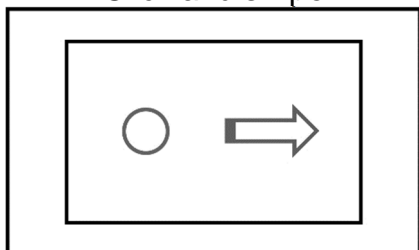

Long click and drag

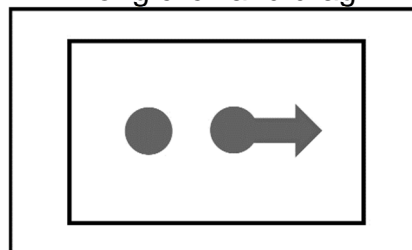


8.1 Switching on the touchscreen

Press and hold the  button on the front (to the right of the screen) for at least 3 seconds.

Switch on the autoclave at the rocker switch below the front cover.





8.2 Switching off the touchscreen

Press the  button on the front (to the right of the screen) for at least 3 seconds. A message will then appear on the screen. Select “Shutdown” to switch off the screen.



8.3 Standby Mode for the touchscreen


Briefly press the  button next to the screen. The background light of the screen then switches off. Press the  button again to reactivate the touchscreen.

9 Instructions for Daily Use



Please make sure that the autoclave is earthed.

9.1 Preparations before use

Switch on the autoclave at the green mains switch. Switch on the touchscreen directly via the screen. To do so, press the  button on the front of the screen (to the right of the screen) for at least 3 seconds. After this the autoclave touchscreen switches on and displays the login window. The autoclave is ready for operation.



Before the first use, the touchscreen displays the notification “Please fill the water tank with distilled water”. As soon as the water is filled, the notification disappears and the autoclave is ready for operation.

9.2 Replenishing with distilled water

The inlet to the freshwater tank is located on the cover of the autoclave. Open the tank cover and full the tank with distilled water.



Only distilled or demineralised water is suitable for operating the autoclave.

Make sure that you do not overfill the tank. Do not fill with more than 4 litres of water maximum. If the tank has been overfilled, the surplus water can be drained via the wastewater hose on the front of the autoclave. If the notification “Please fill the water tank with distilled water” appears on the screen during sterilisation, the sterilisation process must not be terminated. Before the next sterilisation process is started, however, the tank must be replenished with distilled water.



Before filling the water tank with distilled water, please switch on the autoclave and make sure that the autoclave is horizontal with 1° inclination to the rear.



Only fill the fresh water tank with demineralised or distilled water. A water quality sensor continuously monitors the quality of the filled water. If the water used to full the tank is too impure, a warning message will appear on the integrated LCD screen of the autoclave and touchscreen.



Neither the manufacturer nor dealers are liable for damage resulting from the use of contaminated water.

9.3 Opening the autoclave door

To open the door, click on the lock icon at the top right of the Main Menu. The door then opens automatically.




9.4 Steps to start a sterilisation process


First check if the autoclave is ready for operation before starting the program. If no warning or error message is displayed on the touchscreen, the autoclave is ready for operation.

9.4.1 Selecting the sterilisation program on the screen

After logging in, navigate to the Sterilisation tab. You will see 4 programs in the centre. Scroll down in the list to view all available programs.

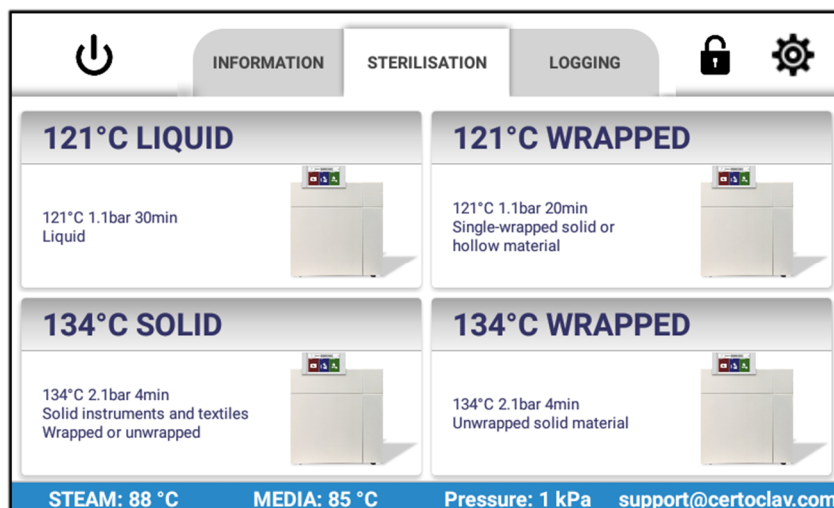

Press on a program to select it. Confirm that you wish to start the program.

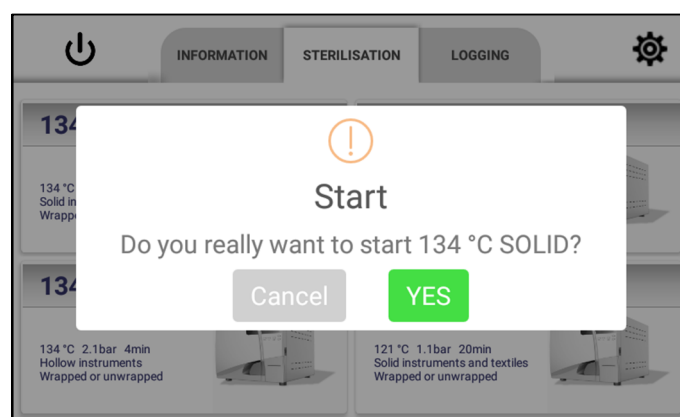


If activated in the settings, step-by-step videos will be displayed before starting the program.

9.4.2 CertoTrace – Documentation and load management

CertoTrace is a software package enabling simple and professional documentation of the entire sterilisation process and tracing of all sterilised products.

If CertoTrace is activated, the CertoTrace Assistant will be executed directly after selecting a program.

Scan the barcode of bottles, instruments or other objects that have been placed in the autoclave, in order to be able to trace their sterilisation with CertoTrace.

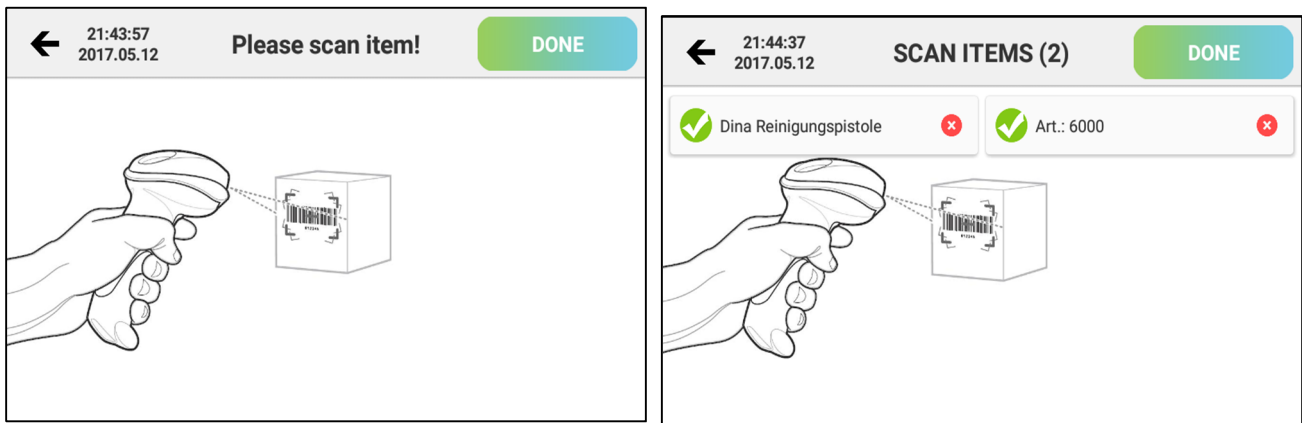


Once an object has been scanned in, it will appear in a list of the CertoTrace Scan Menu. The object and the barcode are noted in the protocol of the sterilisation process.


After all objects have been read in and placed in the sterilisation chamber, press on the Next button.

To terminate the CertoTrace Assistant, press the Back button at the top left of the screen.

If a barcode does not exist for an object, use the label designer to generate and print out a barcode for the object.

To open the label designer, navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Press on “Sterilization” and then “Open label printer designer”.

9.4.3 Step-by-step video assistant

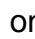To activate the step-by-step video assistant, navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Select the Sterilisation entry on the left of the Settings window. Press on the Step-by-step video assistant to activate or deactivate the step-by-step video assistant.

The step-by-step video assistant guides the user through all steps necessary to operate the autoclave via video instructions.



9.4.4 Loading the autoclave



Please make sure that the total volume of the loads is less than 70 % of the volume of the autoclave chamber.



Please leave at least 10 mm free space between the load and sterilisation chamber wall.

To check the effectiveness of the sterilisation processes independently, place the indicators or bags together with the load into the sterilisation chamber. If you execute the B&D or Helix program, place the indicator in the centre of the sterilisation chamber. A reliable test result is attained in this way.

9.4.5 Closing the autoclave door

To close the autoclave door, press the door gently against the boiler. The vacuum pumps then suction the door so it is automatically locked electrically.




If the temperature in the sterilisation chamber is too high, the door of the autoclave can only be closed with difficulty. Wait 2 minutes before closing the door.

9.4.6 Material tests and automatic repetitions

Material tests can be activated with checkboxes in the settings. To activate or deactivate Material Test Mode, navigate to the Settings Menu by clicking on the ⚙ symbol at the top right of the screen. Press on “Sterilization” and then “Material testing mode”.

If Material Test Mode has been activated, the autoclave asks how often the program is to be executed automatically in succession before the program start.

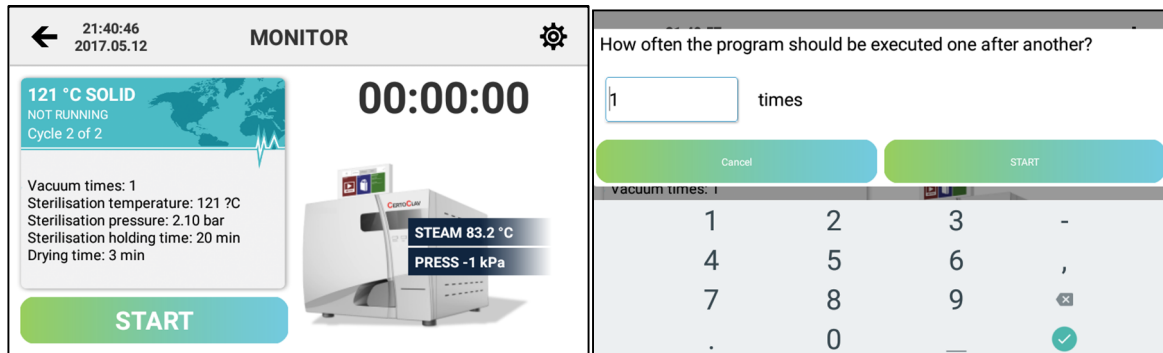


You can enter a number between 1 and 1000. The program will execute the desired number of cycles fully automatically in succession. The status of the program cycle can be called up at any time via a browser at www.certocloud.com or offline on the CertoClav server at <http://certocloud>. If you have activated email or SMS notifications, you can call up the current status of the material test by SMS or email. The email or SMS notifications can be configured in the settings. You can find the notifications in the Settings under the Notifications menu item.

9.4.7 Status screen

The Status screen becomes visible as soon as a program is running.



Information on the current cycle is listed on the left of the screen. The current steam temperature and data on the pressure are on the right of the screen. If the liquid program has been started, the temperature of the liquid will also be visible.

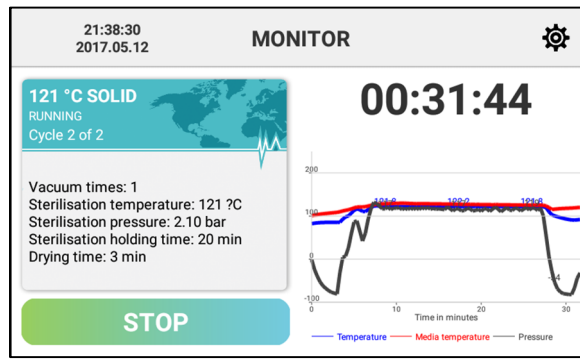

To show a graph of the temperature and pressure, hold and drag the autoclave symbol to the left. The graphic recording will then become visible.

9.5 Unloading the autoclave

As soon as the sterilisation cycle has finished, “Program finished” is displayed on the screen and you will hear three beep sounds. Return to the Main Menu and press the lock icon to open the door. You can then open the door and remove the sterile load.




Risk of scalding! Hot steam or hot water can escape from the chamber when opening the door.

Do not open the door until the pressure in the sterilisation chamber is in the range from -5kPa to 5kPa.

9.6 Emptying the used water tank

If “Empty the used water tank” is displayed on the screen, the used water tank is full. Please empty the used water via the wastewater hose supplied. The left port is for used water, the right port is the drain for distilled fresh water.

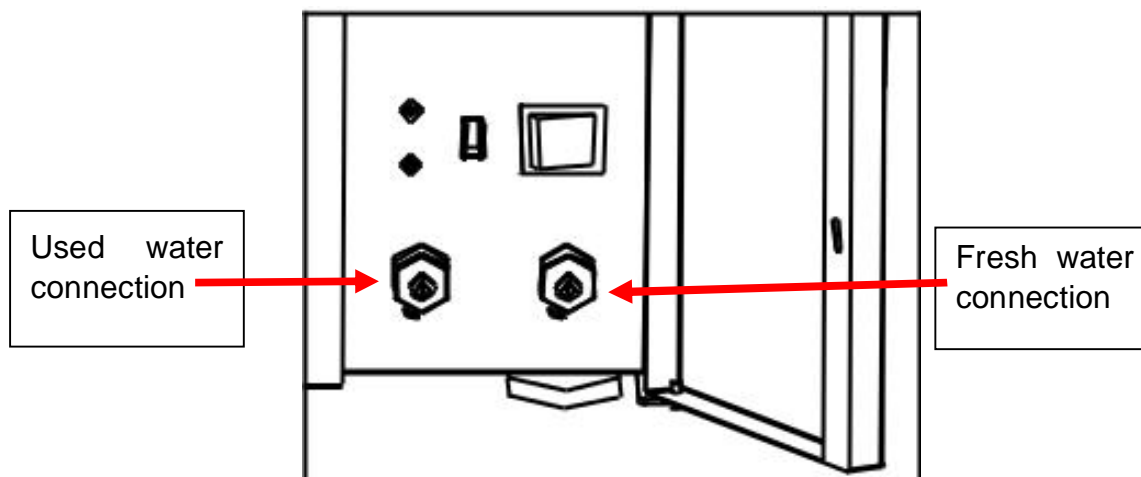


Used water may be hot.

9.7 PC software

CertoCloud is a browser-based management software package for your autoclave. The software is available at www.certocloud.com for PC, smartphone and tablets.

CertoCloud is also available as a laboratory-internal private cloud. The CertoServer must be connected to your laboratory network for this. The CertoServer does not require an Internet connection.

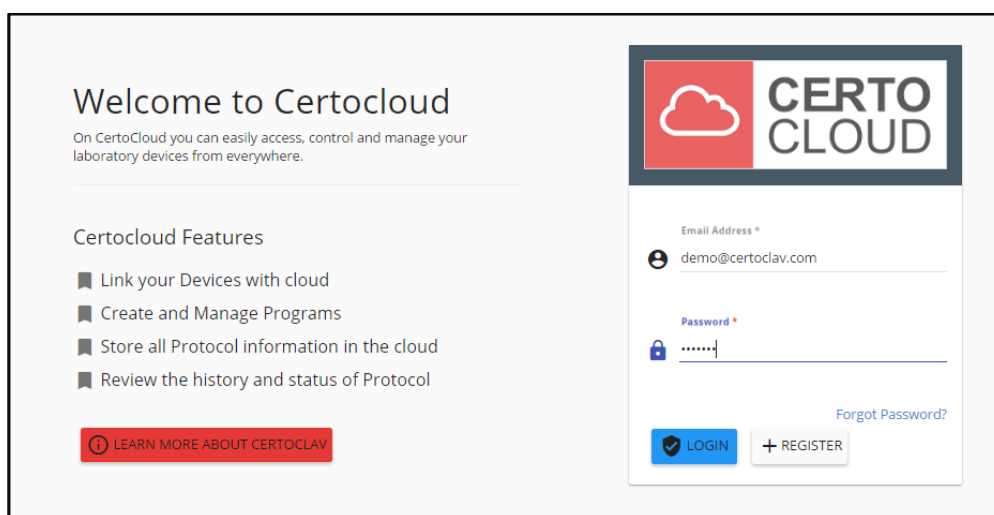

9.7.1 Functions of the CertoCloud

Email Address *

max.muster@mail.com

Password *


.....

[Forgot Password?](#)

LOGIN **+ REGISTER NOW**

Creation of user accounts directly at the autoclave or web-based.

Autoclave Lab 2



Serial
220002576

Model
**Certoclav
22 Vac
Pro**

Firmware
DEFAULT

Tracing real-time data.

If An Error Occured, Send Notification

SMS 0123456789 ✕


EMAIL max.muster@mail.com ✕

Administration of SMS and email notifications. For instance, if maintenance is due.

List of Protocols

Cycle ↑	Program Name	Start
76	cloud generated	7/22/ PM
75	121°C SOLID	7/22/ PM
74	Neues Programm	7/22/ PM
74	121°C SOLID	7/22/ PM

Downloading of all sterilisation protocols as signed PDF, Excel or ZIP file.

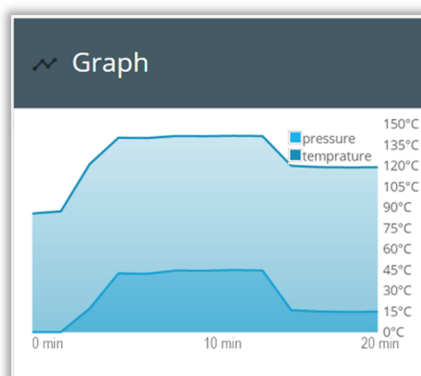

Viewing of all protocols as graphic view, list or detailed view.

Protocol Verification

Protocol ID *

0 / 24

Device Key *

0 / 17

Security Code *


0 / 40

VERIFY

Validity verification of exported protocols.

9.7.2 Display of sterilisation protocols via web browser


To manage the autoclave and view protocols, please use a browser to visit www.certocloud.com or <http://certocloud>. A typical protocol is shown below.

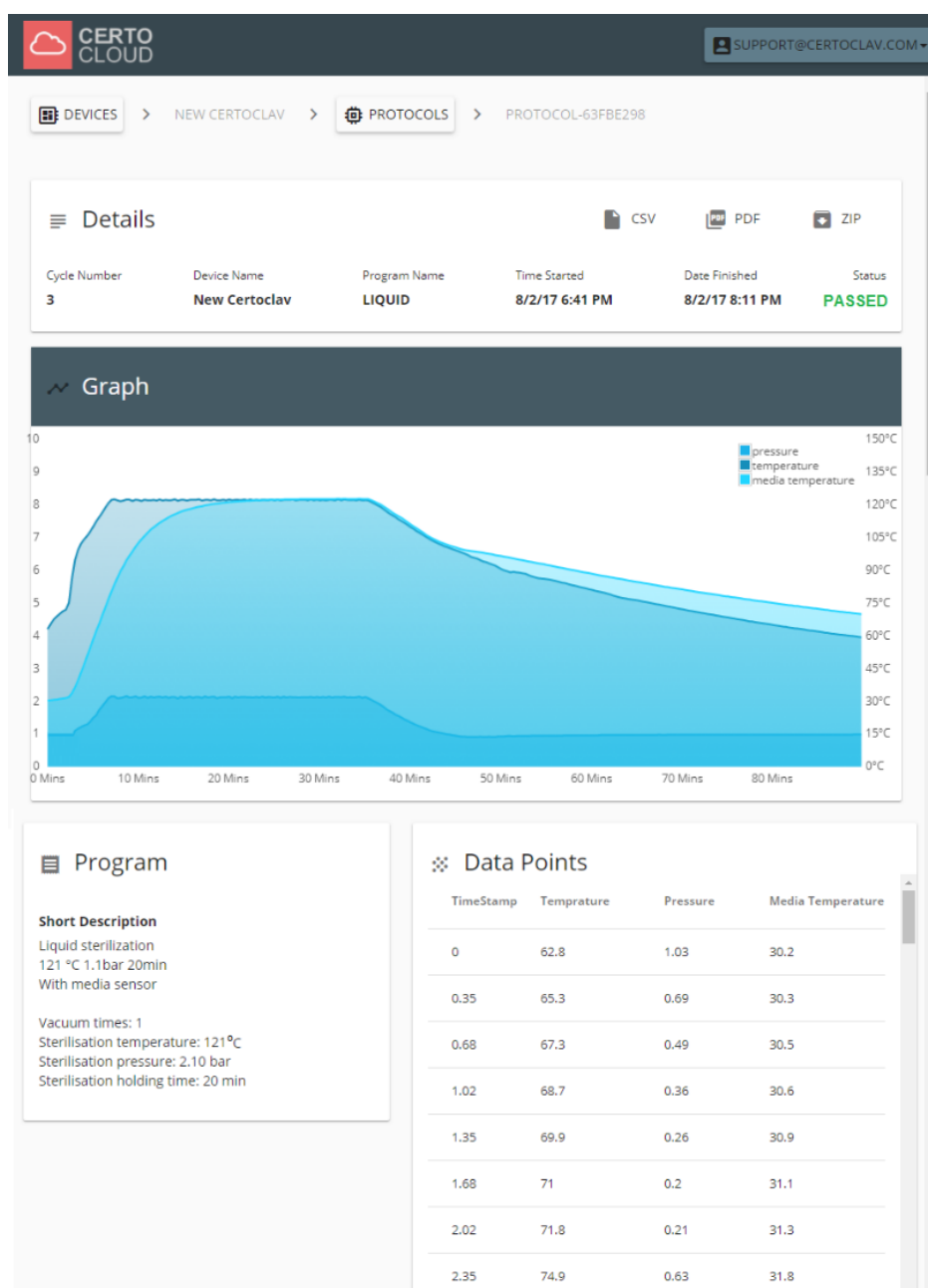


10 Control Panel of the Autoclave

Icons, tabs and text fields can be edited by pressing with your finger.


10.1 Login Menu

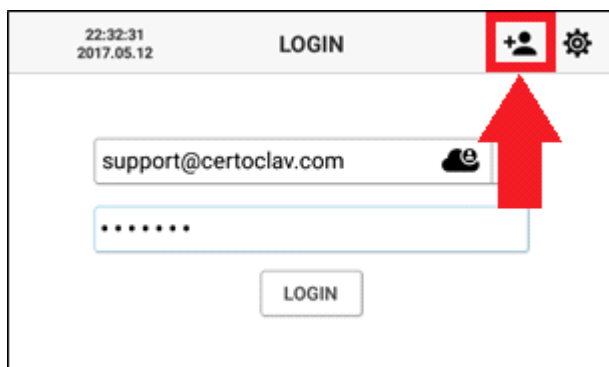

The Login window appears after switching on the touchscreen. To create a new user, press on the plus symbol at the top right of the screen.

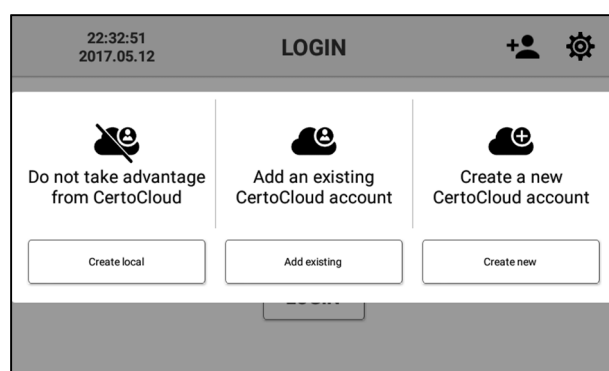

To create a local account without server connection, press on “Do not take advantage from CertoCloud”.
To add an existing CertoCloud user account with server connection, press on “Add an existing CertoCloud account”.
If you wish to create a new CertoCloud user account with server connection, press on “Create a new CertoCloud account”.

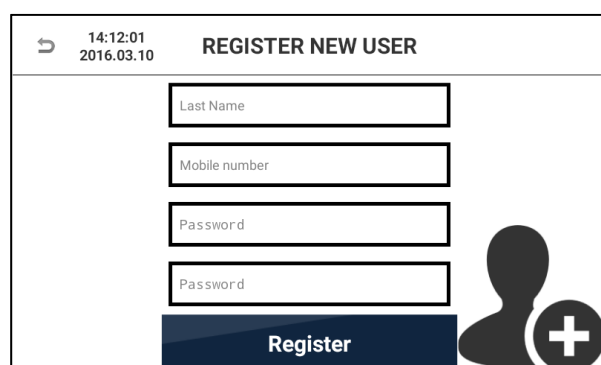

Enter your user data to create a new CertoCloud user account. Enter your phone number including country code. This is needed if you wish to receive status notifications from the autoclave by phone. Enter a password of your choice with at least 4 characters and repeat entry. To cancel the registration, press on the Back button in the top left corner of the screen. Press on “Register” to complete the registration.


To add an existing CertoCloud user account, enter your email address and password. Then press on “Add account”. To cancel the registration, press on the Back button in the top left corner of the screen.

To enable you to log into a CertoCloud user account recently created, first select your email address from the list and then enter your password. Make sure that Online Mode is activated if you wish to use a user account with server connection.

When you log into your CertoCloud user account for the first time, you will be prompted for an activation code. You should have received the activation code by SMS or email after you created your CertoCloud solution. Enter this code in the text field and press on “Activate account”. Your CertoCloud user account is now activated and you can now log into the autoclave with this.


10.2 Support Menu

Navigate to the Information tab to view the Information Menu. Here you can watch training videos, contact the CertoSupport directly from the touchscreen or order spare parts, accessories or service for your autoclave.

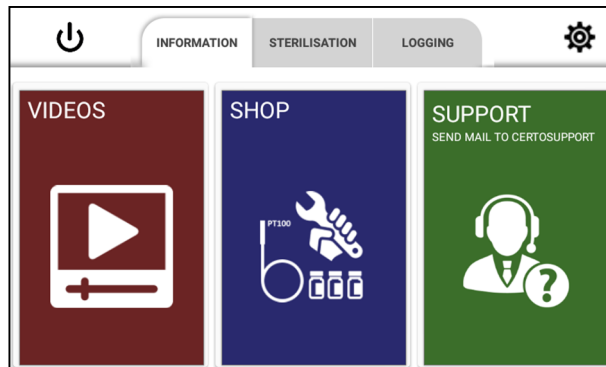


10.2.1 Tutorial videos

To scroll through the list of all available videos, press on the video section in the Information tab, Select the required video to view it in full-screen mode.

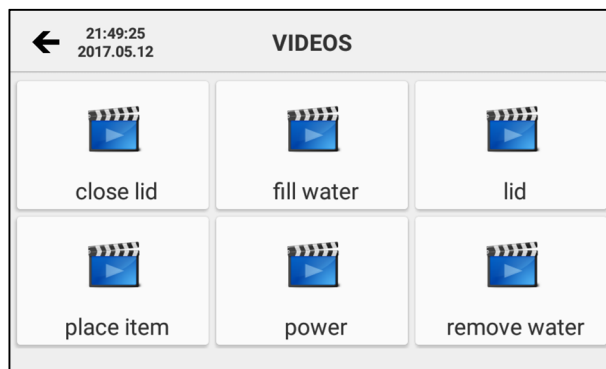


10.2.2 Send email to CertoSupport

An Internet connection must be available to send an email to CertoSupport from the touchscreen. To connect the touchscreen to the Internet, you can either connect a network cable to the Ethernet port or establish a WLAN connection to your WLAN router. Navigate to the Information tab and press Support. In the dialog window that follows select the option “Send an Email to CertoSupport”.

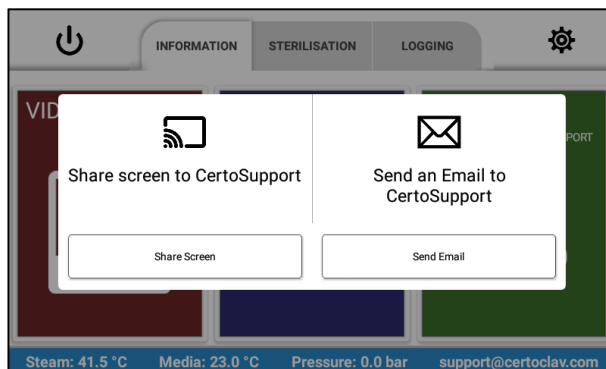


Now select the desired subject from the dropdown list and then enter your message in the text field. Press on “Send” to send the email. We will respond to your question within 24 hours via the email address of your autoclave account.

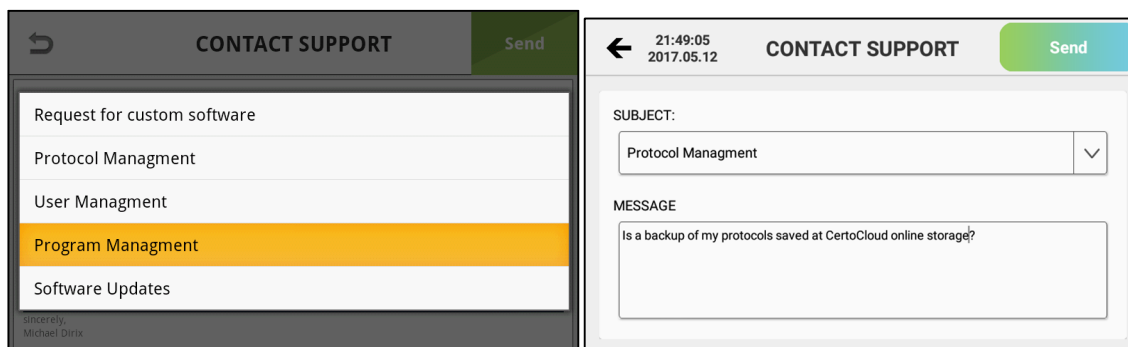

10.2.3 TeamViewer, remote support and share screen


To share the screen via TeamViewer, navigate to the Information tab and press Support. In the dialog window that follows select the option “Share screen to CertoSupport”. The TeamViewer request is now accepted if a schedule with the CertoSupport has been agreed beforehand via email or phone.

To arrange a time window for the remote support, navigate to the Information tab and press Support. In the dialog window that follows select the option “Send an Email to CertoSupport”.

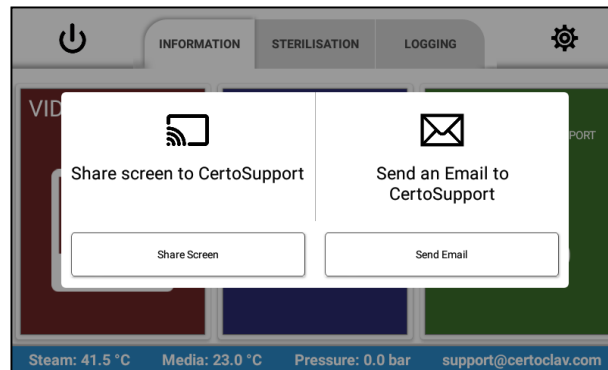


Request a schedule for the remote support. We will usually respond to your request within 12 hours.


As soon as you start a TeamViewer session, a window appears, giving the opportunity to describe the problem roughly once again. Then press on “Start”.

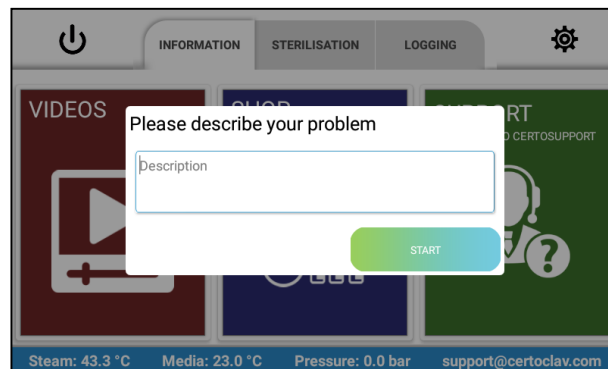

After pressing on Start, wait until CertoSupport has joined the TeamViewer session.

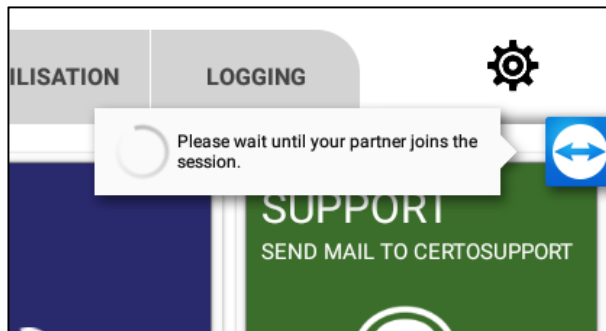

As soon as CertoSupport has joined the session, we will ask your permission to split the screen. Press on “Allow”.

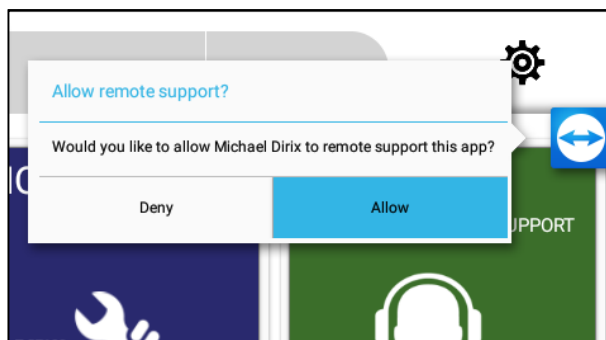


The CertoSupport service technician is now able to service your autoclave. Please wait until the technician has made the required settings to your device. To cancel the screen transmission, drag the TeamViewer icon to the X symbol at the bottom of the screen.

10.3 Protocols Menu


10.3.1 Viewing protocols

To view sterilisation protocols, navigate to the Logging tab. A list of all recorded protocols is displayed on the left half of the screen. The cloud symbol in the protocol list indicates whether the protocols have been backed up in the CertoCloud or on your local CertoServer. Press on the dropdown menu above the protocol list to sort the protocols according to define criteria. To display the details and graphic view, select the protocol by pressing with your finger on the list element. A graph of the protocol will now be shown on the right of the window. To display the protocol as a list, press on the list symbol above the graph.

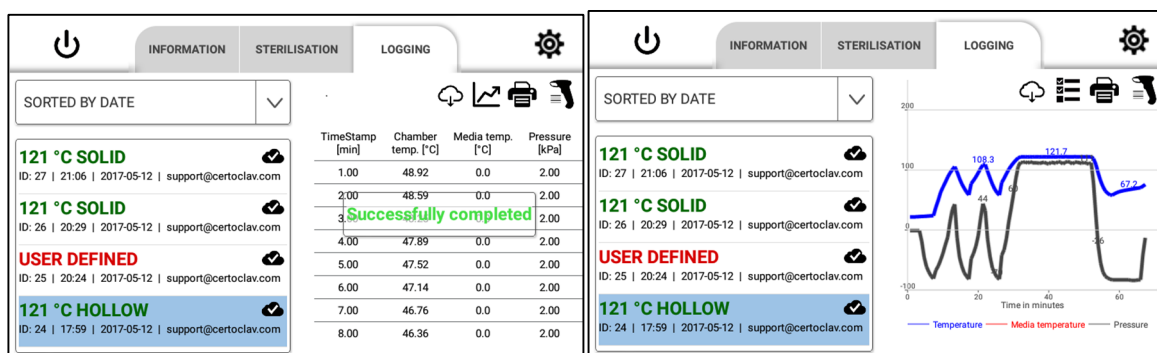

10.3.2 Printing protocols

Navigate to the Logging tab. Select the protocol you wish to print out on the left of the screen. Press on the pressure symbol at the top right of the screen. Now select the option "Print the protocol".

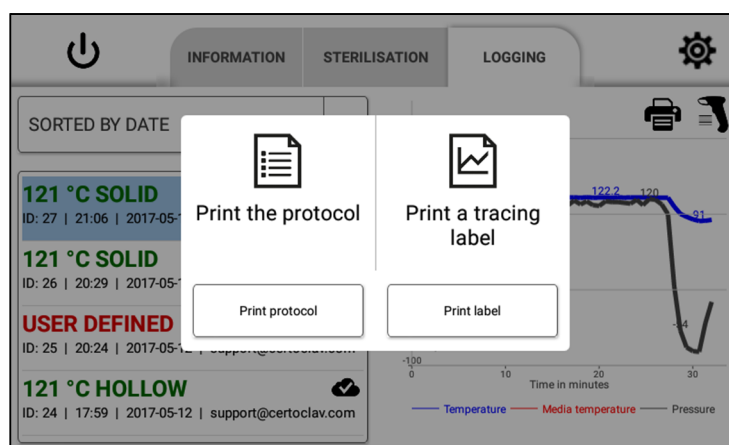

10.3.3 Example of a protocol printout

Optional

```

Header line 1
Header line 2
Header line 3


Protocol      CertoClav Vacuum Pro Series
Model:        VAC-PRO-45L
S/N.:         22ABCDEFGHJ
Program:      134°C HOLLOW
Program description:
    3 x venting
    134 °C   2.1bar   30min
    Hollow instruments
    wrapped or unwrapped
Media:        NO
Cycle no:     971
Start time:   12.01.15 17:42:00
End time:     12.01.15 18:16:00
Result:       PASSED
hh:mm:ss      Temp °C   Press [bar]
17:42:00      27.5      1.00
17:42:20      27.5      0.50
17:42:40      27.6      0.18
17:43:00      80.4      0.80
[...]
18:15:00      65.9      0.18
18:15:20      65.4      0.18
18:15:40      68.0      0.50
18:16:00      70.2      1.00
Program finished successfully
END

Signed by _____
Confirmed by _____


```

10.3.4 Print barcode

Navigate to the Logging tab and select the protocol last recorded on the left of the screen. Press on the pressure symbol at the top right of the screen. Now select the option “Print a tracing label”.

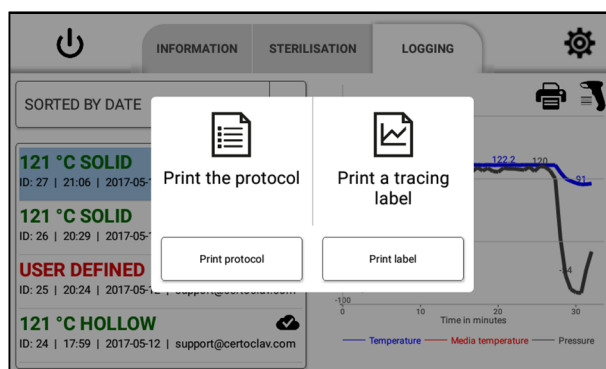


Now enter the shelf-life of the sterilised products.

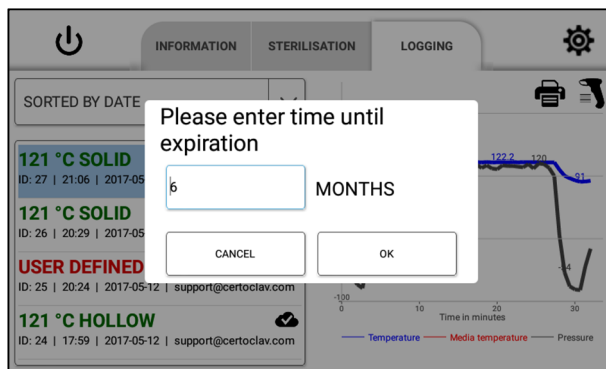


Then press on OK to print the self-adhesive label with shelf-life date and barcode.

10.3.5 Tracing marked products

Navigate to the Logging tab and press on the Scan icon.

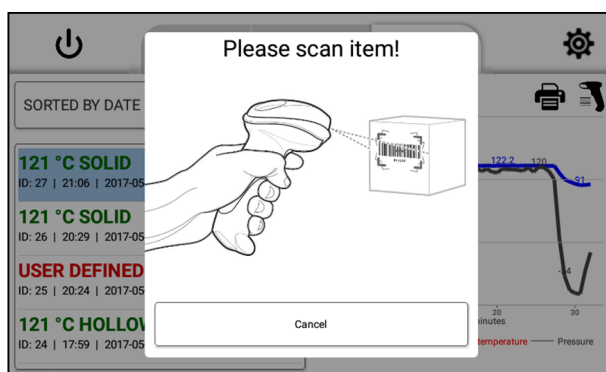


A dialog window appears. Scan the barcode of the sterilised product with the barcode reader. The protocol for the sterilised product now appears on the screen. A graphic view and list view of the protocol can now be seen on the right of the screen.

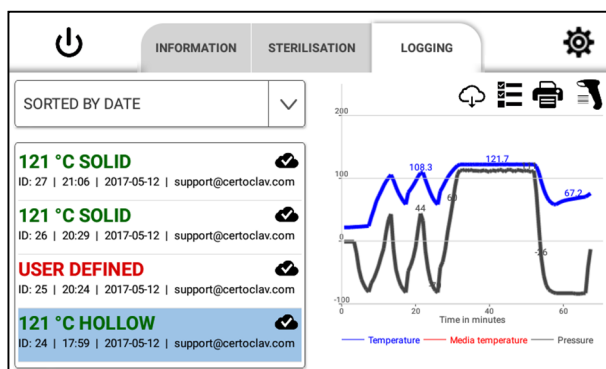

11 Settings


11.1 User administration

Press on the  symbol on the right of the screen to navigate to the settings. Select the User Administration list entry in the left window. The User Administration will now be shown on the right of the window. Press on Edit to change your user data and password or on Delete to delete your user account.




11.2 Network settings

11.2.1 Connecting to a WLAN network


Navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Then press on "Network". If no checkmark is set at "WiFi On/Off", press on the empty checkbox to activate WiFi.

After this press on "Manage WiFi connections" to display a list of available WLAN networks. Select the relevant WLAN network and enter any required access code and further parameters.

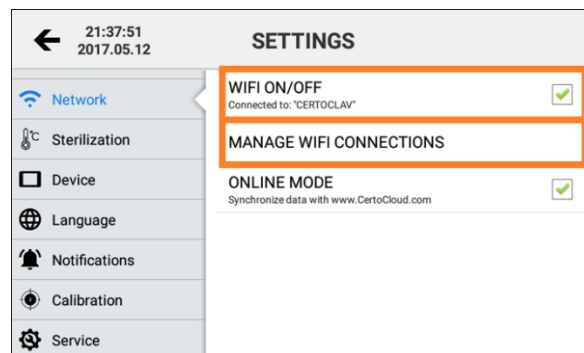


11.2.2 Connecting to a LAN network


The autoclave can be connected to a LAN via an Ethernet cable. Connect the network cable to the Ethernet port of the interface box.

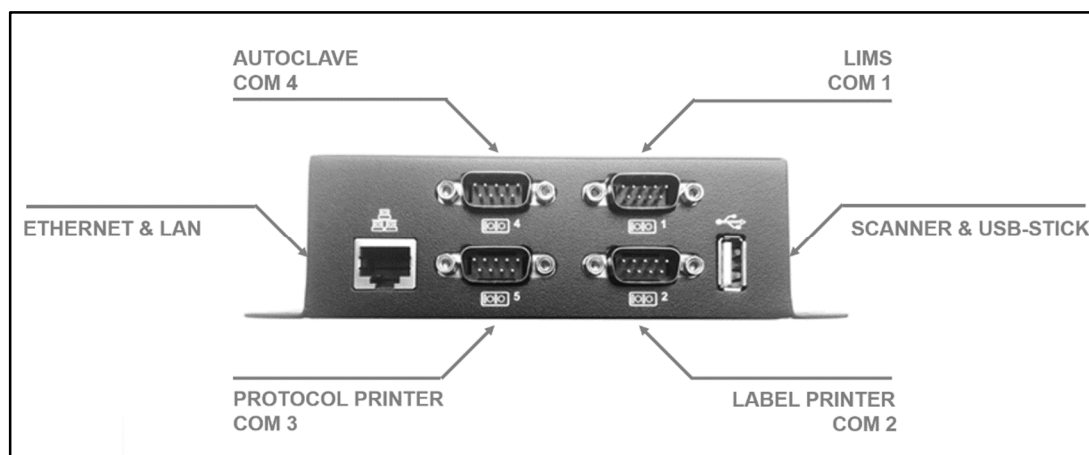

11.2.3 Network communication

Online Mode needs to be activated if the autoclave data is to be synchronised with the CertoCloud or the local CertoServer. If the CertoCloud is used for synchronisation, email and SMS notifications will also be available.

Navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Then press on “Network”. Now select Online Mode to activate or deactivate the data synchronisation.

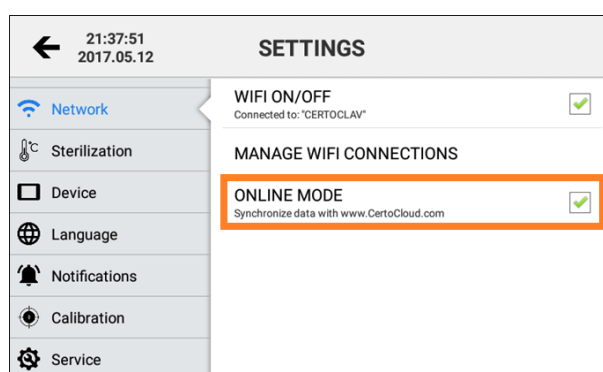


Note: If Online Mode is deactivated, your protocols will only be saved locally on the autoclave. As soon as you reactivate Online Mode, all protocols saved locally up until then are synchronised with the CertoCloud and are thus also available online in your backup memory.

11.2.4 Connecting to the CertoServer


If you do not have an Internet connection but wish to use the CertoCloud, we recommend connecting the local CertoServer to the autoclave.

Connect the CertoServer to the LAN and switch in the server. Connect the autoclave to the same network via Ethernet cable or WLAN.


Navigate to the Settings Menu by clicking on the ☰ symbol at the top right of the screen. Press on “Network” and then on “choose server”.

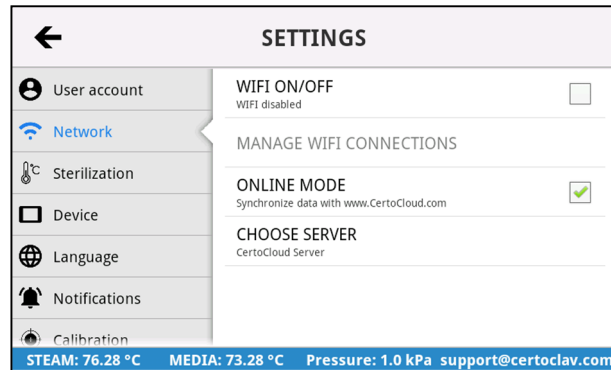


Select the “Add a server from local network” option.

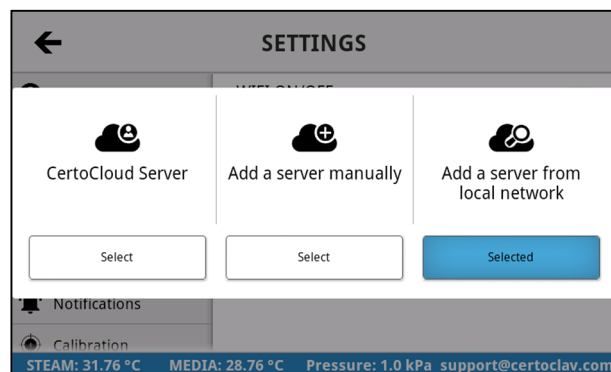

The autoclave now searches for the CertoServer in the local network.

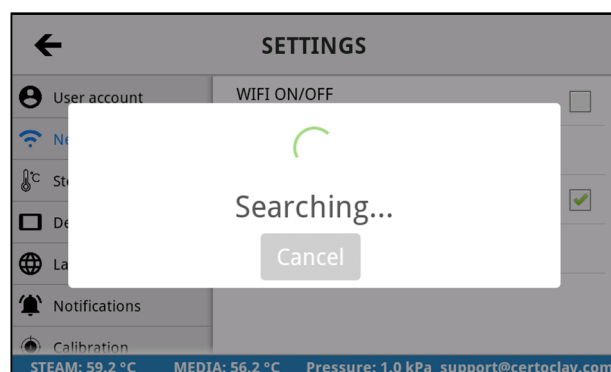


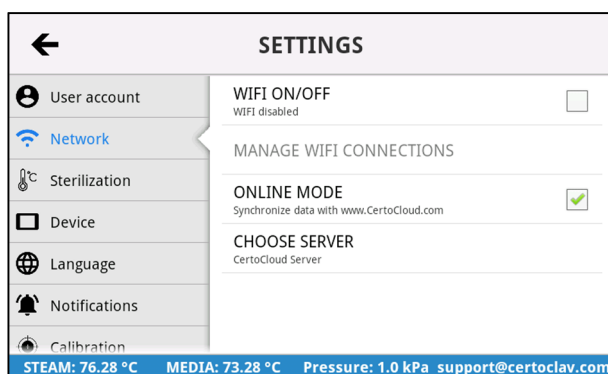
After the CertoServer has been found, confirm with “Save”.

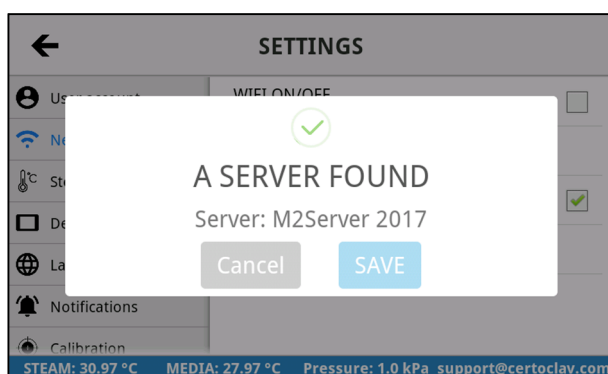


11.2.5 Connecting to the CertoCloud server

If the autoclave is connected to the Internet, you can alternatively use the free CertoCloud for documentation and configuration of the autoclave.

To activate communication with the CertoCloud, navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Press on “Network” and then on “choose server”.




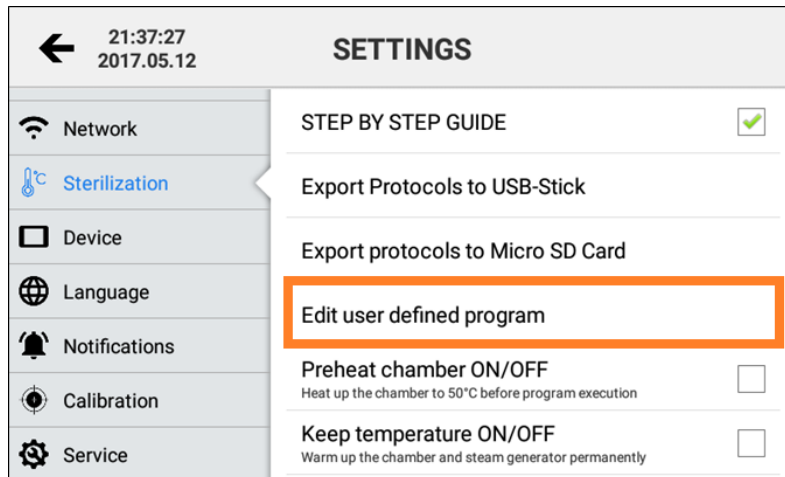
Select the “CertoCloud Server” option.

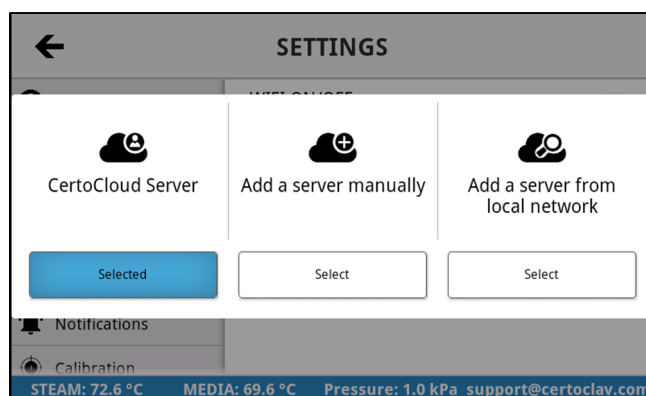

11.3 Sterilisation settings

11.3.1 Customising sterilisation programs


To customise a sterilisation program, navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Press on “Sterilization” and then “Edit user defined program”.



Now define the program parameters individually corresponds to your requirements:

- Number of prevacuum pumps between 1 and 10
- Sterilisation temperature between 105 °C and 134 °C
- Sterilisation time between 4 and 60 minutes
- Drying time between 1 and 25 minutes

Then press on “Apply” to save the program parameters or on “Cancel” to return to the Settings Menu and cancel the process.

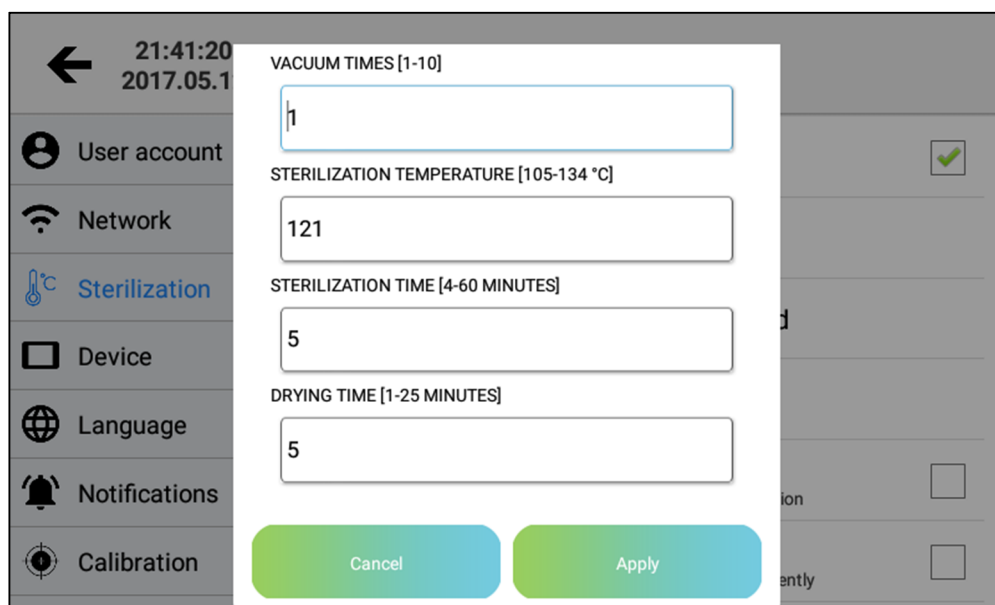

11.3.2 Material tests and automatic repetitions

If Material Test Mode is activated, you will be asked how often a program is to be executed before starting it for the first time.

How often the program should be executed one after another?


times

Cancel START

vacuum times: 1

1	2	3	-
4	5	6	,
7	8	9	✕
.	0	—	✓

This has the advantage that longer material tests can be performed fully automatically without user interaction.

To activate or deactivate Material Test Mode, navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Press on “Sterilization” and then “Material testing mode”.

← SETTINGS

User account

Network

🌡️ Sterilization

📺 Device

🌐 Language

🔔 Notifications

🎯 Calibration

Edit user defined program

Preheat chamber ON/OFF ☐
Heat up the chamber to 50°C before program execution

Keep temperature ON/OFF ☐
Warm up the chamber and steam generator permanently

Material testing mode ☐
On program start, show a dialog to enter how often to run the program automatically


Open label printer designer

CertoTrace load documentation ☐
Enable autoclave load documentation. Barcodescanner and labelprinter required.

11.3.3 CertoTrace — Your support during material tests

CertoTrace is a professional documentation software package for the entire sterilisation process. CertoTrace records which products have been sterilised, when and by whom. The tracing of sterilised products and management of shelf-lives is thus integrated in the work process.


To activate CertoTrace, navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen. Press on “Sterilization” and then “CertoTrace load documentation”. After activation, CertoTrace is automatically integrated in the workflow when starting a program.

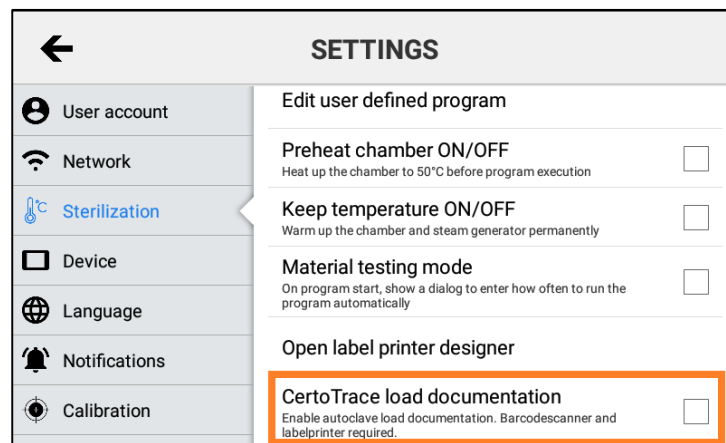

11.3.4 Label designer


The CertoClav label printer prints shelf-life data and a barcode for tracing the product on self-adhesive labels.

You can use the integrated label designer to identify articles with barcodes. To open the label designer, navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen. Press on “Sterilization” and then “Open label printer designer”.

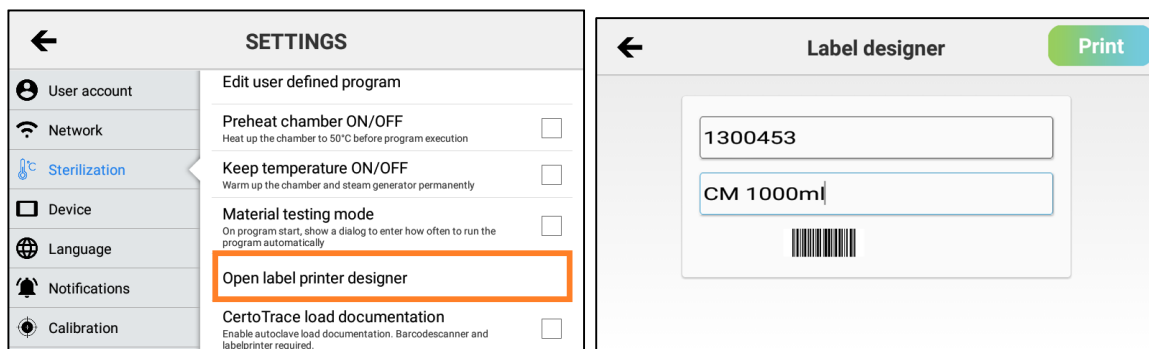


11.3.5 Real-time protocolling to USB


You can use a USB stick to save the sterilisation protocols in real-time during the sterilisation process. For this, please connect a USB stick to the front USB port of the autoclave.

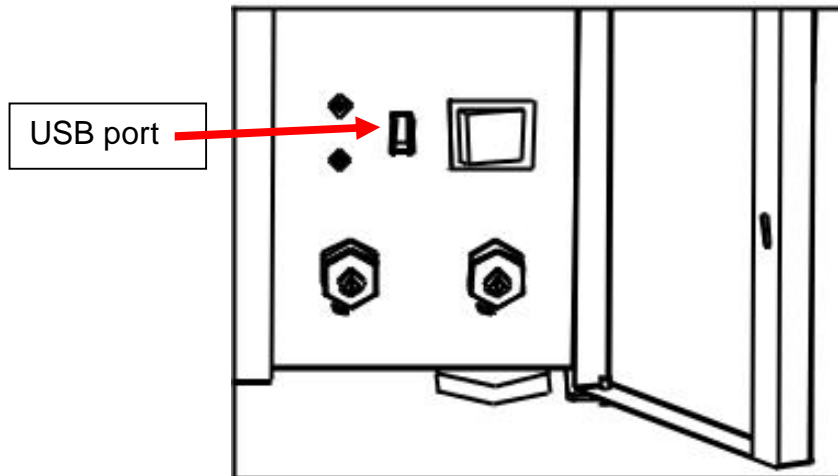


11.3.6 Exporting all protocols to a USB

The autoclave can save any number protocols internally. To export all protocols to a USB stick, plug a USB stick into the interface box.

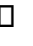Navigate to the Settings Menu by clicking on the  symbol at the top right of the screen. Press on "Sterilization" and then "Export protocols to USB-Stick".

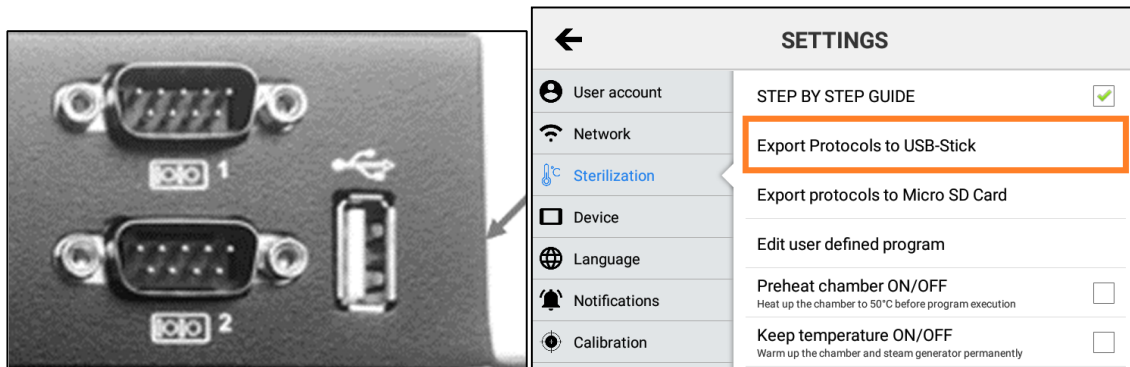


If the protocol export fails, your USB stick has not been recognised. In this case plug the USB stick into the device and restart the touchscreen. Now attempt to export the protocols again.

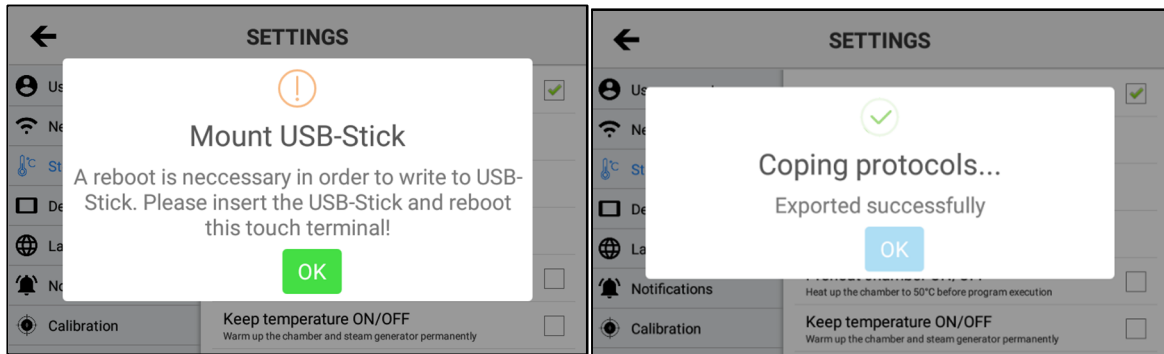


11.3.7 Exporting all protocols to an SD card

The autoclave can save any number of protocols internally. To export all protocols to an SD card, plug a Micro SD card into the opened push-in compartment on the right of the touchscreen.

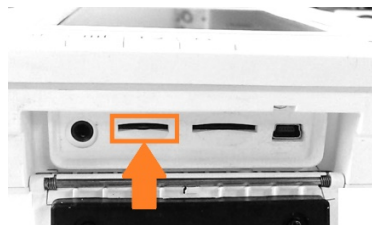


Navigate to the Settings Menu by clicking on the □ symbol at the top right of the screen. Press on “Sterilization” and then “Export protocols to Micro SD Card”.

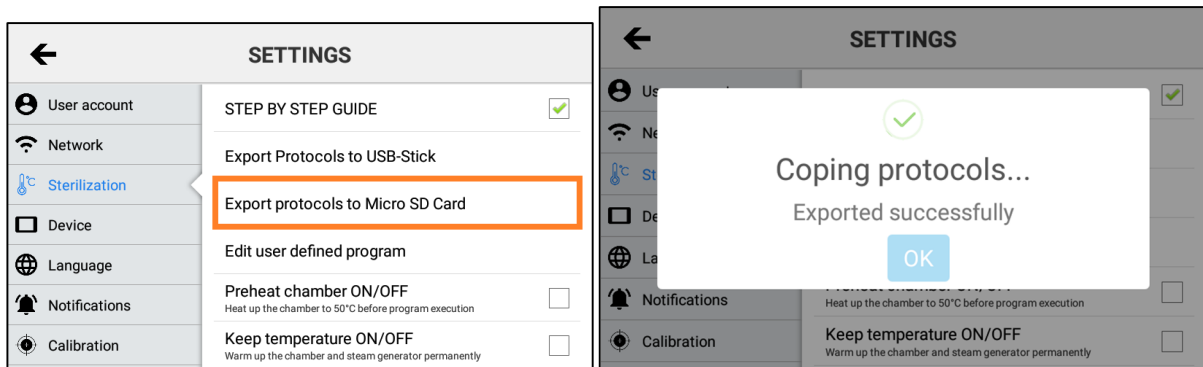

If the protocol export fails, restart the touchscreen with the SD card inserted. Now attempt to export the protocols again.

11.3.8 Exporting all protocols via a network

Log into the autoclave with your CertoCloud or CertoServer user account. Make sure that Online Mode is activated. If the login is successful, all protocols will be synchronised with the CertoCloud. The export has been successful if checkmark icons are displayed in the protocol list.

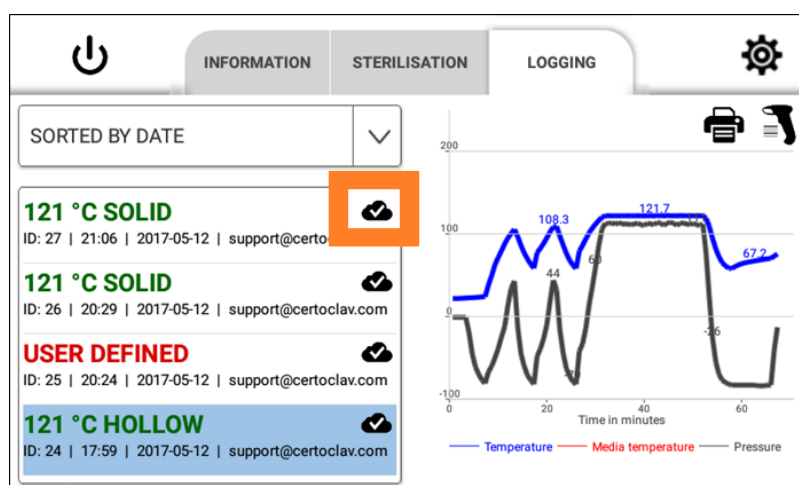


11.4 System settings


11.4.1 Changing the date and time

Use the buttons of the LCB screen to change the date and time of the autoclave. Use the M button to navigate to “Basic Set” and then to “Date” for this. To adapt the required parameters, use the arrow buttons and confirm the correct value with M button.


Program Basic Set Report	DATE:12-04-03 TIME:09:10:08 LANGUAGE:ENG Counter: 12
--------------------------------	---

11.4.2 Installing updates via the Internet

Make sure that autoclave is connected to the Internet. Go to the Settings Menu by clicking on the  symbol at the top right of the screen. Press on “Device” and then on “Check for updates”.

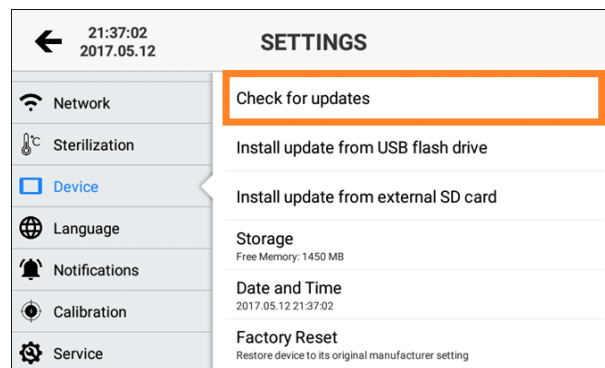

A progress display appears with the download. All updates are automatically installed as soon as they are downloaded. If a new software version is available, an installation assistant will guide you through the installation process.

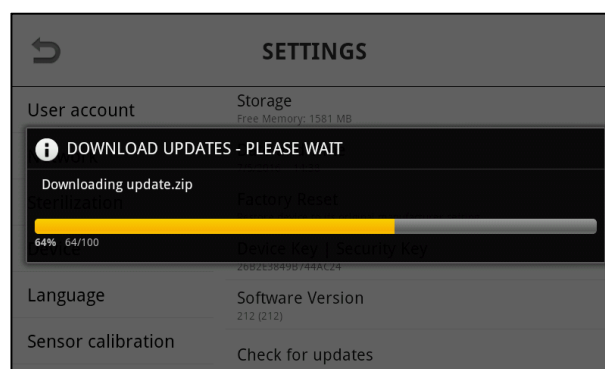


11.4.3 Installing updates via USB

Download the latest update from www.certoclav.com or ask for the latest update for your device by email to support@certoclav.com. Please indicate the serial number of your autoclave for the Support Team. You can find the serial number on the type plate of the autoclave or in the settings. For this, navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen. Press on “Device”.

Copy the “update.zip” to the main directory of a USB stick and plug this USB stick into the extensive cable of the touchscreen.

Navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen. Press on “Device” and then on “Install update from USB flash drive”.

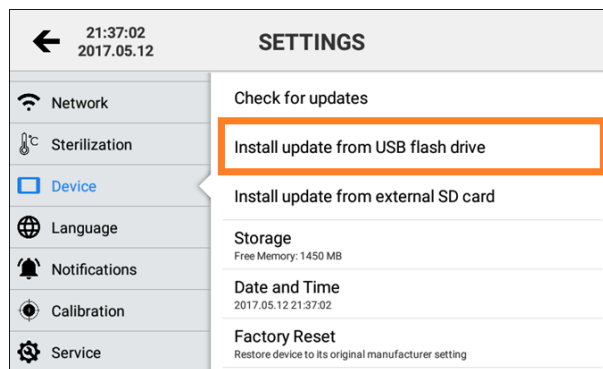


If the USB stick is not recognised, restart the touchscreen and try again. Follow the instructions on the screen to install corresponding updates.

11.4.4 Installing updates via SD card

Download the latest update from www.certoclav.com or ask for the latest update for your device by email to support@certoclav.com. Please indicate the serial number of your autoclave for the Support Team. You can find the serial number on the type plate of the autoclave or in the settings. For this, navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen. Press on “Device”.

Copy the “update.zip” to the main directory of Micro SD card and plug this into the push-in compartment on the right of the touchscreen.

Navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen. Press on “Device” and then on “Install update from external SD card”.

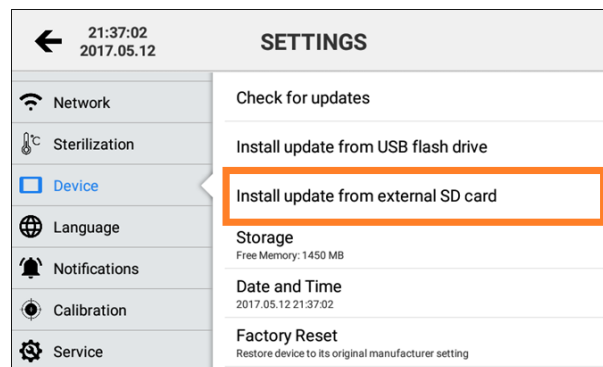

If the SD card is not recognised, restart the touchscreen and try again. Follow the instructions on the screen to install corresponding updates.

11.4.5 Restoring the default settings


Navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen and pressing on “Device”. In the detailed view on the right of the screen scroll to the Restore default settings entry and press on it.



Once you have reset the touchscreen in the default settings, all protocols and user accounts not previously exported and not synchronised with the CertoCloud will be deleted.

11.4.6 Changing the language

Navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen and press on “Language”. Pressing the current language setting in the detailed view on the right of the screen opens a dropdown list with the languages supported. Select a language by clicking on the desired language.

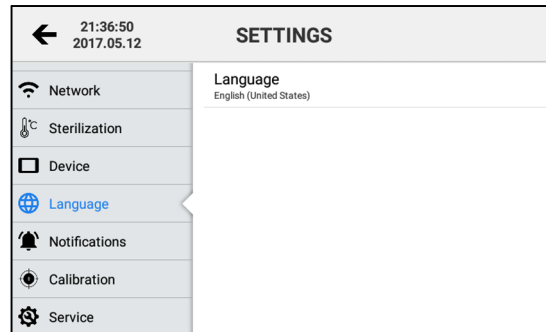

11.5 Configuration of SMS and email notifications


If you still do not have a CertoCloud account, you can create a new CertoCloud account in the Login window.

Note: The autoclave must be connected to the Internet to be able to create the CertoCloud account. Log in with your CertoCloud user account.

To configure SMS and email notifications, navigate to the Settings Menu by clicking on the ☐ symbol at the top right of the screen and pressing on “Notifications”.

If you wish to receive an email if a program is terminated, when maintenance is necessary or once a program has successfully been executed, activate the relevant assigned checkbox.


Enter the email address of your choice, which is to be used for the notification. If you have a CertoCloud account, you can also save a mobile phone number to which notifications via SMS are to be sent. Then press on “Save” to save the changes.

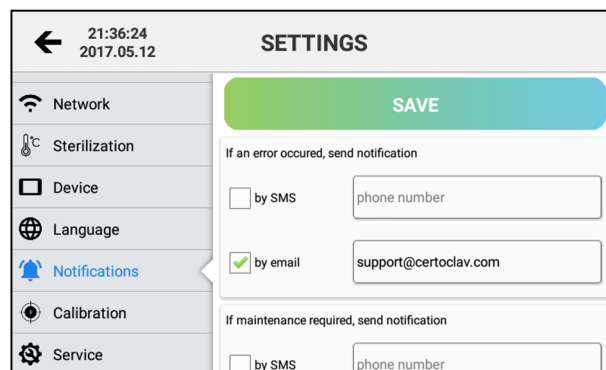

11.6 Sensor calibration




The sensor calibration should be calibrated by trained personnel only.

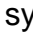To calibrate the sensors, navigate to the Settings Menu by clicking on the  symbol at the top right of the screen and pressing on “Calibration”.

Press on the offset you wish to adjust. The offset parameters must be in the interval [-3°C; 3°C] and [-5 kPa; +5kPa]. Then press on “Apply”.

SETTINGS	
← 21:36:10 2017.05.12	
Network	Offset of steam sensor [°C] <input type="text" value="-1.0"/>
Sterilization	Offset of heater sensor [°C] <input type="text" value="0.0"/>
Device	Offset of steam generator sensor [°C] <input type="text" value="0.0"/>
Language	Offset of media sensor [°C] <input type="text" value="0.0"/>
Notifications	Offset of pressure sensor [kPa] <input type="text" value="4"/>
Calibration	
Service	<input type="button" value="Apply"/>

11.7 GLP and GMP data

To view the GLP and GMP settings, navigate to the Settings Menu by clicking on the  symbol at the top right of the screen and pressing on GLP and GMP data.

11.7.1 Header

The header setting activates the printout of the GLP header. Up to four text lines are available. If this setting is activated, the header will be displayed at the top of each sterilisation protocol and printed out.

11.7.2 Autoclave designation

Designate the autoclave in this setting. An alphanumeric text can be entered. The standard name is empty.

11.7.3 Project designation

Designate your project in this setting. An alphanumeric text of up to 25 characters can be entered. The standard designation is empty.

11.7.4 Advanced GLP and GMP settings

Activate the following checkboxes in the Advanced GLP and GMP settings to configure a printout and digital copy of the sterilisation products.

Setting	Option (Default value is bold)
Print autoclave designation	YES , NO
Print user identification	YES , NO
Print project designation	YES , NO
Print application name	YES , NO
Print signature fields	YES , NO
Print header	YES , NO
Print date	YES , NO
Print autoclave serial number	YES , NO

11.8 Locking settings

To prevent unauthorised changes to the autoclave, sensor parameters, programs and settings, it is only possible to change settings in the password-protected Administrator user account.


12 Error Messages

If an error occurs during a program cycle, the autoclave will sound an acoustic alarm and display an error code on the screen. The autoclave will also automatically release pressure from inside the sterilisation chamber and switch off the heating. This safety system is designed to protect the user. If an alarm occurs, please write down the error coded and switch off the device. Do not open the door. Switch the device on again and wait until the pressure returns to -0.5 kPa ~0.5 kPa.



We recommend rebooting the program in order to see whether the error occurs again.

A repeated long beep sound is audible at each of the following error messages. The user must confirm the error message by pressing the M button for a longer period so that the error memory is deleted.

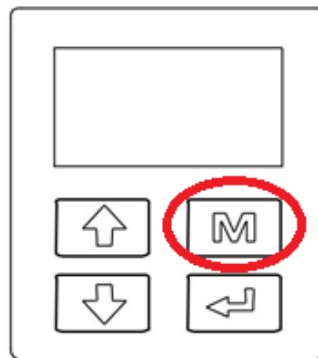

Code	Error	Possible cause	Troubleshooting
E-1	The program was terminated manually by the user.	User has pressed the STOP button during sterilisation.	-
E-2	The communication between the external touchscreen and mainboard is interrupted.	The serial cable (COM 4) between the touchscreen and autoclave has become detached.	Screw the COM4 cable into the RS232 base on the rear of the autoclave again. If the problem is still not resolved, restart the autoclave and touchscreen. If the problem continues to persist, navigate to the Settings to reset the touchscreen.
E1	Sensor error	The sensor of the steam generator is defective.	Contact your dealer.
E2	Sensor error	The sensor of the chamber probe is defective	Contact your dealer.
E3	Sensor error	The sensor of the heating probe is defective	Contact your dealer.

E5	Pressure relief error	The pressure could not be released from the chamber	Contact your dealer.
E6	Door sensor error	The door is open during sterilisation	Contact your dealer.
E9	Temperature control	The temperature could not be retained	Contact your dealer.
E11	Steam generator heating	The steam generator could not be preheated	Contact your dealer.
E12	Chamber heating	The jacket heating could not be preheated	Contact your dealer.
E13	Vacuum pump or leak tightness	The generation of a vacuum during a cycle has failed	Clean the door seal and restart the process. Contact your dealer if the error persists.
E20	Program terminated manually	The user has terminated the program manually during a current cycle	
E22	Vacuum pump or leak tightness	Vacuum test has failed	Clean the door seal and restart the process. Contact your dealer if the error persists.
E23	Vacuum test was terminated unexpectedly	The temperature in the chamber has become too hot during the program cycle	Deactivate the “Keep warm” or “Retain temperature” function in the sterilisation settings and wait until the chamber has cooled down.
E24	The pressure is less than 30 kPa at an temperature of 120°C.	The temperature-pressure ratio does not correspond to that of saturated steam.	Contact your dealer.
E25	Door locking failed	The door could not be locked. The door is possibly set too narrow.	Contact your dealer.
E26	Door opening failed		Contact your dealer.
E27	The vacuum test cannot be started	The temperature in the chamber is too hot to start the vacuum test	Deactivate the “Keep warm” or “Retain temperature” function in the sterilisation settings and wait until the chamber has cooled down.

13 Maintenance



Please use original spare parts of the manufacturer only. Please contact your dealer if you have any questions regarding maintenance and repair.

Make sure that the autoclave is disconnected from the power supply during maintenance and repair.

Make sure that autoclave is at room temperature.

The following components need to be checked or replaced regularly:

- Sealing ring
- Safety valve

Inspection plan:

Component	Inspection by	Cycle	Maintenance
Door	Trained personnel	Every 2 years	See service instructions
Sealing ring	Trained personnel	Annually	See service instructions
Fuse	Trained personnel	If the fuse has tripped.	Operate the fuse lever on the rear of the autoclave.
Safety valve	Trained personnel	Annually	See service instructions

14 Transport and Storage




Always keep the transport crate in which the autoclave was delivered.

14.1 Preparations

Allow the autoclave to cool down and pull the mains plug out of the socket.

14.2 Emptying the water containers

Please empty all internal water tanks: Connect the wastewater hose to the wastewater port for the fresh and used water. Activate the drain by pulling on the connection port. Press the connection port back towards the autoclave to stop the flow.

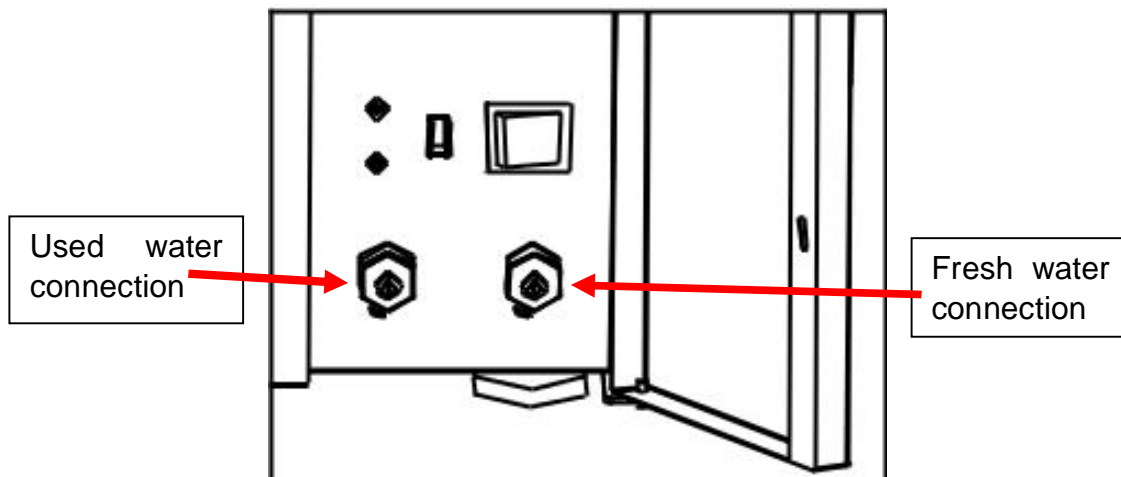

14.3 Transport conditions

Please contact your dealer.



Do not drop the autoclave during transport

14.4 Storage conditions


The autoclave should be stored in a dry place at a room temperature between 5°C to maximum 40°C. The relative air humidity must not be higher than 80%. Make sure that the store is well ventilated and does not contain any corrosive gases.

15 Technical Data

Supply voltage	230VAC, 50Hz, 2300VA
Maximum operating pressure	0.21 □ 0.23 MPa
Maximum operating temperature	135 □ 137 °C
Capacity of water tank	4 l
Device dimensions (W*H*D)	640*560*840 mm
Chamber dimensions	Ø 319, T 617 mm
Chamber material	Stainless steel 304L
Net weight	120 kg
Ambient temperature	5 - 40 °C
Relative air humidity	≤ 80 %
Air pressure	> 70 kPa
Noise level	< 70 dB

16 Appendix

16.1 Appendix 1: Structural diagram

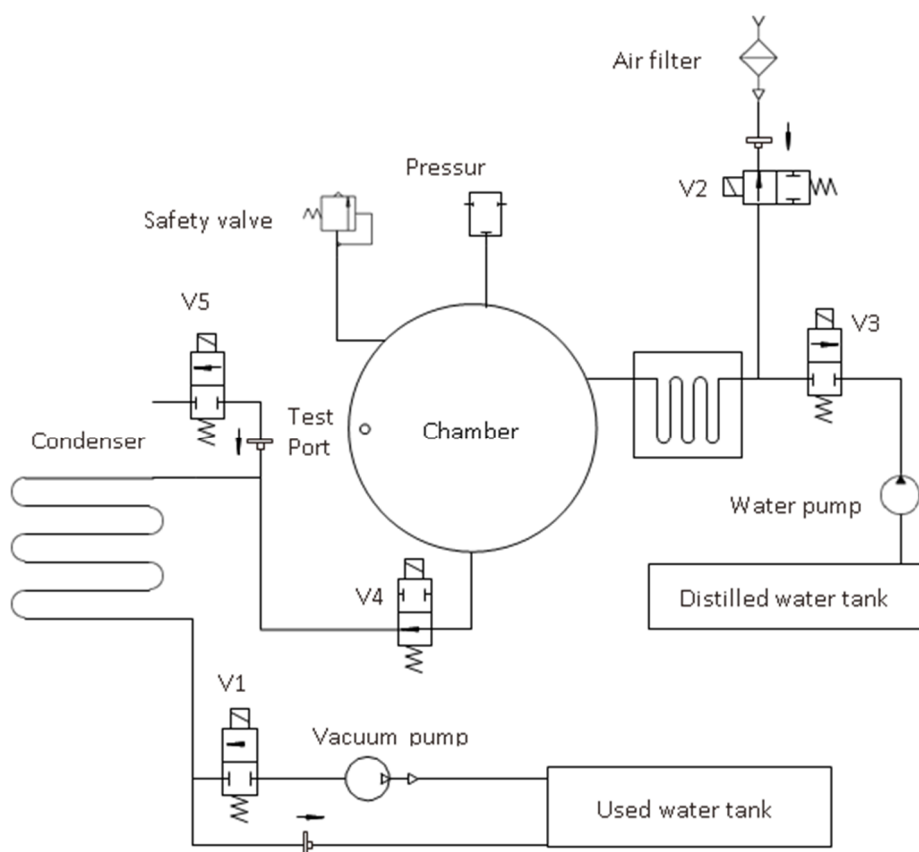

V1: Vacuum pump valve


V2: Air filter valve

V3: Pump valve

V4: Water release valve

V5: Vacuum pump start valve

16.2 Appendix 2: Circuit diagram

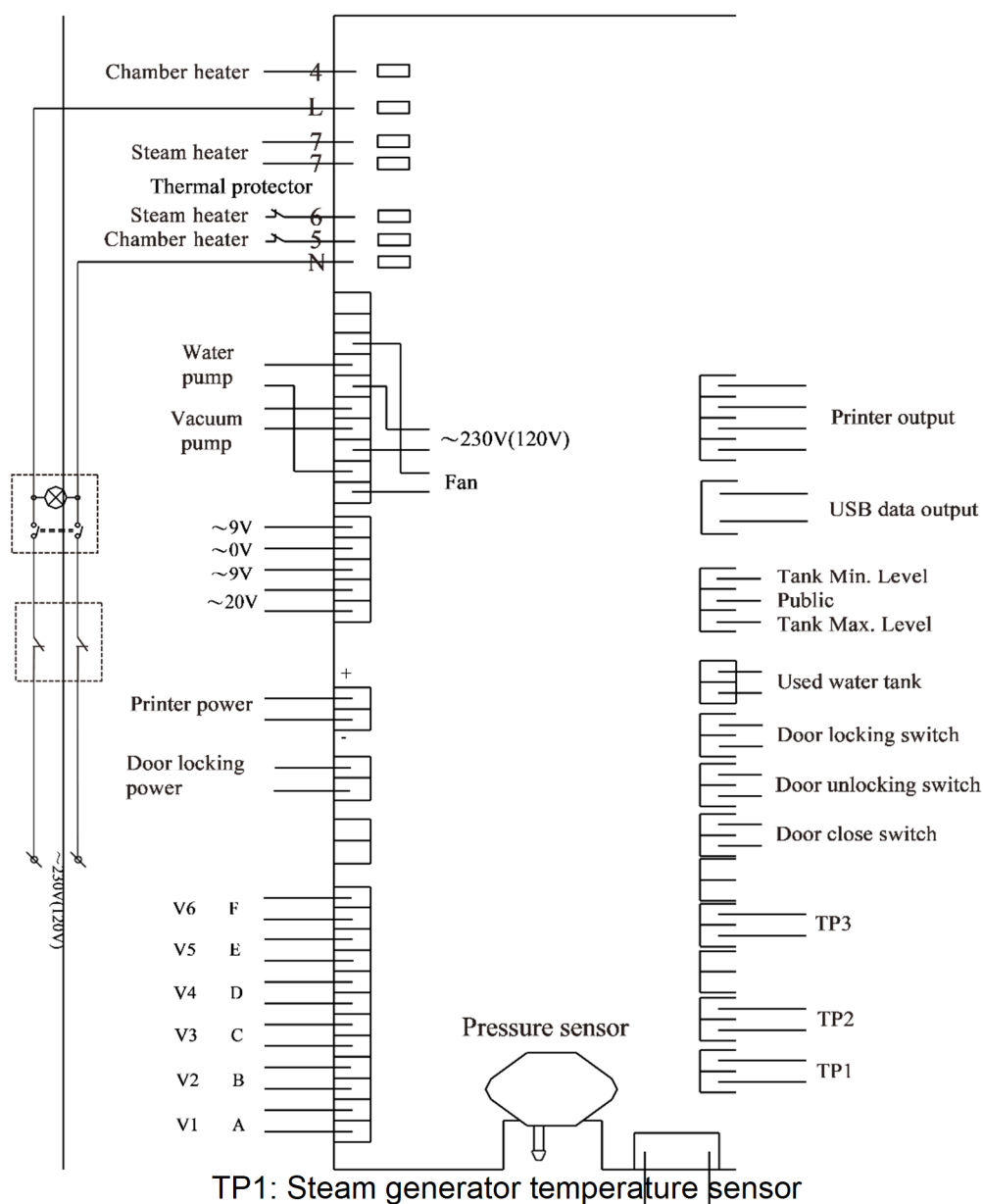

TP1: Steam generator temperature sensor

TP2: Inner temperature sensor 1

TP3: Temperature sensor of chamber wall

V1: Vacuum pump valve

V2: Air filter valve

V3: Pump valve

V4: Water release valve

V5: Vacuum pump start valve

V6: Vacuum pump fan

16.3 Appendix 3: CE Declaration of Conformity for Autoclave

EU/EG Declaration of Conformity

The manufacturer

Ningbo Ican Machines Co., Ltd.
No.77,Yunlin East Road, Gulin Town, Haishu District, Ningbo, China

hereby declares, under its sole responsibility, that the product described below is

Designation: Steam Sterilizer
Property: vessel, piping, assembly, agitator, compressor, according to
EN-2014/68/EU
Model/Type: CB-45T

All relevant harmonization regulations of the following product directives of the European Union fulfilled: (select all applicable product guidelines)

	guideline	Description of the
<input checked="" type="checkbox"/>	2014/68/EU	Directive on the availability of pressure equipment on the market
<input checked="" type="checkbox"/>	2014/30/EU	Electromagnetic Compatibility Directive
<input checked="" type="checkbox"/>	2014/35/EU	Directive on the making available on the market of electrical equipment for use within certain voltage limits

Harmonised standards and technical specifications:

- EN 13480, AD2000, EN ISO 12100:2010, EN764-1:2015+A1L2016, EN 764-5:2014, EN 764-7:2002+AC:2006

Specific information concerning the Pressure Equipment Directive 2014/68/EU

Max. permissible pressure PS: 2.8 bar
Max. permissible temperature TS: 141 °C
Volume: 45 L
Fluid: Steam
Fluid group: Water
Test pressure PT: 3.5 bar

Conformity assessment procedure applied:
Category: II Module: E1

EU Certificate of Conformity

No. 0B170816.NIMDW08

Notified body EU identification number 1282
Ente Certificazione Macchine Srl
Via Ca' Bella, 243 – Loc. Castello di Serravalle – 40053 Valsamoggia (BO) – ITALY

Specific data according to EMC directive 2014/30/EU:

The product has been placed on the market according to article § 14 a), internal production control according to Annex II of the EMC Directive.

Spain, 29.05.2019
Place / Date (day/month/year)


Management Representative

Page: 1 of 1

Template: Declaration of Conformity DG MS NSP EMV
ATEX

Template ID:

creator: customer

Rev.no.: 1

Date: 19.05.2017

16.4 Appendix 4: CE Declaration of Conformity for Safety Valve

NUOVA GENERAL INSTRUMENTS s.r.l. <small>LOC. CAMPASSO - 29010 PIANELLO VALTIDONE (PC) Italy - Tel. +39 0523994629 - Fax +39 0523997219 http://www.nuovageneral.it - E-mail: info@nuovageneral.it - P.IVA 01337040339 - C.F. 01317650065 REA-PC 117793 - Reg. Imp. PC 01317650065 - Capitale Sociale € 100.000,00 i.v.</small>		N. from 018023901 to 018026900	
Tipo: D7/A OT (Typ.): Dis. N° 041-1 Rev. 20/05/08			
DICHIARAZIONE DI CONFORMITÀ Allegato IV D. E. 2014/68/UE DECLARATION OF CONFORMITY Annex IV, E. D. 2014/68/EU KONFORMITÄTSEKHLARUNG Anlage IV Richtlinie 2014/68/UE DECLARATION DE CONFORMITÉ Annexe IV D. E. 2014/68/UE		VALVOLA DI SICUREZZA -- SAFETY VALVE SICHERHEITSVENTIL -- SOUPAPE DE SURETE	
PROGETTO, COSTRUZIONE E COLLAUDO CONFORME ALLA DIRETTIVA 2014/68/UE - 15/05/14 CATEGORIA IV Modulo H1 DESIGN, MANUFACTURE AND TEST ACCORDING TO 2014/68/EU DIRECTIVE - 15/05/14 CLASS IV Module H1 PROJEKT, KONSTRUKTION UND ABNAHMEPRÜFUNG GEMÄß EG-RICHTLINIE 2014/68/EU - 15/05/14 KATEGORIE IV Formblatt H1 PROJET, FABRICATION ET VERIFICATION CONFORME A LA DIRECTIVE 2014/68/UE-15/05/14 CATEGORIE IV Module H1			
Garanzia qualità totale Nr.: Garantie gesamte Qualität Nr.:		Full quality assurance system Nr.: Garantie qualité totale Nr.: Module H1: DGR-0036-QS-1180-15	
Norme e spec. utilizzate: (Codes or spec.ref.): (Angewandte normen und spezifikationen): (Reglements et spec. utilise): AD-Merkblatt A2 - VdTUV-Merkblatt Sicherheitsventil 100 / Racc. "E" ed.1979 / ISO 4126 - 1 / AS1271-2003 / EN-378 / EN13136			
Organismo notificato: (Notified body): (Meldestelle): (Organisme agree):		TÜV SÜD Industrie Service GmbH Westendstraße 199 D 80686 Munich 0036	
CE			
- Marchio del costruttore: (Manufacturer's Mark): (Herstellerzeichen): (Marque du constructeur)			
G			
- P.N.: 60 (Nenndruck): - h / do: 0,28571			
- D.N.: (Nominal Diameter): (Anschlussgewinde): (D.N.):			
- Entrata: (In): (Eintritt): (Entre):			
G.1/4" ISO 228			
- Uscita: (Out): (Austritt): (Sortie):			
- Area: (Area): (Offnung): (Surface):			
0,385 cm²			
- Limiti di temperatura di esercizio dei materiali utilizzati a pressione atmosferica: (Operating temperature limits for materials used at atmospheric pressure): (Betriebstemperaturgrenze der unter Luftdruck verwendeten Materialien): (Limites de température de fonctionnement des matériaux utilisés sous pression atmosphérique):			
-20 / +200 °C			
- Coefficiente di efflusso ridotto: (Reduced flow coefficient): (Reduzierte Ausflussziffer): (Coefficient d'écoulement réduit)			
Kd 0,90			
- Pressione di taratura: (Setting Pressure): (Abgleichungsdruck): (Pression de tarage):			
2,6 bar			
- Sovrappressione: (Overpressure): (Über-Unterdruck): (Surpression):			
10%			
- Scarto di chiusura: (Closing variation): (Unterdruck): (Ecart de fermeture):			
15%			
TIPO FLUIDO: (FLUID TYPE): (MEDIUM): (Type de fluide): ARIA-AIR-LUFT			
- Scarico in atmosfera: (Discharge in Atmosphere): (Abfluß gegen Atmosphäre): (Decharge en atmosphere):			
105,87 kg/h 81,9 Nm³ / h 1364,92 lt/min.			
TIPO FLUIDO: (FLUID TYPE): (MEDIUM): (Type de fluide): VAPORE SATURO-STEAM-DAMPF			
- Scarico in atmosfera: (Discharge in Atmosphere): (Abfluß gegen Atmosphäre): (Decharge en atmosphere):			
64,12 kg/h 0 Nm³ / h 0 lt/min.			
TIPO FLUIDO: (FLUID TYPE): (MEDIUM): (Type de fluide):			
- Scarico in atmosfera: (Discharge in Atmosphere): (Abfluß gegen Atmosphäre): (Decharge en atmosphere):			
kg/h Nm³ / h lt/min.			
1- UNI EN12164 CW614N R360 3- UNI EN12164 CW614N R360 5- UNI EN12164 CW614N 8- UNI EN12164 CW614N 9- UNI EN 10270-1 SM 10 VITON (-20 / +200 °C)			
Anno costruzione: (Year of manufacture): (Baujahr): (Année de fabrication): 2018			
We declare that this valve has passed successfully the test, and the hydraulic test of the body to 1,5 x P.N. Hiermit bestätigen wir, dass das Ventil die Abschlusskontrolle einschließlich der hydraulischen Gehäuseprüfung mit einem Druck von 1,5 x Nenndruck bestanden hat. On declare que la soupape a été soumise avec résultat satisfaisant à la control final y-compris l'épreuve hydraulique du corps a 1,5xP.N.			
EN 10204 3.1 Cod: S.V. Code:			
Date: 02/01/2018			

Nr.Ord.: E-MAIL

S.V. Code:

Date: 02/01/2018

NUOVA GENERAL INSTRUMENTS SRL
 Alberto Patrizia (CEO)
 Valerio

Orders

Dealer verification

Instruction manuals

Technical Support



Certoclav Sterilizer GmbH
Peintner Straße 10
A-4060 Leonding



Tel: +43 732 674 278

E-Mail: support@certoclav.com

Website: www.certoclav.com