
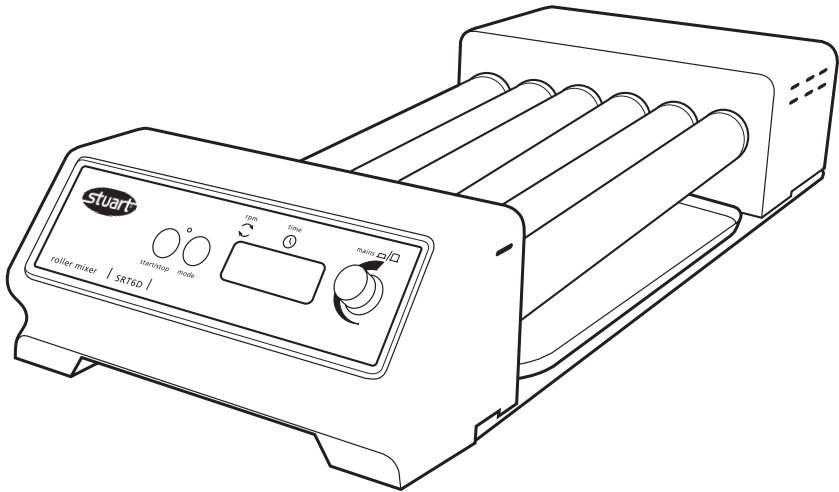
# Roller Mixers

SRT6 SRT6D SRT9 SRT9D


Instructions for use  
Mode d'emploi  
Istruzioni per l'uso  
Instrucciones de uso  
Bedienungsanleitung
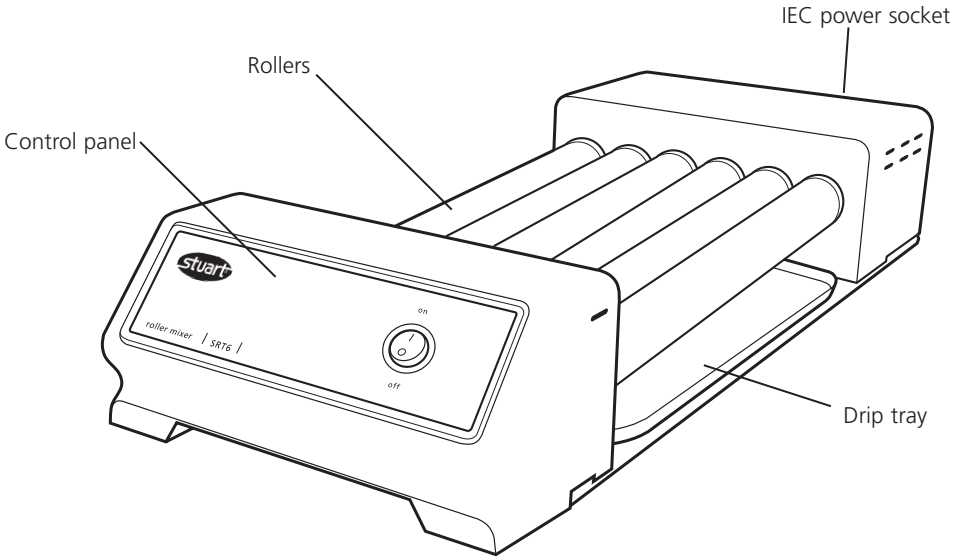
# English

**SRT6 & SRT9**

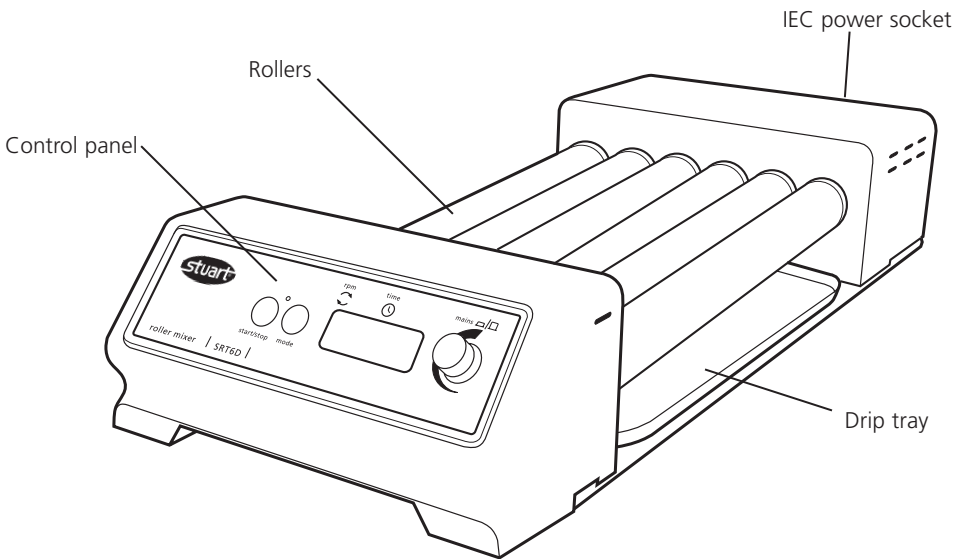**SRT6D & SRT9D**

# Roller Mixers

SRT6 SRT6D SRT9 SRT9D

**Thank you for purchasing this piece of Stuart equipment. To get the best performance from the equipment, and for your own safety, please read these instructions carefully before use.**

This equipment is designed to operate under the following conditions:-

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +4°C to +60°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuation not exceeding 10%
- ❖ Overvoltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664

If the equipment is not used in the manner described in this manual the protection provided by the equipment may be impaired.

## Electrical Installation

**THIS EQUIPMENT MUST BE EARTHED**

**Before connection please ensure that the line supply corresponds to that shown on the rating plate. This model requires a supply rated at 220 - 240V, 50 Hz, ~, single phase.**

### Power requirements

Model	Wattage
SRT6 & 9	50W

There is an IEC socket at the rear of the instrument for connection to the mains supply. The unit is supplied with two mains leads fitted with IEC plugs for connection to the instrument. One has a UK 3 pin plug and the other has a 2 pin "Shuko" plug for connection to the mains. Choose the lead appropriate for your electrical installation and discard the other. Should neither lead be suitable, take the lead with the UK plug and replace the plug with a suitable alternative.

See the enclosed instruction sheet for advice on how to carry out this procedure. Should the mains lead require replacement a cable of 1mm<sup>2</sup> of harmonised code H05W-F connected to an IEC 320 plug should be used.

### IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN

The mains lead should be connected to the instrument BEFORE connection to the mains supply.

## General Description

The Stuart Roller Mixers provide a gentle, but highly efficient, rocking and rolling action. On the SRT6 & 9 the rollers rotate at a fixed speed of 33rpm and rise and fall 16mm. The SRT6D & 9D have adjustable speed control between 5 and 60rpm, and an adjustable digital timer that automatically stops rotation after the set time (1 to 9999 minutes, 1 second to 90 minutes or 1 minute to 9 hours).

Most sizes of tube, Bijoux, Universals and other bottles can be accommodated

**SRT6 & 6D:** Have 6 rollers

**SRT9 & 9D:** Have 9 rollers

**Available separately** are stacking blocks to allow up to three units to be stacked vertically, optimizing bench space. Order: SRT/STACK Stacking system for rollers (4 x stack blocks).

## Safety Advice Before Use

- ❖ The unit should be carried using both hands with fingers under each side frame.
- ❖ Never move or carry the unit when in use or connected to the mains electricity supply.
- ❖ Do not mix flammable liquids or use the equipment in hazardous atmospheres.
- ❖ In case of mains interrupting, a fault or mechanical failure, the unit will continue to operate on removal of fault.
- ❖ Mechanical energy can lead to breakage of glass vessels. Use with care.

## Operation SRT6 & 9

Position the apparatus on a firm level surface and connect to the mains electricity supply. Place the samples to be mixed between the rollers. Once all the samples have been set, switch the Roller Mixer ON using the Mains On/Off switch.

## Operation SRT6D & 9D

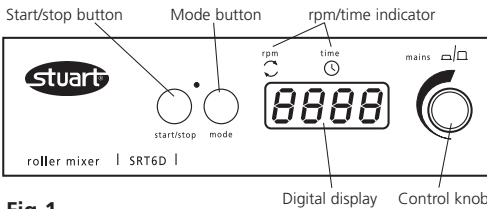
Fig 1

Position the apparatus on a firm level surface and connect to the mains electricity supply. Place the samples to be mixed between the rollers. Once all the samples have been set, switch the Roller Mixer ON by pressing the control knob (see fig 1). The unit will carry out a self test routine and then the display will show the last stored set speed (red dot in the display shows RPM is selected). This can be adjusted by turning the control knob. Once the correct speed is displayed press the start / stop button to begin motion. The speed can be adjusted without halting the unit. The unit can be halted at any time by pressing the start / stop button.

### Setting the timer.

In order to use the timer press the mode button. The red dot on the display moves over to time. The display will show the last stored time in hours and minutes (H is displayed on left hand side of digital display) or seconds and minutes depending on the timer unit selected (see below to change). Select the desired count down time using the control knob (--- indicates continuous, no timer set). When the display shows the correct time in

minutes, press the start / stop button to begin operating. During operation, motion can be paused at any time by pressing the start / stop button. If the button is pressed again, motion will start again and the timer will continue to count down. The current timer or speed can be displayed or adjusted by using the mode button to toggle between the two and altered with the control knob, without halting the unit. When the timer reaches zero, the unit will be automatically halted and an alert will sound. Press the start / stop button to revert back to the saved time and speed values set (see instructions below to adjust saved values) or select new values following instructions given above.

**N.B To disable the timer and revert to continuous rotation, select - - - in the display during timer mode.**

To switch the unit OFF completely, press the control knob in. The display will show OFF.

### Saving speed/time values.

In order to save a commonly used speed/time value, switch the unit ON by pressing the control knob in (see fig 1). Select both a speed value and a time value (either 1 to 9999 minutes, 1 second to 90 minutes or 1 minute to 9 hours or continuous - - -) using the instruction given above. Switch the unit OFF by pressing in the control knob. Wait 5 seconds and switch the unit back ON by pressing the control knob in. The input values are now saved in memory. They can be altered at any time by following the same procedure.

**N.B. When the unit is switched OFF using the control knob, whatever values of speed/time were set prior to turning the unit off will be saved, and will be the values restored when the unit is switched back on.**

### Changing the timer unit

The unit timer is defaulted to minutes only (1 minute to 9999). To change to minutes and hours (1 minute to 9 hours, H is displayed in the left digit of the timer display), or to change to seconds and minutes (1 second to 90 minutes) turn the unit off using the control knob then switch the mains power off at the plug. Hold down both the start/stop button and the mode button whilst simultaneously switching the mains

power back on. The display will temporarily show 88:88 then OFF. Using the control knob turn the unit on, the display will show nn:ss for minutes and seconds, H:nn for hours and minutes and nn for minutes, turn the control knob to select the desired timer unit and turn the unit off by the control knob, OFF will be displayed. When the unit is next used the timer will have been altered to the set unit.

## Maintenance & Servicing

**WARNING:** Ensure the unit is disconnected from the mains electricity supply before attempting maintenance or servicing.

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh or abrasive cleaning agents.

**Any repairs or replacement of parts MUST be undertaken by suitably qualified personnel.**

The following spares and accessories are available from your laboratory dealer.

Description	Catalogue Number
Stacking system	SRT/STACK

For a comprehensive list of parts required by service engineers conducting internal repairs and a service manual, please contact the Technical Service Department of Bibby Scientific Ltd. quoting both the model and serial number.

Only spare parts supplied or specified by Bibby Scientific Ltd. or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed into the instrument.

If in any doubt, please contact the Technical Service Department of Bibby Scientific Ltd. or the point of sale.

Bibby Scientific Ltd.  
Stone, Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0)1785 812121  
Fax: +44 (0)1785 813748  
e-mail [info@bibby-scientific.com](mailto:info@bibby-scientific.com)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)

## Warranty

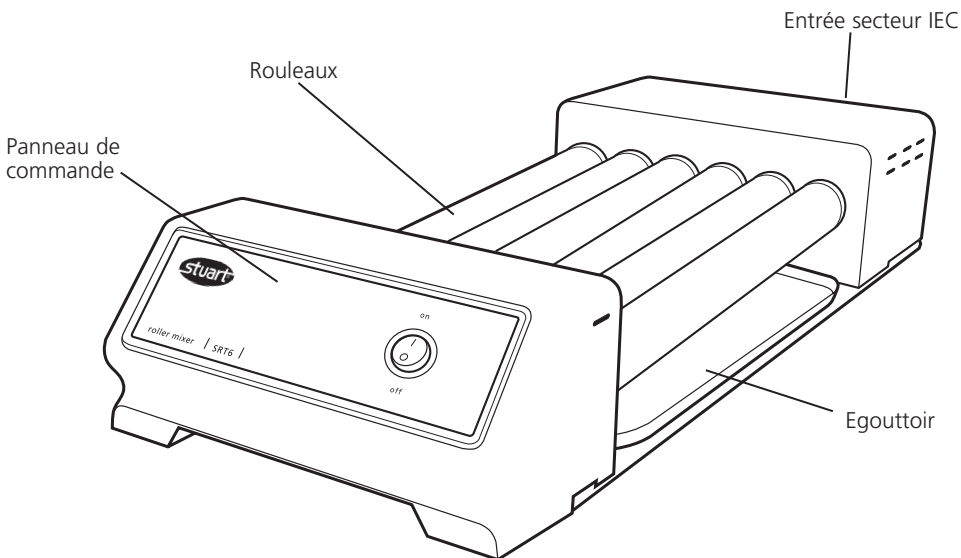
Bibby Scientific Ltd warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of three (3) years. In the event of a justified claim Bibby Scientific will replace any defective component or replace the unit free of charge. This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage caused by incorrect installation, adaptation, modification, fitting of non approved parts or repair by unauthorised personnel.

# Technical Specification


	<b>SRT6</b>	<b>SRT6D</b>
Number of rollers	6	6
Speed	33rpm	5 to 60rpm
Amplitude	16mm	16mm
Maximum load	5kg	5kg
Controls	Analogue	Digital
Timer	No	Yes
Roller size (length x diameter)	340x30mm	340x30mm
Overall dimensions (w x d x h)	565x240x110mm	565x240x110mm
Net weight	5.1kg	5.1kg
Electrical Supply	230V, 50Hz, 50W	230V,50Hz, 50W

	<b>SRT9</b>	<b>SRT9D</b>
Number of rollers	9	9
Speed	33rpm	5 to 60rpm
Amplitude	16mm	16mm
Maximum load	5kg	5kg
Controls	Analogue	Digital
Timer	No	Yes
Roller size (length x diameter)	340x30mm	340x30mm
Overall dimensions (w x d x h)	565x360x110mm	565x360x110mm
Net weight	6.9kg	6.9kg
Electrical Supply	230V, 50Hz, 50W	230V,50Hz, 50W

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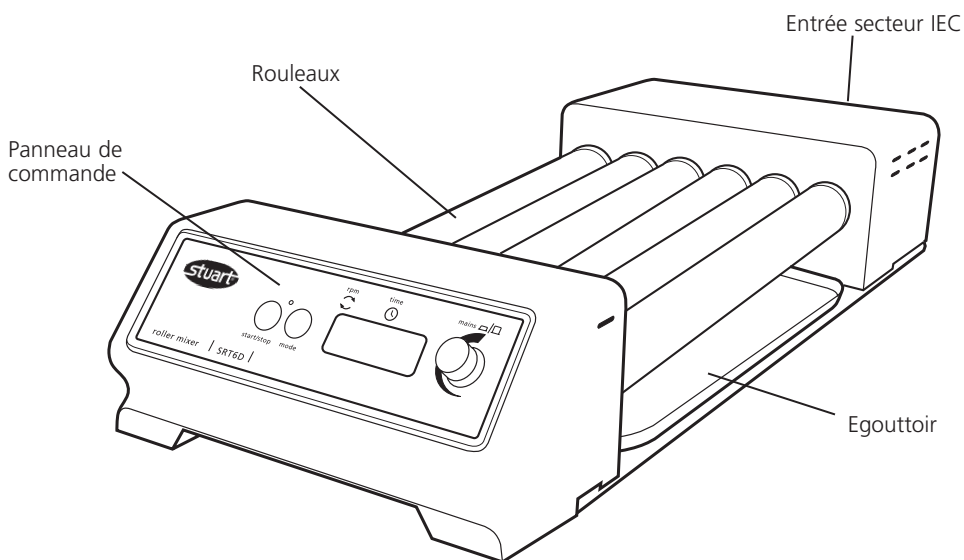These products meet the relevant EC harmonised standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in their vicinity will meet these

standards and we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, please contact the Technical Service Department of Bibby Scientific Ltd.

## Declaration of Conformity

### Stuart Roller Mixers, Models SRT6, SRT6D, SRT9 & SRT9D

These products comply with the requirements of the EU Directives listed below:

**89/336/EEC** Electromagnetic Compatibility Directive amended by 93/68/EEC.  
**73/23/EEC** Low Voltage Directive amended by 93/68/EEC.

Compliance with the requirements of these Directives is claimed by meeting the following standards

EN 61326: 1997 + Amendments (Electrical Equipment for Measurement, Control and Laboratory use)  
EN 61010-1: 2001 (Safety Requirements for Electrical Equipment for Measurement, Control and Laboratory use)  
EN 61010-2-051: 2003 (Particular Requirements for Laboratory Equipment for Mixing and Stirring)

#### Compliance Certificates and Equipment Specification.

The full product specifications, listed in report numbers:

RETS(E)1719/A/1  
RETS(L)1719/A/2  
RETS(L)1719/A/3

from Epsilon Technical Services showing compliance to the above standards, are available on request

#### CE mark affixed '06.

Signed:  (Mr D. E. Hicks)

Date: 21<sup>st</sup> March 2006

Authority: Technical/Development Manager  
Barloworld Scientific Ltd

Barloworld Scientific Ltd - Stone - Staffs - ST15 0SA - UK  
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

# INSPECTION REPORT

**MODEL** \_\_\_\_\_

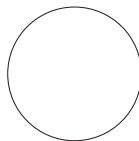
## ELECTRICAL SAFETY

1. Earth continuity
2. Insulation
3. Flash test

## FUNCTIONAL

1. Indicators
2. Roller control
3. Visual acceptance

**QUALITY CONTROL INSPECTOR**

**Bibby Scientific**

### **UK**

Bibby Scientific Ltd.  
Beacon Road, Stone  
Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0)1785 812121  
Fax: +44 (0)1785 810405  
e-mail: sales@bibby-scientific.com  
www.bibby-scientific.com

### **North & South America**

Bibby Scientific US Inc.  
t/a Techne Inc.  
3 Terri Lane, Suite 10  
Burlington, NJ 08016 USA  
Toll Free (in NA): 800-225-9243  
Tel: +1 609 589 2560  
Fax: +1 609-589-2571  
e-mail: labproducts@techneusa.com  
www.techneusa.com

### **Middle East**

Bibby Scientific Middle East Ltd.  
PO Box 27842, Engomi 2433  
Nicosia  
Cyprus  
Tel: +357 22 660 423  
Fax: +357 22 660 424  
e-mail: sales@bibby-scientificme.com

### **France**

Bibby Scientific Limited  
Bâtiment Le Deltaparc Parc Silic PN2  
7 rue du Canal  
BP 55437 VILLEPINTE  
95944 ROISSY Charles de Gaulle  
France  
Tel: +33(0)148 63 78 03  
Fax: +33(0)148 63 78 01  
e-mail: ventes@bibby-scientific.com  
www.bibby-scientific.com

### **Italy**

Bibby Scientific Italia Srl  
Via Alcide de Gasperi 56  
20077 Riozzo di Cerro al Lambro  
Milano Italia  
Tel: +39 (0)2 98230679  
Fax: +39 (0)2 98230211  
e-mail: marketing@bibby-scientific.it  
www.bibby-scientific.it

### **Asia**

Bibby Scientific - Singapore  
Prudential Tower, Level 26  
30 Cecil Street  
Singapore 049712  
Tel: +65 6631 2976  
Fax: +44 (0)1785 810405  
e-mail: info@bibby-scientific.com  
www.bibby-scientific.com