

HI 9811-5


pH/EC/TDS/°C meter

Quick Reference Guide

TAKING MEASUREMENTS

- Connect the HI 1285-5 probe and switch the meter ON.
- Submerge the probe into the sample to be tested.
- Select the desired measurement range (pH, $\mu\text{S}/\text{cm}$, mg/L, $^{\circ}\text{C}$) by pressing the correspondent key.

Note: If the display shows only an "E" on the right side of the LCD, the reading is out of range.

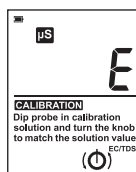

- Wait for a couple of minutes for reaching the thermal equilibrium before taking any measurements.
- After use switch off the instrument, clean and cover the probe with the protective cap.

Note: It is recommended to clean often the probe with HI 700661 Cleaning Solution.

BATTERY REPLACEMENT

When the battery level is so low that it may cause unreliable measurements, the meter turns off.

It is recommended to replace the battery immediately.

Unscrew the three screws on the rear of the meter, remove the battery compartment cover and replace the 9V alkaline battery with a new one while observing its polarity.

Make sure the battery contacts are tight and secure before replacing the cover.

 **HANNA**[®]
instruments
www.hannainst.com