Multi Reax

Betriebsanleitung
Instruction Manual
Mode d'Emploi
Instrucciones de
Empleo
Istruzioni per l'uso

D

DEUTSCH

Seite 3 - 13

E

ENGLISH

page 14 - 24

F

FRANCAISE

page 25 - 35

ES

ESPAÑOL

página 36 - 47

I

ITALIANO

Pagina 48 - 59

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EN 61010-1 + EN 61010-2-051

Thank you for purchasing a Heidolph Instruments product. This item has been designed, made and inspected in compliance with DIN EN ISO 61010 for long-term and flawless operation.

SUMMARY

SUMMARY 14

STANDARD HARDWARE & OPTIONS..... 15

GENERAL..... 15

SAFETY INFORMATION 16

SET-UP 17

1. Machine set-up 17

2. Electric hook-up 17

3. Secure containers or tubes in carousel attachment..... 17

4. Exchange carousel attachment..... 19

OPERATION AND CONTROLS..... 20

A. Continuous shaking..... 20

B. Timer controlled shaking..... 20

C. Repetition..... 21

D. Edit time during shaking operation..... 21

CLEANING & SERVICING..... 21

DISASSEMBLY & STORAGE..... 21

DISPOSAL 22

MALFUNCTION & REMEDY..... 22

SPECIFICATIONS..... 22

WARRANTY, LIABILITY & COPYRIGHT..... 23

FAQ / REPAIR WORK..... 23

CE-DECLARATION OF CONFORMITY..... 24

Important information

Advise about power cord / mains supply

Caution: mandatory action

Caution: fire- and explosion hazard

Advise about maintenance / repair

STANDARD HARDWARE & OPTIONS

	product	quantity	P/N 230/240V 50/60Hz	P/N 115V 50/60Hz
	Multi Reax	1	545-10000-00	545-10000-01
	Instruction Manual	1	01-005-004-28	01-005-004-28
	Power cord	1	14-007-003-81	14-007-003-89

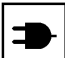
Standard options (shipped with item)

	item	P/N
	carousel attachment for 26 test tubes, 10 - 16mmØ	549-01600-00-0
	carousel attachment for 12 test tubes, 16 - 32mmØ	549-03000-00-0
	rubber ring for carousel attachment for 26 test tubes	23-08-06-01-75-0
	rubber ring for carousel attachment for 12 test tubes	23-08-06-01-76-0

GENERAL

-  Unpack your item carefully. Inspect for damage and report such damage or missing parts to your supplier right away.
-  Read your Instruction Manual carefully. Take time to save time while working with your product. Make sure that every user has read and understood the Instruction Manual.
-  Please store the Instruction Manual in a place easily accessible to every user.

IF ALL ELSE FAILS, READ THESE INSTRUCTIONS !

-  A so-called EURO-plug (DIN 49441 CEE 7/VII 10/ 16 A 250 V) is standard on all of the products.
For the Continental US they feature a US-standard plug (NEMA Pub.No.WDI.1961 ASA C 73.1 . 1961 page 8 15A 125V).
-  For using the item in a country with deviating outlet / plug systems, we recommend to use approved adapters or to have an electrician replace the standard plug with one suiting your needs.
-  As shipped, the item features a protective ground wire. When replacing the original plug, make sure to reconnect this protective ground wire in the new plug !

SAFETY INFORMATION

Please comply with all safety and accident-prevention regulations as in force for laboratory work!

Use extra care when working with flammable substances; refer to safety data sheets.

Use extra care when working in the vicinity of flammable and explosive substances. Motors are of non-sparking type, the item itself however is not explosion-proof.

When connecting your item with your local power supply, please make sure your item is wired for your local voltage; go by the data plate on the item.

Please connect your unit with a protective-ground outlet only.

Turn your power switch OFF whenever the item is not used, or before disconnecting the plug.

Repair work is limited to technicians approved by Heidolph Instruments.

Your item needs a solid stand.

Lab bench needs to be of rigid design, and features an anti-skid surface coat.

Before starting the item, make sure all vessels are attached safely (must not move while shaking).

SET-UP

1. Machine set-up

Please locate the shaker on a stable, horizontal surface. For reasons of safety, keep the area around the machine clear of other items.

2. Electric hook-up

Use the power cord from your hardware bag and connect it with the plug connector in the item's rear panel.

3. Secure containers or tubes in carousel attachment

Just plug containers or tubes into retaining pockets against preload of a rubber ring. This rubber ring is going to safely secure the tubes in the retainer pocket.

To ease installation, lift rubber ring by hand before introducing containers.



Please make sure the rubber ring is undamaged. If not, containers cannot be seated properly. Change if necessary.



Refer to recommendations below when securing containers:

Locate containers concentric with retainer plate all the time !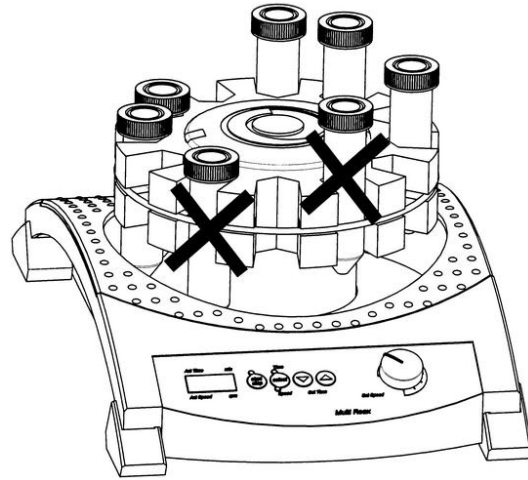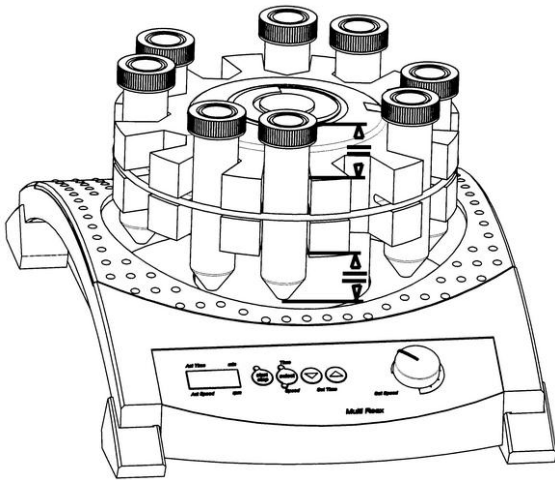

Never install different sizes of containers in neighbored pockets. If not, the smaller diameter container will not be secured properly !

When using different sizes of containers, form groups of items with same diameter and leave one pocket between groups void.

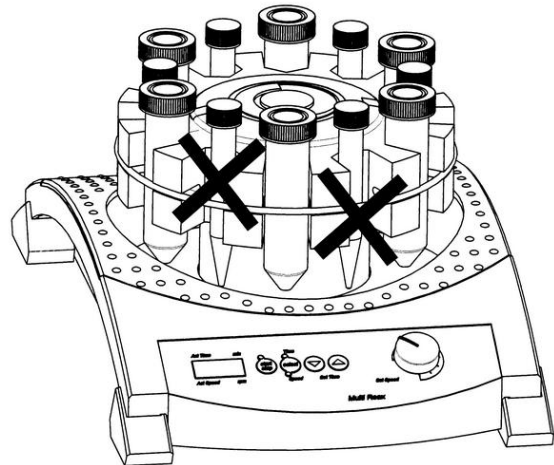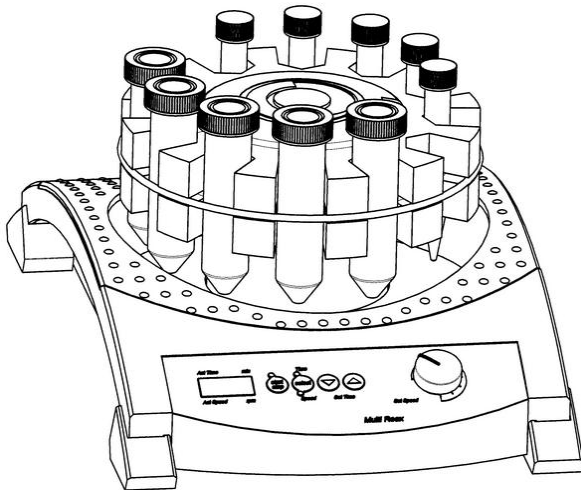Before starting shaking action, make sure all vessels are secured in place properly.

Before starting shaking action, make sure, all containers are oscillating free and cannot hit the machine housing.

The carousel attachment for 26 test tubes accepts round containers of 10 - 16 mm dia. and a max. length of 160 mm only; the carousel attachment for 12 test tubes holds round containers of 16 - 32 mm dia. and a max. length of 120 mm only.

As a function of load applied on the shaker holders, certain points of resonance may be found in a speed range from 1,200 to 1,500 rpm. If such resonance is found (strange vibrations of the shaker), please increase or decrease speed by a minor amount as to get out of the critical frequency range.

4. Exchange carousel attachment


The item arrives with the carousel attachment for 26 test tubes installed. You may replace it with the carousel attachment for 12 test tubes at ease. Remove all containers before installing the retainers. Remove carousel attachment by pulling with both hands, plug new carousel attachment in place.

Check for location of the two "noses". Carefully slip carousel attachment over these "noses". Be sure not to damage the part. These "noses" secure the carousel on the shaker.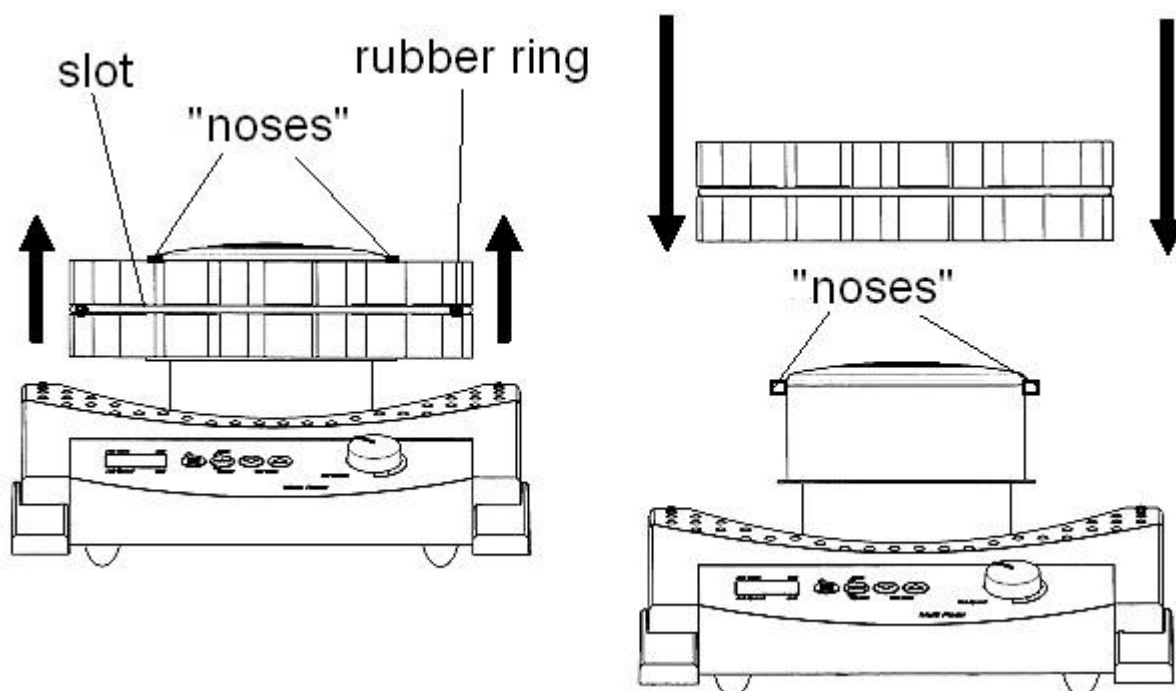

Check rubber ring for proper seat. This ring secures your containers in place. Seat ring in slot of carousel on condition.

OPERATION AND CONTROLS

-  Before connecting power cord with mains outlet, make sure that:
 - your item is wired for your local voltage and frequency (data plate on item).
 - master switch is set to "0" and all controls are in "min." position (turn CCW completely); this way you avoid spilling fluids by too intense shaking action.


 Carefully close your vessels and select appropriate shaking intensity (if vessels remain open) to avoid splashes and spillage.

-  We recommend to start with low shaking speed and gradually increase frequency to avoid accidental shaking at high speed.

 When handling hazardous fluids, make reference to applicable safety information.

The control panel features the following controls (from right to left):

- A 2-pole master switch (toggle-type, lighted green on the right side of the unit)
- B setting knob, shaking frequency
- C Set Time key \uparrow : time up
- D Set Time key \downarrow : time down
- E Select key: display selector: shaking frequency (Act Speed) / time (Act Time)
- F LED Time: time appears in display
- G LED Speed: rotational speed (shaking frequency) appears in display
- H Start / Stop key: to START / STOP shaking action
- I LED Start /Stop: lighting, while item is running
- J 4-digit display

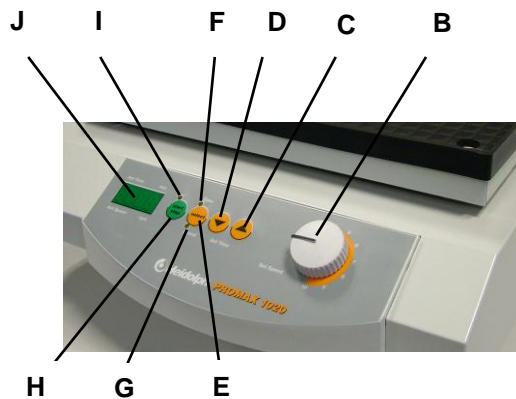A. Continuous shaking

- Turn item ON with master switch (A)
- Set time with \uparrow and \downarrow keys to 000 (depress \uparrow and \downarrow together).
- Start continuous operation by hitting the Start / Stop key. Start and Speed LEDs turn on, shaking frequency is displayed (4-digit display).
- Set shaking frequency with knob (B)
- Hitting Start / Stop key will stop shaking.

B. Timer controlled shaking

- Set time between 0 and 999 minutes with Set Time \uparrow or Set Time \downarrow keys.
Short hits on the keys change time setting slowly, whereas depressing keys is going to

change time rapidly. Depressing the Set Time ↑ and Set Time ↓ keys resets the timer to 000.

- Depressing the Start / Stop key (H) starts shaking at the frequency selected with knob (B) and for the time interval set before. Start LED turns ON.
- Time lapse appears in a 3-digit display. Time LED turns ON.
- Current shaking action may be discontinued with the Start / Stop key (H). Actual time won't change. Hitting this Start / Stop key continues shaking at the moment it had been discontinued.
- In case of a power supply failure or interruption (item turned OFF), this factor is lost.
- Hitting the Select key during timer-controlled operation, you may change display from time to shaking frequency and vice versa. Speed LED turns ON. Timer LED flashes to indicate timer operation mode.
- Timer LED also flashes, when timer-controlled operation was discontinued with the Start / Stop key.

C. Repetition

- Repeating last timer controlled operation does not require new time setting, as long as you did not touch the master switch.

D. Edit time during shaking operation

- Time setting may be changed during shaking, while timer is running and even when timer controlled operation was discontinued.
- Memory timer will keep initial value in its memory.

CLEANING & SERVICING

Cleaning: wipe housing clean with a damp cloth (add some sort of mild liquid soap).

Note

To avoid damage to the surface finish, avoid using chlorine bleach, chlorine-based detergents, abrasive substances, ammonia, rags or cleaning agents containing metal particles.

The item is maintenance-free. Repair work is limited to technicians so approved or appointed by Heidolph Instruments. Please call your local Heidolph Instruments Dealer or a Heidolph Instruments Field Representative.

DISASSEMBLY & STORAGE

Disassembly

1. Turn item OFF and disconnect mains plug.
2. Remove all of the hardware arranged around the shaker to ease disassembly.

Forward & Store

1. We recommend to store the item and its components in its original box, or a similar container that offers adequate protection against damage in transit. Tape the box securely.
2. Store the item in a dry place.

Caution

Do not jolt or shake the item during transport.

DISPOSAL

For disposal, please comply with your local or national regulations.
 Split by metal, plastic, etc.
 Packing material to be treated as described above (material split).

MALFUNCTION & REMEDY

Work on electric, electronic and cryogenic components is limited to qualified personnel.

Master switch on shaker / mixer won't light

1. Check power cord
2. Check circuit breakers

Item won't shake (master switch lighted))

1. Timer run-down
2. Thermal motor circuit breaker triggered by motor overload
 Remedy:
 Wait about 20 minutes, decrease load applied on shaker plate.
3. Mechanical parts broken (humming motor noise) or electronic failure (no motor noise).

SPECIFICATIONS

Model	Multi Reax
Motion	Vibrating
Rotation speed (rpm)	150 – 2000
Rotation speed setting	Digital
Orbit (mm)	3
Operating mode	Timer or continuous
Input power (W)	33
Weight (kg)	9,8
Dimensions (wxhxd) (mm)	270 x 410 x 172
Standard equipment	Carousel attachment for 26 test tubes, 10 - 16mmØ and max. length of 160 mm, or Carousel attachment for 12 test tubes, 16 - 32mmØ and max length of 120 mm
Max. load capacity (kg)	1,5
Supply voltage	230/240V / 50/60 Hz or 115V / 50/60 Hz
Overheating protection	Self-resetting
Protection class (DIN 40050)	IP 30
Permissible ambient temp. (°C)	0 – 50
Permissible relative humidity (%)	80
Designed for continuous use	yes

WARRANTY, LIABILITY & COPYRIGHT

Warranty

Heidolph Instruments warrants that the present product shall be free from defects in material (except wear parts) and workmanship for 3 years from the date shipped off the manufacturer's warehouse.

Damage in transit is excluded from this warranty.

To file for such warranty service, contact Heidolph Instruments (phone ++49-9122-9920-68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation are not covered by this warranty.

Alterations to the present warranty need Heidolph Instruments' consent in writing.

Exclusion Clause

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

Copyright

Copyright in pictures and wording of the present Instruction Manual is held by Heidolph Instruments.

FAQ / REPAIR WORK

If any **aspect** of installation, operation or maintenance remains unanswered in the present Manual, please contact the following address:

For repair services please call Heidolph Instruments (phone: +49 - 9122 - 9920-68) or your local, authorized Heidolph Instruments Dealer.

Note

You will receive approval for sending your defective item to the following address:

Heidolph Instruments GmbH & Co. KG
Lab Equipment Sales
Walpersdorfer Str. 12
D-91126 Schwabach / Germany
phone: ++49-9122-9920-68
Fax: ++49-9122-9920-65
E-Mail: Sales@Heidolph.de

Note

If you are based in the United States of America, please contact Heidolph US:

Heidolph Instruments, LLC
Lab Equipment Sales
2615 River Rd.
Cinnaminson, NJ 08077
Phone: 856-829-6160
Fax: 856-829-7639
E-Mail: heidolph@snip.net

Safety Information

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact *substance*
- take proper *protective* measure to ensure the safety of our receiving and service personnel
- *mark* the pack IAW Hazardous Materials Act

CE-DECLARATION OF CONFORMITY

We herewith declare that the present product complies with the following standards and harmonized documents:

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