20000003968

Vortex 2_032018

IKA

IKA Vortex 2

| | | |
|--------------------------|----|----|
| Betriebsanleitung | DE | 03 |
| Ursprungssprache | | |
| Operating instructions | EN | 08 |
| Mode d'emploi | FR | 13 |
| Руководство пользователя | RU | 18 |
| 使用说明 | ZH | 23 |
| 取扱説明書 | JA | 28 |
| 사용 설명서 | KO | 33 |

| | | |
|---------------------------|----|----|
| Indicaciones de seguridad | ES | 38 |
| Veiligheidsinstructies | NL | 39 |
| Norme di sicurezza | IT | 40 |
| Säkerhetsanvisningar | SV | 41 |
| Sikkerhedshenvisninger | DA | 42 |
| Sikkerhetsanvisninger | NO | 43 |
| Turvallisuusohjeet | FI | 44 |
| Normas de segurança | PT | 45 |
| Wskazówki bezpieczeństwa | PL | 46 |
| Bezpečnostní pokyny | CS | 48 |

| | | |
|---------------------------|----|----|
| Biztonsági utasítások | HU | 49 |
| Varnostna opozorila | SL | 50 |
| Bezpečnostné pokyny | SK | 51 |
| Ohutusjuhised | ET | 53 |
| Drošības norādes | LV | 54 |
| Nurodymai dėl saugumo | LT | 55 |
| Инструкции за безопасност | BG | 56 |
| Indicații de siguranță | RO | 57 |
| Υποδείξεις ασφάλειας | EL | 59 |

IKA-Werke, Germany
Reg. No. 004343

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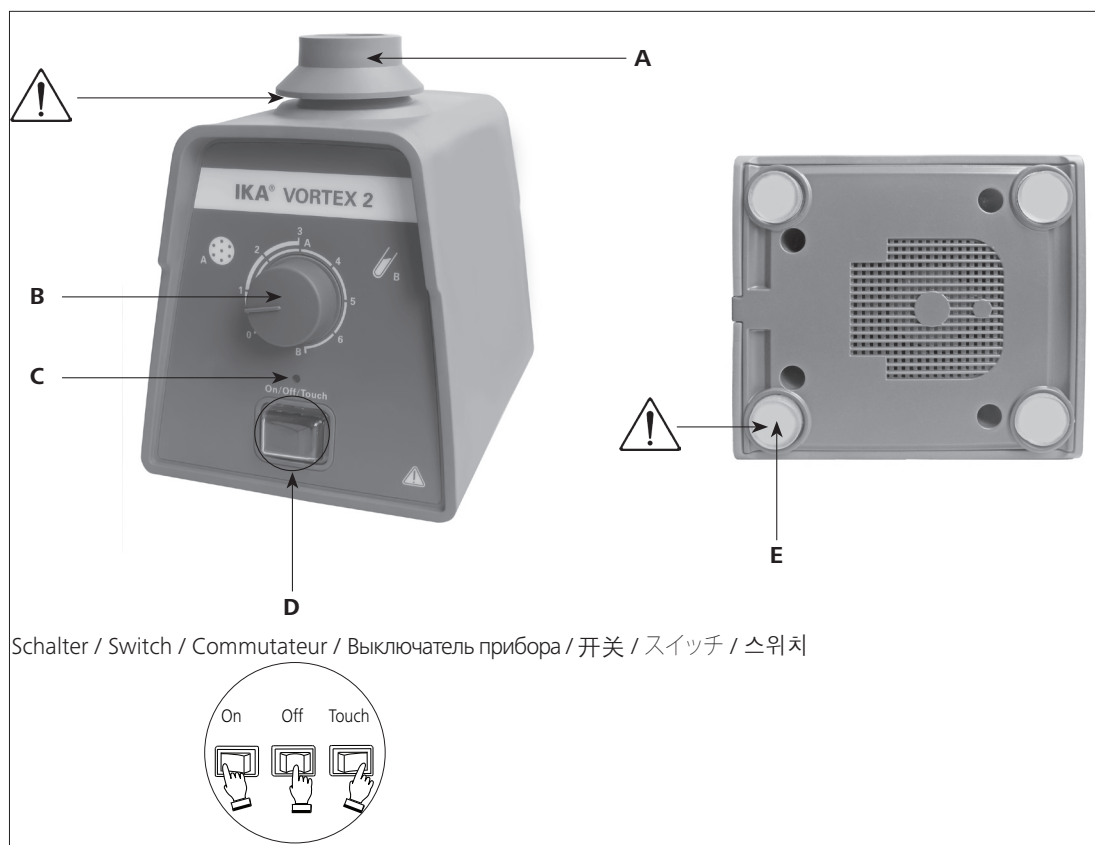
Fig. 1

| Pos. Bezeichnung | Item Designation | Pos. Désignation | Пункт описание |
|------------------------------------|--|-------------------------------------|--|
| A Standardaufsatz VG 3.1 | A Standard attachment VG 3.1 | A Support standard VG 3.1 | A Стандартное крепление VG 3.1 |
| B Drehknopf | B Rotating knob | B Bouton rotatif | B ручку вращения |
| C Kontrollleuchte | C Control light | C Lumière côntrol | C Индикаторная лампа |
| D Schalter | D Switch | D Commutateur | D Выключатель прибора |
| E Gerätefuß | E Base | E Pied | E база устройства |

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Contents

| | Page |
|--------------------------------|------|
| EC-Declaration of conformity | 8 |
| Explication of warning symbols | 8 |
| Safety instructions | 9 |
| Correct use | 10 |
| Unpacking | 10 |
| Commissioning | 10 |
| Accessories | 11 |
| Maintenance and Cleaning | 12 |
| Warranty | 12 |
| Technical data | 12 |

EC-Declaration of conformity

We declare under our sole responsibility that this product is in compliance with the regulations 2006/95/EC, 2004/108/EC and 2011/65/EU and conforms to the standards or standardized documents: EN 61010-1, EN 61010-2-051 and EN 61326-1. A copy of the complete EU Declaration of Conformity can be requested at sales@ika.com.

Explication of warning symbols

Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Indicates practices which, if not avoided, can result in equipment damage.

Safety instructions

- **Read the operating instructions completely before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Because the options for combining products, attachments, shaking vessels, experiment and medium are nearly endless, user safety cannot be ensured simply with design requirements on the part of the product.

DANGER

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise there is a risk of:

- **splashing liquids;**
- **projectile parts;**
- **body parts, hair, clothing and jewellery getting caught.**

- Keep hand/arm vibration to a minimum (Directive 2002/44/EC) when operated manually:
 - Keep the sample container as vertical as possible.
 - Only exert the minimum pressure necessary for the mixing process.
 - Use attachments for multiple samples when there are a large number of samples.
- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.

DANGER

Please pay attention to the dangerous parts of the equipment in Fig. 1.

- The feet of the appliance must be clean and undamaged.
- If the shaking movement of the appliance increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.
- Make certain that the unit is set at the lowest speed before commissioning; otherwise, the unit will begin running at the speed set in last operation. Gradually increase the speed.
- Pay attention when setting the speed to any imbalance of the attachments, shaking vessels and possible spraying of the medium to be mixed.
- Ensure that the shaker does not start to move.

DANGER

Reduce the speed if:

- **the medium splashes out of the vessel because the speed is too high;**
- **the appliance is not running smoothly;**
- **the appliance begins to move around because of dynamic forces;**
- **an error occurs.**

Do not touch rotating parts during operation!

- Firmly secure the accessories and vessels in place, otherwise shaking vessels could be damaged or projected out.
- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
- Position one single shaking vessel in the centre and several shak-

- ing vessels so that they are evenly spread out.
- The appliance is not suitable for hand-held operation.
- Sharp edged vessels cause wear debris on the attachments.

DANGER

Beware of the risk of:

- **flammable materials;**
- **glass breakage as a result of mechanical shaking power.**

- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- **Do not** operate the appliance in explosive atmospheres, with hazardous substances or under water.
- Process pathogenic materials only in closed vessels under a suitable fume hood. Please contact **IKA** application support if you have any question.
- Safe operation is only guaranteed with the accessories described in the "**Accessories**" chapter.
- Always disconnect the plug before fitting accessories.
- The appliance starts up again automatically following a cut in the power supply.
- The instrument can only be disconnected from the mains supply by pulling out the mains plug or the connector plug.
- The socket for the mains cord must be easily accessible.
- The appliance may heat up when in use.
- Never cover the ventilation slots or cooling fins on the motor or on the instrument.

WARNING

This user manual needs to be consulted in all cases where it is used.

- Socket must be earthed (protective ground contact).
- The voltage stated on the nameplate must correspond to the mains voltage.
- Protect the appliance and accessories from bumps and impacts.
- Please observe the permitted speed for the attachments. Never set higher speed.
- During operation please always keep the appliance be covered with the top attachment. Do not allow liquids to get into the appliance when changing attachments, if liquids accidentally splash into the appliance, stop using it immediately and send it back to **IKA** for repair.
- The machine must only be opened by trained specialists, even during repair. The instrument must be unplugged from the power supply before opening. Live parts inside the instrument may still be live for some time after unplugging from the power supply.

WARNING

Covering or parts that are capable of being removed from the unit without accessory equipment have to be reattached to the unit for safe operation in order to prevent, for example, the ingress of fluids, foreign matter, etc.

Correct use

Use

For mixing liquids:

- Touch mode for single test tubes
- Continuous mode for single or several vessels

Intended use: Tabletop device

Range of use (Only for indoor use)

- Laboratories
- Pharmacies
- Schools
- Universities

The device is suitable for use in residential areas and all other areas.

The safety of the user cannot be guaranteed:

- If the instrument is operated with accessories that are not supplied or recommended by the manufacturer;
- If the instrument is operated improperly or contrary to the manufacturer's specifications;
- If the instrument or the printed circuit board are modified by third parties.

Unpacking

Unpack

- Please unpack the device carefully;
- In the case of any damage a fact report must be sent immediately (post, rail or forwarder).

Delivery scope

- Shaker **Vortex 2**
- Standard attachment VG 3.1
- Operating instructions
- A warranty card

Commissioning

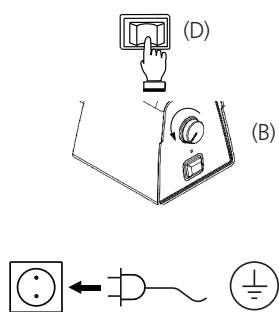
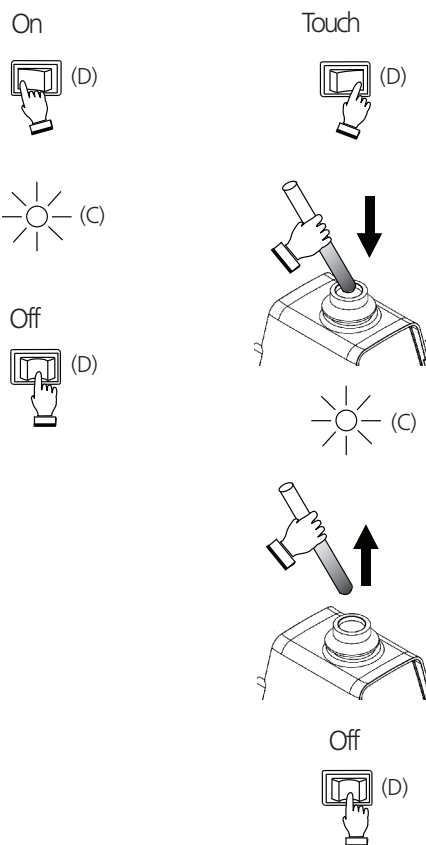
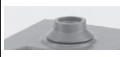
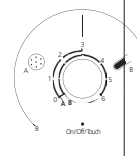
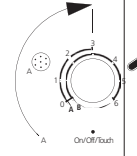
Fig. 2


Accessories

Possible uses and permitted speed ranges of attachments:


| Designation | Description | Id.-No. | Touch Mode | Continuous Mode | Range of speed Max. | Max. load (kg) |
|--|--|-----------|------------|-----------------|---|----------------|
|  VG 3.1 Standard attachment | For test tubes and small vessels up to diameter of 30 mm | 33 412 00 | × | × |  | 0.4 |
|  VG 3.2 One-hand attachment | 88 mm-plate attachment with rubber pad | 33 423 00 | × | × | | |
|  VG 3.3 Universal attachment | 150 mm-plate attachment with rubber pad | 33 424 00 | - | × |  | |

See more accessories on www.ika.com.

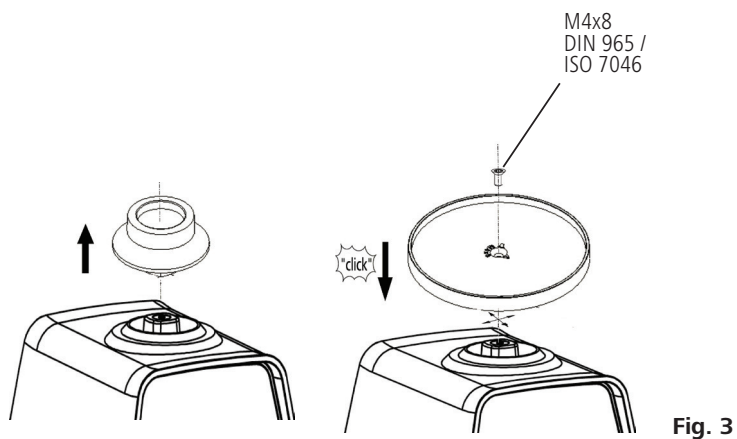Changing attachments

Using inserts

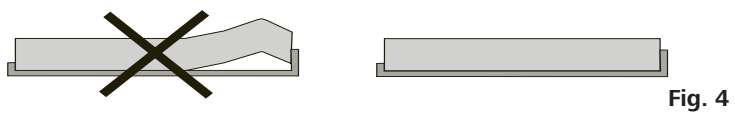Maintenance and Cleaning

The instrument is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Cleaning

For cleaning disconnect the main plug!

Use only cleaning agents which have been approved by **IKA** to clean **IKA** instruments.

Dirt Cleaning agent

| | |
|------------------------|--|
| Dyes | isopropyl alcohol |
| Construction materials | water containing tenside/isopropyl alcohol |
| Cosmetics | water containing tenside/isopropyl alcohol |
| Foodstuffs | water containing tenside |
| Fuels | water containing tenside |

For materials which are not listed, please request information from **IKA** application support.

Wear protective gloves during cleaning the instruments.

Electrical instruments may not be placed in the cleansing agent for the purpose of cleaning.

Do not allow moisture to get into the instrument when cleaning.

Before using another than the recommended method for cleaning

or decontamination, the user must ascertain with **IKA** that this method does not destroy the instrument.

The rubber of the attachments' pads (VG 3.2 and VG 3.3) is a natural product, so talcum can discharge. Clean the rubber pad with glycerin.

Spare parts order

When ordering spare parts, please give:

- Machine type;
- Manufacturing number, see nameplate;
- Item number and designation of the spare part, see www.ika.com.

Repair

Please send instrument in for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.

For this you should request the "Decontamination Clearance Certificate" from **IKA**, or use the download printout of it from the **IKA** website www.ika.com.

If you require servicing, return the instrument in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

Technical data

| | | |
|--|-------------|--|
| Operating voltage / Design frequency | V/Hz | 220...240 ±10% / 50/60 |
| or | V/Hz | 115±10% / 60 |
| or | V/Hz | 100±10% / 50/60 |
| Input power | W | 60 |
| Speed | rpm | (750 - 2750) ± 250, at 50 Hz, infinitely adjustable (750 - 3300) ± 250, at 60 Hz, infinitely adjustable |
| Speed adjustment | | analog (infinitely) rotating knob in front of the device |
| Speed display | | Scale 0 - 6 |
| Drive | | Shaded pole motor |
| Agitation stroke | mm | 4 |
| Shaking motion | | horizontal, approximately circular |
| Perm. duration of operation | % | 100 |
| Perm. ambient temperature | °C | +5 to +40 |
| Perm. relative humidity | % | 80 |
| Protection type (acc. to DIN EN 60529) | | IP 21 (IP 21 can only be achieved when fitted with top attachment) |
| Protection class | | I (protective earth) |
| Overvoltage category | | II |
| Contamination level | | 2 |
| Operation at a terrestrial altitude | m | max. 2000 |
| Dimensions total (W x D x H) (without attachment) | mm | 120 x 138 x 140 |
| Weight without supported load | kg | 3.9 |
| Max. supported load (incl. attachment) | kg | 0.4 |

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IKA - Werke GmbH & Co.KG