40 268 00

**IKA®**

## IKA® ULTRA -TURRA® Tube Drive control

CE

ULTR\_c\_042012

Reg.-No. 4343/01

Betriebsanleitung Ursprungssprache	DF	4
Operating instructions	EN	16
Mode d'emploi	FR	28
Indicaciones de seguridad	ES	40
Veiligheidsinstructies	NL	41
Norme di sicurezza	IT	42
Säkerhetsanvisningar	SV	43
Sikkerhedsanvisninger	DA	44
Sikkerhetsanvisninger	NO	45
Turvallisuusohjeet	FI	46
Instruções de serviço	PT	47
Wskazówki bezpieczeństwa	PL	48
Bezpečnostní pokyny	CS	49
Biztonsági utasítások	HU	50
Varnostna opozorila	SL	51
Bezpečnostné pokyny	SK	52
Ohutusjuhised	ET	53
Drošības norādes	LV	54
Nurodymai dėl saugumo	LT	55
Указания за безопасност	BG	56
Instrucțiuni de securitate	RO	57
Υποδείξεις ασφαλείας	EL	58

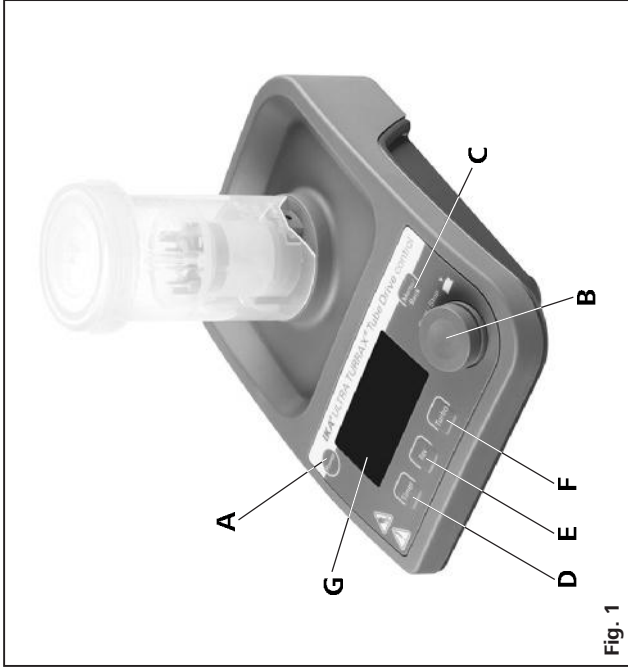
Fig. 1

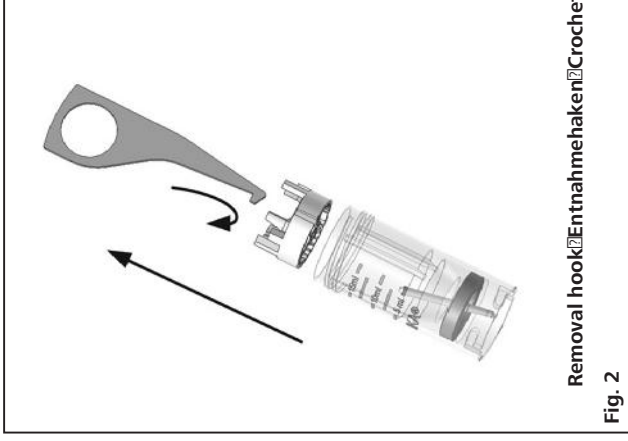
Fig. 2

Pos.	Bezeichnung	Item	Designation	Pos.	Désignation
A	Taste "Power"	A	"Power" button	A	Bouton "Power"
B	Druck- und Drehknopf	B	Rotating Pressing knob	B	Bouton Rotatif Pressage
C	Taste "Menu/Back"	C	"Menu/Back" button	C	Bouton "Menu/Back"
D	Taste "Timer"	D	"Timer" button	D	Bouton "Timer"
E	Taste "Rev"	E	"Rev" button	E	Bouton "Rev"
F	Taste "Turbo"	F	"Turbo" button	F	Bouton "Turbo"
G	Anzeige	G	Display	G	Écran

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## EC-Declaration of conformity

We declare under our sole responsibility that this product conforms to the regulations 89/336EC, 98/79EC and 2006/95EC and conforms with the standards or standardized documents DIN EN IEC 61010-1, DIN EN IEC 61326-1, DIN EN ISO 12100-1, -2 and EN 60204-1.

## Explanation of symbols

General hazard

This symbol identifies information **that is of vital importance for safeguarding your health and safety**. Disregarding this information can lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**. Disregarding this information can result in damage to the instrument or to system components.

This symbol indicates information **which is important for ensuring that the operations of the appliance are performed efficiently and for using the instrument**. Failure to observe this information can result in inaccurate results.

Risk of biohazard infection.

Risk of hand injury.

### 3 Safety instructions

For your protection:

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- The instrument must be supervised at all times when in operation.

**The instrument must only be operated with a tube attached with no tube there is a risk of injury from the rotating locating pin.**

- **Hearing protection must be worn when working with BMT-20 or BMT-50 tubes.**

**Wear your personal protective equipment in accordance with the hazard category of the medium to be processed.**

Otherwise there is a risk of:

- splashing liquids
- projectile parts
- body parts, hair, clothing and jewellery getting caught.

- **Set up the instrument in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.**

- The feet of the instrument must be clean and undamaged.

- The instrument is not suitable for manual operation.

- The instrument may heat up when in use.

- Check the instrument and accessories beforehand for damage each time when you use them. Do not use damaged components.

- Ensure that the cover on the tube is screwed on tightly.

- Ensure that the tube is firmly attached to the bayonet lock connector on the drive unit prior to operating the instrument.

- The tube must only be attached and removed while the motor is stationary.

The IKA® tubes must always be closed when the instrument is in operation. Switch off the instrument immediately if any material leaks from the tube. Clean the instrument.

- **Always open the tube carefully after use as the media in the tube may heat up due to transfer of energy during operation; leading to pressurization of the container: risk of material spraying, protective equipment must be worn.**

**When handling infectious or toxic sample materials, do NOT touch the sample, mixture or waste with your hands. Wear gloves and lab coat and goggles if necessary.**

- **The temperature of the material must not exceed 40 °C.**

- Only use tubes approved by IKA®.

- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.

- Beware of the risk of flammable materials.

- **Do not operate the instrument in explosive atmospheres, with hazardous substances or under water.**

- The instrument doesn't start up again automatically following a cut in the power supply.

- Safe operation is only guaranteed with the accessories described in the 'Accessories' chapter.

**For protection of the instrument**

- The voltage stated on the name plate must correspond to the mains voltage.

- The instrument must only be operated with the original plug-in power supply unit.

- **Protect the instrument and accessories from bumps and impacts.**

- The instrument may only be opened by experts.

## 4 Unpacking

### • Unpack

- Please unpack the device carefully
- In the case of any damage a fact report must be set immediately (for post, rail or forwarder)

### • Delivery scope

- **Tube Drive control**
- Power supply unit with exchangeable adapters\_
- Removal hook
- Operating instructions
- A warranty card
- A decontamination clearance certificate.

Power supply unit: with exchangeable adapters\_

### Power supply

**Exchangeable adapter  
Europe, Switzerland**

**Exchangeable adapter  
USA, China**

**Exchangeable adapter  
England**

**Exchangeable adapter  
Australia**

**Optionally in the storage suite-case:**

- 1 x ST 20
- 1 x DT 20
- 1 x BMT 20 S
- 1 x BMT 20 G

## 5 Correct use

### • Use

	Application	Volume	Mode
▶ with Tube ST-20...	Mixing Stirring	2 - 15 ml	A, B, P1-P9, Turbo, Rev
▶ with Tube ST-50...		15 - 50 ml	
▶ with Tube DT-20...	Dispensing Stirring	5 - 15 ml	A, B, P1-P9, Turbo, Rev
▶ with Tube DT-50...		15 - 50 ml	
▶ with Tube BMT-20...	Crushing Mixing wet/dry	2 - 15 ml	A, B, P1-P9, Turbo, Rev
▶ with Tube BMT-50...		15 - 50 ml	

### • Range of use (indoor use only)

- Laboratories
- Pharmacies
- Schools
- Universities

This device is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network which also supplies residential areas.

The safety of the user cannot be guaranteed:

- if the instrument is operated with accessories that are not supplied or recommended by the manufacturer
- if the instrument is operated improperly contrary to the manufacturer's specifications
- if the instrument is operated improperly contrary to the manufacturer's specifications.

## 6 Useful facts

Changes in viscosity and volume caused by processing of dispersion media may cause small fluctuations in the speed of the device.

## 7 Interfaces and outputs

### Remote control

The device can be operated in "Remote" mode via USB interface using labworldsoft® laboratory software. The USB interface is located on the back side on the device and can be connected to a PC using the USB cable.

**Note:** Please note the system requirements as well as the operating instructions and help section included within the software.

### USB Interface

The Universal Serial Bus (USB) is a serial bus system used to connect the device with the PC. Devices that support USB can be connected to each other whilst they are running (hot pluggable) and provide automatic recognition of the connected devices and their properties. Use the USB interface in conjunction with labworldsoft® for operation in "Remote" mode.

### Installation

Connect the device to the PC using the USB data cable. The device will then transmit information to the Windows operating system to tell which device drivers are required. Windows will then either:

- Load the driver.
- Install the driver automatically if not already installed.
- Prompt the user to perform a manual installation.

Navigate to <http://www.ika.net/ika/lws/download/stmcdc.inf>

### Command syntax and format

The following applies to the command set:

- Commands are generally sent from the computer (Master) to the device (Slave).
- The device sends only at the computer's request. Even fault indications cannot be sent spontaneously from the device to the computer (automation system).
- Commands are transmitted in capital letters.
- Commands and parameters including successive parameters are separated by at least one space (Code: hex 0x20).
- Each individual command (incl. parameters and data) and each response are terminated with Blank CR Blank LF (Code: hex 0x20 hex 0x0d hex 0x20 hex 0x0a) and have a maximum length of 80 characters.
- The decimal separator in a number is a dot (Code: hex 0x2E).

The above details correspond as far as possible to the recommendations of the NAMUR working party (NAMUR recommendations for the design of electrical plug connections for analogue and digital signal transmission on individual items of laboratory control equipment, rev. 2.9).

The NAMUR commands and the additional specific IKA® commands commissioning serve only as low level commands for communication between the tube driver and the PC. With a suitable terminal or communications program these commands can be transmitted directly to the tube driver. The IKA® software package, labworldsoft, provides a convenient tool for controlling tube driver and collecting data under MS Windows, and includes graphical entry features, for motor speed ramps for example.

The following table summarizes the (NAMUR) commands understood by the tube drive control up to Version 2.9.


Abbreviations used:

m	4	Numbering parameter (integer)
m	4	Speed

NAMUR Commands	Function	Remark
IN_NAME	Request designation. Start the Remote control function.	Device keypad not available except "power" button is pressed, after this command is sent to device. "Remote" symbol is displayed.
RESET	Switch to normal operation	Device keypad available again, and clear the "Remote" symbol.
IN_PV_X X=4	Read actual speed value	
IN_SP_X X=4	Read target value input	
OUT_SP_X X=4	Set target speed value, input: rpm	
START_X X=4	Switch on appliance remote function. Start value function with setting.	
IN_SOFTWARE	Request software ID, number, date and version	

## 8 Commissioning

Pay attention to the ambient conditions listed in the "Technical Data".

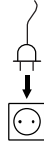If these conditions are met, the unit is ready for operation when the mains plug has been plugged in.

Fit the tube carefully and ensure it is properly attached.

If above procedures are not followed, safe operation can not be guaranteed and/or the equipment may be damaged.


## 9 Switching on

- ⇨ Press "Power" button (A) to switch on the device, see Fig. 1.
- ⇨ The start-up screen will be displayed for several seconds. This screen shows the unit designation and the software version.

- ⇨ Then, the system check screen will be displayed for several seconds.


Symbols are shown on the screen, including the working mode setting, the target speed, the maximum rotation speed (rpm) and timer maximum value etc. of the device.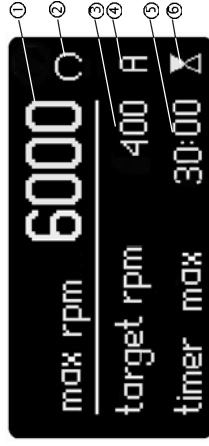

Symbols Explanation

- ① Maximum rotation speed of the device (If "Turbo" function is activated, the maximum rotation speed could reach 8000 rpm).
- ② Rotation symbol (The symbol will change into  $\cup$  or  $\cup$  if "Rev" function is activated).
- ③ Rotation speed setting.
- ④ Operation mode.
- ⑤ "Timer" maximum value of the device.
- ⑥ "Timer" symbol.

⇒ After the system check screen, the working screen will appear and the device is ready for operation.

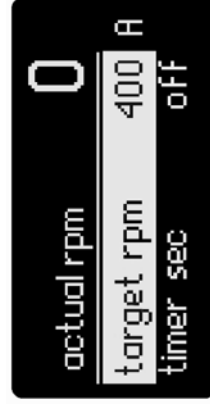**Note:** Figure above shows specified values corresponding to the default settings at delivery.

## 10 Factory setting

The device is supplied with following factory setting:

Menu Settings:		
mode	A	rpm
max rpm	6000	minute
timer max	30	
language	English	
Working Setting:		
target rpm	400	rpm
timer sec	off	
rev sec	off	

### 10.1 Resetting parameters to factory setting

To reset the parameters to the factory settings, proceed as following:

- **Pull out main plug from the device.**
- Hold the controller buttons "Menu/Back" and "Timer" simultaneously and re-insert main plug in power outlet.
- The unit is now reset to the factory settings.

## 11 Setting menu parameters

### 11.1 Operation mode

The device can be operated in mode A, B and P1 to P9. The mode setting will be stored automatically when operation is over, or the device is switched off. User may change the mode setting in the main menu if necessary.

#### • Operation mode A

In this mode, when current operation is over or the device is switched off, all working settings will not be stored.

#### • Operation mode B

In this mode, when current operation is over or the device is switched off, all menu and working settings will be stored.

#### • Operation mode P1 - P9

In mode P1-P9, the user can set and save required parameters, (target speed, timer value and rev value) in program. The parameter settings can only be changed in the menu program. These settings can also be used directly when necessary.

If the timer value in the program setting is more than 3 minutes, the speed upper limit is 4000 rpm.

If the timer is not activated in the program and speed is higher than 4000 rpm, the speed will also decrease to 4000 rpm after running 3 minutes.

#### • Setting operation mode between A and B

- ☞ Press the "Menu/Back" button (C) in the main menu.
- ☞ Turn the "Rotating/Pressing" knob (B) to select the "mode" item in the main menu.
- ☞ Press the "Rotating/Pressing" knob (B).
- ☞ Turn the "Rotating/Pressing" knob (B) to select mode A or B in submenu.
- ☞ Press the "Rotating/Pressing" knob (B) again to complete the process.
- ☞ Press the "Menu/Back" button (C) to return to the working screen.

#### • Setting operation mode between P1 and P9

- ☞ Press the "Menu/Back" button (C) in the main menu.
- ☞ Turn the "Rotating/Pressing" knob (B) to select the "mode" item in the main menu.
- ☞ Press the "Rotating/Pressing" knob (B).
- ☞ Turn the "Rotating/Pressing" knob (B) to select the "mode" item in mode P1 and P9 in submenu.
- ☞ Press the "Rotating/Pressing" knob (B) and enter desired menu program.
- ☞ Turn the "Rotating/Pressing" knob (B) to select desired program item.
- ☞ Press the "Rotating/Pressing" knob (B) .
- ☞ Turn the "Rotating/Pressing" knob (B) to change the value.
- ☞ Press the "Rotating/Pressing" knob (B) again to complete the setting.
- ☞ Press the "Menu/Back" button (C) to complete the process, display will return to working screen.

**Note:** The target speed and timer value in the program should be less than "rpm max" and "timer max" setting in menu. Otherwise, the device will run according to the setting in the menu.

### 11.2 Operation language

The language can be set to: English, German, Chinese, French, Spanish, Italian and Korean. The language setting will be stored automatically when operation is over or the device is switched off. User can change the language setting in the main menu if necessary.

#### • Setting operation language

- ☞ Press the "Menu/Back" button (C) in the main menu.
- ☞ Turn the "Rotating/Pressing" knob (B) to select the "language" item in the main menu.
- ☞ Press the "Rotating/Pressing" knob (B) to confirm "language".
- ☞ Turn the "Rotating/Pressing" knob (B) to select a language.

- Press the "Rotating/Pressing" knob (B) again to complete the setting.
- Press the "Menu/Back" button (C) to return to the working screen.

### 11.3 Maximum rotation speed (max rpm)

The default setting "max rpm" is 6000 rpm in the device. User can change the setting to the range of 400 to 6000 rpm in the main menu.

The "max rpm" setting will be stored automatically when operation is over or the device is switched off.

#### • Setting "max rpm"

- Press the "Menu/Back" button (C) in the main menu.
- Turn the "Rotating/Pressing" knob (B) to select the "max rpm" item in the main menu.
- Press the "Rotating/Pressing" knob (B).
- ⇒ A mark ▶ will appear to indicate the activated item.
- Turn the "Rotating/Pressing" knob (B) to adjust the value.
- Press the "Rotating/Pressing" knob (B) again to complete the process.
- Press the "Menu/Back" button (C) to return to the working screen.

**Note:** The speed „max rpm“ can be adjusted with the "Rotating/Pressing" knob (B) in a range between 400 to 6000 rpm. This is achieved in magnitudes of 10 rpm. In remote control mode via computer, adjusting magnitude is 1 rpm.

### 11.4 Timer maximum value (timer max)

The default "timer max" is 30 minutes in the device. User can change the setting in 10 second intervals and up to 30 minutes in the main menu.

The "timer max" setting will be stored automatically when operation is over or the device is switched off.

#### • Set "timer max"

- Press the "Menu/Back" button (C) in the main menu.
- Turn the "Rotating/Pressing" knob (B) to select the "timer max" in the main menu.
- Press the "Rotating/Pressing" knob (B)
- ⇒ A mark ▶ will appear to indicate the activated item.
- Turn the "Rotating/Pressing" knob (B) to change the value.
- Press the "Rotating/Pressing" knob (B) again to complete the process.
- Press the "Menu/Back" button (C) to return to the working screen.

**Note:** The timer „timer max“ can be adjusted with the "Rotating/Pressing" knob (B) in a range from 10 seconds to 30 minutes. This is achieved in magnitudes of 10 seconds. In remote control mode via computer, adjusting magnitude is 1 second.

## 12 Setting working parameters

### 12.1 Target speed (target rpm)

Set "target rpm" within "rpm max" range in the working screen by rotating the "Rotating/Pressing" knob (B). Adjustment-magnitude is 10 rpm. In remote control mode via computer, adjustment magnitude is 1 rpm.

**Note:** For tube protection, if the "Timer" setting is more than 3 minutes, the speed will be limited to less than 4000 rpm. Or the speed can be set to any value within "rpm max" range which is set in the main menu.  
If the "Timer" isn't activated and the speed value is over 4000 rpm, the speed will be reduced to 4000 rpm after running 3 minutes. User can change the speed by rotating "Rotating/Pressing" knob (B) during running in mode A and B.

#### • Setting "target rpm" in working screen

- ☞ Turn the "Rotating/Pressing" knob (B) to change the value.
- ☞ Press "Rotating/Pressing" knob (B) to start working.
- ⇒ Rotation symbol (C, D, E) indicates running.

### 12.2 Timer function

Set "Timer" within "timer max" range in the working screen by rotating the "Rotating/Pressing" knob (B). Adjustment magnitude is 10 seconds. In remote control mode via computer, adjustment magnitude is 1 second.

If the "Timer" is deactivated when the device is started, the device will automatically count from 0 to 30 minutes and then the unit will stop running.

**Note:** If the "Timer" button (D) is pressed in mode B after processing has stopped by pressing "Start/Stop" knob (B), the "Timer" function will be activated automatically and device will store last running time as "timer" value in next operation.

- Setting "Timer" value (timer sec) in working screen
- ☞ Press "Timer" button (D) to activate the timer setting in the working screen.

- ☞ Turn the "Rotating/Pressing" knob (B) to change the value.
- ☞ Press "Timer" button (D) again to complete the setting.
- ⇒ Current values are stored.

### 12.3 Reverse (Rev) function

When setting "Rev" value with the "Rotating/Pressing" knob (B), the value can be set from 10 to 60 seconds where adjustment magnitude is 1 second.

**Note:** If both "Rev" function and "Timer" function are activated, the "Rev" value should be less than "Timer" value. Otherwise, the reverse function will not work.

If the reverse function is activated, press "Rev" button (E) during running, the reverse function will be deactivated. Press "Rev" button (E) again, the reverse function will be reactivated and the screen will indicate "Rev" value setting.

#### • Setting "Rev" timer (rev sec) in working screen

- ☞ Press "Rev" button (E) to activate the reverse setting in the working screen.
- ☞ Turn the "Rotating/Pressing" knob (B) to change the value.
- ☞ Press "Rev" button (E) again to complete the setting.
- ⇒ Current values are stored.

### 12.4 Turbo Function

The rotation speed can increase to 8000 rpm in several seconds by pressing down "Turbo" button (F) during operating. For tube protection, Turbo function can only last 60 seconds at most if you press down "Turbo" button (F). Several seconds later, Turbo function can be activated again after symbol has disappeared.

- Activate and hold "Turbo" function in working screen:
  - Press down (Turbo) button (F) over 2 seconds when the device is working, to activate Turbo function.
  - ⇒ The Turbo mark  appears and changes alternately to indicate that the Turbo function is activated and rotation speed increase to 8000 rpm in a short time.
  - Release "Turbo" button (F).
  - ⇒ Turbo function will be stopped and rotation speed will return to original setting. A symbol  appears to indicate that the Turbo function is stopped. Before this symbol disappears, turbo function can not be activated again.

Only use cleansing agents which have been recommended by IKA®.

#### Cleaning agent

- isopropyl alcohol
- water containing tenside/isopropyl alcohol
- water containing tenside/isopropyl alcohol
- water containing tenside
- water containing tenside

- Dyes
- Construction materials
- Cosmetics
- Foodstuffs
- Fuels

For materials with are not listed, please request information from IKA® application support.

Wear protective gloves during cleaning the instruments.

Electrical instruments may not be placed in the cleansing agent for the purpose of cleaning.

**Do not allow moisture to get into the instrument when cleaning.**  
**Before using another than the recommended method for cleaning or decontamination, the user must ascertain that this method does not destroy the instrument.**

#### Spare parts order

- When ordering spare parts, please give:
  - machine type
  - manufacturing number, see type plate
  - item and designation of the spare part, see [www.ika.com](http://www.ika.com), spare parts diagram and spare parts list.

#### Repair

In case of repair the device has to be cleaned and free from any materials which may constitute a health hazard.

For repair, use the form supplied or printed copy on the IKA® website: [www.ika.com](http://www.ika.com).

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

### 13 Error message

It is possible for the media to become lodged between the cutters of the DT-tubes. The unit will automatically change the direction of rotation up to three times in an attempt to free the cutters. The unit will switch off automatically if this is not possible.

If the device becomes blocked on three occasions within a certain period of time the following error message will occur:

→ Display Err 1

→ An audible signal will be given.

**Corrective action:** Switch off the device, check contents, switch on the device.

### 14 Maintenance

The drive unit is maintenance-free but not immune from wear of carbon brushes of the motor wear down over the time.

For cleaning disconnect the main plug.

## 15 Accessories

- **ST-20** Tube with stirring device 20 ml
- **ST-20-M** Tube with stirring device 20 ml, with pierceable membrane
- **ST-20-M-gamma** Tube with stirring device 20 ml,  $\gamma$ -sterilized, with pierceable membrane
- **ST-50** Tube with stirring device 50 ml
- **ST-50-M** Tube with stirring device 50 ml, with pierceable membrane
- **ST-50-M-gamma** Tube with stirring device 50 ml,  $\gamma$ -sterilized, with pierceable membrane
- **DT-20** Tube with dispersing device 20 ml
- **DT-20-M** Tube with dispersing device 20 ml, with pierceable membrane
- **DT-20-M-gamma** Tube with dispersing device 20 ml,  $\gamma$ -sterilized, with pierceable membrane
- **DT-50** Tube with dispersing device 50 ml
- **DT-50-M** Tube with dispersing device 50 ml, with pierceable membrane
- **DT-50-M-gamma** Tube with dispersing device 50 ml,  $\gamma$ -sterilized, with pierceable membrane
- **BMT-20-S** Tube for grinding with stainless steel balls 20 ml,
- **BMT-20-S-M** Tube for grinding with stainless steel balls 20 ml, with pierceable membrane
- **BMT-20-S-M-gamma** Tube for grinding with stainless steel balls 20 ml,  $\gamma$ -sterilized, with pierceable membrane
- **BMT-50-S** Tube for grinding with stainless steel balls 50 ml,
- **BMT-50-S-M** Tube for grinding with stainless steel balls 50 ml, with pierceable membrane
- **BMT-50-S-M-gamma** Tube for grinding with stainless steel balls 50 ml,  $\gamma$ -sterilized, with pierceable membrane

- **BMT-20-G**
- **BMT-20-G-M**

Tube for grinding with glass balls 20 ml  
Tube for grinding with glass balls 20 ml, with pierceable membrane

- **BMT-50-G**
- **BMT-50-G-M**

Tube for grinding with glass balls 50 ml  
Tube for grinding with glass balls 50 ml, with pierceable membrane

- **TC-20**
- **TC-20-M**

Tube cover 20 ml  
Tube cover 20 ml with pierceable membrane

- **TC-50**
- **TC-50-M**

Tube cover 50 ml  
Tube cover 50 ml with pierceable membrane

- **Glass balls**
- **Stainless steel balls**
- **labworldsoft®**

## 16 Warranty

In accordance with **KA**'s warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

## 17 Technical data

### Power supply unit

Input	<b>VAC</b>	100 - 240
	<b>A</b>	0,6
	<b>Hz</b>	50-60
Output	<b>VDC</b>	24
		24W, LPS (limited power source)

Protection class

2 (double insulated) 

### Tube Drive

Operating voltage	<b>VDC</b>	24
	<b>mA</b>	1000
	<b>W</b>	20

Power consumption, normal operation

**W** 2

Power consumption, standby operation

**W** 17

Power output

**W** EC fiat motor

Drive

400 - 6000

Speed range

infinitely adjustable, adjustment magnitude 10 rpm

Turbo speed

Mode A

Mode B

Mode P1-P9

Speed setting

rotating/pressing knob

Speed display

digital

Timer

10 seconds-30 minutes

Reverse timer

10 - 60 Seconds

Display-timer

digital

Perm. ambient temperature

**°C** +5 to +40

Perm. rel. humidity

**%** 80

Protection to DIN EN 60529

IP 20

Contamination level

2

Overvoltage category

II

Operation at a terrestrial altitude level

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**IKA® - Werke GmbH & Co.KG**