



CL-20

INSTRUCTION MANUAL





Description

CL-20 is a fluorescence ring light illuminator designed to use with the SZM and SZR series.

Specifications

Light source:	8 W, fluorescence ring light
Mains power:	230 V

Instructions

Attach the ring in Fig. 1 to the objective holder (the inferior part of the SZM or SZR microscope head). Fix the ring light illuminator around it, using the side screws (see Fig. 2). Connect the mains cable to a 230 V power source and switch the illuminator on and off using the switch located close to the power cable on the top.



Fig. 1

Clean with a soft dry cloth. Do not use solvents.

Note: during use, the lamp may become warm. This is absolutely normal, and does not cause malfunction.



Fig. 2 The final setup



MAD Apparecchiature scientifiche

Via Rigla 32, Ponteranica (BG) - ITALY

Tel.: +39 035 571392 (6 linee) Telefax: +39 035 571435

Optika Mikroskop ve Laboratuvar Cihazlari Ticaret Ltd. Şti.

Yürekli Adam Sokak N0:20 Kat:6 Daire:14 Kavacık BEYKOZ/İSTANBUL

Tel: 090 216 413 4100 Fax: 090 216 413 1617

MAD Iberica

C/ La Lluna, 11 - 08001- Barcelona - ESPAÑA

Tel.: +34 93 3248682 Fax: +34 93 3248683