

In 1973 NTC services introduced their first automatic gas cylinder change-over unit for use with CO<sub>2</sub> incubators in Tissue Culture, IVF and Microbiology laboratories. Since that date our Pneuchange units have become the best selling gas bottle change-over equipment for these applications.

This equipment is recommended and sold by most of Europe's incubator manufacturers as a result of which there are currently in excess of ten thousand of these units in round the clock use.

NTC has become a major player in the gas control and handling market and now presents it's range, manufactured to the highest quality standards and guaranteed for 12 months.

Our range of gas control equipment is used by:

Chemists  
Biochemists  
Physicists  
Biologists  
Biotechnologists  
Microbiologists  
Embryologists  
Pathologists  
Engineers.

# Laboratory Gas Control Equipment



## Over 40 Years Successfully Serving Laboratories With Innovative Gas Control Equipment



# PC SERIES

## Pneuchange Mk3



Pneuchange is the unit of choice in thousands of laboratories for ensuring the continuity of CO<sub>2</sub> and other inert gases supplies to CO<sub>2</sub> incubators and other equipment where loss of gas would be inconvenient, costly or dangerous. These compact units can be bench or wall mounted using the keyholes on the back of the unit. Full installation instructions are included.

The unique pneumatic logic construction of Pneuchange means that it can be placed anywhere between the gas bottles and the equipment being served, no electrical supplies are involved. These units are suitable for use with all non-corrosive, non flammable gases.

The inlet ports are connected with suitable tubing to the gas bottle regulators. Two pressure gauges indicate the inlet pressures and two coloured indicators show which cylinder is active. On exhaustion of the active bottle the unit switches automatically to the second bottle. The empty cylinder can be replaced at leisure and without fear of loss of gas from the disconnected regulator or inlet port. An audible battery operated change-over alarm is fitted. Max inlet 100 PSI. Min. inlet 30 PSI.

Standard port fittings are 3/8 BSP male and are supplied with (6mm) hose tails. Fittings for 6mm. OD tube as used on some European incubators and other hoesail sizes are also available.

### How to order:

NTC PC15	Outlet pressure range 0-15 PSI
NTC PC30	Outlet pressure range 0-30 PSI
NTC PC60	Outlet pressure range 0-60 PSI
NTC PC100	Outlet pressure range 0-100 PSI

Add VFC suffix for volt free contacts option to allow connectivity to remote alarm system instead of integral buzzer.

## PC Mk3 Variants

### Pneuchange Trio. PC-Trio



This unit has three outlets into which special outlet pressure modules (OPMs) can be plugged. The modules consist of a pressure regulator and gauge. Using these OPMs, up to three incubators can be served at the **same or different** pressures.

PC-Trio is supplied without OPM's. Additional modules can be supplied at the same time or later. The OPM ports are self-sealing and on removal of the OPM will completely isolate the incubator from the gas supply.

### How to order:

NTC PCTRIO	PC-Trio base unit
------------	-------------------


### Output Pressure Modules (OPM):

NTC OPM 35	OPM with 0-35 PSI gauge.
NTC OPM 100	OPM with 0-100 PSI gauge.



## N.B Pneuchange units are not suitable for Gas Chromatography



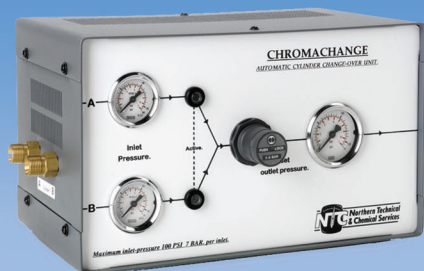

# Chromachange CC100

Chromachange was developed to meet the needs of Chromatographers and other laboratory workers requiring a reliable, compact device for maintaining continuity of gas supplies to GCs, ICPs and other gas dependent equipment.

Unlike large industrial change-over systems, CHROMACHANGE is a self-contained, non-electrical automatic device, requiring no resetting after cylinder change. The specially developed valve system requires inlet gas pressures of 100 PSI (7 BAR) and hence the gas cylinders must be fitted with reliable gas regulators. Change-over takes place when the gas cylinder empties and the outlet pressure falls to 70% of the original value. It is recommended that inlet pressures should be maintained at 100 PSI. The outlet pressure regulator and gauge allows the pressure to your equipment to be set at any value up to that of the inlet pressure.

The status of the feed cylinders is indicated on the inlet pressure gauges and the currently active (on-line) cylinder is indicated by coloured indicators. Gas ports are 3/8" BSP male (as fitted on UK cylinder regulators).

Chromachange is suitable for all non-corrosive, non-flammable gases only.

All units can be bench or wall mounted. Preset, fixed output cylinder regulators are available and are recommended for use with this equipment. A range of optional equipment is available including battery operated change-over alarms and remote indicator panels. Our technical department is always ready to discuss and advise on your particular requirements

How to order:

NTC CC100      Chromachange

## Gas Cylinder Regulators

We can supply a full range of gas cylinder regulators for all laboratory gases. Our prices are some of the most competitive on the market and many models can be supplied from stock.

All regulators are made to the highest international quality standards and are guaranteed for 12 months.





# WolfLabs

**Pricing on any accessories shown can be found by keying the part number into the search box on our website.**

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

**[www.wolflabs.co.uk](http://www.wolflabs.co.uk)**

**Tel : 01759 301142**

**Fax : 01759 301143**

**[sales@wolflabs.co.uk](mailto:sales@wolflabs.co.uk)**

Please contact us if this literature doesn't answer all your questions.