# **Instruction Manual**

HI300N - HI310N


HI301N - HI311N

HI302N - HI312N

HI303N - HI304N

HI322N - HI324N

# **Magnetic Stirrers**

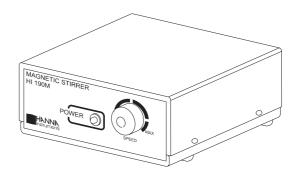



Dear Customer,

Thank you for choosing a Hanna product.

Please read this instruction manual carefully before using the meter. If you need additional technical information, do not hesitate to e-mail us at tech@hannainst.com.

These meters are in compliance with <€ directives.

#### **WARRANTY**

All Hanna Instruments products are warranted for two years against defects in workmanship and materials when used for their intended purpose and maintained according to the instructions. Damages due to accidents, misuse, tampering or lack of prescribed maintenance are not covered. This warranty is limited to repair or replacement free of charge. If service is required, contact the dealer from whom you purchased the instrument. If under warranty, report the model number, date of purchase, serial number and the nature of the failure. If the repair is not covered by the warranty, you will be notified of the charge for repair or replacement. If the instrument is to be returned to Hanna Instruments, obtain a Return Goods Authorization from the Customer Service Department first and then send it with shipment cost prepaid. When shipping any instrument, make sure it is properly packaged for complete protection.

#### **TABLE OF CONTENTS**

| Preliminary Examination                 | 3  |
|---|----|
| General Description                     |    |
| Application Guide                       | 5  |
| Electrical Connections                  | 6  |
| Accessories                             | 6  |
| Functional Description & Specifications | 7  |
| HI 300N & HI 310N                       |    |
| HI 301N & HI 311N                       | 10 |
| HI 302N & HI312N                        | 12 |
| HI 303N                                 | 14 |
| HI 304N                                 | 16 |
| HI 322N & HI 324N                       | 18 |
| Operational Guide for                   |    |
| HI 300N & HI 310N                       | 20 |
| HI 301N, HI 303N & HI 311N              | 22 |
| HI 302N, HI 304N & HI 312N              |    |
| HI 322N & HI 324N                       | 26 |
|   |    |

#### PRELIMINARY EXAMINATION

Remove the instrument from the packing material and examine it to make sure that no damage has occurred during shipping. If there is any damage, notify your Dealer.

Each stirrer is supplied complete with a magnetic stir bar.

**Note:** Save all packing material until you are sure that the instrument functions correctly. All defective items must be returned in the original packing with the supplied accessories.

#### **GENERAL DESCRIPTION**

The Hanna series of magnetic stirrers are designed with simplicity of use for stirring and mixing solutions in a wide range of laboratory applications. Several models are available in an extensive range covering:

- general purpose stirrers
- dual-speed range
- auto-reverse cycle stirring
- · rpm indication
- timer control

All models employ a motorized magnet to rotate a stirring bar to create a forced vortex rotational flow in the solution with a constant angular speed which is adjustable with an external speed knob.

Unlike conventional stirrers, all Hanna stirrers use the concept of pulse-driven drive to give the full power required to effectively stir the fluid regardless of the speed.

This eliminates the need to start the motor at high speed and the user is now able to control the speed precisely. Each stirrer is equipped with an opto-sensor for speed monitoring, and a speed limiter circuit that prevents hazardous operation when the load is suddenly removed.

The power required to effectively stir a fluid is affected by the torque load on the stirrer which is dependent on the capacity, type and viscosity of the fluid. The speed of the stirrer decreases as the torque load increases and vice versa, unless the power delivered changes accordingly.

For Newtonian fluids like water, the viscosity remains constant during stirring. General purpose, normal stirrers are ideal for this kind of fluids.

For some fluids like oil and clay, the viscosity changes during the stirring process, making it necessary to adjust the power to maintain a constant speed during the stirring.

Hanna Auto-Speed stirrers (HI 31X series) are designed with automatic feedback control to maintain constant speed when the torque load changes and are therefore ideal for stirring non-Newtonian fluids where the viscosity changes during stirring.

They are also ideal for mixing applications where one solution is added to another which changes the capacity and thereby the torque load on the stirrer.


The Auto-Speed stirrers maintain a constant stirring speed during the mixing up to the full capacity of 5 liters. They can also be used with Newtonian fluids up to a capacity of 5 liters.

Hanna Instruments reserves the right to modify the design, construction and appearance of its products without advance notice.

#### **APPLICATION GUIDE**

| APPLICATION  | MODEL                                    |
|--|--|
| Newtonian fluids up to 2.5 liters  | HI 300N<br>HI 301N<br>HI 302N<br>HI 304N |
| Non-Newtonian fluids<br>up to 5 liters whose<br>viscosity changes<br>during stirring or<br>Newtonian fluids up<br>to 5 liters or mixing<br>operation where load<br>changes during stirring | HI 310N<br>HI 311N<br>HI 312N            |
| Repetitive stirring in<br>Newtonian fluids up<br>to 2.5 liters   | HI 303N                                  |
| Repetitive stirring in multiple samples with the same stirring speed   | HI 304N                                  |
| Routine stirring application with a predetermined stirring time  | HI 322N                                  |
| Repetitive stirring in multiple samples with the same stirring speed and predetermined stirring time   | HI 324N                                  |

#### **ELECTRICAL CONNECTIONS**

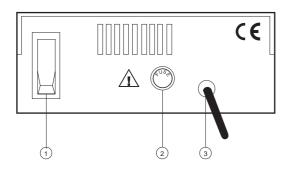

- 1. ON/OFF switch for power supply
- 2. Fuse holder for 400 mA/250 V fuse
- 3. AC power cord.

Unplug the stirrers from the power supply before replacing the fuse.

#### **ACCESSORIES**

HI 731320 Magnetic stir bar (10 pcs)

# **FUNCTIONAL DESCRIPTION AND SPECIFICATIONS**

See the following pages for specifications of each model.

The capacity is referred to stirring water with a viscosity of 0.001 N·s/m<sup>2</sup> at 25°C.

#### Specifications for all models

Speed ON/OFF and continuously Control variable with potentiometer

knob

**Drive** Direct drive D.C. Motor with

magnet

4 W **Power** 

Consumption

**Torque** 85 g-cm Delivered

Power 115 or 230 Vac, 50/60 Hz Supply depending on models

400 mA/250 V; 20x5 mm **Fuse** 

Stirring Bar Teflon® coated,

L 50 x dia 7 mm

#### Recommendations for Users

Before using these products, make sure that they are entirely suitable for the environment in which they are used.

Operation of these instruments in residential area could cause unacceptable interferences to radio and TV equipments, requiring the operator to take all necessary steps to correct interferences

Any variation introduced by the user to the supplied equipment may degrade the instruments' EMC performance.


Unplug the stirrers from the power supply before replacing the fuse.


To avoid damages or burns, do not place these stirrers in microwave ovens.

### <u>HI 300N & HI 310N</u> GENERAL PURPOSE STIRRERS

These general purpose stirrers are suitable for daily stirring applications in the laboratory.

**HI 300N** normal stirrer and **HI 310N** autospeed stirrer are covered with a corrosion resistant stainless steel top.

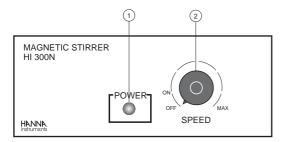

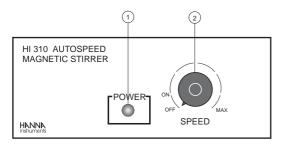



- 1. **Power indicator** (LED lights up when the stirrer is powered)
- 2. **ON/OFF/SPEED knob**, to turn the stirrer ON and OFF, and set the stirring speed (continuously variable up to the maximum)

|  | HI300N   | HI310N                    |
|--|--|---------------------------|
| Maximum<br>Stirring<br>Capacity                      | 2.5 Liters<br>(0.66 gallons)   | 5 Liters<br>(1.3 gallons) |
| Speed Range<br>Min<br>Max                            | 100 rpm<br>800 to 1000 rpm   |                           |
| Auto-Speed   | notapplicable  | standard                  |
| Power Supply HI 300N-1 HI 300N-2 HI 310N-1 HI 310N-2 | 110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz<br>110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz |                           |
| Installation<br>Category                             | П  |                           |
| Cover<br>Material                                    | AISI 316 stainless steel   |                           |
| Environment  | 0 to 50°C (32 to 122°F);<br>RH max 95%   |                           |
| Dimensions   | 180 x 180 x 70 mm<br>(7.1 x 7.1 x 2.8")  |                           |
| Weight   | 1.4 kg (3.   | 1 lb.)                    |

### HI 301N & HI 311N TWO-SPEED STIRRERS

**HI 301N** normal stirrer and **HI 311N** autospeed stirrer, are ideal for long duration stirring which require slow speed and accurate speed control (for example in electrolysis applications). The selectable HIGH/LOW range allows to set the speed precisely.

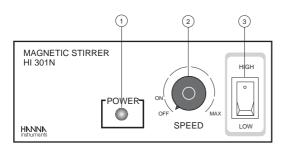

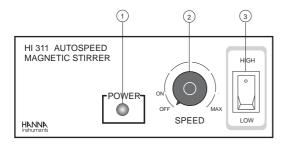


- 1. **Power indicator** (LED lights up when the stirrer is powered)
- ON/OFF/SPEED knob, to turn the stirrer ON and OFF, and set the stirring speed (continuously variable up to the maximum)
- 3. **HIGH/LOW switch** (for selecting low or high speed range)

|  | HI301N   | HI311N                    |
|--|--|---------------------------|
| Maximum<br>Stirring<br>Capacity                      | 2.5 Liters<br>(0.66 gallons)   | 5 Liters<br>(1.3 gallons) |
| Speed Range<br>Low<br>High                           | 100 to 500 rpm<br>100 to 800/1000 rpm  |                           |
| Auto-Speed   | not applicable   | standard                  |
| Power Supply HI 301N-1 HI 301N-2 HI 311N-1 HI 311N-2 | 110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz<br>110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz |                           |
| Installation<br>Category                             | П  |                           |
| Cover<br>Material                                    | AISI 316 stainless steel   |                           |
| Environment  | 0 to 50°C (32 to 122°F);<br>RH max 95% RH  |                           |
| Dimensions   | 180 x 180 x 70 mm<br>(7.1 x 7.1 x 2.8")  |                           |
| Weight   | 1.4 kg (3.1 lb.)   |                           |

#### HI 302N & HI 312N **AUTO-REVERSE STIRRERS**

HI 302N normal stirrer and HI 312N autospeed stirrer allow to stir the fluid in both directions and are ideal for powder mixing applications to ensure a homogeneous final product and complete dissolution of hard-to-wet powders. The reverse stirring period is adjustable from 30 seconds to 3 minutes. When set to 1 minute, the fluid will rotate in one direction for minute and then reverse its rotation for another minute at the desired speed in a repeated cycle until the stirrer is switched off.

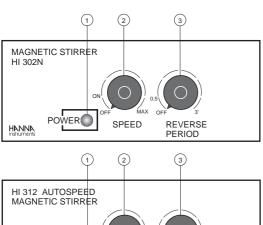

|                                   | 3)                                 |
|-----------------------------------|------------------------------------|
| HI 312 AUTOSPEE<br>MAGNETIC STIRR |                                    |
| HANNA POWER                       | ON 0.5 0.5 3' SPEED REVERSE PERIOD |


- 1. Power indicator (LED lights up when the stirrer is powered)
- 2. **ON/OFF/SPEED knob**, to turn the stirrer ON and OFF, and set the stirring speed (continuously variable up to the maximum)
- 3. REVERSE PERIOD knob with the ON/ OFF and continuously variable selection of inversion cycle time in minutes, from 0.5 to 3 minutes

|  | HI 302N  | HI312N                    |
|--|--|---------------------------|
| Maximum<br>Stirring<br>Capacity                      | 2.5 Liters<br>(0.66 gallons)   | 5 Liters<br>(1.3 gallons) |
| Speed Range<br>Min.<br>Max.                          | 100 rpm<br>800 to 1000 rpm   |                           |
| Reverse<br>Interval                                  | from 30 seconds<br>to 3 minutes  |                           |
| Auto-Speed   | not applicable   | standard                  |
| Power Supply HI 302N-1 HI 302N-2 HI 312N-1 HI 312N-2 | 110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz<br>110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz |                           |
| Installation<br>Category                             | П  |                           |
| Cover<br>Material                                    | AISI 316 stainless steel   |                           |
| Environment  | 0 to 50°C (32 to 122°F)<br>95% RH max.   |                           |
| Dimensions   | 180 x 180 x 70 mm<br>(7.1 x 7.1 x 2.8")  |                           |
| Weight   | 1.4 kg (3.1 lb.)   |                           |

### <u>HI 303N</u> <u>TWO-SPEED STIRRER WITH TACHO-</u> <u>METER</u>

 $\mbox{\bf HI~303N}$  stirrer incorporates an LCD tachometer.

This model allows precise stirring repeatability by using two separate ranges. In each range, the speed can be fine tuned for exacting accuracy.

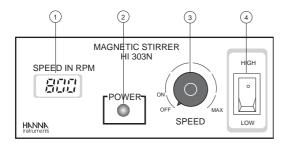


- 1. rpm LCD displays the rpm speed
- 2. **Power indicator** (LED lights up when the stirrer is powered)
- 3. **ON/OFF/SPEED knob**, to turn the stirrer ON and OFF, and set the stirring speed (continuously variable up to the maximum)
- 4. **HIGH/LOW switch** (for selecting low speed range or high speed range)

|                                  | HI 303N  |  |
|----------------------------------|--|--|
| Maximum<br>Stirring<br>Capacity  | 2.5 Liters<br>(0.66 gallons)                   |  |
| Speed Range<br>Low<br>High       | 100 to 500 rpm<br>100 to 1000 rpm              |  |
| Tachometer                       | 4-digit LCD                                    |  |
| Power Supply HI 303N-1 HI 303N-2 | 110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz |  |
| Installation<br>Category         | П  |  |
| Cover<br>Material                | AISI 316 Stainless Steel                       |  |
| Environment                      | 0 to 50°C (32 to 122°F);<br>RH max 95%         |  |
| Dimensions                       | 180 x 180 x 70 mm<br>(7.1 x 7.1 x 2.8")        |  |
| Weight                           | 1.4 kg (3.1 lb.)                               |  |

#### <u>HI 304N</u> <u>AUTO-REVERSE STIRRER WITH TACHO-</u> <u>METER</u>

**HI 304N** is a heavy-duty stirrer with a built-in tachometer. An advanced circuit allows to reverse the direction of the stir at an interval that can be adjusted from 30 seconds up to 3 minutes.

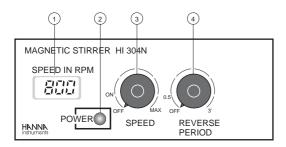



- 1. rpm LCD displays the rpm speed
- 2. **Power indicator** (LED lights up when the stirrer is powered)
- ON/OFF/SPEED knob, to turn the stirrer ON and OFF, and set the stirring speed (continuously variable up to the maximum)
- REVERSE PERIOD knob with the ON/OFF and continuously variable selection of inversion cycle time in minutes, from 0.5 to 3 minutes

|                                  | HI 304N  |
|----------------------------------|--|
| Maximum<br>Stirring<br>Capacity  | 2.5 Liters<br>(0.66 gallons)                   |
| Speed Range<br>Min.<br>Max.      | 100 rpm<br>800 to 1000 rpm                     |
| Tachometer                       | 4-digit LCD                                    |
| Power Supply HI 304N-1 HI 304N-2 | 110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz |
| Installation category            | П  |
| Cover<br>Material                | AISI 316 stainless steel                       |
| Environment                      | 0 to 50°C (32 to 122°F);<br>RH max 95% RH      |
| Dimensions                       | 180 x 180 x 70 mm<br>(7.1 x 7.1 x 2.8")        |
| Weight                           | 1.4 kg (3.1 lb.)                               |

#### HI 322N & HI 324N TIMER-CONTROLLED STIRRERS

HI 322N and HI 324N are autospeed stirrers equipped with timer control to start and automatically stop the stirring process after an user-selectable period (from 5 minutes to 2 hours). They are the ideal stirrers for routine applications where the required stirring time is known, and for non-Newtonian fluids whose viscosities change during stirring or Newtonian fluids up to 5 liters capacity.

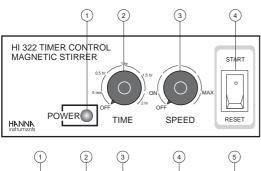

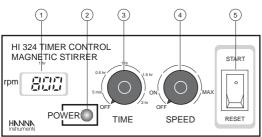

- 1. rpm LCD displays the rpm speed
- 2. **Power indicator** (LED lights up when the stirrer is powered)
- 3. **TIME knob** for setting the stirring time from 5 minutes to 2 hours
- 4. **SPEED knob**, to turn the stirrer ON and OFF, and set the stirring speed (continuously variable up to the maximum)
- 5. **START/RESET switch** for starting and resetting the time control

|  | HI322N   | HI324N                    |
|--|--|---------------------------|
| Maximum<br>Stirring<br>Capacity                      | 5 Liters<br>(1.3 gallons)  | 5 Liters<br>(1.3 gallons) |
| Speed Range<br>Min.<br>Max.                          | 100 rpm<br>800 to 1000 rpm   |                           |
| Timer Range  | user selectable from 5<br>minutes to 2 hours   |                           |
| Tachometer   | notapplicable  | 4-digit<br>LCD            |
| Auto-Speed   | standard   | standard                  |
| Power supply HI 322N-1 HI 322N-2 HI 324N-1 HI 324N-2 | 110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz<br>110/115 Vac, 50/60 Hz<br>220/240 Vac, 50/60 Hz |                           |
| Installation<br>Category                             | П  |                           |
| Cover<br>Material                                    | AISI 316 stainless steel   |                           |
| Environment  | 0 to 50°C (32 to 122°F);<br>RH max 95%   |                           |
| Dimensions   | 180 x 180 x 70 mm<br>(7.1 x 7.1 x 2.8")  |                           |
| Weight   | 1.4 kg (3.1 lb.)   |                           |

# OPERATIONAL GUIDE FOR HI 300N and HI 310N

- Make sure that the SPEED knob (on the front panel) and the ON/OFF switch (on the rear panel) are in OFF position.
- MAGNETIC STIRRER
  HI 190M
  POWER OF SPEED

  A O
- Place a beaker or glass container at the center of the top of the stirrer, and place the stirring bar in the beaker.
- MACNETIC STIRRER
  HI 190M
  POWER OF OUT
- Pour the liquids to be mixed into the beaker.

Do not fill up to the brim and leave sufficient space to prevent spillage during stirring.




- Connect the power cord to the mains and turn the stirrer on.
- Adjust the SPEED knob slowly until the required stirring speed is achieved.

20




 Allow sufficient time for the mixing operation to attain uniform rotational flow.

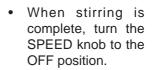




 If no further stirring is required, turn the stirrer off.

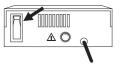

# OPERATIONAL GUIDE FOR HI 301N, HI 303N and HI 311N

 Make sure that the SPEED knob and the ON/OFF switch are in the OFF position.

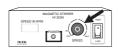




 Place a beaker or glass container at the center of the top of the stirrer and place the stirring bar in the beaker.



 Pour the liquids to be mixed into the beaker. Do not fill up to the brim and leave sufficient space to prevent spillage during stirring.



 Connect the power cord to the mains and turn the stirrer on with the ON/OFF switch.

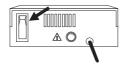

 Select the low or high range with the HIGH/ LOW switch as required.



 Adjust the SPEED knob slowly until the required stirring speed is achieved.




 Allow sufficient time for the mixing operation to attain uniform rotational flow.

The stirring speed is indicated in the rpm LCD (**HI 303N** only).

 When stirring is complete, turn the SPEED knob to the OFF position.





 If no further stirring is required, turn the stirrer off.

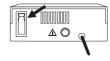

# OPERATIONAL GUIDE FOR HI 302N, H I304N and HI 312N

 Make sure that the SPEED knob and the ON/OFF switch are in the OFF position.

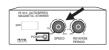

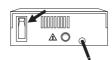


 Place a beaker or glass container at the center of the top of the stirrer and place the stirring bar in the beaker.




 Pour the liquids to be mixed into the beaker. Do not fill up to the brim and leave sufficient space to prevent spillage during stirring.



 Connect the power cord to the mains and turn the stirrer on with the ON/OFF switch.

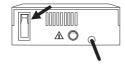

 Set the REVERSE PE-RIOD knob to the desired reverse stirring period from 30 seconds to 3 minutes.

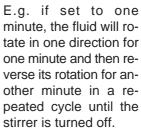




- Adjust the SPEED knob slowly until the required stirring speed is achieved.
- Allow sufficient time for the mixing operation to attain uniform rotational flow.
   The stirring speed is indicated in the rpm LCD for HI 304N.
- When stirring is complete, turn the SPEED knob to the OFF position.




 If no further stirring is required, turn the stirrer off.

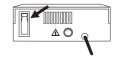

# OPERATIONAL GUIDE FOR HI 322N and HI 324N


- Make sure that the SPEED knob and the ON/OFF switch are in the OFF position and the START/RESET switch is in RESET position.
- Place a beaker or glass container at the center of the top of the stirrer and place the stirring bar in the beaker.
- Pour the liquids to be mixed into the beaker. Do not fill up to the brim and leave sufficient space to prevent spillage during stirring.
- Connect the power cord to the mains and turn the stirrer on with the ON/OFF switch.



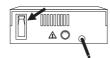






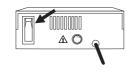

 Set the TIME knob to the desired stirring time from 5 minutes to 2 hours. Then set the START/RESET switch to START. The stirrer will only start stirring when the START switch is in the START position and will end when the selected stirring time has passed.

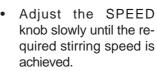






- Allow sufficient time for the mixing operation to attain uniform rotational flow.
   The stirring speed is indicated in the rpm LCD for HI 324N.
- When stirring is complete, turn the SPEED knob to the OFF position.



 If no further stirring is required, turn the stirrer off.

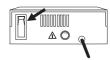

# **SALES AND TECHNICAL SERVICE CONTACTS**

Australia:

Tel. (03) 9769.0666 • Fax (03) 9769.0699

China:

Tel. (10) 88570068 • Fax (10) 88570060

**Egypt:** Tel. & Fax (02) 2758.683

**Germany:** Tel. (07851) 9129-0 • Fax (07851) 9129-99

Greece:

Tel. (210) 823.5192 • Fax (210) 884.0210

Indonesia:

Tel. (21) 4584.2941 • Fax (21) 4584.2942

Japan:

Tel. (03) 3258.9565 • Fax (03) 3258.9567

Korea:

Tel. (02) 2278.5147 • Fax (02) 2264.1729

Malaysia:

Tel. (603) 5638.9940 • Fax (603) 5638.9829

Singapore:

Tel. 6296.7118 • Fax 6291.6906

South Africa:

Tel. (011) 615.6076 • Fax (011) 615.8582

Taiwan:

Tel. 886.2.2739.3014 • Fax 886.2.2739.2983

Thailand:

Tel. 66.2619.0708 • Fax 66.2619.0061

**United Kingdom:** 

Tel. (01525) 850.855 • Fax (01525) 853.668

Tel. (401) 765.7500 • Fax (401) 765.7575

For e-mail contacts and complete list of Sales and Technical offices, please see www.hannainst.com