

# Operating Manual

**CB (E7)**

**CO<sub>2</sub> Incubators**

**CO<sub>2</sub> Incubators with O<sub>2</sub> Control**

**CO<sub>2</sub> Incubators with Active Humidification**

with sterilizable NDIR sensor system for CO<sub>2</sub>  
and microprocessor program controller MB2

Model	Model version	Voltage	Equipment	Art. no.
CB 170	CB170-230V	200-230 V		9040-0131, 9140-0131
CB 170	CB170-230V-O	200-230 V	O <sub>2</sub> control 0-20 vol.-%	9040-0132, 9140-0132
CB 170	CB170-230V-F	200-230 V	Active humidification	9040-0133, 9140-0133
CB 170	CB170-230V-FO	200-230 V	Active humidification O <sub>2</sub> control 0-20 vol.-%	9040-0134, 9140-0134
CB 170-UL	CB170-120V	100-120 V		9040-0139, 9140-0139
CB 170-UL	CB170-120V-O	100-120 V	O <sub>2</sub> control 0-20 vol.-%	9040-0140, 9140-0140
CB 170-UL	CB170-120V-F	100-120 V	Active humidification	9040-0141, 9140-0141
CB 170-UL	CB170-120V-FO	100-120 V	Active humidification O <sub>2</sub> control 0-20 vol.-%	9040-0142, 9140-0142

## **BINDER GmbH**

- Address: Post office box 102, 78502 Tuttlingen, Germany ► Phone: +49 7462 2005 0
- Fax: +49 7462 2005 100 ► Internet: <http://www.binder-world.com>
- E-mail: [info@binder-world.com](mailto:info@binder-world.com) ► Service Hotline: +49 7462 2005 555
- Service Fax: +49 7462 2005 93 555 ► Service E-Mail: [service@binder-world.com](mailto:service@binder-world.com)
- Service Hotline USA: +1 866 885 9794 or +1 631 224 4340 x3
- Service Hotline Asia Pacific: +852 390 705 04 or +852 390 705 03
- Service Hotline Russia and CIS: +7 495 988 15 16

## Contents

<b>1. SAFETY.....</b>	<b>7</b>
1.1 Legal considerations .....	7
1.2 Structure of the safety instructions.....	7
1.2.1 Signal word panel .....	7
1.2.2 Safety alert symbol .....	8
1.2.3 Pictograms .....	8
1.2.4 Word message panel structure .....	9
1.3 Localization / position of safety labels on the chamber .....	10
1.4 Type plate .....	11
1.5 General safety instructions on installing and operating the CO <sub>2</sub> incubator .....	12
1.6 Precautions when working with gases .....	13
1.7 Precautions when handling gas cylinders.....	15
1.8 Intended use .....	16
1.9 Measures to prevent accidents .....	17
1.10 Resistance of the humidity sensor against harmful substances .....	18
<b>2. CHAMBER DESCRIPTION .....</b>	<b>19</b>
2.1 Chamber overview .....	20
2.2 Inner chamber .....	21
2.3 Control panel on the rear of the chamber .....	22
2.4 Instrument panel .....	23
<b>3. COMPLETENESS OF DELIVERY, TRANSPORTATION, STORAGE, AND INSTALLATION.....</b>	<b>23</b>
3.1 Unpacking, and checking equipment and completeness of delivery .....	23
3.2 Guidelines for safe lifting and transportation.....	24
3.3 Storage.....	24
3.4 Location of installation and ambient conditions .....	25
<b>4. INSTALLATION AND CONNECTIONS .....</b>	<b>28</b>
4.1 Shelves.....	28
4.2 Permadry™ water pan .....	28
4.3 Connecting the O <sub>2</sub> sensor (chamber with O <sub>2</sub> control) .....	29
4.4 Gas connections .....	30
4.4.1 Connection of the CO <sub>2</sub> gas cylinder.....	31
4.4.2 Connection of the O <sub>2</sub> gas cylinder (chamber with O <sub>2</sub> control with optional alternative control range 10 up to 95 vol.-% O <sub>2</sub> ) .....	32
4.4.3 Connection of the N <sub>2</sub> gas cylinder (chamber with O <sub>2</sub> control).....	33
4.4.4 Connecting the gas hose to the chamber rear (for CO <sub>2</sub> , O <sub>2</sub> , N <sub>2</sub> ) .....	34
4.4.5 Gas cylinder connection kits (option).....	35
4.5 Water supply for the chamber with active humidification .....	36
4.5.1 Types of suitable water quality .....	36
4.5.2 Water supply installation .....	37
4.5.3 BINDER Pure Aqua Service (option) .....	37
4.6 Electrical connection .....	38
4.7 Handling and aligning the divided inner door, gas proof (optional equipment).....	39
<b>5. FUNCTIONAL OVERVIEW OF THE MB2 CHAMBER CONTROLLER.....</b>	<b>40</b>
5.1 Operating functions in normal display .....	41
5.2 Display views: Normal display, program display, chart-recorder display.....	42
5.3 Controller icons overview .....	43
5.4 Operating modes.....	45
5.5 Controller menu structure.....	46
5.5.1 Main menu .....	47
5.5.2 "Settings" submenu.....	48
5.5.3 "Service" submenu.....	48

5.6	Principle of controller entries.....	49
5.7	Performance during and after power failures.....	49
5.8	Performance when opening the door.....	50
<b>6.</b>	<b>START UP .....</b>	<b>50</b>
6.1	Turning on the chamber .....	50
6.2	Controller settings upon start up .....	51
6.3	Entering the altitude of the installation site .....	52
6.4	Factory settings.....	52
6.5	Equilibration time.....	53
6.6	Operating the DuoDoor™ door lock.....	53
6.7	Required gas supply for the chamber with O <sub>2</sub> control.....	54
6.7.1	Hypoxic control range 0.2 vol.-% up to 20 vol.-% O <sub>2</sub> (regular).....	54
6.7.2	Alternative control range 10 vol.-% up to 95 vol.-% (option) .....	54
6.7.3	Operation without O <sub>2</sub> control.....	54
6.8	Activating / deactivating CO <sub>2</sub> , O <sub>2</sub> and humidity control .....	55
6.9	Humidity control of the Permady™ system.....	56
<b>7.</b>	<b>SET-POINT ENTRY IN “FIXED VALUE” OPERATING MODE .....</b>	<b>57</b>
7.1	Set-point entry through the “Setpoints” menu .....	58
7.2	Direct setpoint entry via Normal display.....	59
7.3	Special controller functions via operation lines .....	60
7.4	Safety instructions for setting high gas concentrations.....	61
<b>8.</b>	<b>TIMER PROGRAM: STOPWATCH FUNCTION .....</b>	<b>62</b>
8.1	Starting a timer program .....	62
8.1.1	Performance during program delay time .....	62
8.2	Stopping a running timer program .....	63
8.2.1	Pausing a running timer program .....	63
8.2.2	Cancelling a running timer program.....	63
8.3	Performance after the end of the program.....	63
<b>9.</b>	<b>TIME PROGRAMS .....</b>	<b>64</b>
9.1	Starting an existing time program .....	64
9.1.1	Performance during program delay time .....	65
9.2	Stopping a running time program.....	65
9.2.1	Pausing a running time program.....	65
9.2.2	Cancelling a running time program.....	65
9.3	Performance after the end of the program.....	65
9.4	Creating a new time program.....	66
9.5	Program editor: program management.....	66
9.5.1	Deleting a time program .....	67
9.6	Section editor: section management.....	68
9.6.1	Add a new program section .....	69
9.6.2	Copy and insert or replace a program section.....	69
9.6.3	Deleting a program section.....	70
9.7	Value entry for a program section.....	71
9.7.1	Section duration.....	71
9.7.2	Set-point ramp and set-point step.....	72
9.7.3	Special controller functions via operation lines.....	74
9.7.4	Setpoint entry .....	75
9.7.5	Tolerance range.....	75
9.7.6	Repeating one or several sections within a time program.....	76
9.7.7	Saving the time program.....	77
<b>10.</b>	<b>WEEK PROGRAMS .....</b>	<b>78</b>
10.1	Starting an existing week program.....	78
10.2	Cancelling a running week program .....	78
10.3	Creating a new week program .....	79

10.4	Program editor: program management.....	80
10.4.1	Deleting a week program.....	81
10.5	Section editor: section management.....	82
10.5.1	Add a new program section .....	83
10.5.2	Copy and insert or replace a program section.....	83
10.5.3	Deleting a program section .....	84
10.6	Value entry for a program section.....	84
10.6.1	Set-point ramp and set-point step modes.....	84
10.6.2	Weekday .....	85
10.6.3	Start time.....	85
10.6.4	Setpoint entry.....	86
10.6.5	Special controller functions via operation lines.....	86
<b>11.</b>	<b>NOTIFICATION AND ALARM FUNCTIONS.....</b>	<b>87</b>
11.1	Notification and alarm messages overview.....	87
11.1.1	Notifications .....	87
11.1.2	Alarm messages .....	88
11.2	State of alarm.....	90
11.3	Resetting an alarm, list of active alarms .....	90
11.4	Activating / deactivating the audible alarm (alarm buzzer) .....	91
11.5	Tolerance range settings and alarm delay times .....	91
11.6	Zero-voltage relay alarm output.....	93
<b>12.</b>	<b>TEMPERATURE SAFETY DEVICES.....</b>	<b>94</b>
12.1	Over temperature protective device (class 1) .....	94
12.2	Overtemperature safety controller class 3.1 .....	94
12.2.1	Safety controller modes .....	95
12.2.2	Setting the safety controller .....	95
12.2.3	Message and measures in the state of alarm.....	96
12.2.4	Function check.....	96
<b>13.</b>	<b>USER MANAGEMENT.....</b>	<b>97</b>
13.1	Authorization levels and password protection.....	97
13.2	Log in.....	100
13.3	Log out .....	101
13.4	User change.....	101
13.5	Password assignment and password change .....	102
13.5.1	Password change .....	102
13.5.2	Deleting the password for an individual authorization level.....	104
13.5.3	New password assignment for "service" or "admin" authorization level when the password function was deactivated .....	105
13.6	Activation code.....	106
<b>14.</b>	<b>GENERAL CONTROLLER SETTINGS.....</b>	<b>107</b>
14.1	Selecting the controller's menu language .....	107
14.2	Setting date and time .....	107
14.3	Selecting the temperature unit.....	109
14.4	Display configuration.....	109
14.4.1	Adapting the display parameters .....	109
14.4.2	Touchscreen calibration.....	110
14.5	Network and communication.....	111
14.5.1	Serial RS485 interface (available via BINDER INDIVIDUAL Customized Solutions) .....	111
14.5.2	Ethernet .....	112
14.5.2.1	Configuration .....	112
14.5.2.2	Display of MAC address.....	113
14.5.3	Web server.....	113
14.5.4	E-Mail.....	114
14.6	USB menu: Data transfer via USB interface .....	115

<b>15. GENERAL INFORMATION .....</b>	<b>116</b>
15.1 Service contact page.....	116
15.2 Current operating parameters.....	116
15.3 Event list.....	117
15.4 Technical chamber information .....	117
<b>16. CHART RECORDER DISPLAY .....</b>	<b>118</b>
16.1 Views.....	118
16.1.1 Show and hide legend .....	118
16.1.2 Switch between legend pages .....	118
16.1.3 Show and hide specific indications .....	119
16.1.4 History display.....	119
16.2 Setting the parameters.....	122
<b>17. HUMIDIFICATION SYSTEM (CHAMBER WITH ACTIVE HUMIDIFICATION) ..</b>	<b>123</b>
17.1 Function of the humidifying system.....	123
<b>18. OPTIONS.....</b>	<b>124</b>
18.1 Communication software APT-COM™ 3 DataControlSystem (option).....	124
18.2 Silicone access ports 30 mm / 1.18 in, closable from both sides with silicon plugs (option) .....	124
18.3 Interior socket 230V (option, available via BINDER INDIVIDUAL Customized Solutions) .....	125
18.4 Analog outputs for temperature and CO <sub>2</sub> (option) .....	127
18.5 Data Logger kits (option).....	127
18.6 Access port for extra-low voltage (option).....	128
18.7 BINDER Gas Supply Service – External bottle changer for CO <sub>2</sub> , N <sub>2</sub> or O <sub>2</sub> (option) .....	129
18.8 Stands .....	129
18.8.1 Stacking stand (option) .....	129
18.8.2 Stacking adapter for direct thermal decoupled stacking (option) .....	129
18.8.3 Base on castors (option).....	130
<b>19. REFERENCE MEASUREMENTS .....</b>	<b>130</b>
19.1 CO <sub>2</sub> reference measuring .....	130
19.1.1 Measuring CO <sub>2</sub> concentration indirectly via the pH of the cell medium.....	130
19.1.2 Measuring CO <sub>2</sub> directly via chemical indicator tubes .....	131
19.1.3 Measuring CO <sub>2</sub> directly with an electronic infrared measuring device .....	131
19.2 Temperature reference measurement .....	131
<b>20. AVOIDING MICROBIAL CONTAMINATION.....</b>	<b>132</b>
20.1 Cells and media .....	132
20.2 Laboratory conditions / equipment around the incubator.....	132
20.3 Working and behavior in the lab .....	132
20.4 Chamber design and equipment of the CO <sub>2</sub> incubator .....	133
20.5 Handling the CO <sub>2</sub> incubator .....	134
<b>21. CLEANING, DECONTAMINATION / DISINFECTION, AND STERILIZATION ..</b>	<b>135</b>
21.1 Cleaning .....	135
21.2 Decontamination / chemical disinfection of the chamber.....	137
21.3 Hot-air sterilization at 180 °C / 356 °F.....	138
21.3.1 Overview .....	138
21.3.2 Preparation for a hot-air sterilization.....	139
21.3.3 Starting the sterilization cycle and running the hot-air sterilization.....	140
21.3.4 Prematurely terminating the sterilization cycle .....	141
21.3.4.1 Prematurely terminating the sterilization cycle after less than 4 hours: Ineffective sterilization .....	142
21.3.4.2 Prematurely terminating the sterilization cycle after more than 4 hours, i.e., during the cooling-down phase: successful sterilization .....	142
21.3.5 Completing the entire sterilization cycle .....	143



<b>22. MAINTENANCE AND SERVICE .....</b>	<b>144</b>
22.1 Maintenance intervals, service.....	144
22.2 Checking the humidity system fan .....	145
22.3 Gas inlet fine filter .....	145
22.4 Replacing the CO <sub>2</sub> sensor.....	145
22.5 Sending the chamber back to BINDER GmbH .....	146
<b>23. DISPOSAL .....</b>	<b>147</b>
23.1 Disposal of the transport packing.....	147
23.1.1 Outer chamber packing .....	147
23.1.2 Packing inside the chamber and equipment.....	147
23.2 Decommissioning.....	148
23.3 Disposal of the chamber in the Federal Republic of Germany .....	148
23.4 Disposal of the chamber in the member states of the EU except for the Federal Republic of Germany.....	149
23.5 Disposal of the chamber in non-member states of the EU .....	150
<b>24. TROUBLESHOOTING .....</b>	<b>151</b>
24.1 General.....	151
24.2 Heating .....	151
24.3 Gas cylinder pressure too low.....	152
24.4 Gas concentration .....	154
24.5 Sterilization.....	156
24.6 Humidity (chamber without active humidification).....	156
24.7 Humidity (chamber with active humidification).....	156
24.8 Controller.....	157
24.9 Open door .....	157
<b>25. TECHNICAL DESCRIPTION.....</b>	<b>158</b>
25.1 Factory calibration and adjustment .....	158
25.2 Over current protection .....	158
25.3 Definition of usable volume .....	159
25.4 CB technical data .....	159
25.5 Equipment and Options (extract) .....	161
25.6 Accessories and spare parts (extract) .....	163
25.7 Important conversion data for non-SI units .....	164
25.8 Conversion table for gas inlet pressures, bar – psi.....	164
25.9 Dimensions .....	165
<b>26. CERTIFICATES AND DECLARATIONS OF CONFORMITY.....</b>	<b>166</b>
26.1 EU Declaration of conformity .....	166
<b>27. PRODUCT REGISTRATION .....</b>	<b>168</b>
<b>28. CONTAMINATION CLEARANCE CERTIFICATE .....</b>	<b>169</b>
28.1 For chambers located outside the USA and Canada .....	169
28.2 For chambers in the USA and Canada.....	172

**Dear customer,**

For the correct operation of the CO<sub>2</sub> incubator CB, it is important that you read this operating manual completely and carefully and observe all instructions as indicated. Failure to read, understand and follow the instructions may result in personal injury. It can also lead to damage to the chamber and/or poor equipment performance

## 1. Safety

This operating manual is part of the components of delivery. Always keep it handy for reference. The device should only be operated by laboratory personnel especially trained for this purpose and familiar with all precautionary measures required for working in a laboratory. Observe the national regulations on minimum age of laboratory personnel. To avoid injuries and damage observe the safety instructions of the operating manual.

	 <b>WARNING</b>
	<p><b>Failure to observe the safety instructions.</b>  <b>Serious injuries and chamber damage.</b></p> <ul style="list-style-type: none"> <li>➤ Observe the safety instructions in this operating manual.</li> <li>➤ Carefully read the complete operating instructions of the chamber.</li> </ul>

### 1.1 Legal considerations

This operating manual is for informational purposes only. It contains information for installing, start-up, operation and maintenance of the product. Note: the contents and the product described are subject to change without notice.

Understanding and observing the instructions in this operating manual are prerequisites for hazard-free use and safety during operation and maintenance. In no event shall BINDER be held liable for any damages, direct or incidental arising out of or related to the use of this manual.

This operating manual cannot cover all conceivable applications. If you would like additional information, or if special problems arise that are not sufficiently addressed in this manual, please ask your dealer or contact us directly by phone at the number located on page one of this manual

Furthermore, we emphasize that the contents of this operating manual are not part of an earlier or existing agreement, description, or legal relationship, nor do they modify such a relationship. All obligations on the part of BINDER derive from the respective purchase contract, which also contains the entire and exclusively valid statement of warranty administration. The statements in this manual neither augment nor restrict the contractual warranty provisions.

### 1.2 Structure of the safety instructions

In this operating manual, the following safety definitions and symbols indicate dangerous situations following the harmonization of ISO 3864-2 and ANSI Z535.6.

#### 1.2.1 Signal word panel

Depending on the probability of serious consequences, potential dangers are identified with a signal word, the corresponding safety color, and if appropriate, the safety alert symbol.

 <b>DANGER</b>
<p>Indicates an imminently hazardous situation that, if not avoided, will result in death or serious (irreversible) injury.</p>





## WARNING

Indicates a potentially hazardous situation which, if not avoided, could result in death or serious (irreversible) injury.



## CAUTION

Indicates a potentially hazardous situation which, if not avoided, may result in moderate or minor (reversible) injury.

## CAUTION

Indicates a potentially hazardous situation which, if not avoided, may result in damage to the product and/or its functions or of a property in its proximity.
















### 1.2.2 Safety alert symbol














Use of the safety alert symbol indicates a **risk of injury**.

Observe all measures that are marked with the safety alert symbol in order to avoid death or injury.

### 1.2.3 Pictograms

Warning signs			
			
Electrical hazard	Hot surface	Explosive atmosphere	Stability hazard
			
Lifting hazard	Gas cylinders	Suffocation hazard	CO <sub>2</sub> suffocation and poisoning hazard
			
Explosive substances	Fire promoting agents	Harmful substances	Risk of corrosion and / or chemical burns
			
Biohazard	Pollution Hazard	Magnetic field	



Mandatory action signs			
			
Mandatory regulation	Read operating instructions	Disconnect the power plug	Lift with several persons
			
Environment protection	Wear protective gloves	Wear safety goggles	
Prohibition signs			
			
Do NOT touch	Do NOT spray with water	Do NOT climb	
	Information to be observed in order to ensure optimum function of the product.		

#### 1.2.4 Word message panel structure

**Type / cause of hazard.**





**Possible consequences.**

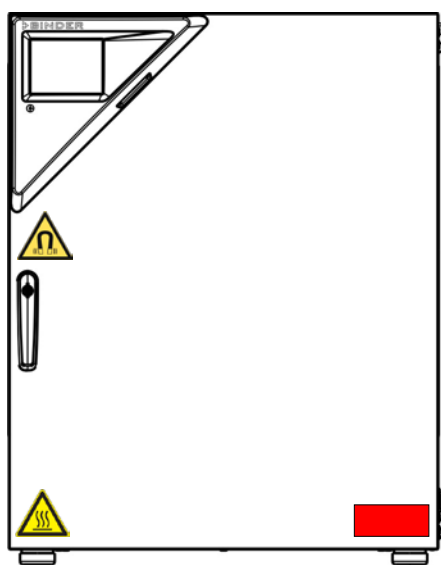
- Instruction how to avoid the hazard: prohibition
- Instruction how to avoid the hazard: mandatory action

Observe all other notes and information not necessarily emphasized in the same way, in order to avoid disruptions that could result in direct or indirect injury or property damage.


### 1.3 Localization / position of safety labels on the chamber

The following labels are located on the chamber:

Pictograms (Warning signs)	Service label
 <p>Hot surface</p> <ul style="list-style-type: none"> <li>on the outer chamber door</li> </ul>	
 <p>Risk of injury</p> <ul style="list-style-type: none"> <li>on the outer door: CB-UL only</li> <li>above the access ports (option)</li> </ul>	
 <p>Magnetic field</p> <ul style="list-style-type: none"> <li>on the outer door above the door handle</li> </ul>	



CO<sub>2</sub> incubator CB

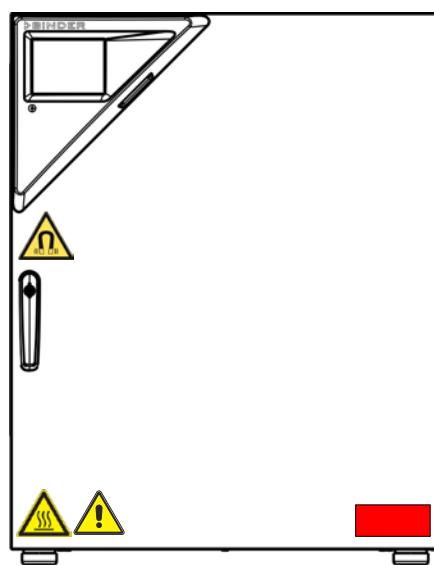

CO<sub>2</sub> incubator CB-UL

Figure 1: Position of labels on the chamber



Keep safety labels complete and legible.

Replace safety labels that are no longer legible. Contact BINDER service for these replacements.

## 1.4 Type plate

Position of type plate: left chamber side (seen from front), at the bottom in the middle

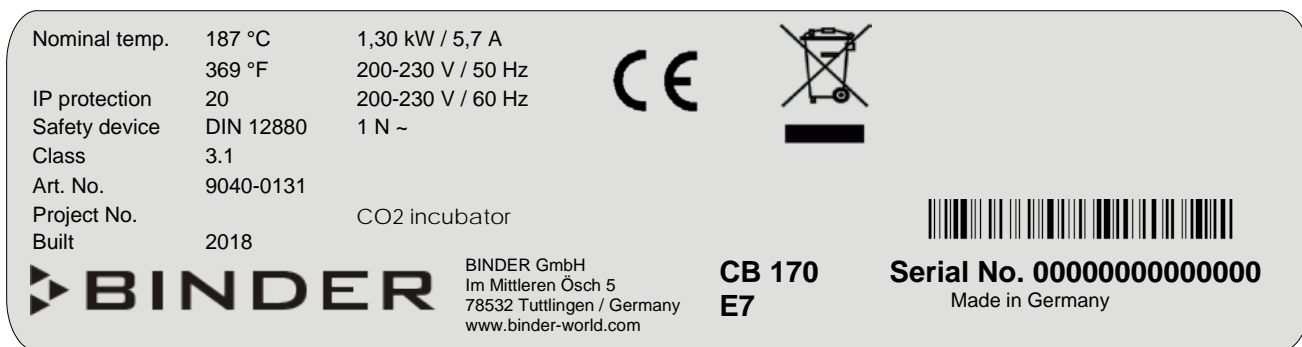




Figure 2: Type plate (example of CB 170 regular chamber)

Indications of the type plate (example)		Information
BINDER		Manufacturer: BINDER GmbH
CB 170		Model designation
CO2 incubator		Device name
Serial No.	0000000000000000	Serial no. of the chamber
Built	2018	Year of construction
Nominal temperature	187 °C 369 °F	Nominal temperature
IP protection	20	IP type of protection acc. to standard EN 60529
Temp. safety device	DIN 12880	Temperature safety device acc. to standard DIN 12880
Class	3.1	Class of temperature safety device
Art. No.	9040-0131	Art. No. of the chamber
Project No.	---	Optional: Special application acc. to project no.
1,30 kW		Nominal power
200-230 V / 50 Hz		Nominal voltage range +/-10% at the indicated power frequency
200-230 V / 60 Hz		
1 N ~		Current type
5,7 A		Nominal current


Symbol on the type plate	Information
	CE conformity marking
	Electrical and electronic equipment manufactured / placed on the market in the EC after 13 August 2005 and to be disposed of in separate collection according to Directive 2012/19/EU on waste electrical and electronic equipment (WEEE).

## 1.5 General safety instructions on installing and operating the CO<sub>2</sub> incubator



With regard to operating the chamber and to the installation location, please observe the guideline BGI/GUV-I 850-0 on safe working in laboratories (formerly BGR/GUV-R 120 or ZH 1/119 laboratory guidelines issued by the employers' liability insurance association) (for Germany).

BINDER GmbH is only responsible for the safety features of the chamber provided skilled electricians or qualified personnel authorized by BINDER perform all maintenance and repair, and if components relating to chamber safety are replaced in the event of failure with original spare parts.



To operate the chamber, use only original BINDER accessories or accessories from third-party suppliers authorized by BINDER. The user is responsible for any risk caused by using unauthorized accessories.

	<b>CAUTION</b>
	<p><b>Danger of overheating.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT install the chamber in unventilated recesses.</li> <li>➤ Ensure sufficient ventilation for dispersal of the heat.</li> </ul>

Do not operate the chamber in hazardous locations.



	 <b>DANGER</b>
	<p><b>Explosion hazard.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT operate the chamber in potentially explosive areas.</li> <li>➤ KEEP explosive dust or air-solvent mixtures AWAY from the chamber.</li> </ul>

The chamber does not dispose of any measures of explosion protection.

	 <b>DANGER</b>
	<p><b>Explosion hazard.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT introduce any substance into the chamber which is combustible or explosive at working temperature.</li> <li>⊘ NO explosive dust or air-solvent mixture in the inner chamber.</li> </ul>



Any solvent contained in the charging material must not be explosive or inflammable. I.e., irrespective of the solvent concentration in the steam room, NO explosive mixture with air must form. The temperature inside the chamber must lie below the flash point or below the sublimation point of the charging material. Familiarize yourself with the physical and chemical properties of the charging material, as well as the contained moisture constituent and its behavior with the addition of heat energy and humidity (chamber with active humidification).




Familiarize yourself with any potential health risks caused by the charging material, the contained moisture constituent or by reaction products which may arise during the temperature process. Take adequate measures to exclude such risks prior to putting the chamber into operation.

	 <b>DANGER</b>
	<p><b>Electrical hazard.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>⊘ The chamber must NOT become wet during operation or maintenance.</li> </ul>

The chambers were produced in accordance with VDE regulations and were routinely tested in accordance to VDE 0411-1 (IEC 61010-1).

During and after a sterilization the temperature of the inner surfaces almost equals the set-point.



	 <b>CAUTION</b>
	<p><b>The inner doors, the inner door and glass door handles, and the inner chamber will become hot during a sterilization.</b></p> <p><b>Danger of burning.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT touch the inner doors, the inner door and glass door handles, and the inner surfaces after a sterilization.</li> </ul>

 	 <b>WARNING</b>
	<p><b>Stability hazard.</b></p> <p><b>Danger of injury.</b></p> <p><b>Damage to the chamber and the charging material.</b></p> <p><b>Housing cover breakaway.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT climb on the lower housing cover.</li> <li>⊘ Do NOT load the lower housing cover with heavy objects while the chamber door is open.</li> </ul>

## 1.6 Precautions when working with gases




### Notes on handling carbon dioxide (CO<sub>2</sub>)

Carbon dioxide (CO<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Vent out any CO<sub>2</sub> gas that may escape via good room ventilation or a suitable connection to an exhaust system. We recommend installing a CO<sub>2</sub> warning system


	 <b>WARNING</b>
	<p><b>High concentration of CO<sub>2</sub> (&gt; 4 Vol.-%).</b></p> <p><b>Risk of death by suffocation.</b></p> <p><b>Danger of poisoning.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling CO<sub>2</sub>.</li> </ul>

### Chamber with O<sub>2</sub> control: Notes on handling oxygen (O<sub>2</sub>)

Oxygen (O<sub>2</sub>) is colorless and almost odorless and therefore practically imperceptible. It promotes burns, which can proceed explosively. There is a fire hazard for flammable oxygenated materials, e.g. clothes and hair. O<sub>2</sub> is heavier than air and may accumulate in low-lying areas.



 	 <b>WARNING</b>
	<p><b>High concentration of O<sub>2</sub> (&gt; 21 % O<sub>2</sub>).</b></p> <p><b>Fire and explosion hazard through contact of combustible materials with O<sub>2</sub>.</b></p> <p><b>Risk of burns and other injuries.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling O<sub>2</sub>.</li> </ul>

Take appropriate measures to prevent oxygen enrichment and fire and explosion hazards in areas where oxygen enrichment is possible.


	<p>General information for safe handling of oxygen:</p> <ul style="list-style-type: none"> <li>• Make sure training of personnel on hazards of oxygen enrichment and necessary safety measures.</li> <li>• Make sure adequate labeling of all oxygen equipment and facilities.</li> <li>• Make sure gas tightness of all gas connections by checking them for leaks (e.g. with leak spray or diluted soap solution).</li> <li>• Close the main valve of the source of oxygen after work when not using the chamber.</li> <li>• Never lubricate O<sub>2</sub> equipment with oil or fat. Use only materials and spare parts which are approved for use with oxygen.</li> <li>• Regularly inspect fire extinguishers for proper condition.</li> <li>• Set up emergency showers where oxygen enrichment is possible.</li> <li>• Strictest smoking ban and no ignition sources in areas where oxygen enrichment is possible.</li> <li>• Make sure good ventilation of areas where oxygen enrichment is possible (location of the chamber and/or O<sub>2</sub> cylinders).</li> <li>• Persons who may have been in a possibly oxygen-enriched atmosphere must keep away from ignition sources (flames, cigarettes, etc.) and ventilate their clothes at least 15 minutes.</li> <li>• Always keep emergency routes free.</li> </ul>
---	---

#### **Chamber with O<sub>2</sub> control: Notes on handling nitrogen (N<sub>2</sub>)**



Nitrogen (N<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Any N<sub>2</sub> gas that may escape must be safely led out via good room ventilation or a suitable connection to an exhaust system.

	<div data-bbox="858 1265 1136 1332">  <b>WARNING</b> </div> <p><b>High concentration of N<sub>2</sub>.</b>  <b>Risk of death by suffocation.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling N<sub>2</sub>.</li> </ul>
---	--



## 1.7 Precautions when handling gas cylinders


	<p>General information for safe handling of gas cylinders:</p> <ul style="list-style-type: none"> <li>• Store and use gas cylinders only in well-ventilated locations.</li> <li>• Open the gas cylinder valve slowly to avoid pressure surges.</li> <li>• Secure gas cylinders during storage and use against falling (chaining).</li> <li>• Transport gas cylinders with a cylinder cart, do not carry, roll, or throw them.</li> <li>• Always close the valve even with apparently empty cylinders; screw on the cap when not in use. Return gas cylinders with the valve closed.</li> <li>• Do not open gas cylinders by force. Mark them when damaged.</li> <li>• Protect gas cylinders against fire, e.g. do not store together with flammable liquids.</li> <li>• Observe relevant regulations for dealing with gas cylinders.</li> </ul>
---	---

Secure the gas cylinders against falling and other mechanical damage.

	<div data-bbox="842 757 1125 824">  <b>WARNING</b> </div> <p><b>Safety valve tearing off.</b>  <b>Sudden release of the stored pressure energy.</b>  <b>Risk of injury.</b></p> <ul style="list-style-type: none"> <li>➤ Secure gas cylinders against falling (chaining).</li> <li>➤ Transport gas cylinders with a cylinder cart.</li> </ul>
---	--

The valve of the gas cylinder **always** must be closed before screwing on or unscrewing the gas hose.

	<div data-bbox="842 1104 1125 1171">  <b>WARNING</b> </div> <p><b>Opening the cylinder valve when the cylinder is not connected.</b>  <b>Sudden release of the stored pressure energy.</b>  <b>Risk of injury.</b></p> <ul style="list-style-type: none"> <li>➤ Close the gas cylinder valve before connecting or removing the gas hose.</li> </ul>
---	--

	<p>After connecting the gas cylinder, check all gas connections for leaks (e.g. with leak spray or diluted soap solution).</p>
---	--






## 1.8 Intended use


Series CB incubators are suitable for the cultivation of mammal cells under typical conditions of approx. 37 °C / 98.6°F. The chamber permits setting defined pH conditions by common NaHCO<sub>3</sub> buffer systems of commercial cell media by keeping an exact CO<sub>2</sub> atmosphere inside. The chambers guarantee high humidity inside to avoid osmolality increasing caused by the evaporation of the cell media.

With the chamber with O<sub>2</sub> control, a variable oxygen atmosphere can additionally influence the growth of the cells.


The chambers are suitable for exact conditioning of harmless materials. Any possible solvent any solvent must not be explosive and flammable. Components of the charging material must NOT form an explosive mixture with air. The operating temperature must lie below the flash point or below the sublimation point of the charging material. Any component of the charging material must NOT be able to release toxic gases.


 	<div style="background-color: red; color: white; padding: 5px; text-align: center;">  <b>DANGER</b> </div> <p><b>Explosion or implosion hazard.</b></p> <p><b>Danger of poisoning.</b></p> <p><b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT introduce any substance combustible or explosive at working temperature into the chamber, in particular no energy sources such as batteries or lithium-ion batteries.</li> <li>Ø NO explosive dust or air-solvent mixture in the inner chamber.</li> <li>Ø Do NOT introduce any substance which could lead to release of toxic gases.</li> </ul>
---	---


**Other applications are not approved.**

	<p>Following the instructions in this operating manual and conducting regular maintenance work (chap. 22.1) are part of the intended use.</p>
---	---

**The chambers are not classified as medical devices as defined by the Medical Device Directive 93/42/EEC.**

	<p>Due to the special demands of the Medical Device Directive (MDD), these chambers are not qualified for sterilization of medical devices as defined by the directive 93/42/EEG.</p>
---	---

	<p><b>WARNING:</b> If customer should use a BINDER chamber running in non-supervised continuous operation, we strongly recommend in case of inclusion of irrecoverable specimen or samples to split such specimen or samples and store them in at least two chambers, if this is feasible.</p>
---	--

	<p>The charging material shall not contain any corrosive ingredients that may damage the machine components made of stainless steel, aluminum, and copper. Such ingredients include in particular acids and halides. Any corrosive damage caused by such ingredients is excluded from liability by BINDER GmbH.</p>
---	---

In case of foreseeable use of the chamber there is no risk for the user through the integration of the chamber into systems or by special environmental or operating conditions in the sense of EN 61010-1:2010. For this, the intended use of the chamber and all its connections must be observed.

## 1.9 Measures to prevent accidents

The manufacturer took the following measures to prevent dangers:

- **Indications on the type plate**  
See operating manual chap. 1.4.
- **Operating manual**  
An operating manual is available for each chamber.
- **Overtemperature monitoring**  
The chamber is equipped with a temperature display, which can be read from outside.  
The chamber is equipped with an additional safety controller (temperature safety device class 3.1 acc. to DIN 12880:2007). Visual and audible (buzzer) signals indicate temperature exceeding.
- **Safety, measurement, and control equipment**  
The safety, measuring, and control equipment is easily accessible.
- **Electrostatic charge**  
The interior parts are grounded.
- **Non-ionizing radiation**  
Non-ionizing radiation is not intentionally produced, but released only for technical reasons by electrical equipment (e.g. power cables). The machine is equipped with a permanent magnet located behind the inner panel of the outer chamber door. If persons with active implants (e.g. pacemakers, defibrillators) keep a safe distance (distance of field source to implant) of 30 cm, an influence of these implants can be excluded with high probability.
- **Protection against touchable surfaces**  
Tested according to EN ISO 13732-1:2008.
- **Floors**  
See operating manual chap. 3.4 for correct installation
- **Cleaning**  
See operating manual chap. 21.

## 1.10 Resistance of the humidity sensor against harmful substances

The following list of harmful substances refers only to the humidity sensor and does not include any other materials incorporated in the chamber or prohibited substances in relation to explosion protection.

Some gases - especially clean gases - do not have any influence on the humidity sensor. Others do have a very small influence, whereas others may influence the sensor to a larger extent.

- The following gases do not influence the sensor and the humidity measurement: Argon (Ar), carbon dioxide (CO<sub>2</sub>), helium (He), hydrogen (H<sub>2</sub>), neon (Ne), nitrogen (N<sub>2</sub>), nitrous oxide (N<sub>2</sub>O), oxygen (O<sub>2</sub>)
- The following gases do not or to a minor extent influence the sensor and the humidity measurement: Butane (C<sub>4</sub>H<sub>10</sub>), ethane (C<sub>2</sub>H<sub>6</sub>), methane (CH<sub>4</sub>), natural gas propane (C<sub>3</sub>H<sub>8</sub>)
- The following gases do not, or to a minor extent influence the sensor and the humidity measurement, provided that the indicated loads are not exceeded:

		Maximum work place threshold limit value		Tolerated concentration with permanent load	
Substance	Formula	ppm	mg/m <sup>3</sup>	ppm	mg/m <sup>3</sup>
Ammonia	NH <sub>3</sub>	20	14	5500	4000
Acetone	CH <sub>3</sub> COCH <sub>3</sub>	500	1200	3300	8000
Benzene		300	1200		150000
Chlorine	Cl <sub>2</sub>	0.5	1.5	0.7	2
Acetic acid	CH <sub>3</sub> COOH	10	25	800	2000
Ethyl acetate	CH <sub>3</sub> COOC <sub>2</sub> H <sub>5</sub>	400	1400	4000	15000
Ethanol	C <sub>2</sub> H <sub>5</sub> OH	500	960	3500	6000
Ethylene glycol	HOCH <sub>2</sub> CH <sub>2</sub> OH	10	26	1200	3000
Formaldehyde	HCHO	0.3	0.37	2400	3000
Isopropanol	(CH <sub>3</sub> ) <sub>2</sub> CHOH	200	500	4800	12000
Methanol	CH <sub>3</sub> OH	200	260	3500	6000
Methyl ethyl ketone	C <sub>2</sub> H <sub>5</sub> COCH <sub>3</sub>	200	590	3300	8000
Ozone	O <sub>3</sub>	0.1	0.2	0.5	1
Hydrochloric acid	HCl	2	3	300	500
Hydrogen sulphide	H <sub>2</sub> S	10	15	350	500
Nitrogen oxides	NO <sub>x</sub>	5	9	5	9
Sulphur dioxide	SO <sub>2</sub>	5	13	5	13
Toluol	C <sub>6</sub> H <sub>5</sub> CH <sub>3</sub>	100	380	1300	5000
Xylene	C <sub>6</sub> H <sub>4</sub> (CH <sub>3</sub> ) <sub>2</sub>	100	440	1300	5000

These values are to be considered as approximate values. The sensor resistance largely depends on the temperature and humidity conditions during the time of exposure to harmful substances. Avoid simultaneous condensation. Tolerated error of measurement: ± 2 %r.H. The maximum work place threshold limit value is one that can be regarded as harmless for humans.

- Vapors of oil and fat are dangerous for the sensor because they may condensate at the sensor and thus prevent its function (insulating layer). For similar reasons it is not possible to measure smoke gases.

## 2. Chamber description

The CO<sub>2</sub> incubators CB are equipped with a multifunctional microprocessor display controller for temperature, CO<sub>2</sub>, and O<sub>2</sub> (chamber with O<sub>2</sub> control) levels and humidity (chamber with active humidification) and a digital display accurate to one-tenth of a degree resp. 0.1 vol.-%. With its comprehensive program control functions, the display program controller MB2 permits the high precision performance of temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control), and humidity (chamber with active humidification) cycles.

**Material:** The inner chamber, the pre-heating chamber and the inside of the doors are all made of stainless steel V2A (German material no. 1.4301, US equivalent AISI 304). The inner surfaces are smooth and therefore easy to clean. The inner chamber is deep-drawn from one piece, polished (suitable for pharmaceutical applications) and has no welds or inaccessible corners. The hinges and the seal of the inner glass door are glued from the outside to aid cleaning of the inner chamber. When operating the chamber at high temperatures (sterilization), the impact of the oxygen in the air may cause discoloration of the metallic surfaces (yellowish-brown or blue) by natural oxidation processes. These colorations are harmless and will in no way impair the function or quality of the chamber. The perforated shelves are also made of stainless steel. You can insert a maximum of 3 (CB 56), 6 (CB 170), resp. 8 (CB 220) shelves. The housing is RAL 7035 powder-coated. All corners and edges are also completely coated.

**Door lock:** The DuoDoor™ door lock offers two to open the outer door and glass door independently or together. The outer door is regularly equipped with a door lock with keys.

**Sterilization:** The chamber's heating system permits hot-air auto-sterilization at a setpoint of 187.5 °C / 369.5°F. Thus, a temperature of 180 °C / 356°F is maintained for at least 30 minutes on all internal surfaces, resulting in sterilization of the entire inner chamber. Therefore, this procedure meets all international guidelines regarding hot air sterilization, e.g. AAMI ST63, DIN 58947, European Pharmacopoeia.

**Temperature safety device** Thanks to the regular safety controller (temperature safety device class 3.1 acc. to DIN 12880:2007), the set temperature is maintained in case of failure.

**CO<sub>2</sub> system:** A highly precise, drift-free CO<sub>2</sub> infrared measuring system in combination with the permanent mixture of CO<sub>2</sub> gas through a special proprietary gas mixing head developed by BINDER allows precise and constant CO<sub>2</sub> concentrations for long periods. This creates optimum growth conditions for cultures. The gas enters the chamber via a fine filter (aseptic filter) with a high filtration efficiency that also filters the smallest particles.

Fast reaction times, maximum accuracy and selectivity characterize the CO<sub>2</sub> measuring procedure of the CB incubator series. The accuracy of the CO<sub>2</sub> measuring system is based on an infrared measuring cell with NDIR (non-dispersive infrared) sensor, which continuously regulates to a reference value. Therefore, disturbance variables and aging phenomena in the measuring system are almost completely eliminated, so that this measuring system, in contrast to other measuring procedures, remains practically drift-free between calibrations and is entirely selective for CO<sub>2</sub>. The sensor is built into the chamber and can be sterilized.

The accuracy of the indicated values of CO<sub>2</sub> depends on the ambient air pressure (approx. 0.08 vol.-% per 10 mbar / 0.15 psi). To compensate for this effect in the CO<sub>2</sub> measurement, the altitude of the installation site is entered into the controller.

**O<sub>2</sub> control (chamber with O<sub>2</sub> control):** The CO<sub>2</sub> incubator is available with O<sub>2</sub> control in addition to CO<sub>2</sub> control. There are two different control ranges:

- Regular equipment: Hypoxic control range **0.2 to 20 vol. % O<sub>2</sub>**. Only N<sub>2</sub> can be connected to reduce O<sub>2</sub> concentration; it is not possible to connect O<sub>2</sub> gas bottles to increase O<sub>2</sub> concentration. Control in the low O<sub>2</sub> range is very precise, in particular in the range below 1 vol. % O<sub>2</sub>.
- Alternative control range **10 to 95 vol. % O<sub>2</sub>** (option). Although the high control range is intended in particular for hyperoxic applications (> 21 vol. % O<sub>2</sub>), it is also suitable for slightly hypoxic applications between 10 and 20 vol. % O<sub>2</sub>.

The O<sub>2</sub> sensor is a semiconductor gas sensor with ZrO<sub>2</sub> ceramic.

**Humidity control (chamber with active humidification):** The chamber is available with humidity control and a microprocessor-controlled humidification system.

Freshwater is supplied by manually filling a freshwater bag, which is placed behind the chamber door in a recess.

A resistance humidifying system humidifies the air. For this purpose, use deionized (demineralized) water. The option BINDER Pure Aqua Service allows using the chamber with any degree of water hardness.

**Controller:** The efficient program controller is equipped with a multitude of operating functions, in addition to recorder and alarm functions. Programming of test cycles is easily accomplished via the modern MB2 touch screen controller and is also possible directly with a computer via Intranet in connection with the BINDER communication software APT-COM™ 3 DataControlSystem (option, chap. 18.1). The chambers are equipped with an Ethernet interface for computer communication. The APT-COM™ 3 software permits networking up to 30 chambers and connecting them to a PC for controlling and programming, as well as recording and representing temperature, CO<sub>2</sub> and O<sub>2</sub> and humidity data. For further options, see chap. 25.5.

Temperature range:	4 °C / 7.2 °F above ambient temperature up to +60 °C / 140 °F
Temperature range (chamber with O <sub>2</sub> control):	6 °C / 10.8 °F above ambient temperature up to +60 °C / 140 °F
CO <sub>2</sub> range:	0 vol.-% up to 20 vol.-%
O <sub>2</sub> range (chamber with O <sub>2</sub> control):	0.2 vol.-% up to 20 vol.-% (hypoxic control range) or 10 vol.-% up to 95 vol.-% (optional alternative control range)
Humidity range (chamber with active humidification):	60 % r.h. up to 93 % r.h.

## 2.1 Chamber overview

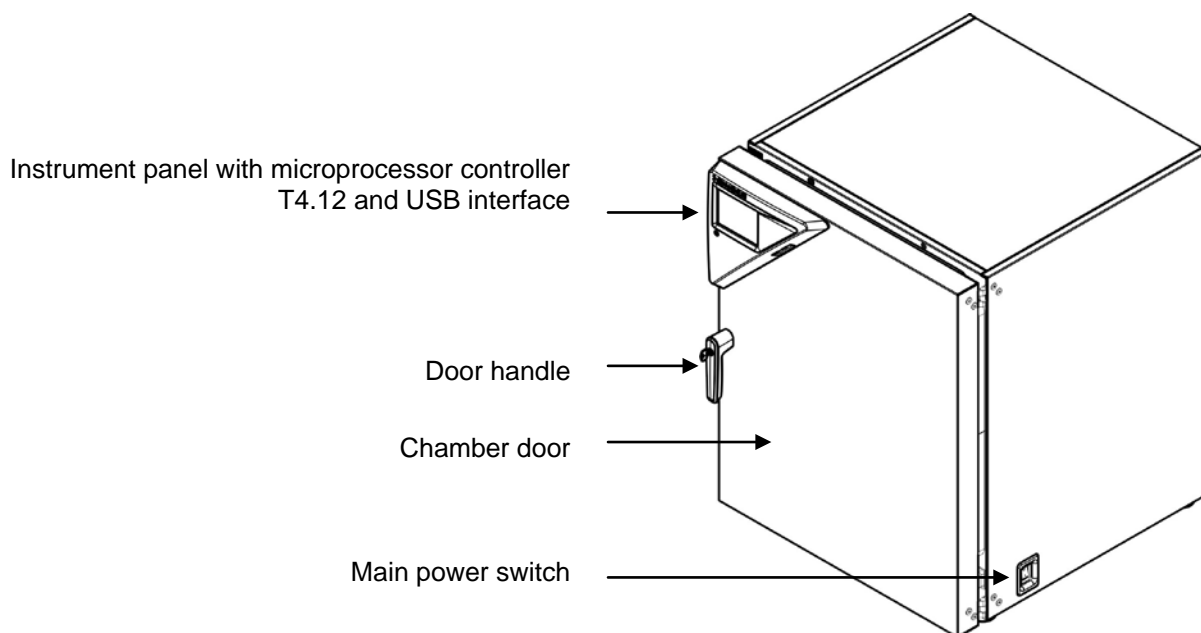


Figure 3: CO<sub>2</sub> incubator CB (example: model CB 170)

## 2.2 Inner chamber

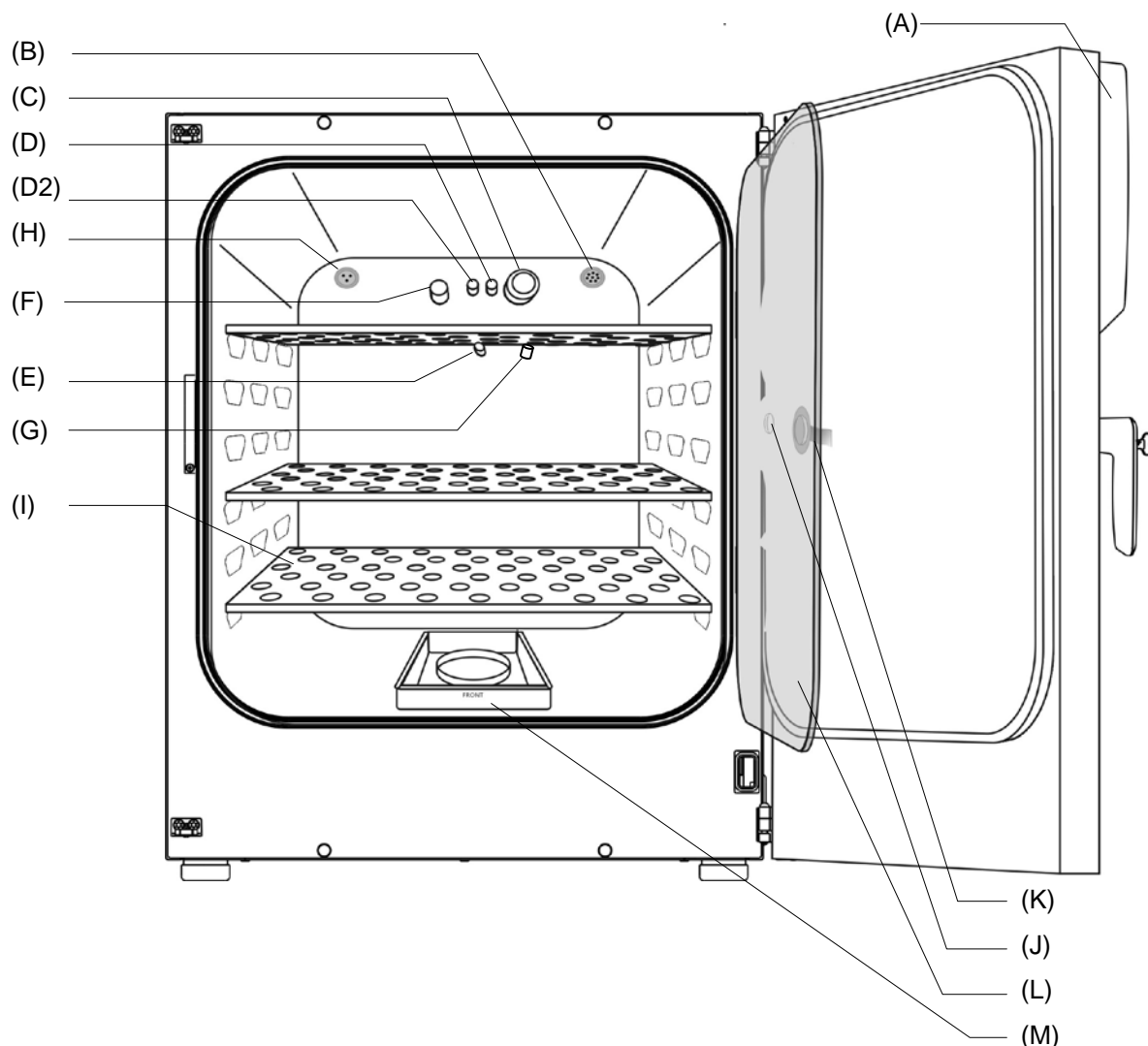


Figure 4: CB 170 with O<sub>2</sub> control and options

- (A) Instrument panel with microprocessor program controller MB2, indicating temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification)
- (B) Connection socket for extra-low voltage supply (option, chap. 18.6)
- (C) CO<sub>2</sub> sensor
- (D) Gas mixing head CO<sub>2</sub>
- (D2) Additional gas mixing head O<sub>2</sub>/ N<sub>2</sub> (chamber with O<sub>2</sub> control)
- (E) Pt 100 temperature probe
- (F) O<sub>2</sub> sensor (chamber with O<sub>2</sub> control)
- (G) Humidity sensor (chamber with active humidification)
- (H) Internal socket 230V (max. 3 A) (option, available via BINDER INDIVIDUAL Customized Solutions, chap. 18.3)
- (I) Perforated shelves, made of stainless steel
- (J) Glass door handle
- (K) Measuring access port
- (L) Inner glass doors
- (M) Permadry™ water pans

## 2.3 Control panel on the rear of the chamber

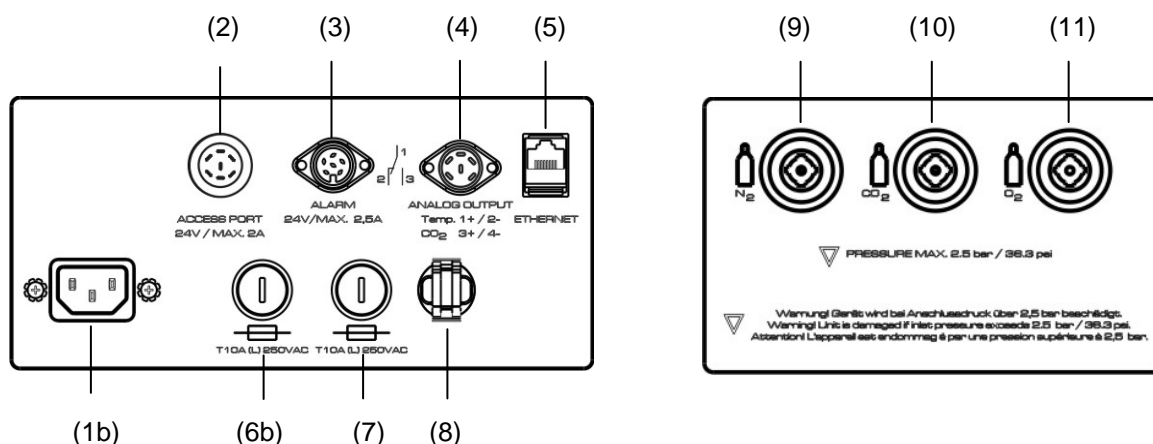


Figure 5: Rear control panel CB with O<sub>2</sub> control and options


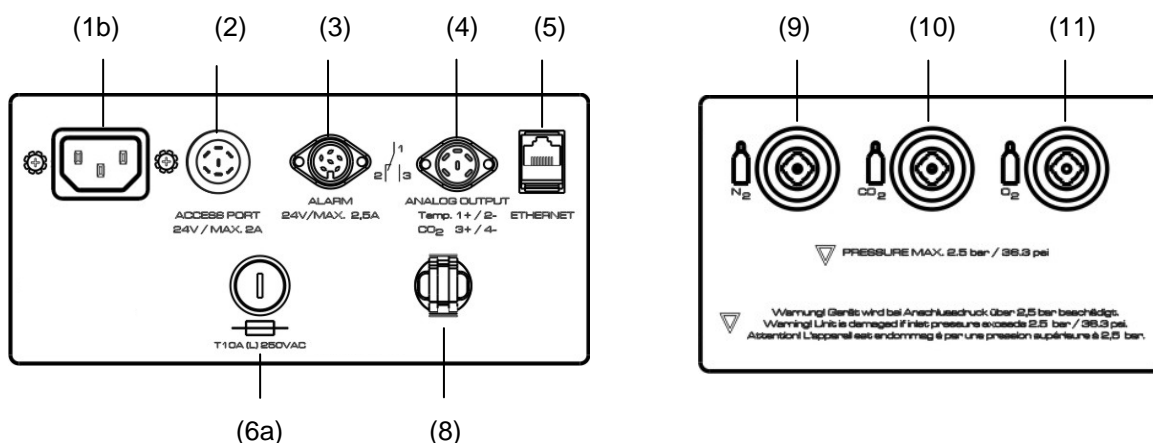


Figure 6: Rear control panel CB-UL with O<sub>2</sub> control and options

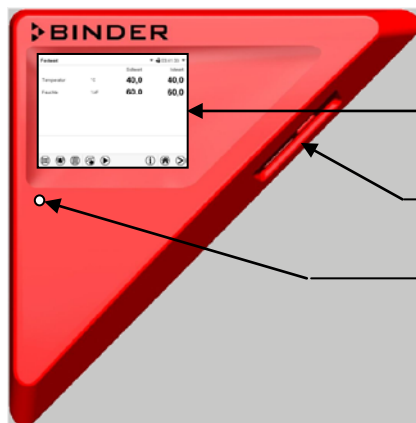- (1a) Socket for IEC connector plug / power cable 100-120 V AC
- (1b) Socket for IEC connector plug / power cable 230 V AC
- (2) External socket for extra-low voltage supply (option for CB 170 / CB 220, chap. 18.6)
- (3) DIN-socket for zero-voltage relay alarm outputs
- (4) DIN socket for analog outputs 4-20 mA (option, chap. 18.4)
- (5) Ethernet interface for computer communication
- (6a) Miniature fuse T12,5 A (L) 250 V AV for 100-120 V chamber
- (6b) Miniature fuse T10 A (L) 250 V AC for 200-230 V chamber
- (7) Miniature fuse T10 A (L) 250 V AC for 200-230 V chamber
- (8) Strain relief for power cable
- (9) Quick acting closure socket for N<sub>2</sub> (chamber with O<sub>2</sub> control)
- (10) Quick acting closure socket for CO<sub>2</sub>
- (11) Quick acting closure socket for O<sub>2</sub> (chamber with O<sub>2</sub> control with optional alternative control range 10 up to 95 vol.-% O<sub>2</sub>)



## 2.4 Instrument panel



MB2 controller display  
5,7" with touchscreen

USB interface

Pilot lamp

Figure 7: Instrument panel with MB2 program controller and USB interface

## 3. Completeness of delivery, transportation, storage, and installation

### 3.1 Unpacking, and checking equipment and completeness of delivery

After unpacking, please check the chamber and its optional accessories, if any, based on the delivery receipt for completeness and for transportation damage. Inform the carrier immediately if transportation damage has occurred.

The final tests of the manufacturer may have caused traces of the shelves on the inner surfaces. This has no impact on the function and performance of the chamber.

Please remove any transportation protection devices and adhesives in/on the chamber and on the doors and remove the operating manuals and accessory equipment.



Remove any protective lamination sheet on the inner metal surfaces prior to commissioning.



**CAUTION**

**Sliding or tilting of the chamber.**

**Damage to the chamber.**

**Risk of injury by lifting heavy loads.**

- Do NOT lift or transport the chamber using the door handle, the door or the lower housing.
- Lift the chamber from the pallet at the four lower corners with the aid of four people.

If you need to return the chamber, please use the original packing and observe the guidelines for safe lifting and transportation (chap. 3.2).

For disposal of the transport packing, see chap. 23.1.





#### Note on second-hand chambers (Ex-Demo-Units):

Second-hand chambers are chambers that were used for a short time for tests or exhibitions. They are thoroughly tested before resale. BINDER ensures that the chamber is technically sound and will work flawlessly.

Second-hand chambers are marked with a sticker on the chamber door. Please remove the sticker before commissioning the chamber.

### 3.2 Guidelines for safe lifting and transportation

After operation, please observe the guidelines for temporary decommissioning (chap. 23.2). Empty the Permady™ water pan before moving the chamber. In case of any spilling of the contents, shut down the chamber and dry it out carefully and completely.

  	<div style="background-color: yellow; text-align: center; padding: 5px;"> <b>CAUTION</b></div> <p><b>Sliding or tilting of the chamber.</b>  <b>Damage to the chamber.</b>  <b>Risk of injury by lifting heavy loads.</b></p> <ul style="list-style-type: none"> <li>➤ Transport the chamber in its original packaging only.</li> <li>➤ For moving or shipping, secure the chamber with transport straps.</li> <li>⊘ Do NOT lift or transport the chamber using the door handle, the door or the lower housing.</li> <li>➤ Lift the chamber at the four lower corners with the aid of 4 people and place it on a rolling pallet.</li> <li>➤ Move the chamber to the desired location and lift it from the rolling pallet with the aid of four people.</li> </ul>
---	---


- Permissible ambient temperature range during transport: 10 °C / 14 °F to +60 °C / 140 °F.

You can order transport packing for moving or shipping purposes from BINDER service.

#### Permissible ambient temperature range during transport:

- If the humidifying system has NOT been emptied: +3 °C / 37.4 °F to +60 °C / 140 °F.
- After BINDER Service has emptied the humidifying system: -10 °C / 14 °F to +60 °C / 140 °F.

With temperatures below +3 °C / 37.4 °F, water must be completely removed from the humidifying system.

	<div style="background-color: yellow; text-align: center; padding: 5px;"><b>CAUTION</b></div> <p><b>Transport below +3 °C / 37.4 °F with filled steam humidifying system.</b>  <b>Freezing in the steam generator.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>➤ Contact BINDER Service before any transportation below +3 °C / 37.4 °F.</li> </ul>
---	---


### 3.3 Storage

Intermediate storage of the chamber is possible in a closed and dry room. Observe the guidelines for temporary decommissioning (chap. 23.2).


#### Permissible ambient temperature range during storage:

- the humidifying system has NOT been emptied: +3 °C / 37.4 °F to +60 °C / 140 °F.
- After BINDER Service has emptied the humidifying system: -10 °C / 14 °F to +60 °C / 140 °F

With temperatures below +3 °C / 37.4 °F, water must be completely removed from the humidifying system.

	<div style="background-color: yellow; text-align: center; padding: 5px;"><b>CAUTION</b></div> <p><b>Storage below +3 °C / 37.4 °F with filled steam humidifying system.</b>  <b>Freezing in the steam generator.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>➤ Contact BINDER Service before any transportation below +3 °C / 37.4 °F.</li> </ul>
---	---

**Permissible ambient humidity:** max. 70 % r.H., non-condensing.

	CAUTION
	<p><b>Condensation by excess humidity.</b>  <b>Danger of corrosion on the housing after operating at humidity values &gt; 70 % r.H. for a long period.</b></p> <ul style="list-style-type: none"> <li>➤ Let the chamber dry for several days before shut-down: <ul style="list-style-type: none"> <li>• Empty the Permadry water pan.</li> <li>• <i>Chamber with active humidification:</i> Turn off humidity control (chap. 6.8).</li> <li>• Set the temperature set point to 60 °C / 140 °F for approx. 2 hours.</li> <li>• Only then, shut down the chamber at the main power switch (1) and empty the freshwater bag.</li> </ul> </li> </ul>

When after storage in a cold location you transfer the chamber to its warmer installation site, condensation may form. Before start-up, wait at least one hour until the chamber has attained ambient temperature and is completely dry.

*Chamber with active humidification:* In case of a prolonged temporal decommissioning: Leave the chamber door open or remove the optional access port plugs.

### 3.4 Location of installation and ambient conditions

#### Notes on the location of installation

Set up the chamber on a flat, even surface, free from vibration and in a well-ventilated, dry location. The chambers are designed for setting up inside a building (indoor use).

Freestanding chamber are suitable for installation on tables or on the optionally available stand (height 200 mm / 0.5 ft). Note: The site of installation must be capable of supporting the chamber's weight (see technical data, chap.25.4).


Align the chamber using a spirit level to ensure even covering of the cell-cultures with the medium. For this purpose, manually adjust the four chamber feet.

The chambers can be stacked on top of each other (two chambers maximum). For safe stacking that is easy to maintain, use the original BINDER stacking stand (chap. 18.8.1) or the stacking adapter (chap. 18.8.2).



To completely separate the chamber from the power supply, you must disconnect the power plug. Install the chamber in a way that the power plug is easily accessible and can be easily pulled in case of danger.

For the user there is no risk of temporary overvoltages in the sense of EN 61010-1:2010.

In order to avoid contamination, never place the chamber directly on the floor.


	CAUTION
	<p><b>Danger of overheating.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure sufficient ventilation for dispersal of the heat.</li> </ul>


Do not install or operate the chamber in potentially explosive areas.

	 <b>DANGER</b>
	<p><b>Explosion hazard.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT operate the chamber in potentially explosive areas.</li> <li>➤ KEEP explosive dust or air-solvent mixtures AWAY from the vicinity of the chamber.</li> </ul>


## Ambient conditions

- Permissible ambient temperature range for operation: +18 °C / 64.4 °F to +30 °C / 86 °F  
At elevated ambient temperature values, fluctuations in temperature and humidity may occur.
- Ideal ambient temperature: at least 7 °C / 12.6 °F below the intended working temperature. E.g., working temperature 37 °C / 98.6 °F = ambient temperature 30 °C / 86 °F and lower. In the event of working temperatures of less than 7 °C / 12.6 °F above the ambient temperature, the setpoint can be exceeded.

	The ambient temperature should not be substantially higher than the indicated ambient temperature of 22 +/-3 °C / 71.6 +/-5.4 °F to which the specified technical data relates. For other ambient conditions, deviations from the indicated data are possible.
---	--



	Avoid direct solar radiation on the chamber.
---	--

- Permissible ambient humidity: 70 % r.H. max., non-condensing.
- Installation height: max. 2000 m / 6561.7 ft above sea level.
- Wall distances: rear 100 mm / 3.94 in, sides 50 mm / 1.97 in.

	To completely separate the chamber from the power supply, you must disconnect the power plug. Install the chamber in a way that the power plug is easily accessible and can be easily pulled in case of danger.
---	---

## Notes on handling carbon dioxide (CO<sub>2</sub>)

Carbon dioxide (CO<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Vent out any CO<sub>2</sub> gas that may escape via good room ventilation or a suitable connection to an exhaust system. We recommend installing a CO<sub>2</sub> warning system.

	 <b>WARNING</b>
	<p><b>High concentration of CO<sub>2</sub> (&gt; 4 Vol.-%).</b></p> <p><b>Danger of death by suffocation.</b></p> <p><b>Danger of poisoning.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling CO<sub>2</sub>.</li> </ul>

Observe the **occupational exposure limit OEL** for CO<sub>2</sub> set by the national authorities (formerly maximum permitted workplace concentration). Check compliance when operating all chambers located in the room.

- OEL for Germany: 5000 ml/m<sup>3</sup> (ppm) = 0,5 Vol.-%
- CO<sub>2</sub> lost from CB 170 at 5 vol.-% CO<sub>2</sub> with each opening the door: about 14.8 g, i.e. 0.00758 cubic meters / 0.268 cubic feet (under normal pressure)
- CO<sub>2</sub> lost during 12h at 5 vol.-% without opening the door: approx. < 2 g, i.e. 0.001 cubic meter / 0.035 cubic feet (under normal pressure 1013 mbar / 14.7 psi)

### An example of how to evaluate laboratory volume and air change rate:

**Question:** Is an air change rate of 1/h sufficient for a lab with a volume of 100 cubic meters / 3,531.5 cubic feet with 10 incubators CB 170, opened 4 times per hour?

**Calculation:** CO<sub>2</sub> concentration = CO<sub>2</sub> lost by opening the door, multiplied by 10 incubators, multiplied by opening the door 4 times per hour, divided by lab volume

0.00758 cubic meters x 10 x 4 div. 100 cubic meters = 0.003032, i.e. 0.303 % or 3032 ppm.




0.2676852 cubic feet x 10 x 4 div. 3,531.467 cubic feet = 0.003032, i.e. 0.303 % or 3032 ppm.

**Result:** The maximum permissible value of 5000 ppm is not exceeded under these operation conditions.


Even when CO<sub>2</sub> or systems operated with CO<sub>2</sub> are handled carefully and appropriately, a residual risk remains, which can lead to life-threatening situations under certain circumstances. Therefore we strongly recommend continuous monitoring of CO<sub>2</sub> concentration in the ambient air of the CO<sub>2</sub> incubator. It must be ensured permanently that the maximum permissible occupational exposure limit OEL for CO<sub>2</sub> (0.5 vol - % CO<sub>2</sub> for Germany) is not exceeded.

### Chamber with O<sub>2</sub> control: Notes on handling oxygen (O<sub>2</sub>)

Oxygen (O<sub>2</sub>) is colorless and almost odorless and therefore practically imperceptible. It promotes burns, which can proceed explosively. There is a fire hazard for flammable oxygenated materials, e.g. clothes and hair. O<sub>2</sub> is heavier than air and may accumulate in low-lying areas.



 	<div style="background-color: orange; text-align: center; padding: 5px;">  <b>WARNING</b> </div> <p><b>High concentration of O<sub>2</sub> (&gt; 21 % O<sub>2</sub>).</b>  <b>Fire and explosion hazard through contact of combustible materials with O<sub>2</sub>.</b>  <b>Risk of burns and other injuries.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling O<sub>2</sub>.</li> </ul>
---	--

Take appropriate measures to prevent oxygen enrichment and fire and explosion hazards in areas where oxygen enrichment is possible.

	<p>General information for safe handling of oxygen:</p> <ul style="list-style-type: none"> <li>• Make sure training of personnel on hazards of oxygen enrichment and necessary safety measures.</li> <li>• Make sure adequate labeling of all oxygen equipment and facilities.</li> <li>• Make sure gas tightness of all gas connections by checking them for leaks (e.g. with leak spray or diluted soap solution).</li> <li>• Close the main valve of the source of oxygen after work when not using the chamber.</li> <li>• Never lubricate O<sub>2</sub> equipment with oil or fat. Use only materials and spare parts which are approved for use with oxygen.</li> <li>• Regularly inspect fire extinguishers for proper condition.</li> <li>• Set up emergency showers where oxygen enrichment is possible.</li> <li>• Strictest smoking ban and no ignition sources in areas where oxygen enrichment is possible.</li> <li>• Make sure good ventilation of areas where oxygen enrichment is possible (location of the chamber and/or O<sub>2</sub> cylinders).</li> <li>• Persons who may have been in a possibly oxygen-enriched atmosphere must keep away from ignition sources (flames, cigarettes, etc.) and ventilate their clothes at least 15 minutes.</li> <li>• Always keep emergency routes free.</li> </ul>
---	---

## Chamber with O<sub>2</sub> control: Notes on handling nitrogen (N<sub>2</sub>)

Nitrogen (N<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Any N<sub>2</sub> gas that may escape must be safely led out via good room ventilation or a suitable connection to an exhaust system.

	 <b>WARNING</b>
	<p><b>High concentration of N<sub>2</sub>.</b>  <b>Risk of death by suffocation.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling N<sub>2</sub>.</li> </ul>

## 4. Installation and connections

### 4.1 Shelves

You can put the shelves in different positions at the line of channel beads in the inner chamber. Hold the shelf straight and then insert it so it will go smoothly inside the chamber.

#### Permitted shelf loads:

Maximum load of one single shelf: 10 kg / 22 lb  
 Maximum total load of all shelves: 30 kg / 66 lb

### 4.2 Permadry™ water pan

The Permadry™ system developed by BINDER is an effective and easy to handle system that ensures high humidity inside the chamber without any condensation forming on the inner surfaces. The Permadry™ water pan consists of two pans in which the outer one is heated and the inner one cooled. With the slight difference of temperature caused by that cooling, the central pan is the specific point for condensation of the surplus humidity. Therefore, all other inner surfaces remain dry.



Figure 8: Permadry™ water pan

- Put the Permadry™ water pan on the bottom of the inner chamber in a way that both notches lock into place.
- The front side of the CB 170 / CB 220 Permadry™ water pan is marked "FRONT".



Figure 9: Letters "FRONT" indicating the front of the CB 170 / CB 220 Permadry™ water pan



- Please make sure that the Permadyr™ water pan has firm contact to the inner chamber bottom and rests tightly on it (notches locked into place).
- Fill only the outer pan with distilled, sterilized water up to the filling level marking on the edge of the inner pan.

Maximum filling quantity of the outer pan: CB 56: approx. 0.7 liters, CB 170: approx. 1.3 liters, CB 260: approx. 2.0 liters.



Figure 10: Filling height line of the outer basin CB 170 / CB 220

- We recommend cleaning and refilling the pans 2 to 3 times a week. For evacuation, remove the Permadyr™ water pan.
- We recommend using distilled, sterile water to achieve optimum growth results. Any corrosive damage that may arise following the use of water of different quality or by additives is excluded from the liability agreement.
- If required, you can add microbiologically inhibiting substances such as copper chips, copper sulfate or ethylene diamine tetra-vinegar acid (EDTA) in a concentration of 1 to 5 mmol/l.

Empty the Permadyr™ water pan before moving the chamber. In case of the contents spilling, immediately shut down the chamber and dry it carefully and completely.

### 4.3 Connecting the O<sub>2</sub> sensor (chamber with O<sub>2</sub> control)

The O<sub>2</sub> sensor is supplied with the chamber in a separate package.



Connect or remove the O<sub>2</sub> sensor only when the chamber is turned off.

Open the door of the inner chamber and plug the O<sub>2</sub> sensor (F) into the left connection socket located in the upper part of the rear of the inner chamber. Pay attention to the correct positioning of the pins.



Figure 11: O<sub>2</sub> sensor

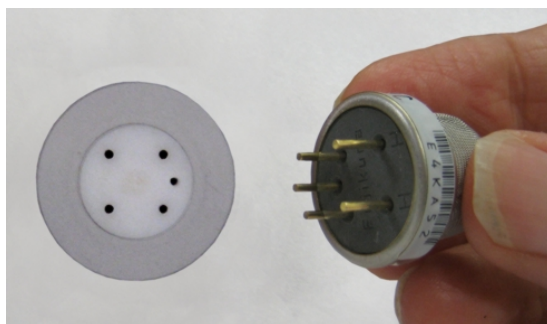



Figure 12: Connecting the O<sub>2</sub> sensor





The O<sub>2</sub> sensor must be plugged in during a hot-air sterilization.





## 4.4 Gas connections


	<p>General information for safe handling of gas cylinders:</p> <ul style="list-style-type: none"> <li>• Store and use gas cylinders only in well ventilated areas.</li> <li>• Open the gas cylinder valve slowly to avoid pressure surges.</li> <li>• Secure gas cylinders during storage and use against falling (chaining).</li> <li>• Transport gas cylinders with a cylinder cart, do not carry, roll, or throw them.</li> <li>• Always close the valve even with apparently empty cylinders; screw on the cap when not in use. Return gas cylinders with the valve closed.</li> <li>• Do not open gas cylinders by force. Mark them when damaged.</li> <li>• Protect gas cylinders against fire, e.g. do not store together with flammable liquids.</li> <li>• Observe relevant regulations for dealing with gas cylinders.</li> </ul>
---	---

Secure the gas cylinder against falling and other mechanical damage.

	<div data-bbox="842 757 1125 831" style="background-color: orange; padding: 5px; text-align: center;">  <b>WARNING</b> </div> <p><b>Safety valve tearing off.</b>  <b>Sudden release of the stored pressure energy.</b>  <b>Risk of injury.</b></p> <ul style="list-style-type: none"> <li>➤ Secure gas cylinders against falling (chaining).</li> <li>➤ Transport gas cylinders with a cylinder cart.</li> </ul>
---	--



The valve of the gas cylinder **always** must be closed before screwing on or unscrewing the gas hose.


	<div data-bbox="842 1108 1125 1182" style="background-color: orange; padding: 5px; text-align: center;">  <b>WARNING</b> </div> <p><b>Opening the cylinder valve when the cylinder is not connected.</b>  <b>Sudden release of the stored pressure energy.</b>  <b>Risk of injury.</b></p> <ul style="list-style-type: none"> <li>➤ Close the gas cylinder valve before connecting or removing the gas hose.</li> </ul>
---	--


	<p>After connecting the gas cylinder, check all gas connections for leaks (e.g. with leak spray or diluted soap solution).</p>
---	--

#### 4.4.1 Connection of the CO<sub>2</sub> gas cylinder

Carbon dioxide (CO<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Vent out any CO<sub>2</sub> gas that may escape via good room ventilation or a suitable connection to an exhaust system. We recommend installing a CO<sub>2</sub> warning system.


	 <b>WARNING</b>
<p><b>High concentration of CO<sub>2</sub> (&gt; 4 Vol.-%).</b>  <b>Risk of death by suffocation.</b>  <b>Danger of poisoning.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling CO<sub>2</sub>.</li> </ul>	

	<p>The CO<sub>2</sub> gas necessary for operation must have a technical grade of 99.5 %.</p>
---	--

	<p>The gas connections must be established by qualified personnel who are trained in handling the respective gases and familiar with the required safety measures.</p>
---	--


The following steps are required:

##### Ensuring the correct CO<sub>2</sub> output pressure

	<p>A gas supply pressure above 2.5 bar / 36 <i>psi</i> will result in chamber damage.</p>
---	---

Use a pressure reducer and make sure to avoid any excessive outlet pressure when connecting the gas hose to the chamber.

The real outlet pressure of gas cylinders, sets of gas cylinders or central gas supplies am on the second manometer must **not** exceed 2.5 bar / 36 *psi*.

	<b>CAUTION</b>
<p><b>Excessive outlet pressure &gt; 2.5 bar / 36 <i>psi</i>.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>Ø The outlet pressure must NOT exceed the indicated value of 2.5 bar / 36 <i>psi</i>.</li> <li>➤ Before connecting, check the outlet pressure on the pressure reducer of the cylinder.</li> <li>➤ Adjust the outlet pressure to 2.0 bar / 29 <i>psi</i> above the ambient pressure.</li> </ul>	

Observe the correct outlet pressure also when replacing the gas cylinders.


##### Establishing the connection to the chamber

Connect the supplied gas hose (internal diameter 6 mm / 0.24 *inches*) to the pressure reducer of the gas cylinders or central gas supply and secure the connection with the supplied hose clamp.

Connect the pre-assembled hose nozzle of the gas hose to the quick acting closure socket (10) DN 6 on the chamber rear, as described in chap. 4.4.4.


##### Leak test

After connecting the gas cylinder, check all gas connections for leaks (e.g. with leak spray or diluted soap solution).




	The recovery times of the gas concentrations inside the chamber after opening the door are indicated in the technical data (chap. 25.4) and refer to a connection pressure of 2.0 bar / 29 <i>psi</i> . Decreasing supply pressure will result in longer recovery times.
---	--


Conversion table for gas inlet pressures, bar – psi, see chap. 25.8.


#### 4.4.2 Connection of the O<sub>2</sub> gas cylinder (chamber with O<sub>2</sub> control with optional alternative control range 10 up to 95 vol.-% O<sub>2</sub>)

	Note: Do not connect the O <sub>2</sub> cylinder or disconnect the gas supply (by pulling off the gas hose) when operating at setpoints below 19 vol.-% O <sub>2</sub> .
---	--

Oxygen (O<sub>2</sub>) is colorless and almost odorless and therefore practically imperceptible. It promotes burns, which can proceed explosively. There is a fire hazard for flammable oxygenated materials, e.g. clothes and hair. O<sub>2</sub> is heavier than air and may accumulate in low-lying areas.


 	 <b>WARNUNG</b>
	<p><b>High concentration of O<sub>2</sub> (&gt; 21 % O<sub>2</sub>).</b>  <b>Fire and explosion hazard through contact of combustible materials with O<sub>2</sub>.</b>  <b>Risk of burns and other injuries.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling O<sub>2</sub>.</li> </ul>

	The O <sub>2</sub> gas necessary for operation must have a technical grade of 99.5 %.
---	---

	The gas connections must be established by qualified personnel who are trained in handling the respective gases and familiar with the required safety measures.
---	---


The following steps are required:

##### Ensuring the correct O<sub>2</sub> output pressure

	A gas supply pressure above 2.5 bar / 36 <i>psi</i> will result in chamber damage.
---	--

Use a pressure reducer and make sure to avoid any excessive outlet pressure when connecting the gas hose to the chamber.

The real outlet pressure of gas cylinders, sets of gas cylinders or central gas supplies am on the second manometer must **not** exceed 2.5 bar / 36 *psi*.

	<b>CAUTION</b>
	<p><b>Excessive outlet pressure &gt; 2.5 bar / 36 <i>psi</i>.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>Ø The outlet pressure must NOT exceed the indicated value of 2.5 bar / 36 <i>psi</i>.</li> <li>➤ Before connecting, check the outlet pressure on the pressure reducer of the cylinder.</li> <li>➤ Adjust the outlet pressure to 2.0 bar / 29 <i>psi</i> above the ambient pressure.</li> </ul>

Observe the correct outlet pressure also when replacing the gas cylinders.


## Establishing the connection to the chamber

Connect the supplied gas hose (internal diameter 6 mm / 0.24 inches) to the pressure reducer of the gas cylinders or central gas supply and secure the connection with the supplied hose clamp.

Connect the pre-assembled hose nozzle of the gas hose to the quick acting closure socket (11) DN 6 on the chamber rear, as described in chap. 4.4.4.

### Leak test



After connecting the gas cylinder, check all gas connections for leaks (e.g. with leak spray or diluted soap solution).


	The recovery times of the gas concentrations inside the chamber after opening the door are indicated in the technical data (chap. 25.4) and refer to a connection pressure of 2.0 bar / 29 psi. Decreasing supply pressure will result in longer recovery times.
---	--


Conversion table for gas inlet pressures, bar – psi, see chap. 25.8.

### 4.4.3 Connection of the N<sub>2</sub> gas cylinder (chamber with O<sub>2</sub> control)

Nitrogen (N<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Any N<sub>2</sub> gas that may escape must be safely led out via good room ventilation or a suitable connection to an exhaust system.


	 <b>WARNING</b>
	<p><b>High concentration of N<sub>2</sub>.</b>  <b>Risk of death by suffocation.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling N<sub>2</sub>.</li> </ul>

	The N <sub>2</sub> gas necessary for operation must have a technical grade of 99.5 %.
---	---

	The gas connections must be established by qualified personnel who are trained in handling the respective gases and familiar with the required safety measures.
---	---


The following steps are required:

### Ensuring the correct N<sub>2</sub> output pressure

	A gas supply pressure above 2.5 bar / 36 psi will result in chamber damage.
---	---

Use a pressure reducer and make sure to avoid any excessive outlet pressure when connecting the gas hose to the chamber.

The real outlet pressure of gas cylinders, sets of gas cylinders or central gas supplies am on the second manometer must **not** exceed 2.5 bar / 36 psi.

	<b>CAUTION</b>
	<p><b>Excessive outlet pressure &gt; 2.5 bar / 36 psi.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>Ø The outlet pressure must NOT exceed the indicated value of 2.5 bar / 36 psi.</li> <li>➤ Before connecting, check the outlet pressure on the pressure reducer of the cylinder.</li> <li>➤ Adjust the outlet pressure to 2.0 bar / 29 psi above the ambient pressure.</li> </ul>

Observe the correct outlet pressure also when replacing the gas cylinders.


#### Establishing the connection to the chamber

Connect the supplied gas hose (internal diameter 6 mm / 0.24 inches) to the pressure reducer of the gas cylinders or central gas supply and secure the connection with the supplied hose clamp.

Connect the pre-assembled hose nozzle of the gas hose to the quick acting closure socket (9) DN 6 on the chamber rear, as described in chap. 4.4.4.

#### Leak test

After connecting the gas cylinder, check all gas connections for leaks (e.g. with leak spray or diluted soap solution).



	<p>The recovery times of the gas concentrations inside the chamber after opening the door are indicated in the technical data (chap. 25.4) and refer to a connection pressure of 2.0 bar / 29 psi. Decreasing supply pressure will result in longer recovery times.</p>
---	---

Conversion table for gas inlet pressures, bar – psi, see chap. 25.8.


#### 4.4.4 Connecting the gas hose to the chamber rear (for CO<sub>2</sub>, O<sub>2</sub>, N<sub>2</sub>)

The procedure of connecting the gas hose to the chamber rear is the same for any gas connection. All quick acting closure sockets (CO<sub>2</sub>, N<sub>2</sub> with O<sub>2</sub> control, and O<sub>2</sub> with O<sub>2</sub> control and optional alternative control range 10 up to 95 vol.-% O<sub>2</sub>) are degreased and supplied with a FKM gasket.

**Note for chambers with O<sub>2</sub> control:** The quick acting closure socket and the hose nozzle for the O<sub>2</sub> connection must be degreased.

	 <b>WARNING</b>
	<p><b>Fire and explosion hazard through contact of fat with O<sub>2</sub>.</b>  <b>Risk of burns and other injuries.</b></p> <ul style="list-style-type: none"> <li>Ø All connection parts for the O<sub>2</sub> connection must be degreased.</li> </ul>

The gas hose, which will be used to establish the connection to a gas cylinder, is already attached to the hose nozzle and secured by a hose clamp. Plug the hose nozzle into the corresponding quick acting closure socket (a) located at the rear of the chamber. This quick acting closure socket is closed by a rubber cover (b).

	<p>Only use the supplied hose nozzle to connect to the quick acting closure socket. Otherwise, the quick acting closure socket may leak, and/or it may become impossible to connect the original hose nozzle. In this case, please contact BINDER Service.</p>
---	--

Remove the rubber cover (b) by pulling it off.


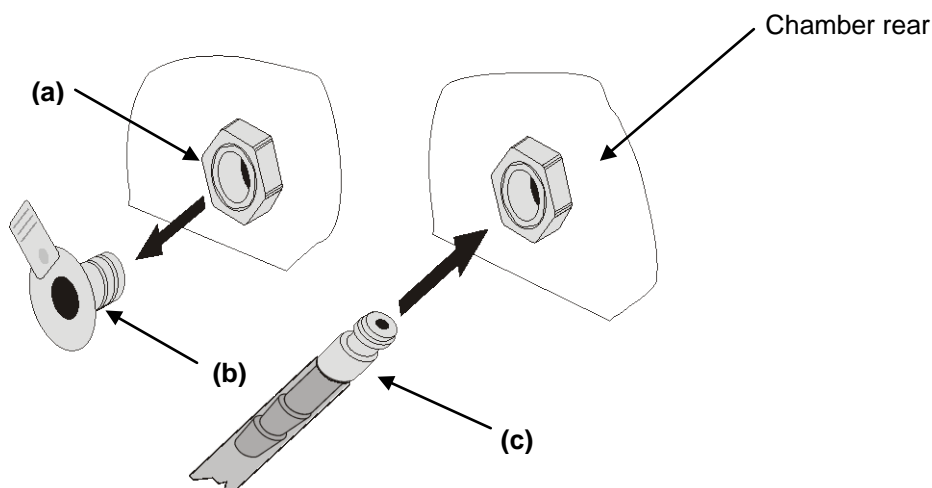


Figure 13: Connecting the hose lead to the gas cylinder

Now fit the hose nozzle (c) in the quick acting closure socket. To remove the connection, pull the hose nozzle off the quick acting closure socket.

#### 4.4.5 Gas cylinder connection kits (option)

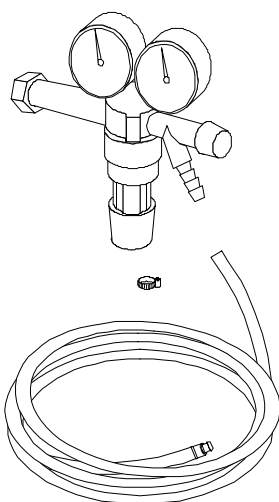

Gas cylinder connection kits are available for CO<sub>2</sub> (Art. No. 8012-0014), O<sub>2</sub> (Art. No. 8012-0015) and N<sub>2</sub> (Art. No. 8012-0016).

The connection kit includes the following parts for connecting a gas cylinder to the CO<sub>2</sub> incubator:

- Pressure reducer with manometers for cylinder pressure (high pressure gauge) and outlet pressure (low pressure gauge)
- 5 m pressure hose with pre-assembled hose nozzle for quick acting closure socket
- 1 hose clamp to connect the gas hose to the pressure reducer

Figure 14: Gas cylinder connection kit



Instructions 7001-0249 included with the connection kit describe connecting and setting the gas cylinder connection kit.



### CAUTION

**Excessive outlet pressure > 2.5 bar / 36 psi.**

**Damage to the chamber.**

- ⊘ The outlet pressure must NOT exceed the indicated value of 2.5 bar / 36 psi.
- Before connecting, check the outlet pressure on the pressure reducer of the cylinder.
- Adjust the outlet pressure to 2.0 bar / 29 psi above the ambient pressure.




The gas connections must be established by qualified personnel who are trained in handling the respective gases and familiar with the required safety measures.

## 4.5 Water supply for the chamber with active humidification

Freshwater is supplied by manually filling the freshwater bag. It is not possible to connect the chamber directly to a water pipe.

The freshwater bag has a volume of 0.9 liters / 0.32 cu.ft. and is placed behind the chamber door in a recess.

	To guarantee humidification during 24 hours even at high humidity set-points, we recommend filling the freshwater bag daily at the end of the day.
---	--


When the freshwater can is empty, the message "Freshwater supply" will be displayed on the controller, the buzzer sounds, and the humidification module turns off. After acknowledging the alarm, the humidification module tries to fill up and start operating.


Water consumption at 37 °C / 98.6 °F and 93 % r.h.:


- Without opening the door: very low.
- When opening the door 10 times a day: approx. 70 ml per day.

### 4.5.1 Types of suitable water quality

- Deionized water from a water treatment installation already existing at the customer's site. Conductivity from 1 µS /cm up to a maximum of 20 µS/cm. (Water, which is in equilibrium with the CO<sub>2</sub> in the air, and has a conductivity below 1 µS/cm (ultrapure water), may cause acid corrosion due to its low pH).
- Tap water that has been treated by the optional water treatment system BINDER Pure Aqua Service (disposable system). A reusable measuring equipment to assess the water quality is included (chap. 4.5.3).

	<p>BINDER GmbH is NOT responsible for the water quality at the user's site.</p> <p>Any problems and malfunctions that might arise following use of water of deviating quality is excluded from liability by BINDER GmbH.</p> <p>The warranty becomes void in the event of use of water of deviating quality.</p>
---	--

	<p style="text-align: center;"><b>CAUTION</b></p> <p><b>Calcification of the humidifying system.</b>  <b>Damage to the chamber.</b></p> <p>➤ Operate the chamber with deionized (demineralized) water only.</p>
---	---

	Water intake temperature NOT below +5 °C / 41 °F and not exceeding 40 °C / 104 °F.
---	--



#### 4.5.2 Water supply installation

Please proceed in the following order:

1. Fill the freshwater bag with sterilized water.  
Observe water quality and temperature (chap. 4.5.1).  
As long as no tube is connected to the outlet of the freshwater bag, the bag outlet must be closed
2. Place the freshwater bag with the opening to the left in the recess behind the chamber door.
3. Establish the connection of the tube to the freshwater bag
4. Lead the tube along the pump
5. For operation with humidity, the freshwater bag outlet must be open

#### 4.5.3 BINDER Pure Aqua Service (option)

The optional BINDER water treatment system (disposable system) is available to treat tap water. The lifetime depends on water quality and the amount of treated water used. The measuring equipment to assess the water quality is reusable.




For detailed information on operating the water treatment system BINDER Pure Aqua Service and its function, please refer to the operating manual supplied with BINDER Pure Aqua Service.

## 4.6 Electrical connection


The chambers are supplied ready for connection. They come with an IEC connector plug.

Model	Art. No. (x = 0 or 1)	Power plug	Voltage +/-10 %	Power frequency	Chamber fuse
CB 170	9x40- 0131 9x40- 0132 9x40- 0133 9x40- 0134	Grounded plug IEC 7/7	200-230 V (1N~)	50/60 Hz	10 A
CB 170-UL	9x40- 0139 9x40- 0140 9x40- 0141 9x40- 0142	NEMA 5-20P	100-120 V (1N~)	50/60 Hz	16 A

- The domestic socket must also provide a protective conductor. Make sure that the connection of the protective conductor of the domestic installations to the chamber's protective conductor meets the latest technology. The protective conductors of the socket and plug must be compatible!
- Prior to connection and start-up, check the power supply voltage. Compare the values to the specified data located on the chamber's type plate (left chamber side, bottom right-hand, see chap. 1.4). We recommend the use of a residual current circuit breaker.
- When connecting, please observe the regulations specified by the local electricity supply company as well as the VDE directives (for Germany).
- Pollution degree (acc. to IEC 61010-1): 2
- Over-voltage category (acc. to IEC 61010-1): II

	<b>CAUTION</b>
	<p><b>Danger of incorrect power supply voltage.</b>  <b>Damage to the equipment.</b></p> <ul style="list-style-type: none"> <li>➤ Check the power supply voltage before connection and start-up.</li> <li>➤ Compare the power supply voltage with the data indicated on the type plate.</li> </ul>

See also electrical data (chap. 25.4).

	<p>To completely separate the chamber from the power supply, you must disconnect the power plug. Install the chamber in a way that the power plug is easily accessible and can be easily pulled in case of danger.</p>
---	--

#### 4.7 Handling and aligning the divided inner door, gas proof (optional equipment)

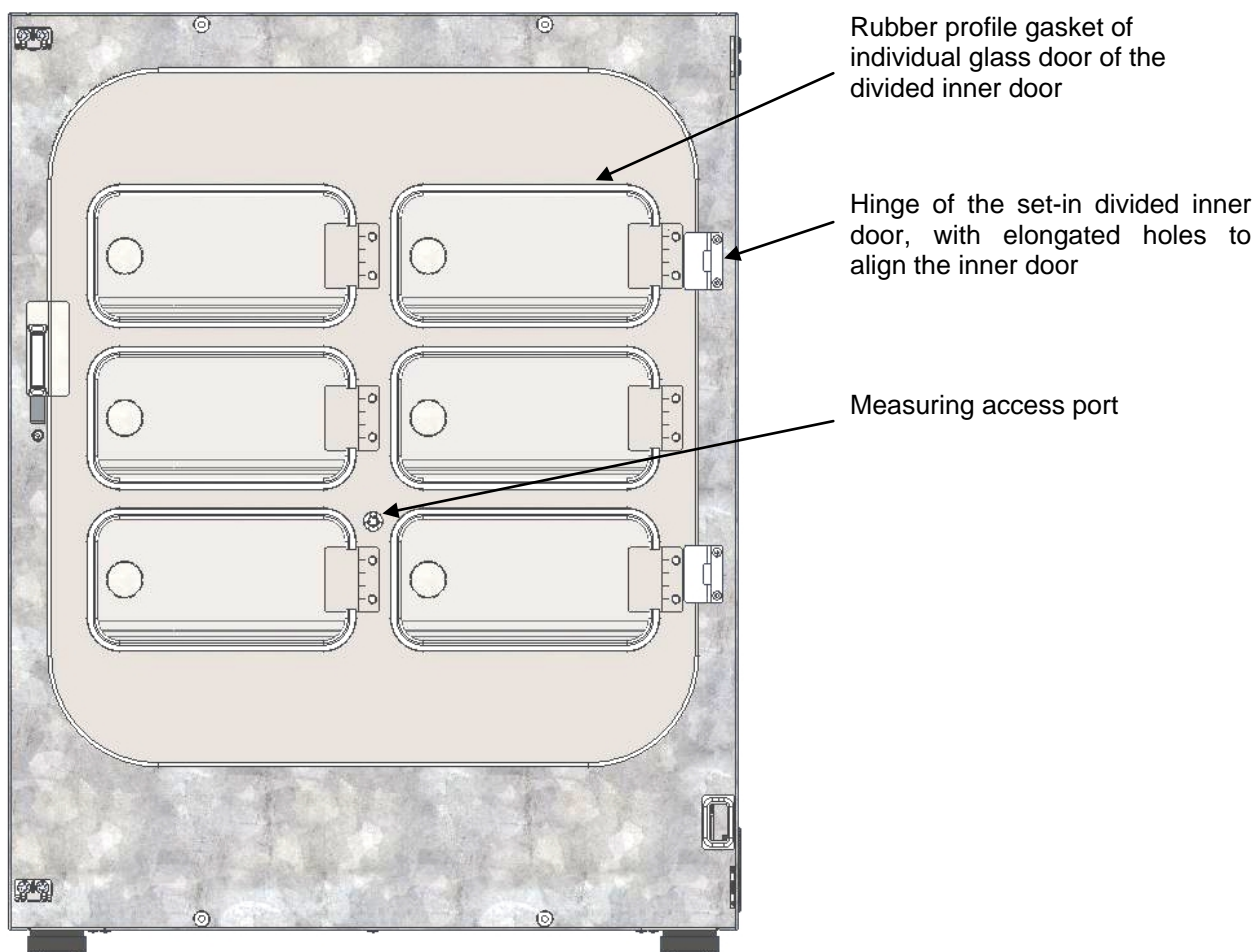


Figure 15: Divided inner door, gas proof (example: CB 170)

- Pull the handles to open each of the individual glass doors. Do not press too hard while closing them to avoid pushing the doors through the gasket into the inner chamber.
- The rubber profile gaskets of the individual glass doors are easy to replace.
- If the divided inner door is not in an aligned position, fasten the hinge screws in the elongated holes of the hinge to align it. Align the door and operate the locking smoothly.

## 5. Functional overview of the MB2 chamber controller

The MB2 chamber controller controls following parameters inside the chamber:

- Temperature in °C
- Carbon dioxide concentration in vol.-%
- Oxygen concentration in vol.-% (chamber with O<sub>2</sub> control)
- Humidity in % r.h. (chamber with active humidification)



Note: In this manual the chambers are described in the maximum equipment variant and including the available options. Therefore, some of the pictured and described features are not available on every chambers

You can enter the desired set point values in fixed value operation mode directly on the display surface or via the setpoint menu. For program operation the controller offers programming week and time programs. In addition there is a timer program available (stopwatch function).

The controller offers various notifications and alarm messages with visual and audible indication and remote alarms via e-mail, an event list (trace file) and the graphical display of the measuring values in the in der chart recorder view. The MB2 program controller permits programming temperature and CO<sub>2</sub> / O<sub>2</sub> and humidity cycles, and specifying special controller functions for each program section. You can enter values or programs directly at the controller or use the APT-COM™ 3 DataControlSystem software (option) specially developed by BINDER.

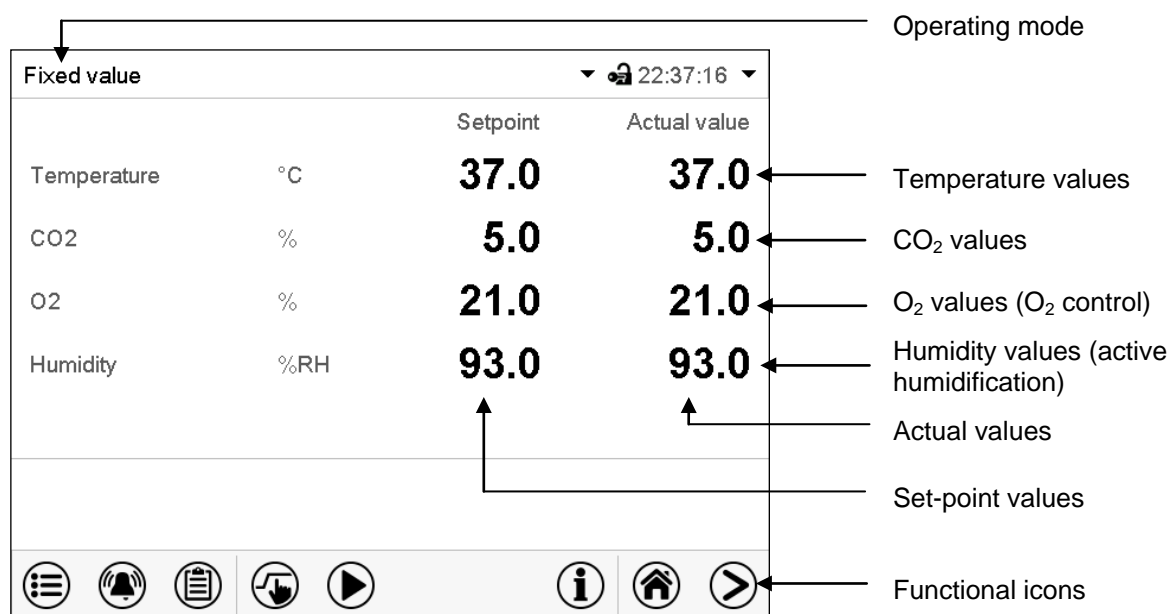


Figure 16: Normal display of the MB2 program controller (sample values, chamber with O<sub>2</sub> control and active humidification)

## 5.1 Operating functions in normal display

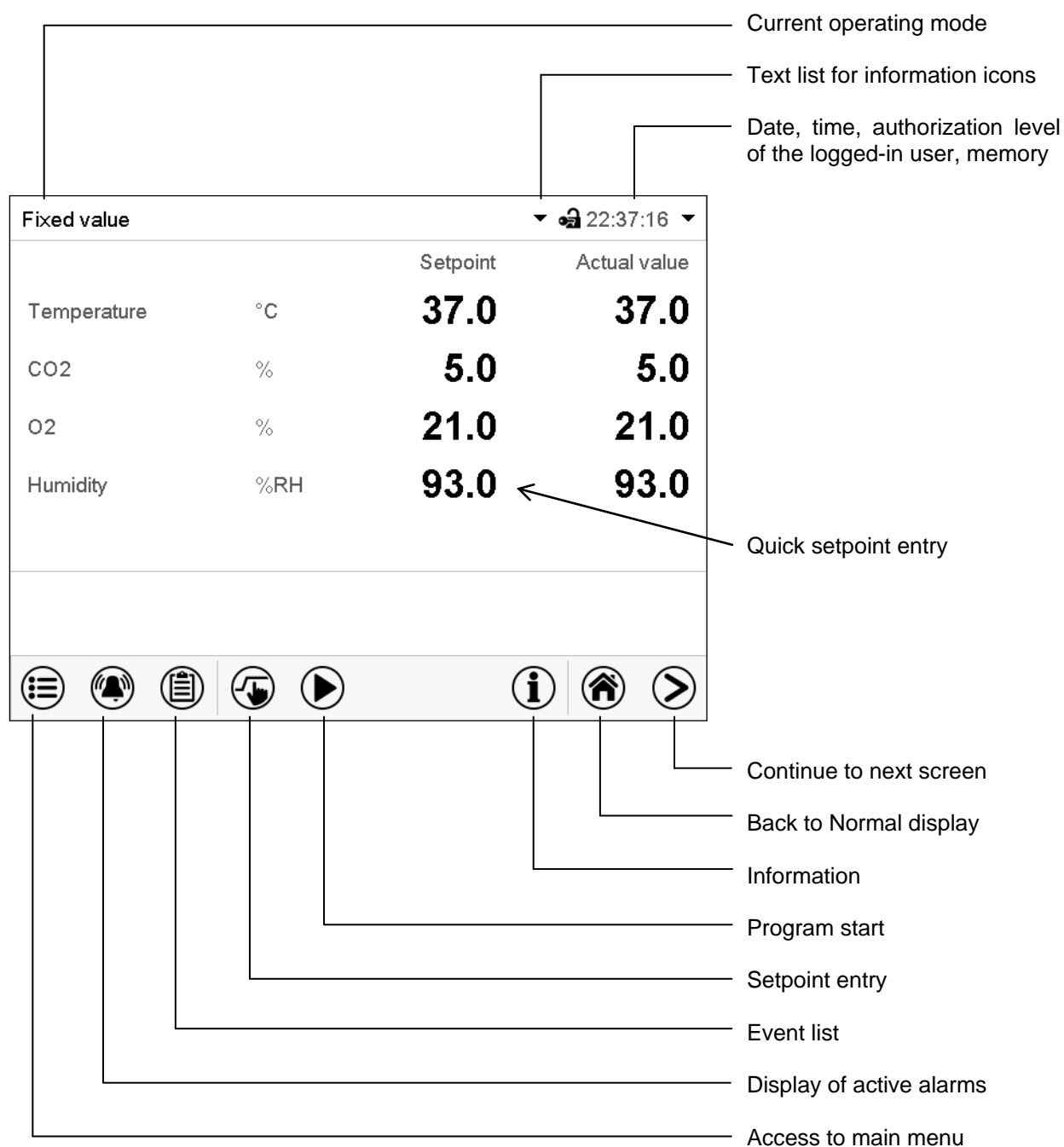


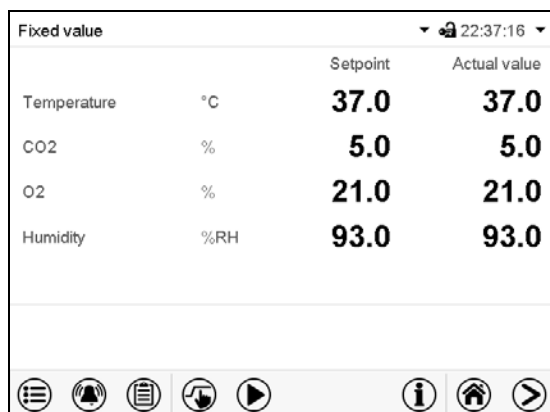



Figure 17: Operating functions of the MB2 controller in normal display (sample values, chamber with O<sub>2</sub> control)

## 5.2 Display views: Normal display, program display, chart-recorder display

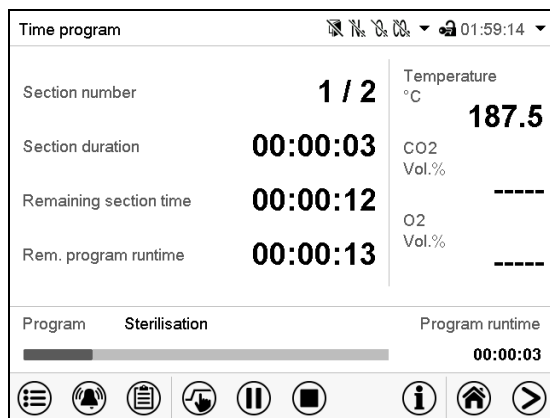	Press the <b>Change view</b> icon to toggle between normal display, program display and chart-recorder display.
	Press the <b>Normal display</b> icon to return from program display and chart recorder display back to Normal display.



Normal display (actual values / setpoint values)



Program display (example: time program)












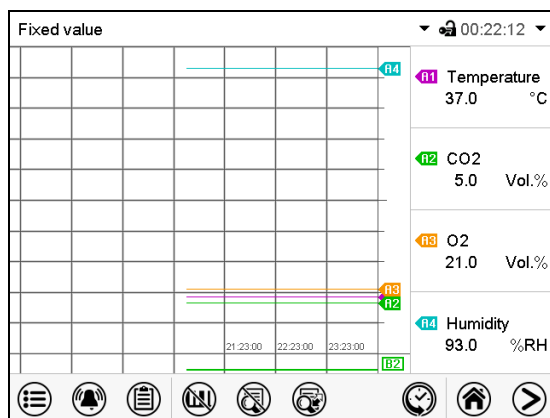

Chart recorder display








(Sample values, chamber with O<sub>2</sub> control and active humidification)

### 5.3 Controller icons overview

#### Navigation icons in Normal display









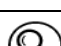


Icon	Signification	Function
	<b>Main menu</b>	Access from Normal display to the main menu
	<b>Alarm</b>	Access from Normal display to the list of active alarms
	<b>Event list</b>	Access from Normal display to the event list
	<b>Setpoint setting</b>	Access from Normal display to the setpoint entry menu: setpoint entry for Fixed value operation, turning on/off CO <sub>2</sub> / O <sub>2</sub> and humidity control, safety controller settings
	<b>Program start</b>	Start a previously entered time or week program, continue a paused time program
	<b>Program pause</b>	Pause a running time program
	<b>Program cancelling</b>	Cancel a running time or week program
	<b>Information</b>	Information on program operation, setpoints, actual values, and the safety controller
	<b>Normal display</b>	Return from program display or chart recorder display to Normal display
	<b>Change view</b>	Toggle between Normal display, program display, and chart recorder display

#### Functional icons in individual menus







Icon	Signification	Function
	<b>Back</b>	Return from each menu to Normal display
	<b>Update</b>	Update the event list and alarm messages
	<b>Confirm</b>	Take over the entries and exit the menu / continue menu sequence.
	<b>Close</b>	Exit the menu / cancel menu sequence. Entries are not taken over. When terminating a menu sequence, an information window appears, which must be confirmed.
	<b>Reset alarm</b>	Acknowledge the alarm and mute the buzzer.
	<b>Change keyboard</b>	Change between uppercase and lower case characters, digits and special characters
	<b>Edit</b>	Edit settings of time and week programs




### Functional icons in the chart recorder display

Icon	Signification	Function
	<b>Show legend</b>	Show legend
	<b>Hide legend</b>	Hide legend
	<b>Switch legend</b>	Switch between legend pages
	<b>Show indications</b>	Show indication "Door open" (B2)
	<b>Hide indications</b>	Hide indication "Door open" (B2)
	<b>History display</b>	Pause chart recorder and change to history display. Data recording continues.
	<b>Curve selection</b>	Go to "Curve selection" submenu in the history display
	<b>Search</b>	Go to "Search" submenu in the history display to select the required instant
	<b>Zoom</b>	Go to "Zoom" submenu in the history display to select the zoom factor
	<b>Show scroll buttons</b>	Show scroll buttons in the history display to scroll to an instant
	<b>Hide scroll buttons</b>	Hide scroll buttons in the history display to scroll to an instant
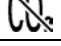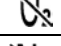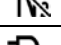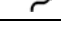
### Information icons referring to chamber conditions

Icon	Text information	Condition
	"Idle mode"	Controller is in Idle mode
	"Door open"	Chamber door is open
	"CO2 pressure alarm off"	CO <sub>2</sub> pressure alarm deactivated
	"O2 pressure alarm off"	O <sub>2</sub> pressure alarm deactivated (chamber with O <sub>2</sub> control)
	"N2 pressure alarm off"	N <sub>2</sub> pressure alarm deactivated (chamber with O <sub>2</sub> control)
	"Interior socket"	Interior socket (option) activated

### Information icon for data processing

Icon	Information
	Waiting icon: Data processing is running. Remaining time to touch the display when calibrating the touchscreen.

## 5.4 Operating modes

The MB2 program controller operates in the following operating modes:

- **Idle mode**

The controller is not functional, i.e., there is no heating, no CO<sub>2</sub> inlet or O<sub>2</sub> inlet (O<sub>2</sub> control) and no humidification (active humidification). The chamber approximates ambient values.

You can activate and deactivate this operating mode with the "Idle mode" control contact in Fixed value operating mode (chap. 7.3), time program operation (chap. 9.7.3) and week program operation (chap. 10.6.5).

- **Fixed value operating mode**

The controller operates as a fixed-point controller, i.e., set-points for temperature, CO<sub>2</sub>, O<sub>2</sub> (O<sub>2</sub> control) and humidity (active humidification) can be defined, which are then maintained until the next manual change (chap. 7.1).

- **Timer program operation**

Stopwatch function: during an entered duration the controller constantly equilibrates to the setpoints entered in Fixed value operation mode.


- **Time program operation**

An entered time program for temperature, CO<sub>2</sub>, O<sub>2</sub> (O<sub>2</sub> control) and humidity (active humidification) is running. The controller offers 25 program memory places with 100 program sections each. The total number of program sections of all programs is unlimited

- **Week program operation**






An entered week program for temperature, CO<sub>2</sub>, O<sub>2</sub> (O<sub>2</sub> control) and humidity (active humidification) is running. The controller offers 5 program memory places with 100 switching points each. The switching points can be distributed over all days of the week.

## 5.5 Controller menu structure

Fixed value		Setpoint	Actual value
Temperature	°C	37.0	37.0
CO <sub>2</sub>	%	5.0	5.0
O <sub>2</sub>	%	21.0	21.0
Humidity	%RH	93.0	93.0
			

Use the **navigation icons** in the screen footer in Normal display to access the desired controller functions.



The available functions depend on the current **authorization level** “Service”, “Admin” or “User” (chap. 13.1). This is selected either during login or can be available without password protection.

	<b>Main menu:</b> program settings, further information, “Service” submenu. The “Settings” submenu allows general configuration of the controller.	chap. 5.5.1
	List of <b>active alarms</b>	chap. 11
	Access to the <b>event list</b>	chap. 15.3
	<b>Setpoint entry</b> for Fixed value operation, turning on/off <b>CO<sub>2</sub> control</b> , <b>O<sub>2</sub> control (chamber with O<sub>2</sub> control)</b> and <b>humidity control (active humidification)</b> , <b>safety controller</b> settings	chap. 7, 6.8, 12.2
	Start/ pause/ cancel an already entered, respectively a running <b>time program</b> or start / cancel an already entered, respectively a running <b>week program</b>	chap. 9.1, 9.2, 10.1









Unless noted otherwise, the figures show the functional range, which is available for the user with “Admin” authorization level.

### 5.5.1 Main menu

The main menu provides access to the general configuration of the controller as well as to program entry and the user administration. Additionally there are support functions like a contact page or the display calibration depending on the available angle.

	Press the <b>Main menu</b> icon to access the main menu from Normal Display.
	Press the <b>Back</b> icon to return from each setting menu to Normal Display.

The main menu provides the following functions and submenus.

Main menu		
 User	User management: login and logout, password management	chap. 13
 Device info	Chamber information	chap. 15.2
 Settings	"Settings" submenu (not visible for user with "User" authorization level)	chap. 14
 Programs	Program entry submenu for time and week programs	chap. 9 and 10
 Service	"Service" submenu	chap. 5.5.3
 Contact	BINDER Service contact page	chap. 15.1
 Calibrate touchscreen	Calibrating the touch screen	chap. 14.4.2
	Back to Normal Display	

#### "Settings" submenu

- Settings of many general controller functions and network settings (chap. 14).
- Available only for users with "Service" and "Admin" authorization level

#### "Service" submenu

- Access to service data, controller reset to factory settings (chap. 5.5.3)
- Available only for users with "Service" and "Admin" authorization level. Full functional range only for BINDER Service (users with "Service" authorization level).











#### "Programs" submenu

- Access to the controller's program functions (chap. 8, 9, 10)

### 5.5.2 “Settings” submenu

The “Settings” submenu is available for users with “Service” or “Admin” authorization level. It serves to enter date and time, select the language for the controller menus and the desired temperature unit and to configure the controller’s communication functions.








Path: [Main menu](#) > [Settings](#)

<div> <div>Main</div> <div>Settings</div> <div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> </div> <div> <div>Chamber</div> <div>Date and time</div> <div>Display</div> <div>Measurement chart</div> <div>Various</div> <div>Serial interfaces</div> <div>Ethernet</div> <div>Web server</div> <div>email</div> <div>Back to main menu</div> </div> </div>		
	Setting the temperature unit, menu language...	chap. 14.1, 14.2
	Setting date and time	chap. 14.2
	Setting the display brightness, continuous operation and screen saver	chap. 14.4
	Settings for the measurement chart: storage interval, storage values, minimum and maximum values	chap. 16.2
	Setting the tolerance range and delay time for tolerance range alarm	chap. 11.4
	Configuration of the RS485 interface (available via BINDER INDIVIDUAL Customized Solutions)	chap. 14.5.1
	Entry of the MAC address and IP address	chap. 14.5.2
	Password protection for web server access	chap. 14.5.3
	Configuration of the e-mail server, assignment of e-mail addresses	chap. 14.5.4
	Back to main menu	

### 5.5.3 “Service” submenu

The “Service” submenu is available for users with “Service” or “Admin” authorization level. When logged-in with “Admin” authorization level the user will find information to tell the BINDER Service in service case.

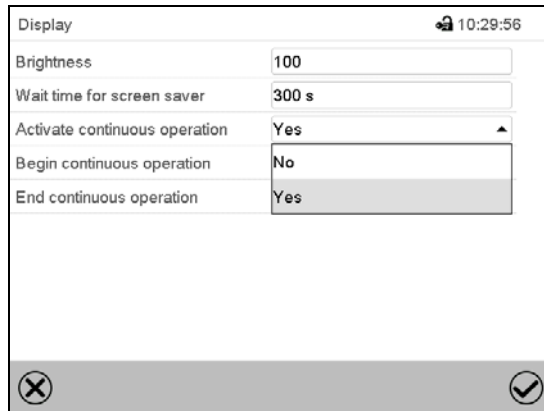
Path: [Main menu](#) > [Service](#)

<div> <div>Main</div> <div>Service</div> <div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> <div>  </div> </div> <div> <div>Service data</div> <div>Counter</div> <div>ST code</div> <div>Factory settings</div> <div>Back to main menu</div> </div> </div>		
	Serial number of the chamber, setup version of the controller software	chap. 14.2
	No function	
	Information for BINDER Service	
	Reset to factory settings	
	Back to main menu	

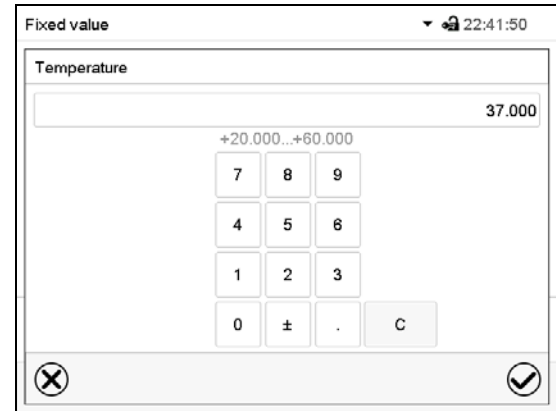
(view with “Admin” authorization level)

## 5.6 Principle of controller entries

In the selection and entry menus there are icons displayed in the footers which you can use to take over the entry or cancel it.





Selection menu (example)



Entry menu (example)

After completing the settings there are the following possibilities:

	Press the <b>Confirm</b> icon to take over the entries and exit the menu or continue the menu sequence.
	Press the <b>Close</b> icon to exit the menu or cancel the menu sequence without taking over the entries. When terminating a menu sequence, an information window appears, which must be confirmed.

## 5.7 Performance during and after power failures

During a power failure, all controller functions are shut down. The gas inlet valves are closed so that no gas can escape into the ambient air. The zero-voltage relay alarm output (3) (chap. 11.6) is switched to alarm position for the whole duration of the power failure.

After the power returns, all functions return to the same status the chamber had before power failure. . The controller continues to function in the original operating mode it was in previously before the power failure occurred.

- Performance after power failure in Idle mode  
Control is deactivated
- Performance after power failure in Fixed value operation mode  
All functions return to the same status the chamber had before power failure. The set-points are immediately resumed.
- Performance after power failure during time program operation  
The program is resumed at the point where the interruption occurred with the latest set-points reached during the program run.
- Performance after power failure during week program operation  
The week program continues with the values corresponding to the current time.
- Performance after power failure during sterilization operation  
The sterilization process is cancelled and the controller changes to Fixed value operation mode with the setpoints already entered. All setpoint values remain in memory.

Power failure and power return are noted in the event list (chap. 15.3).

If during power failure the inner chamber temperature or the CO<sub>2</sub> / O<sub>2</sub> concentration has fallen below the alarm thresholds, confirm the alarms (e.g., tolerance range, safety controller). See chap. 11.3.

## 5.8 Performance when opening the door

When you open the door, CO<sub>2</sub> control and O<sub>2</sub> control (chamber with O<sub>2</sub> control) immediately turns off.

After 60 seconds from opening the door, the heating and humidification (chamber with active humidification) turn off.

After closing the door, heating, CO<sub>2</sub> control, O<sub>2</sub> control (chamber with O<sub>2</sub> control), humidification (chamber with active humidification) turn on again.

## 6. Start up

### 6.1 Turning on the chamber

After connecting the supply lines (chap. 4), turn on the chamber by its main power switch. The lit pilot lamp shows the chamber is ready for operation.

When the main power switch is turned on and yet the controller display is dark, the display is in stand-by mode. Press on the touchscreen to activate it.

Check that the CO<sub>2</sub> control is activated (chap. 6.8).

*Chamber with active humidification:*

Fill the freshwater bag (chap. 4.5). Check that the humidifying system is activated (chap. 6.8).

After the first turning on of the chamber or after an interruption of the power supply the relative humidity will increase after a delay of about 20 minutes. During this period, the relative humidity can drop considerably.

*Chamber with O<sub>2</sub> control:*

Check that the O<sub>2</sub> control is activated (chap. 6.8).

Warming chambers may release odors in the first few days after commissioning. This is not a quality defect. To reduce odors quickly we recommend heating up the chamber to its nominal temperature for one day and in a well-ventilated location.



**WARNING:** If customer should use a BINDER chamber running in non-supervised continuous operation, we strongly recommend in case of inclusion of irrecoverable specimen or samples to split such specimen or samples and store them in at least two chambers, if this is feasible.



## 6.2 Controller settings upon start up

The window „Language selection“ enables the **language selection**, in case that it's activated in the „Start-up“ menu. Afterwards occurs a request of the **time zone** and the **temperature unit**.

Language selection

German

English

✓

Start-up

Temperature unit: Degrees Celsius

Time zone: UTC+1h (CET)

Daylight saving time switch: Automatic

▼ Start of daylight saving time

▼ End of daylight saving time

Language query after restart: Yes


✗ ✓

The controller will function in the **operating mode**, which was active before the last shut-down. It controls temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification) in fixed value operating mode to the last entered values and in the program mode to the set points achieved beforehand.

### Locked operation

Provided that the user administration has been activated by the assignment of passwords for the different authorization types, the **controller operation** is first locked after turning on the unit, recognizable by the closed lock icon in the header.

Fixed value		Setpoint	Actual value
Temperature	°C	37.0	37.0
CO <sub>2</sub>	%	5.0	5.0
O <sub>2</sub>	%	21.0	21.0
Humidity	%RH	93.0	93.0



☰ 🔔 📄 🔄 ▶

ℹ 🏠 ➡

In the locked view the controller provides all display functions. No setting functions are available.

The setpoints are shaded (light grey) in normal display. Changing them by direct entry in the fixed value operating mode is not possible. The functional icons for setpoint entry and program start in the footer are without function.

After turning on the unit, user log-in is required to operate the controller (chap. 13.2)

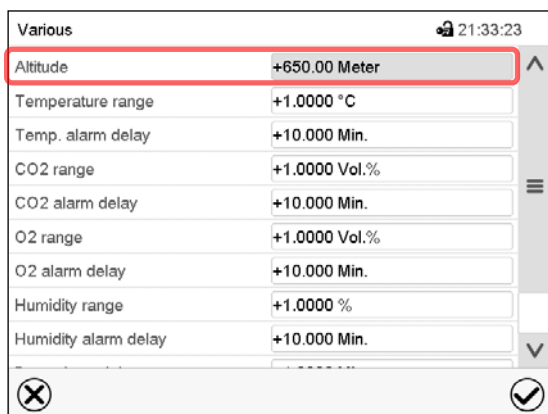
### Operation without user log-in / without password-protection

If the password function has been deactivated, after turning on the unit without user log-in there are those controller functions available, which correspond to the highest authorization level without a password protection. There is no lock icon in the header.


### 6.3 Entering the altitude of the installation site

After first turning on the chamber, it is required to enter the altitude above sea level of the installation site into the controller. This entry serves to correct the calculation of CO<sub>2</sub> concentration in vol.-% from the measurement of partial pressure. The setting will remain stored after shutting the power off.

Path: **Main menu > Settings > Various**



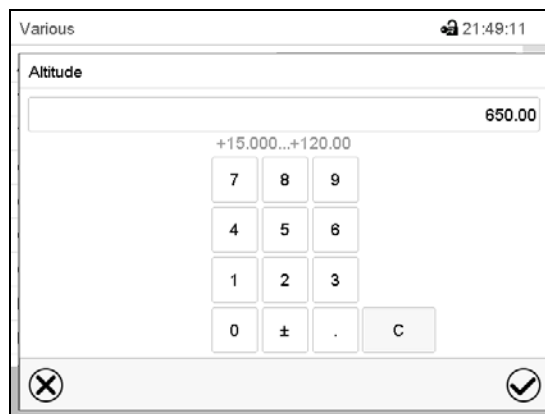Submenu "Various".  
Select the field "Altitude"



"Altitude" entry menu. Enter the desired altitude above sea level in meters and confirm entry with **Confirm** icon.  
Factory setting: 650 m

### 6.4 Factory settings

The chamber is supplied with the following basic preset parameters:

- |   |                     |
|---|---------------------|
| • Temperature setpoint  | 37 °C / 98.6 °F     |
| • CO <sub>2</sub> setpoint                                      | 5 vol.-%            |
| • O <sub>2</sub> setpoint (chamber with O <sub>2</sub> control) | 21 vol.-%           |
| • Humidity setpoint (chamber with active humidification)        | 93 % r.h.           |
| • Safety controller class 3.1                                   | Offset 2 °C         |
| • Sterilization temperature                                     | 187.5 °C / 369.5 °F |
| • Audible alarm signal (buzzer)                                 | activated           |
| • Altitude above sea level                                      | 650 m               |

The set temperature determines the target working temperature in the inner chamber, i.e. set value 37 °C / 98.6 °F = target working temperature 37 °C / 98.6 °F. The same is valid for the CO<sub>2</sub> and O<sub>2</sub> concentration (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification). For the hot-air sterilization, the set value is 187.5 °C / 369.5 °F and cannot be changed.



As long as there is a difference between the actual and set value shown in the display, adequate operation of the chamber is not guaranteed.

## 6.5 Equilibration time

### Temperature

- Equilibration time is max. 4 hours.

### CO<sub>2</sub>

- After turning in the chamber, after approx. 5 minutes the CO<sub>2</sub> concentration equilibrates automatically to the pre-set value of 5 vol.-% CO<sub>2</sub>.
- Equilibration time is max. 0.5 hours.

### O<sub>2</sub> (chamber with O<sub>2</sub> control)

- After turning in the chamber, the ambient oxygen concentration of approx. 20.9 vol.-% is displayed after a delay of 10 minutes. During the first 10 minutes, O<sub>2</sub> control is not defined. Then O<sub>2</sub> is regulated to the displayed setpoint value of 21 vol.-%.
- For high O<sub>2</sub> setpoint values (with alternative control range option), equilibration time is max. 1.5 hours.

### Humidity (without active humidification)

- With humidity control effected via the Permadyr<sup>®</sup> system equilibration time is max. 15 hours

### Humidity (chamber with active humidification)

- Equilibration time is max. 1 hour.

## 6.6 Operating the DuoDoor™ door lock

The DuoDoor™ door lock offers two different ways to open the doors, dependent on the glass door lock position.

### Opening outer door and glass door independently

When closing the glass door, close the glass door lock.

Then you can close and open the outer door using its handle. The glass door will remain closed until you manually open the glass door lock.

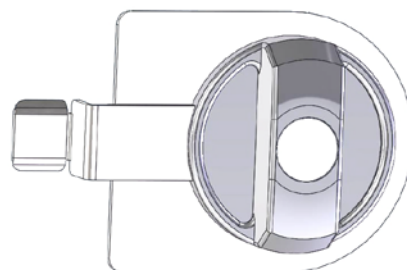


Figure 18: Closed position

### Opening outer door and glass together

When closing the glass door leave the glass door lock in open position. Then close the outer door. The glass door is pressed against the gasket so that both doors are securely closed.

When opening the outer door, a magnetic device also opens the glass door.

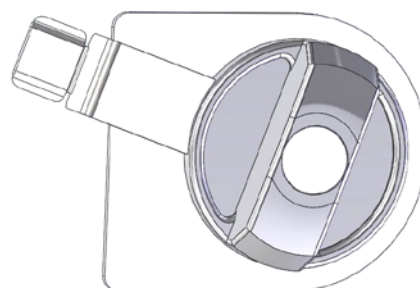


Figure 19: Open position

## 6.7 Required gas supply for the chamber with O<sub>2</sub> control

In the chamber with O<sub>2</sub> control, cell growth can be additionally influenced by a variable oxygen content. Depending on the required range of O<sub>2</sub> concentration, the type of gas which is required may vary. To this end, two equipment variants are available:

### 6.7.1 Hypoxic control range 0.2 vol.-% up to 20 vol.-% O<sub>2</sub> (regular)

In addition to the CO<sub>2</sub> connection, the chamber provides a gas connection for N<sub>2</sub>.

As compared to the ambient air (O<sub>2</sub> concentration of 20.9 vol.-%) only the N<sub>2</sub> concentration needs to be increased, no oxygen is required for the O<sub>2</sub> control. Therefore, it is not necessary and possible to connect an O<sub>2</sub> cylinder. Even with activated O<sub>2</sub> control no O<sub>2</sub> pressure alarm will be emitted.

### 6.7.2 Alternative control range 10 vol.-% up to 95 vol.-% (option)

In addition to the CO<sub>2</sub> connection, the chamber provides gas connections for O<sub>2</sub> and N<sub>2</sub>.

- **Operation in the hyperoxic range (> 22 vol.-% O<sub>2</sub>)**

As compared to the ambient air (O<sub>2</sub> concentration of 20.9 vol.-%) only the O<sub>2</sub> concentration needs to be increased, no nitrogen is required for the O<sub>2</sub> control. Therefore, it is not necessary to connect a N<sub>2</sub> cylinder.

You can deactivate the N<sub>2</sub> pressure alarm in the controller (chap. 7.3). Then also with activated O<sub>2</sub> control no N<sub>2</sub> pressure alarm will be emitted.

- **Alternating operation in the hyperoxic *and* hypoxic range or in the range from 20 vol.-% O<sub>2</sub> up to 22 vol.-% O<sub>2</sub>**

N<sub>2</sub> and O<sub>2</sub> are required for O<sub>2</sub> control. It is necessary to connect all gas supplies.

You can deactivate the O<sub>2</sub> and / or N<sub>2</sub> pressure alarm in the controller (chap. 7.3). Then also with activated O<sub>2</sub> control no O<sub>2</sub> and / or N<sub>2</sub> pressure alarm will be emitted.


### 6.7.3 Operation without O<sub>2</sub> control

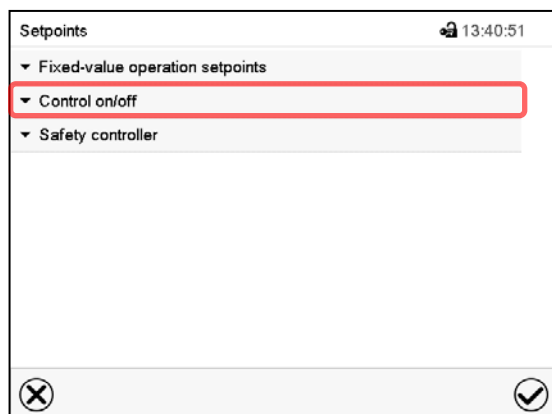
You can deactivate the O<sub>2</sub> control in the controller (chap. 6.8). With this setting, no O<sub>2</sub> and N<sub>2</sub> pressure alarms and no O<sub>2</sub> tolerance range alarm will be issued. The zero-voltage relay alarm contact (chap. 11.6) will also not be triggered.

Since neither nitrogen nor oxygen are required, it is not necessary to connect an O<sub>2</sub> or N<sub>2</sub> cylinder.


## 6.8 Activating / deactivating CO<sub>2</sub>, O<sub>2</sub> and humidity control

If CO<sub>2</sub> control and/or O<sub>2</sub> control (chamber with O<sub>2</sub> control) and/or humidity control (chamber with active humidification) shall not be used, it is possible to deactivate them in the controller.

 Press the **Setpoint setting** icon to access the “Setpoint” setting menu from Normal display.



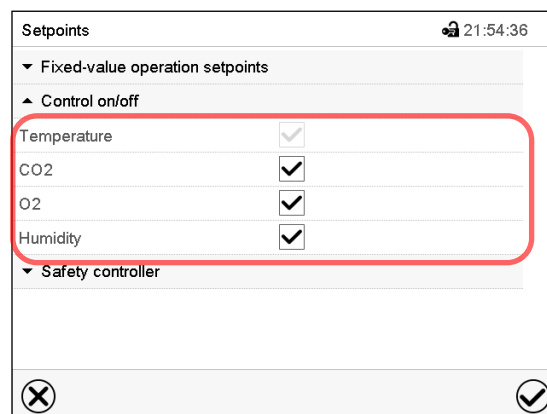“Setpoints” menu.  
Select “Control on/off”.



You can turn on or off CO<sub>2</sub> control and O<sub>2</sub> control (chamber with O<sub>2</sub> control) and humidity control (active humidification)

Mark / unmark the checkbox to activate / deactivate the desired function and press the **Confirm** icon.



When operating the chamber with deactivated CO<sub>2</sub> control, no CO<sub>2</sub> pressure alarm and no CO<sub>2</sub> tolerance range alarm will be emitted.



*Chamber with O<sub>2</sub> control:* When operating the chamber with deactivated O<sub>2</sub> control, no O<sub>2</sub> and N<sub>2</sub> pressure alarms and no O<sub>2</sub> tolerance range alarm will be emitted.



*Chamber with active humidification:* When operating the chamber with deactivated humidity control no humidity tolerance range alarm will be emitted. The controller continues to show the actual humidity value but no longer the humidity set-point.

When operating the chamber with deactivated humidity control, humidification can be realized using the Permadyr™ water pan (chap. 6.9).

## 6.9 Humidity control of the Permadyr™ system

With the chamber with active humidification, humidity setting via the Permadyr™ System is not required.

The Permadyr™ system with its 2-pan water system guarantees a maximum humidity of up to 95 % r.H. in the inner chamber, which remains condensation-free. This performance assumes an average ambient temperature of 22 +/- 3 °C / 71.6 +/- 5.4 °F and a working temperature in the inner chamber of 37 °C / 98.6 °F.

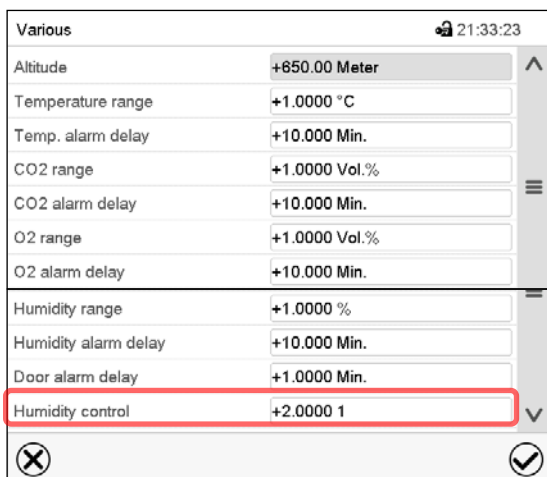
You can increase or decrease humidity slightly via the heating of the water pan if required.

Humidity is infinitely adjustable in whole numbers from -5 (low) to +5 (high).


Factory setting: "0" (medium).

### Setting:

Path: [Main menu](#) > [Settings](#) > [Various](#)



Various		21:33:23
Altitude	+850.00 Meter	^
Temperature range	+1.0000 °C	
Temp. alarm delay	+10.000 Min.	
CO2 range	+1.0000 Vol.%	
CO2 alarm delay	+10.000 Min.	
O2 range	+1.0000 Vol.%	
O2 alarm delay	+10.000 Min.	
Humidity range	+1.0000 %	
Humidity alarm delay	+10.000 Min.	
Door alarm delay	+1.0000 Min.	
Humidity control	+2.0000 1	v

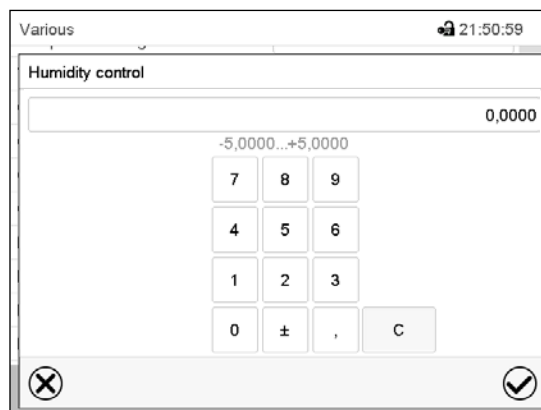Humidity control		21:50:59
-5,0000...+5,0000		0,0000
7	8	9
4	5	6
1	2	3
0	±	,
		C

"Humidity control" entry menu. Enter the desired humidity and confirm entry with **Confirm** icon.

Submenu "Various".

Select the field "Humidity control"




If the ambient temperature deviates by more than +/- 5 °C from the values recommended by the manufacturer, the conditions for maximum air humidity with condensation-free inner chamber are no longer guaranteed. Contact BINDER Service for assistance.


## 7. Set-point entry in “Fixed value” operating mode


In Fixed value operating mode you can enter set-points for temperature, CO<sub>2</sub> and O<sub>2</sub> (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification), and the switching-state of up to 16 operation lines.

All settings made in Fixed value operating mode remain valid until the next manual change. They are saved also when turning off the chamber or in case of toggling to Idle Mode or Program Mode.

	Setting ranges	Control ranges
<b>Temperature</b>	20 °C / 68 °F up to 60 °C / 140 °F	4 °C / 7.2 °F above ambient temperature up to 60 °C / 140 °F
<b>Temperature</b> (chamber with O <sub>2</sub> control)		6 °C / 10.8 °F above ambient temperature up to 60 °C / 140 °F
<b>CO<sub>2</sub></b>	0 vol.-% up to 20 vol.-%	0 vol.-% up to 20 vol.-%
<b>O<sub>2</sub></b> (chamber with O <sub>2</sub> control, with regular hypoxic control range)	0.2 vol.-% up to 21 vol.-%	0.2 vol.-% up to 20 vol.-%
<b>O<sub>2</sub></b> (chamber with O <sub>2</sub> control, with optional alternative control range)	5 vol.-% up to 95 vol.-%	10 vol.-% up to 95 vol.-%
<b>Humidity</b> (chamber with active humidification)	60 % r.h. up to 93 % r.h.	60 % r.h. up to 93 % r.h.

	<p>With set-point type “<b>Limit</b>”, adapt the safety controller (chap. 12.2) always when you changed the temperature set-point. Set the safety controller set-point by approx. 2 °C above the controller temperature set-point.</p> <p>Recommended setting: Set-point type “<b>Offset</b>” with safety controller set-point 2 °C.</p>
---	--

	<p>When setting a lower temperature setpoint, in order to save time, we recommend cooling down the chamber by turning it off and opening both chamber doors.</p>
---	--

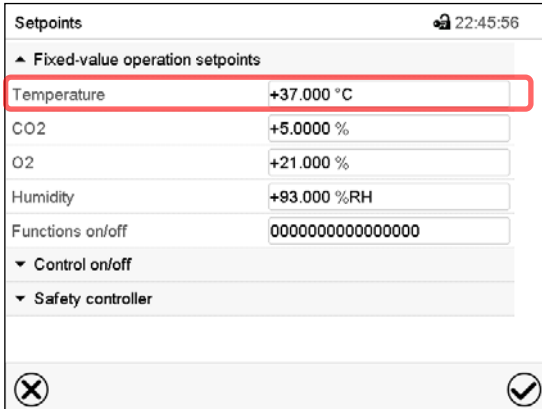
	<p>When setting a lower CO<sub>2</sub>, O<sub>2</sub> or humidity setpoint, the gas or humidity must be able to escape first. Open both chamber doors for this purpose. Observe the precautions when working with gases (chap. 1.6).</p>
---	--



## 7.1 Set-point entry through the “Setpoints” menu



Press the **Setpoint setting** icon to access the “Setpoint” setting menu from Normal display.



“Setpoints” menu.

Select “Fixed value operation setpoints” to access the individual parameters.

- Select the field “Temperature” and enter the desired temperature setpoint.  
Setting range: 20 °C up to 60 °C. Confirm entry with **Confirm** icon.
- Select the field “CO<sub>2</sub>” and enter the desired CO<sub>2</sub> setpoint.  
Setting range: 0 vol.-% up to 20 vol.-%.  
Confirm entry with **Confirm** icon.

*Chamber with O<sub>2</sub> control:*

- Select the field “O<sub>2</sub>” and enter the desired O<sub>2</sub> setpoint.  
Setting range with regular hypoxic control range: 0.2 vol.-% up to 20 vol.-%.  
Setting range with alternative control range: 5 vol.-% up to 95 vol.-%.  
Confirm entry with **Confirm** icon.

*Chamber with active humidification:*

- Select the field “Humidity” and enter the desired humidity setpoint.  
Setting range: 60 % r.h. up to 93 % r.h. Confirm entry with **Confirm** icon.



When entering a value outside the setting range, the message: “Value outside of limits! (Min: xxx, Max: xxx)” appears (xxx is a wildcard for the limits of the respective parameter). Press the **Confirm** icon and repeat the entry with a correct value.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

## 7.2 Direct setpoint entry via Normal display

Alternatively you can also enter the setpoints directly via Normal display.

Fixed value		22:37:16	
		Setpoint	Actual value
Temperature	°C	37.0	37.0
CO2	%	5.0	5.0
O2	%	21.0	21.0
Humidity	%RH	93.0	93.0

Normal display. Select the setpoint you want to change.



Fixed value		22:41:50	
Temperature			
			37.000
+20.000...+60.000			
7	8	9	
4	5	6	
1	2	3	
0	±	.	C

Example: "Temperature" entry menu. Enter the desired setpoint and confirm entry with **Confirm** icon.

### 7.3 Special controller functions via operation lines



Press the **Setpoint setting** icon to access the “Setpoint” setting menu from Normal display.

You can define the switching state of up to 16 operation lines (control contacts). They are used to activate / deactivate special controller functions.

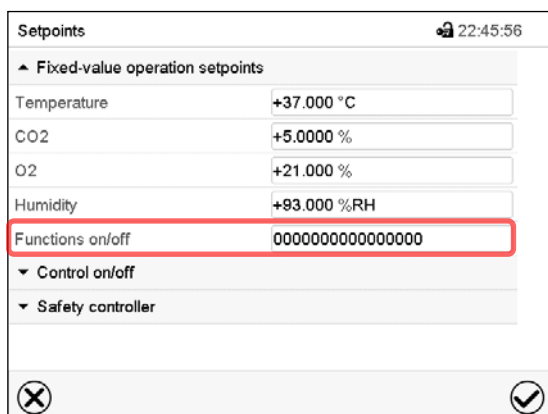
*Chamber with O<sub>2</sub> control:* Depending on the required range of O<sub>2</sub> concentration, the type of gas which is required may vary (chap. 6.6). Therefore, the pressure alarms for O<sub>2</sub> and N<sub>2</sub> can be deactivated individually with active O<sub>2</sub> control.

For chambers equipped with the optional interior socket (chap. 18.3) you can turn on and off the voltage of the interior socket via the controller.

- Operation line “Idle mode” activates / deactivates the operating mode “Idle mode”.
- Operation line “O<sub>2</sub> pressure alarm off” activates / deactivates the O<sub>2</sub> pressure alarm (chamber with O<sub>2</sub> control)
- Operation line “N<sub>2</sub> pressure alarm off” activates / deactivates the N<sub>2</sub> pressure alarm (chamber with O<sub>2</sub> control)
- Operation line “Interior socket” activates / deactivates interior socket (option, available via BINDER INDIVIDUAL Customized Solutions)


The other operation lines are without function.

Use the “Setpoints” menu to configure the operation lines.



“Setpoints” menu.

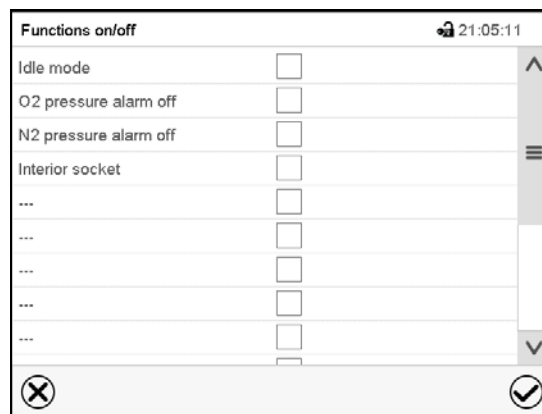Select the field “Functions on/off”.

“Functions on/off” entry menu (view with chamber with O<sub>2</sub> control with optional interior socket).

Mark / unmark the checkbox to activate / deactivate the desired function and press the **Confirm** icon

Activated operation line: switching status “1” (On)

Deactivated operation line: switching status “0” (Off)

The operation lines count from right to left.

#### Example:



Activated operation line “Idle mode” = 000000000000000001

Deactivated operation line “Idle mode” = 000000000000000000

## 7.4 Safety instructions for setting high gas concentrations

### Notes on handling carbon dioxide (CO<sub>2</sub>)




Carbon dioxide (CO<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Vent out any CO<sub>2</sub> gas that may escape via good room ventilation or a suitable connection to an exhaust system. We recommend installing a CO<sub>2</sub> warning system.

	 <b>WARNING</b>
	<p><b>High concentration of CO<sub>2</sub> (&gt; 4 Vol.-%).</b>  <b>Danger of death by suffocation.</b>  <b>Danger of poisoning.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling CO<sub>2</sub>.</li> </ul>

If CO<sub>2</sub> is released, leave the area und inform the security service or fire department.

### Chamber with O<sub>2</sub> control: Notes on handling oxygen (O<sub>2</sub>)

Oxygen (O<sub>2</sub>) is colorless and almost odorless and therefore practically imperceptible. It promotes burns, which can proceed explosively. There is a fire hazard for flammable oxygenated materials, e.g. clothes and hair. O<sub>2</sub> is heavier than air and may accumulate in low-lying areas.



 	 <b>WARNING</b>
	<p><b>High concentration of O<sub>2</sub> (&gt; 21 % O<sub>2</sub>).</b>  <b>Fire and explosion hazard through contact of combustible materials with O<sub>2</sub>.</b>  <b>Risk of burns and other injuries.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses</li> <li>➤ Ensure technical ventilation measures</li> <li>➤ Observe the relevant regulations for handling O<sub>2</sub>.</li> </ul>

Take appropriate measures to prevent oxygen enrichment and fire and explosion hazards in areas where oxygen enrichment is possible.

Observe the general information for safe handling of oxygen (chap. 1.6).

### Chamber with O<sub>2</sub> control: Notes on handling nitrogen (N<sub>2</sub>)


Nitrogen (N<sub>2</sub>) in high concentrations is hazardous to health. It is colorless and almost odorless and therefore practically imperceptible. Any N<sub>2</sub> gas that may escape must be safely led out via good room ventilation or a suitable connection to an exhaust system.

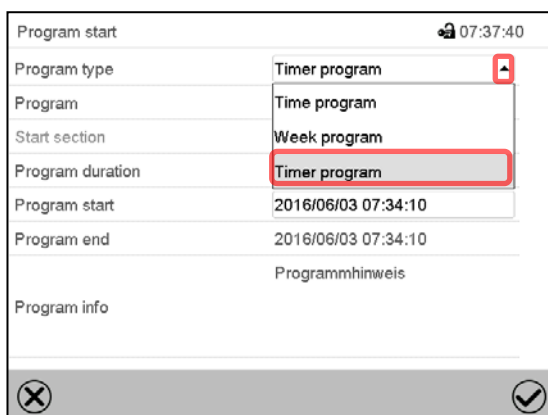
	 <b>WARNING</b>
	<p><b>High concentration of N<sub>2</sub>.</b>  <b>Risk of death by suffocation.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT set up chambers in non-ventilated recesses</li> <li>➤ Ensure technical ventilation measures</li> <li>➤ Observe the relevant regulations for handling N<sub>2</sub>.</li> </ul>

## 8. Timer program: stopwatch function

During an entered duration the controller constantly equilibrates to the setpoints entered in Fixed value operation mode (temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control), humidity (chamber with active humidification), configuration of the operation lines). This duration can be entered as a "Timer program". During the program runtime, any setpoint changes do not become effective; the controller equilibrates to the values which were active during program start.

### 8.1 Starting a timer program

 In Normal display press the **Program start** icon to access the "Program start" menu.



"Program start" menu

- In the field "Program type" select "Timer program".
- Select the field "Program duration" and enter the desired program duration. Press the **Confirm** icon.
- Select the field "Program start" and enter the desired start time of the program in the "Program start" entry menu. Press the **Confirm** icon. The program delay time until program start begins.



Normal display.


Information on the bottom of the screen indicates the currently running program and the time already passed. The grey bar shows how much time of the whole time is elapsed.

#### 8.1.1 Performance during program delay time

During the configured program delay time until program start, the controller equilibrates to the current setpoints of Fixed value operation mode. Modifications of these setpoints are possible but become effective only after the timer program is finished. When the configured moment for program start is reached, the program delay time ends and the program starts running. The controller equilibrates to the values which had been active during program start.



## 8.2 Stopping a running timer program

### 8.2.1 Pausing a running timer program


	Press the <b>Program pause</b> icon to interrupt the program.
---	---

The program is paused. The program runtime stops running down, the time display flashes.

There are the following options:

	Press the <b>Program start</b> icon to continue the program
	Press the <b>Cancelling</b> icon to cancel the program

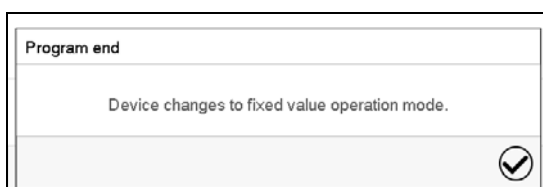
### 8.2.2 Cancelling a running timer program

	Press the <b>Program cancelling</b> icon to cancel the program.
---	---

A confirmation prompt is displayed. Press the **Confirm** icon to confirm that the program shall really be cancelled.

After confirming the message the controller changes to Fixed value operation mode. Temperature and humidity will then equilibrate to the setpoints of Fixed value operation mode.

## 8.3 Performance after the end of the program



After the end of the program the message "Device changes to fixed value operation mode" appears on the screen.


Press the **Confirm** icon.

After confirming the message the controller changes to Fixed value operation mode. Temperature and humidity will then equilibrate to the setpoints of Fixed value operation mode.

## 9. Time programs

The MB2 program controller permits programming time programs with real-time reference. It offers 25 program memory positions with up to 100 program sections each.

For each program section you can enter set-points for temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control), and humidity (chamber with active humidification), section duration, type of set-point transition (ramp or step) and the tolerance ranges.



If the safety controller has been set to "limit" mode, check the setting of the safety controller when changing the temperature set-point, (chap. 12.2).

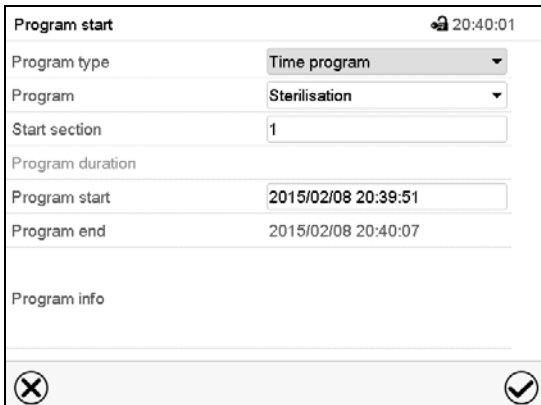
Programming remains saved in case of a power failure or after turning off the unit.

Path: [Main menu](#) > [Programs](#) > [Time program](#)

### 9.1 Starting an existing time program



In Normal display press the **Program start** icon to access the "Program start" menu.



"Program start" menu

- In the field "Program type" select the setting "Time program".
- In the field "Program" select the desired program.
- Select the field "Program start" and enter the desired program start time in the "Program start" entry menu. Press the **Confirm** icon. The program delay time until program start begins.

The program end is adapted automatically depending on the entered program duration.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu. The program starts running.

If instead you press the **Close** icon to exit the menu without taking over the entries, the program will not start.




Normal display. Information on the bottom of the screen indicates the currently running program and the time already passed. The grey bar shows how much time of the whole time is elapsed. If program duration has been set to infinite, the grey bar is not displayed.

### 9.1.1 Performance during program delay time

During the configured program delay time until program start, the controller equilibrates to the current setpoints of Fixed value operation mode. Modifications of these setpoints are effective. When the configured moment for program start is reached, the program delay time ends and the program starts running.



## 9.2 Stopping a running time program

### 9.2.1 Pausing a running time program


	Press the <b>Program pause</b> icon to interrupt the program..
---	--

The program is paused. The program runtime stops running down, the time display flashes.

There are the following options:

	Press the <b>Program start</b> icon to continue the program
	Press the <b>Cancelling</b> icon to cancel the program

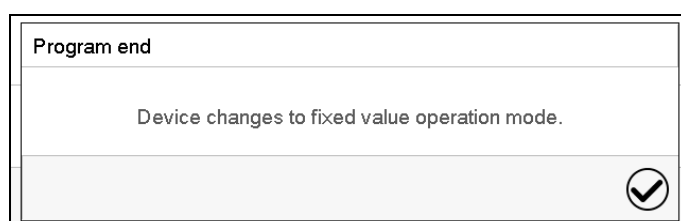
### 9.2.2 Cancelling a running time program

	Press the <b>Program cancelling</b> icon to cancel the program.
---	---

A confirmation prompt is displayed. Press the **Confirm** icon to confirm that the program shall really be cancelled.

After confirming the message, the controller changes to Fixed value operation mode. Temperature and humidity will then equilibrate to the setpoints of Fixed value operation mode.

## 9.3 Performance after the end of the program



After the end of the program the message "Device changes to fixed value operation mode" appears on the screen.

Press the **Confirm** icon.


As long as the message has not been confirmed, the setpoint of the last program section remains effective. Program the last section as desired. If the heating, CO<sub>2</sub> control, O<sub>2</sub> control (chamber with O<sub>2</sub> control) and humidification (chamber with active humidification) shall turn off, activate operation line "Idle mode" in the last program section.

After confirming the message the controller changes to Fixed value operation mode. The setpoints of Fixed value operation mode are then equilibrated.




## 9.4 Creating a new time program

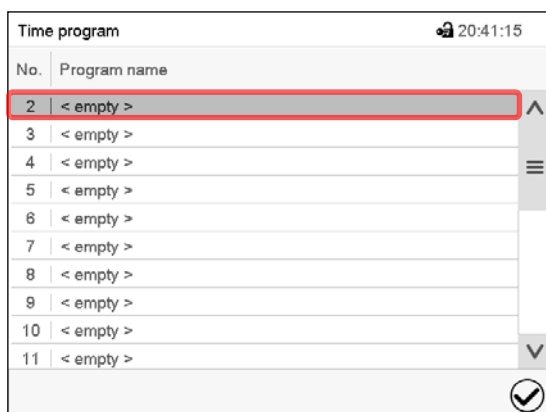Path: [Main menu > Programs > Time program](#)



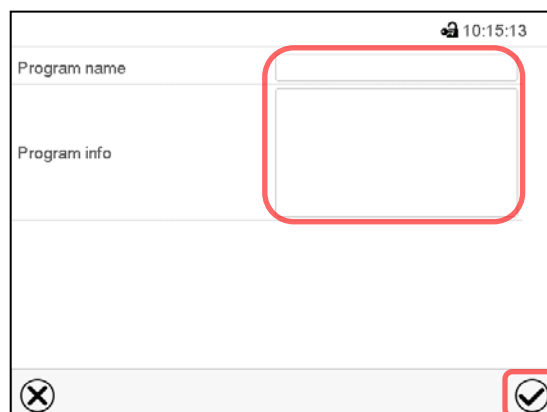
"Time program" menu:  
overview of the existing programs.  
Select an empty program place.




Enter the program name and, if desired,  
additional program information in the  
corresponding fields.  
Press the **Confirm** icon.  
The program view opens (chap. 9.5).

## 9.5 Program editor: program management

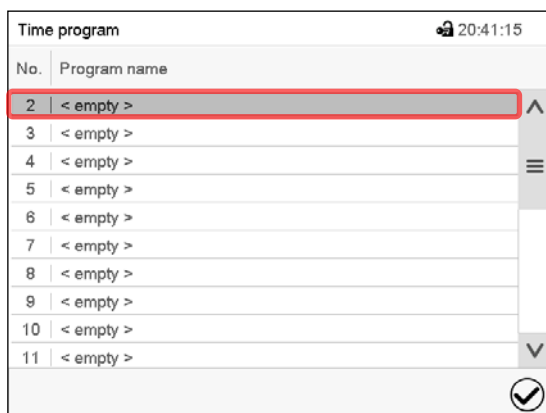Path: [Main menu > Programs > Time program](#)



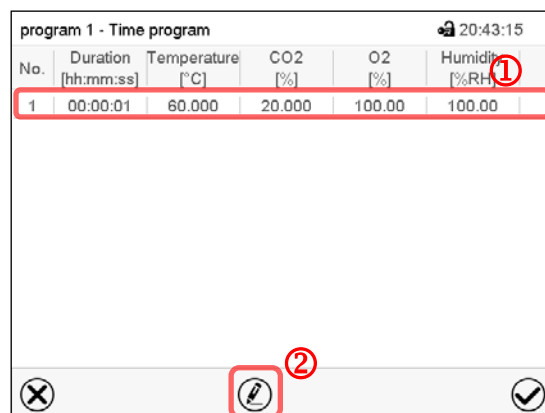
"Time program" menu:  
overview of the existing programs.  
Select an existing program (example: program 3)  
or create a new program (chap. 9.4).  
The program view opens.



Program view (example: program 1).  
If a new program has been created, there is  
just one program section.

There are the following options:

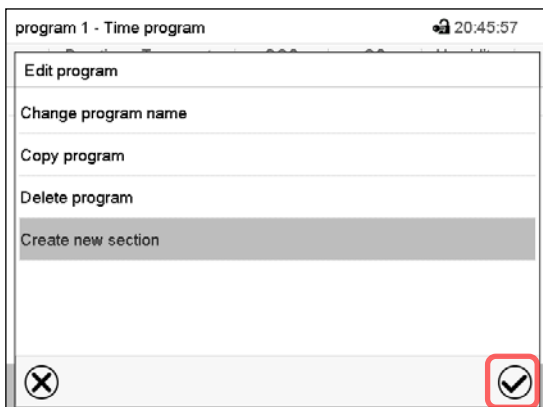- ① Select a program section to open the section editor (chap. 9.6)
- ② Press the **Edit** icon to open the program editor




Program editor: "Edit program" menu

Select the desired function and press the **Confirm** icon.


The program editor offers following options:

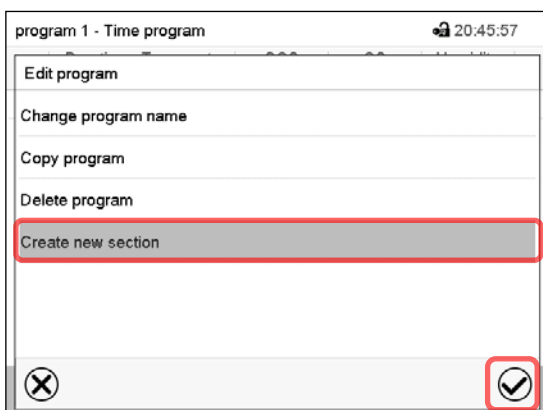- Change the program name
- Copy program
- Replace program: Replacing an new or an existing program with the copied program. This menu point is visible only after a program has been copied.
- Delete program
- Create new section



To add a new section, select "Create new section" and press the **Confirm** icon.

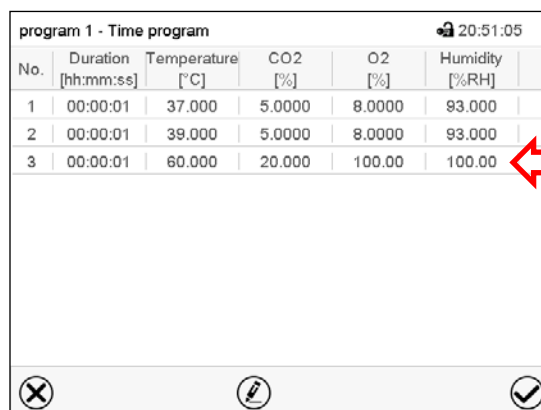The program view opens.

No.	Duration [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]	Humidity [%RH]
1	00:00:01	37.000	5.0000	8.0000	93.000
2	00:00:01	39.000	5.0000	8.0000	93.000
3	00:00:01	60.000	20.000	100.00	100.00

Program view.

A new section is always added at the very bottom (example: section 2).

### 9.5.1 Deleting a time program

Path: [Main menu](#) > [Programs](#) > [Time program](#)

In the "Time program" menu select the program to be deleted. The program view opens.



In the **program view** press the **Edit** icon to open the program editor



In the **program editor** select "Delete program" and press the **Confirm** icon.  
The program is deleted. The controller returns to the program view.

## 9.6 Section editor: section management

Path: [Main menu](#) > [Programs](#) > [Time program](#)

Select the desired program.

program 1 - Time program 20:52:55

No.	Duration [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]	Humidity [%RH]
1	00:00:01	37.000	5.0000	8.0000	93.000
2	00:00:01	39.000	5.0000	8.0000	93.000
3	00:00:01	37.000	5.0000	8.0000	93.000

Program view.

Select the desired program section  
(example: section 1)



program 1 - Section number 1 20:55:07

Duration 00:00:01

Course Ramp

Functions on/off 0000000000000000

Number of repetitions 0

Start section for repetition 1

Temperature +37.000

Tolerance band min. +0.0000

Tolerance band max. +0.0000

CO2 +5.0000

Section view (example: section 1).

There are the following options:

- ① Select a parameter to enter or modify the according value (chap. 9.7)
- ② Press the **Edit** icon to open the program editor



program 3 - Section number 1 11:01:06

Edit section

Copy section

Delete section

Add new section

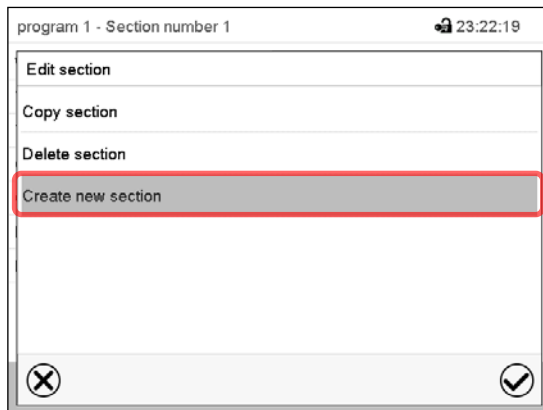
Section editor: "Edit section" menu

Select the desired function and press the **Confirm** icon.

The section editor offers following options:

- Copy section
- Replace section: Replacing an existing section with the copied section. This menu point is visible only after a section has been copied.
- Insert section: Adding the copied section. This menu point is visible only after a section has been copied.
- Delete section
- Add new section


### 9.6.1 Add a new program section



Section editor: "Edit section" menu.


Select "Create new section" and press the **Confirm** icon.

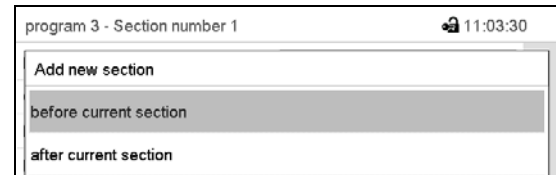Then select whether to insert the new section before or after the current section.




Press the **Confirm** icon. The new section opens.

### 9.6.2 Copy and insert or replace a program section

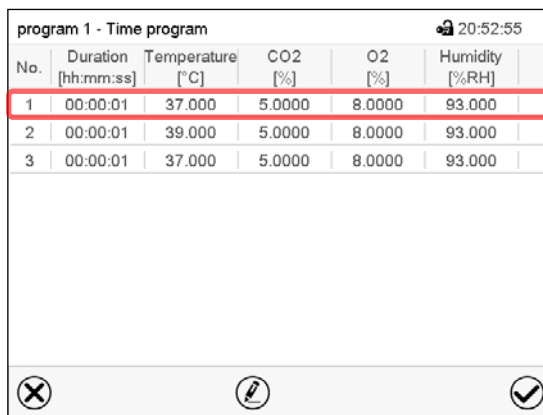


Program view.

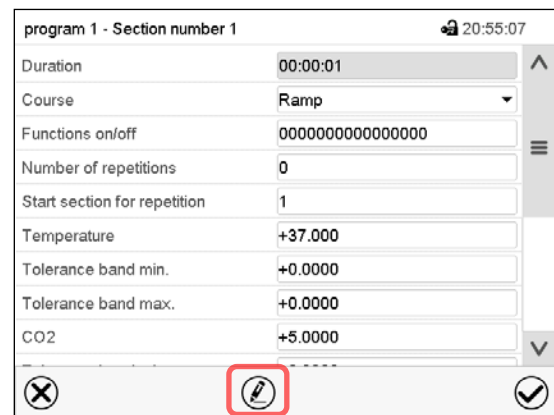Select the program section to be copied (example: section 1)



Section view (example: section 1).

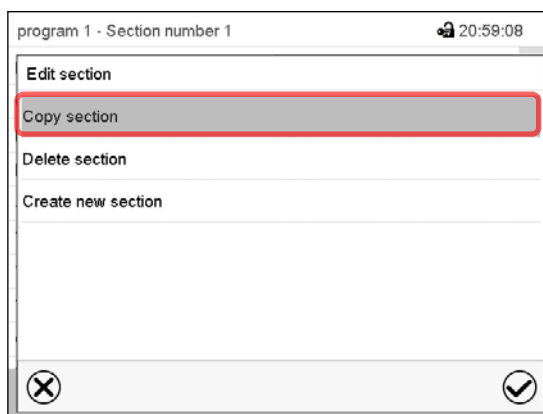Press the **Edit** icon to open the section editor.




Section editor: "Edit section" menu

Select "Copy section" and press the **Confirm** icon.

The current section (example: section 1) is copied.

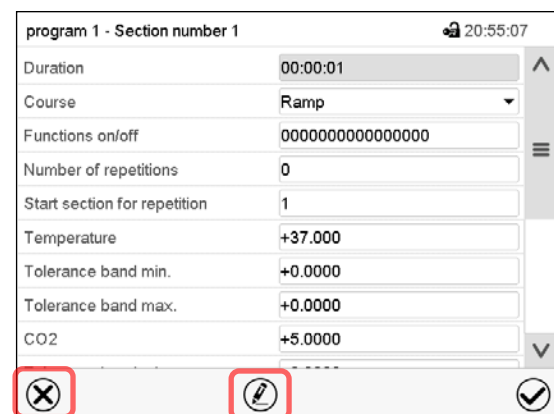The controller returns to the section view.



Section view (example: section 1).

Press the **Close** icon to change to the program view, if you want to select another section to be replaced or before or after which the copied section shall be inserted...



**or**

Press the **Edit** icon to open the section editor if you want the current section to be replaced or the copied section to be inserted before or after it

program 1 - Time program 20:52:55

No.	Duration [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]	Humidity [%RH]
1	00:00:01	37.000	5.0000	8.0000	93.000
2	00:00:01	39.000	5.0000	8.0000	93.000
3	00:00:01	37.000	5.0000	8.0000	93.000

✕ ✎ ✓

Program view.

Select the section to be replaced or before or after which the copied section shall be inserted (example: section 2) and press the **Confirm** icon.

program 1 - Section number 1 20:55:07

Duration 00:00:01

Course Ramp

Functions on/off 0000000000000000

Number of repetitions 0

Start section for repetition 1

Temperature +37.000

Tolerance band min. +0.0000

Tolerance band max. +0.0000

CO2 +5.0000

✕ ✎ ✓

Section view (example: section 1).

Press the **Edit** icon to open the section editor

program 1 - Section number 1 21:00:40

Edit section

Copy section

Replace section

Insert section

Delete section

Create new section

✕ ✓

Section editor: "Edit section" menu

Select "Replace section" to replace the selected section with the copied section

**or**

Select "Insert section" to additionally add the copied section.

In this case select whether to insert it before or after the selected section.

Insert section

before current section

after current section

Press the **Confirm** icon

### 9.6.3 Deleting a program section

In the **program view** select the program section to be deleted. The section view opens.



In the **section view** press the **Edit** icon to open the section editor



In the **section editor** select "Delete section" and press the **Confirm** icon.  
The section is deleted. The controller returns to the section view.



### 9.7.2 Set-point ramp and set-point step

You can define the type of transitions for temperature, CO<sub>2</sub> concentration, O<sub>2</sub> concentration (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification) for each individual program section.

#### **“Ramp” mode: Gradual set-points changes**

The set-point of a given program section functions as the section's start value. During the section's duration, each set-point gradually passes to the set-point of the subsequent program section. The actual value follows the continually changing set-point.

If the last program section is in “ramp” mode and the setpoint shall change within this section, then you must program an additional section (with the shortest possible section duration) to provide the target value of the last program section. Otherwise, the setpoint would remain constant during the section's duration.

Programming in the “ramp” mode allows all kinds of temperature and CO<sub>2</sub> / O<sub>2</sub> and humidity transitions:

- Gradual changes of temperature and CO<sub>2</sub> / O<sub>2</sub> concentration and humidity  
The setpoint changes its value gradually during the entered section duration. The actual value follows the continually moving set-point at any time.
- Program sections with constant temperature and CO<sub>2</sub> / O<sub>2</sub> concentration and humidity  
The setpoints (initial values) of two subsequent program sections are identical; so the temperature and humidity remain constant during the entire duration of the first program section.
- Sudden changes of temperature and CO<sub>2</sub> / O<sub>2</sub> concentration and humidity  
Steps can be programmed in ramp mode as temperature / CO<sub>2</sub> / O<sub>2</sub> / humidity changes (ramps) that occur during a very short interval. If the duration of this transitional program section is very short (minimum entry 1 sec), the temperature, CO<sub>2</sub> / O<sub>2</sub> or humidity change will proceed rapidly within the minimum amount of time.

#### **“Step” mode: Sudden set-points changes**

The set-point of any program section functions as the section's target value. At the start of the program section, heating up, attaining the set gas concentration and humidification occur with the maximum speed to reach the set-point values and then remain constant for the remaining section time. Therefore the set-points remain constant for the section's duration. These changes occur rapidly within the minimum amount of time (minimum entry: 1 second).

Programming in the “step” mode allows only two kinds of temperature and humidity transitions:

- Programming gradual changes of temperature / CO<sub>2</sub> / O<sub>2</sub> / humidity (ramps) is impossible in the “step” mode
- Program sections with constant temperature and humidity  
The setpoints (target values) of two subsequent program sections are identical; so the temperature and humidity remain constant during the entire duration of the first program section.
- Sudden changes of temperature and humidity  
The entered setpoint of the section is reached as fast as possible and then held constant for the remaining section duration.

## Selecting the setting “Ramp” or “Step”

program 3 - Section number 2 11:17:48

Duration

Course

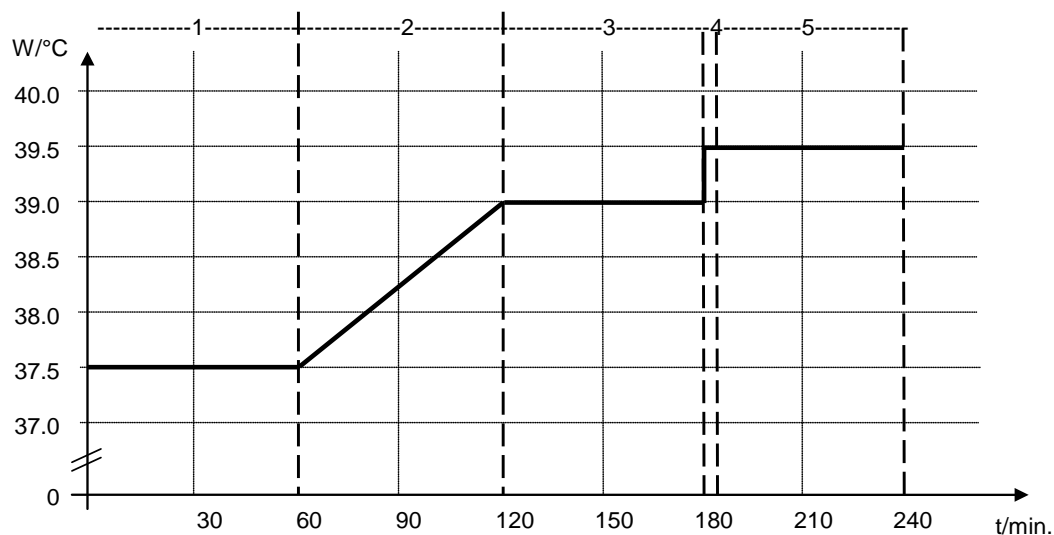
Functions on/off

Number of repetitions

Section view (partial view).

In the field “Course” select the desired setting “Ramp” or “Step”.

## “Ramp” and “Step” mode example (representation of a temperature course)



## Corresponding program table

Section No.	Duration [hh:mm:ss]	Temperature [°C]	CO2 [Vol.-%]	O2 [Vol.-%]	Humidity [% rH]	Ramp or Step
1	01:00:00	37.5	xxxx	xxxx	xxxx	Step
2	01:00:00	37.5	xxxx	xxxx	xxxx	Ramp
3	01:00:00	39.0	xxxx	xxxx	xxxx	Step
4	00:00:10	39.5	xxxx	xxxx	xxxx	Step
5	01:00:00	39.5	xxxx	xxxx	xxxx	Step



### 9.7.3 Special controller functions via operation lines

You can define the switching state of up to 16 operation lines (control contacts). They are used to activate / deactivate special controller functions.

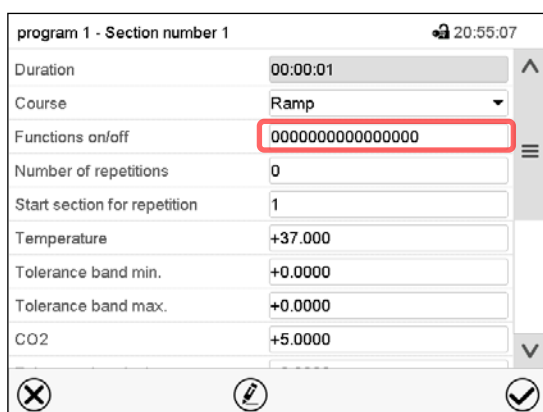
*Chamber with O<sub>2</sub> control:* Depending on the required range of O<sub>2</sub> concentration, the type of gas which is required may vary (chap. 6.6). Therefore, the pressure alarms for O<sub>2</sub> and N<sub>2</sub> can be deactivated individually with active O<sub>2</sub> control.

For chambers equipped with the optional interior socket (chap. 18.3) you can turn on and off the voltage of the interior socket via the controller.

- Operation line "Idle mode" activates / deactivates the operating mode "Idle mode".
- Operation line "O<sub>2</sub> pressure alarm off" activates / deactivates the O<sub>2</sub> pressure alarm (chamber with O<sub>2</sub> control)
- Operation line "N<sub>2</sub> pressure alarm off" activates / deactivates the N<sub>2</sub> pressure alarm (chamber with O<sub>2</sub> control)
- Operation line "Interior socket" activates / deactivates interior socket (option, available via BINDER INDIVIDUAL Customized Solutions)


The other operation lines are without function.

Use the Section editor to configure the operation lines.



Section view.

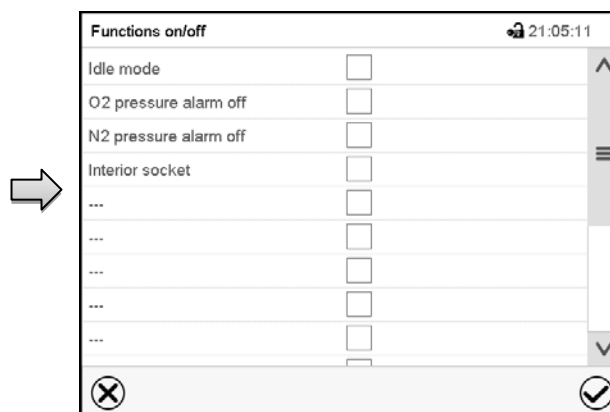Select the field "Functions on/off".



"Functions on/off" entry menu (view with chamber with O<sub>2</sub> control with optional interior socket).

Mark / unmark the checkbox of the desired function to activate / deactivate it and press the Confirm icon.

The controller returns to the section view.

Activated operation line: switching status "1" (On)

Deactivated operation line: switching status "0" (Off)

The operation lines count from right to left.

#### Example:

Activated operation line "Idle mode" = 00000000000000001

Deactivated operation line "Idle mode" = 00000000000000000

### 9.7.4 Setpoint entry

- Select the field “Temperature” and enter the desired temperature setpoint.  
Setting range: 20 °C up to 60 °C.  
Confirm entry with **Confirm** icon. The controller returns to the section view.
- Select the field “CO<sub>2</sub>” and enter the desired CO<sub>2</sub> setpoint.  
Setting range: 0 vol.-% up to 20 vol.-%.  
Confirm entry with **Confirm** icon. The controller returns to the section view.

*Chamber with O<sub>2</sub> control:*

- Select the field “O<sub>2</sub>” and enter the desired O<sub>2</sub> setpoint.  
Setting range with regular hypoxic control range: 0.2 vol.-% up to 20 vol.-%.  
Setting range with alternative control range: 5 vol.-% up to 95 vol.-%.  
Confirm entry with **Confirm** icon. The controller returns to the section view.

*Chamber with active humidification:*

- Select the field “Humidity” and enter the desired humidity setpoint.  
Setting range: 60% r.h. up to 93% r.h.  
Confirm entry with **Confirm** icon. The controller returns to the section view.

### 9.7.5 Tolerance range

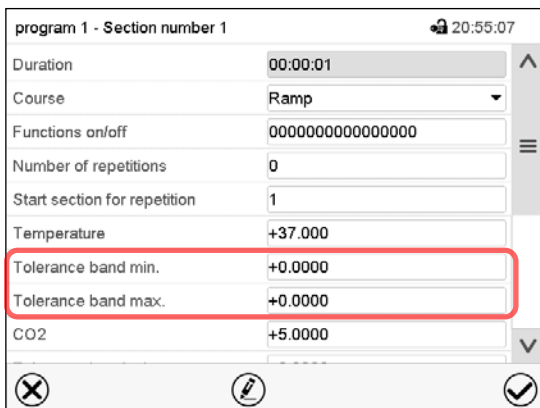
You can specify tolerance ranges for temperature, CO<sub>2</sub> and O<sub>2</sub> (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification) for each program section with different values for the tolerance minimum and maximum. When the actual value exceeds the given threshold, the program is interrupted. This is indicated on the display (see below). When the actual temperature is situated again within the entered tolerance limits, the program automatically continues. Therefore, the duration of the program may be extended due to the programming of tolerances.



Programming of tolerances may extend program duration.

An entry of “-99999” for the tolerance minimum means “minus infinite” and an entry of “999999” for the tolerance maximum means “plus infinite”. Entry of these values will never lead to program interruption. The entry of “0” for the tolerance minimum and/or maximum deactivates the respective tolerance function.

When requesting rapid value transitions, we recommend not programming tolerance values in order to enable the maximum speed for heating-up and CO<sub>2</sub> / O<sub>2</sub> control, and humidification.



Section view, showing the temperature tolerance band

- Select the field “Tolerance band min” and enter the desired lower tolerance band value. Setting range: -99999 to 99999. Confirm entry with **Confirm** icon. The controller returns to the section view.
- Select the field “Tolerance band max” and enter the desired upper tolerance band value. Setting range: -99999 to 99999. Confirm entry with **Confirm** icon. The controller returns to the section view.

Set the tolerance ranges for other parameters accordingly, if desired.

If one of the actual values (temperature, CO<sub>2</sub>, O<sub>2</sub>, humidity) is outside the program tolerance range the whole program course is interrupted. During this program interruption time the controller equilibrates to the set-points of the current section.

The screen header indicates “Program pause (tolerance band)”. The program runtime indication flashes and does not proceed any further.


When the temperature or CO<sub>2</sub> / O<sub>2</sub> or humidity values are back within the entered program tolerance range, the program continues automatically.

### 9.7.6 Repeating one or several sections within a time program

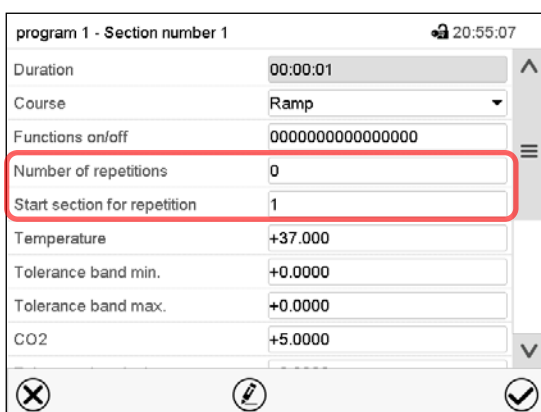You can repeat several subsequent sections together. It is not possible to define the start section the same time also as the target section, therefore you cannot repeat a single individual section.

Enter the desired number of repetitions in the field „Number of repetitions“ and the number of the section to start the repetition cycle with in the field “Start section for repetition” To have sections repeated infinitely, enter the number of repetitions as “-1”.

The selected sections are repeated as many times as selected. Then the program continues.



program 1 - Section number 1		20:55:07
Duration	00:00:01	↑
Course	Ramp	▼
Functions on/off	0000000000000000	≡
Number of repetitions	0	
Start section for repetition	1	
Temperature	+37.000	
Tolerance band min.	+0.0000	
Tolerance band max.	+0.0000	
CO2	+5.0000	▼

Section view, showing the repetition function

- Select the field “Number of repetitions” and enter the desired number of repetitions. Setting range: 1 to 99, and -1 for infinite. Confirm entry with **Confirm** icon. The controller returns to the section view.
- Select the field “Start section for repetition” and enter the section number, at which the repetition should start. Setting range: 1 up to the section before the currently selected section. Confirm entry with **Confirm** icon. The controller returns to the section view.

### 9.7.7 Saving the time program

program 1 - Section number 1 20:55:07

Duration 00:00:01

Course Ramp

Functions on/off 0000000000000000

Number of repetitions 0

Start section for repetition 1

Temperature +37.000

Tolerance band min. +0.0000

Tolerance band max. +0.0000

CO2 +5.0000

✕ ⓘ **✓**

Section view.

After the all desired values of the program section have been configured, press the **Confirm** icon to take over the programming.  
The controller changes to the program view.



program 1 - Time program 20:51:05

No.	Duration [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]	Humidity [%RH]
1	00:00:01	37.000	5.0000	8.0000	93.000
2	00:00:01	39.000	5.0000	8.0000	93.000
3	00:00:01	60.000	20.000	100.00	100.00

✕ ⓘ **✓**

Program view.

Press the **Confirm** icon to take over the programming.

The controller changes to the Normal display.




To save the programming it is absolutely required to press the **Confirm** icon. Otherwise all settings will be lost! There is no confirmation prompt!

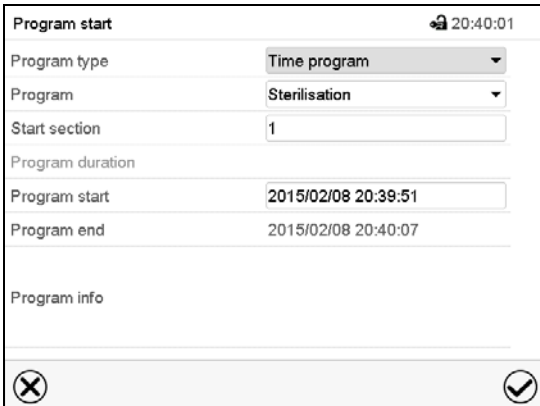
## 10. Week programs

The MB2 program controller permits programming week programs with real-time reference. It offers 5 week program places in total with up to 100 shift points for each week program.

Path: [Main menu > Programs > Week program](#)

### 10.1 Starting an existing week program

 In Normal display press the **Program start** icon to access the “Program start” menu.




“Program start” menu.

- In the field “Program type” select the setting “Week program”.
- In the field “Program” select the desired program.
- There are no further settings available in the “Program start” menu for week programs, as they are needed only for time programs.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu. The program starts running.


If instead you press the **Close** icon to exit the menu without taking over the entries, the program will not start.

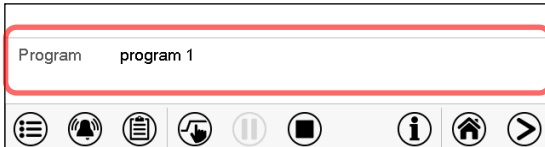After starting the week program, the previously entered week program setpoints are active and will be equilibrated according to the current time.



Information on the bottom of the screen indicates the currently running program.

### 10.2 Cancelling a running week program

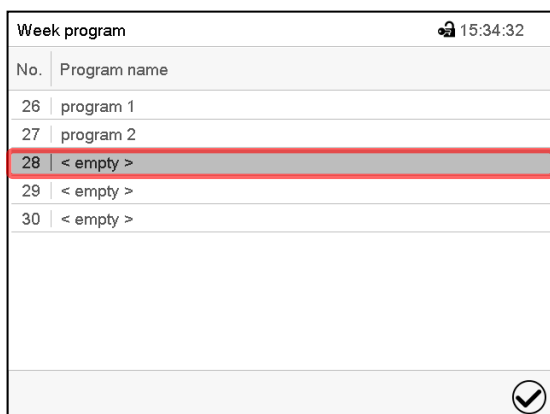
 Press the **Program cancelling** icon to cancel the program.

A confirmation prompt is displayed. Press the **Confirm** icon to confirm that the program shall really be cancelled.

After confirming the message the controller changes to Fixed value operation mode. Temperature and humidity will then equilibrate to the setpoints of Fixed value operation mode.


### 10.3 Creating a new week program

Path: [Main menu](#) > [Programs](#) > [Week program](#)



No.	Program name
26	program 1
27	program 2
28	< empty >
29	< empty >
30	< empty >

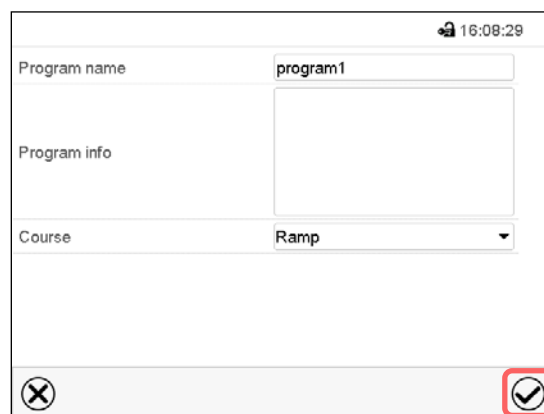"Week program" menu:  
overview of the existing programs.  
Select an empty program place.

Program name:

Program info:

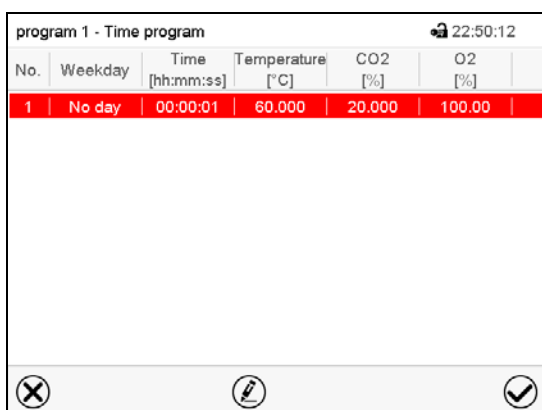
Course: Ramp



Enter the program name and, if desired,  
additional program information in the  
corresponding fields.

Select the set-point course "Ramp" or "Step"  
(chap. 10.6.1).

Press the **Confirm** icon. The program view  
opens.



No.	Weekday	Time [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]
1	No day	00:00:01	60.000	20.000	100.00

Program view.

For the first section no weekday is specified.  
Therefore the section is first marked in red and  
cannot be saved.

## 10.4 Program editor: program management

Path: **Main menu > Programs > Week program**

Week program		08:39:23	
No.	Program name		
26	program 1		
27	program 2		
28	< empty >		
29	< empty >		
30	< empty >		

“Week program” menu:  
overview of the existing programs.  
Select an existing program  
(example: program 1).

program 1 - Time program						23:00:01	
No.	Weekday	Time [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]		
1	Monday	03:00:00	37.000	5.0000	8.0000		
2	Tuesday	00:00:01	39.000	8.0000	100.00		

Program view (example: program 1).  
If a new program has been created, there is just one program section.

There are the following options:

- ① Select a program section to open the section editor (chap. 10.4.1)
- ② Press the **Edit** icon to open the program editor

program 1 - Week program

15:43:30

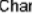
Edit program


Change program name

Copy program

Delete program

Create new section







Program editor: “Edit program” menu.  
Select the desired function and press the **Confirm** icon.

The program editor offers following options:

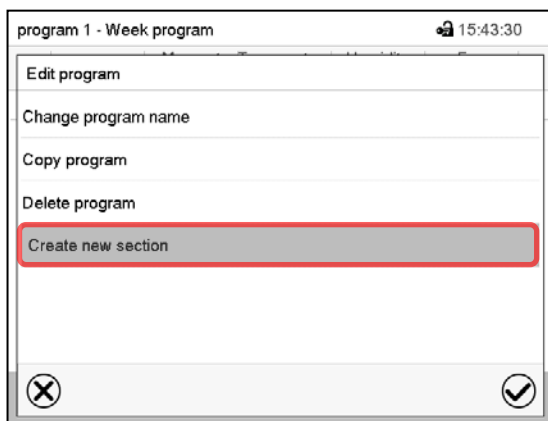- Change program name. This menu also offers to configure the ramp / step mode setting (chap. 10.6.1).
- Copy program
- Replace program: Replacing an new or an existing program with the copied program. This menu point is visible only after a section has been copied.
- Delete program
- Add new section



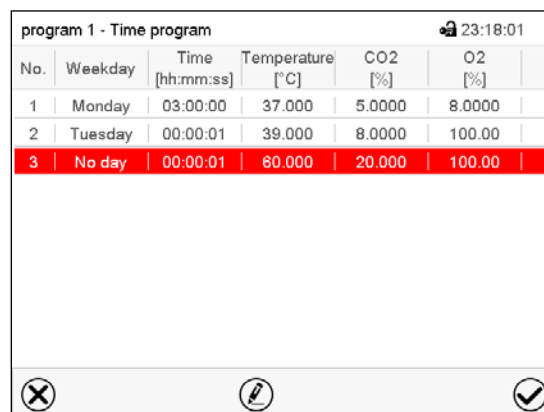To add a new section, select “Create new section” and press the **Confirm** icon.  
The program view opens.

No.	Weekday	Time [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]
1	Monday	03:00:00	37.000	5.0000	8.0000
2	Tuesday	00:00:01	39.000	8.0000	100.00
3	No day	00:00:01	60.000	20.000	100.00

Program view.

With a new section no weekday is specified.  
Therefore the section is first marked in red and cannot be saved.

A new section is always added at the very bottom (example: section 3). When the section start is specified the sections are automatically arranged in the correct chronological order.

#### 10.4.1 Deleting a week program

Path: [Main menu](#) > [Programs](#) > [Week program](#)

In the “Week program” menu select the program to be deleted. The program view opens.



In the **program view** press the **Edit** icon to open the program editor



In the **program editor** select “Delete program” and press the **Confirm** icon.  
The program is deleted. The controller returns to the program view.



## 10.5 Section editor: section management

Path: [Main menu](#) > [Programs](#) > [Week program](#)

Select the desired program.

program 1 - Time program 23:00:01

No.	Weekday	Time [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]
1	Monday	03:00:00	37.000	5.0000	8.0000
2	Tuesday	00:00:01	39.000	8.0000	100.00

✕ ✎ ✓

Program view.

Select the desired program section  
(example: section 1)



program 1 - Section number 1 22:58:41

Weekday: Monday

Time: 03:00:00 ①

Temperature: +37.000 ②

CO2: +5.0000

O2: +8.0000

Humidity: +93.000

Functions on/off: 0000000000000000

✕ ✎ ✓

Section view (example: section 1).

There are the following options:

- ① Select a parameter to enter or modify the according value (chap. 10.6)
- ② Press the **Edit** icon to open the program editor



program 1 - Section number 1 23:22:19

Edit section

Copy section

Delete section

Create new section

✕ ✓

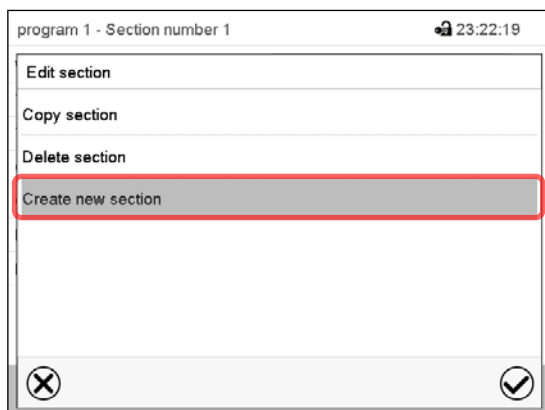
Section editor: "Edit section" menu

Select the desired function and press the **Confirm** icon.

The section editor offers following options:


- Copy section
- Replace section: Replacing an existing section with the copied section. This menu point is visible only after a section has been copied.
- Insert section: Adding the copied section. This menu point is visible only after a section has been copied.
- Delete section
- Create new section

### 10.5.1 Add a new program section



Section editor: "Edit section" menu.

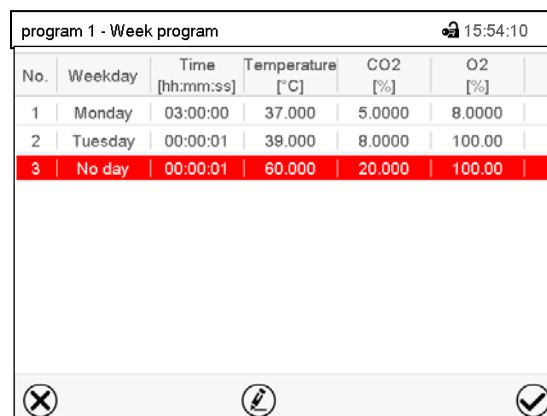Select "Create new section" and press the **Confirm** icon.




No.	Weekday	Time [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]
1	Monday	03:00:00	37.000	5.0000	8.0000
2	Tuesday	00:00:01	39.000	8.0000	100.00
3	No day	00:00:01	60.000	20.000	100.00

Program view.

With a new section no weekday is specified. Therefore the section is first marked in red and cannot be saved.

A new section is always added at the very bottom (example: section 3). When the section start is specified the sections are automatically arranged in the correct chronological order.

### 10.5.2 Copy and insert or replace a program section

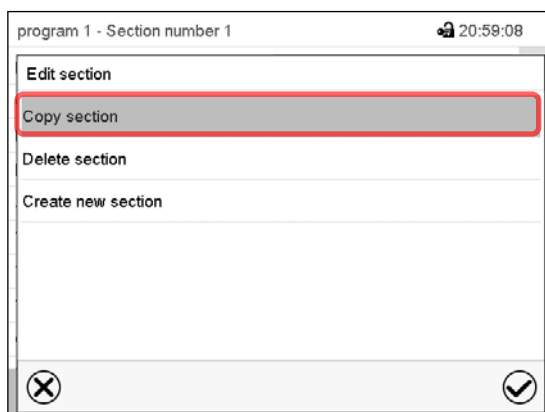


Section editor: "Edit section" menu

Select "Copy section" and press the **Confirm** icon.

The current section (example: section 1) is copied.

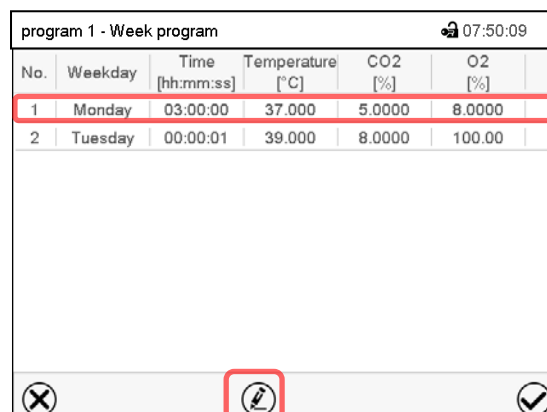The controller returns to the program view.




No.	Weekday	Time [hh:mm:ss]	Temperature [°C]	CO2 [%]	O2 [%]
1	Monday	03:00:00	37.000	5.0000	8.0000
2	Tuesday	00:00:01	39.000	8.0000	100.00

Program view

Select the section to be replaced or before or after which the copied section shall be inserted (example: section 2).

Press the **Confirm** icon

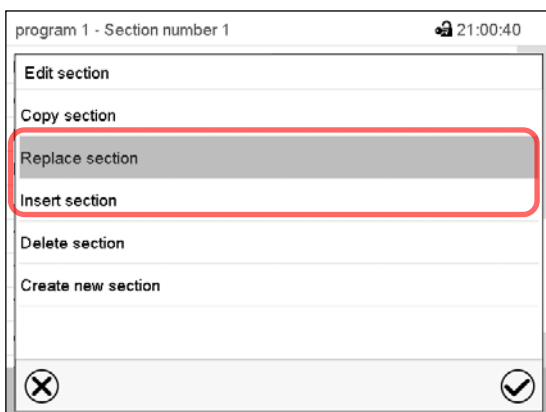The controller returns to the section editor



Section editor: "Edit section" menu

Select "Replace section" to replace the selected section with the copied section

or

Select "Insert section" to additionally add the copied section.

Press the **Confirm** icon.

If you selected "Insert section" the sections are automatically arranged in the correct chronological order.

### 10.5.3 Deleting a program section

In the **program view** select the program section to be deleted. The section view opens.

➡ In the **section view** press the **Edit** icon to open the section editor

➡ In the **section editor** select "Delete section" and press the **Confirm** icon.  
The section is deleted. The controller returns to the section view.

## 10.6 Value entry for a program section

Path: [Main menu > Programs > Week program](#)

Select the desired program and section.

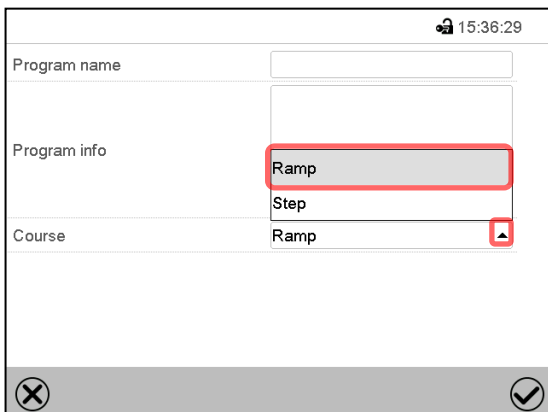
The setting and control ranges for the individual parameters are the same as for "Fixed value" operating mode (chap. 7).

### 10.6.1 Set-point ramp and set-point step modes

The explanation of the settings "Ramp" or "Step" is given in chap. 9.7.2.

You can define the type of transitions for temperature, CO<sub>2</sub> concentration, O<sub>2</sub> concentration (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification) for the entire week program.


Select the desired program and press the **Edit** icon to open the program editor. In the program editor select the "Change program name" function and press the **Confirm** icon.



“Change program name” menu.


In the field “Course” select the desired setting “Ramp” or “Step” and press the **Confirm** icon.

## 10.6.2 Weekday

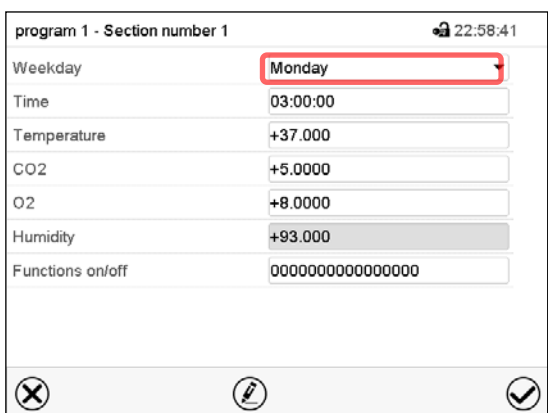


Section view.

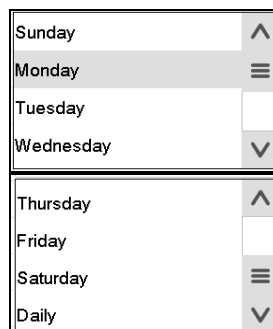In the field “Weekday” select the desired weekday.




With “Daily” selected, this section will run every day at the same time.

## 10.6.3 Start time

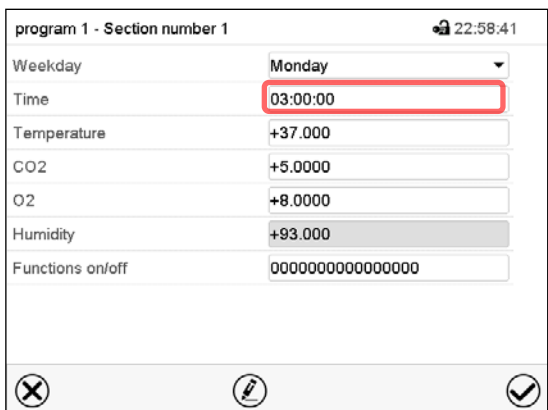

Section view.

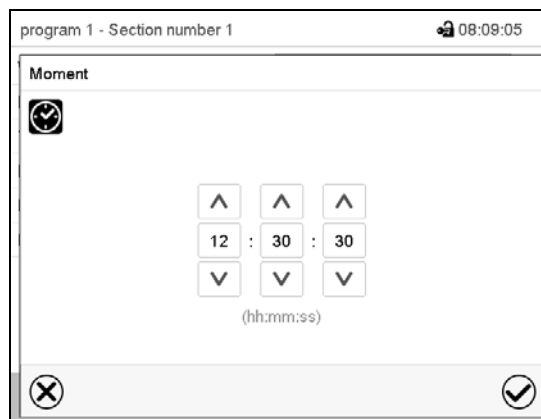Select the field “Moment”.

Entry menu “Moment”.

Select with the arrow keys the desired start moment of the section and press the **Confirm** icon.

#### 10.6.4 Setpoint entry

- Select the field “Temperature” and enter the desired temperature setpoint.  
Setting range: 20 °C up to 60 °C.  
Confirm entry with **Confirm** icon. The controller returns to the section view.
- Select the field “CO<sub>2</sub>” and enter the desired CO<sub>2</sub> setpoint.  
Setting range: 0 vol.-% up to 20 vol.-%.  
Confirm entry with **Confirm** icon. The controller returns to the section view.

*Chamber with O<sub>2</sub> control:*

- Select the field “O<sub>2</sub>” and enter the desired O<sub>2</sub> setpoint.  
Setting range with regular hypoxic control range: 0.2 vol.-% up to 20 vol.-%.  
Setting range with alternative control range: 5 vol.-% up to 95 vol.-%.  
Confirm entry with **Confirm** icon. The controller returns to the section view.

*Chamber with active humidification:*

- Select the field “Humidity” and enter the desired humidity setpoint.  
Setting range: 60% r.h. up to 93% r.h.  
Confirm entry with **Confirm** icon. The controller returns to the section view.

#### 10.6.5 Special controller functions via operation lines

You can define the switching state of up to 16 operation lines (control contacts). They are used to activate / deactivate special controller functions.

*Chamber with O<sub>2</sub> control:* Depending on the required range of O<sub>2</sub> concentration, the type of gas which is required may vary (chap. 6.6). Therefore, the pressure alarms for O<sub>2</sub> and N<sub>2</sub> can be deactivated individually with active O<sub>2</sub> control.

For chambers equipped with the optional interior socket (chap. 18.3) you can turn on and off the voltage of the interior socket via the controller.

- Operation line “Idle mode” activates / deactivates the operating mode “Idle mode”.
- Operation line “O<sub>2</sub> pressure alarm off” activates / deactivates the O<sub>2</sub> pressure alarm (chamber with O<sub>2</sub> control)
- Operation line “N<sub>2</sub> pressure alarm off” activates / deactivates the N<sub>2</sub> pressure alarm (chamber with O<sub>2</sub> control)
- Operation line “Interior socket” activates / deactivates interior socket (option, available via BINDER INDIVIDUAL Customized Solutions)

The other operation lines are without function.

Select the desired program and section. You can set the operation lines in the “Functions on/off” field.

*For details please refer to chap. 9.7.3.*

## 11. Notification and alarm functions

### 11.1 Notification and alarm messages overview

#### 11.1.1 Notifications

Notifications are indicated by **information icons** displayed in the screen header in Normal display

An information icon serves as an indication of a certain condition.







If this condition persists, in some cases an alarm will be triggered after a fix or configurable interval. As long as the condition persists, the information icon therefore continues to be displayed also in state of alarm. If during alarm the conditions ends, e.g., if during a tolerance range alarm the actual value returns to within the tolerance range, the information icon disappears, whereas the alarm will continue until manual acknowledgement.

Press the flash icon next to the information icon to access the corresponding text information.



Normal display showing the text information.

The currently valid information texts are highlighted in black (example: "O2 pressure alarm off")

Condition	Information icon	Text information	Start after condition occurred
The controller is in Idle mode (chap. 5.4).		"Idle mode"	immediately
Chamber door open		"Door open"	immediately
CO <sub>2</sub> pressure alarm deactivated		"CO2 pressure alarm off"	immediately
Chamber with O <sub>2</sub> control: O <sub>2</sub> pressure alarm deactivated		"O2 pressure alarm off"	immediately
Chamber with O <sub>2</sub> control: N <sub>2</sub> pressure alarm deactivated		"N2 pressure alarm off"	immediately
Interior socket (option) activated		"Interior socket"	immediately

Notifications are not shown in the event list.

### 11.1.2 Alarm messages

Condition	Alarm message	Start after condition occurred	Zero-voltage relay alarm output
<p>Having reached the setpoint, the actual temperature value deviates by more than the set tolerance range value and longer than the set delay time from the setpoint</p> <p><b>or</b></p> <p>the temperature doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door.</p>	"Temperature range"	After configurable delay time (chap. 11.5). Factory setting: 10 minutes	With alarm start
<p>Having reached the setpoint, the actual CO<sub>2</sub> concentration deviates by more than the set tolerance range value and longer than the set delay time from the setpoint</p> <p><b>or</b></p> <p>the CO<sub>2</sub> concentration doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door</p>	"CO <sub>2</sub> range"	After configurable delay time (chap. 11.5). Factory setting: 10 minutes	With alarm start
<p><i>Chamber with O<sub>2</sub> control:</i></p> <p>Having reached the setpoint, the actual O<sub>2</sub> concentration deviates by more than the set tolerance range value and longer than the set delay time from the setpoint</p> <p><b>or</b></p> <p>the O<sub>2</sub> concentration doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door</p>	"O <sub>2</sub> range"	After configurable delay time (chap. 11.5). Factory setting: 10 minutes	With alarm start
<p><i>Chamber with active humidification:</i></p> <p>Having reached the setpoint, the actual humidity value deviates by more than the set tolerance range value and longer than the set delay time from the setpoint</p> <p><b>or</b></p> <p>the humidity doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door.</p>	"Humidity range"	After configurable delay time (chap. 11.5). Factory setting: 10 minutes	With alarm start
Chamber door is open for longer than the set delay time.	"Door open"	After configurable delay time (chap. 11.5). Factory setting: 1 minute	----
Power failure	---	----	immediately

Condition	Alarm message	Start after condition occurred	Zero-voltage relay alarm output
Exceeded setpoint of the safety controller class 3.1	"Safety controller	immediately	----
Low CO <sub>2</sub> outlet pressure (< 0.3 bar / 4.4 psi)	"CO <sub>2</sub> pressure"	immediately	----
<i>Chamber with O<sub>2</sub> control:</i> Low O <sub>2</sub> outlet pressure (< 0.3 bar / 4.4 psi)	"O <sub>2</sub> pressure"	immediately	----
<i>Chamber with O<sub>2</sub> control:</i> Low N <sub>2</sub> outlet pressure (< 0.3 bar / 4.4 psi)	"N <sub>2</sub> pressure"	immediately	----
Temperature sensor defective	e.g. "- - - -" or "<-<-<" or ">->->"	immediately	immediately
Safety controller temperature sensor defective	Safety controller sensor	immediately	immediately
CO <sub>2</sub> sensor defective. Contact BINDER Service.	"CO <sub>2</sub> sensor fault"	immediately	immediately
<i>Chamber with O<sub>2</sub> control:</i> O <sub>2</sub> sensor defective. Contact BINDER Service	"O <sub>2</sub> sensor fault"	immediately	immediately
<i>Chamber with active humidification:</i> Humidity sensor defective. Contact BINDER Service	"Humid sensor fault"	immediately	immediately
Hot air sterilization running	"Sterilization"	immediately	----

Alarm messages are displayed in the list of active alarms until acknowledging them. They are also shown in the event list.



When operating the chamber without CO<sub>2</sub> / O<sub>2</sub> / N<sub>2</sub> connection, turn off the corresponding gas control in the "setpoints" menu (chap. 6.8) in order to avoid tolerance range and pressure alarms.

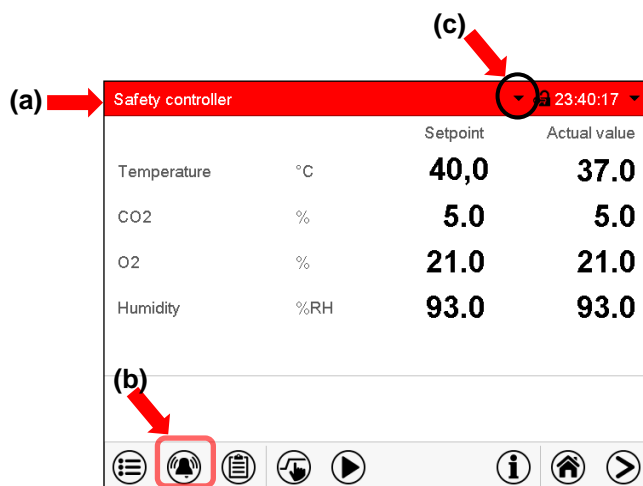


*Chamber with active humidification:* When operating the chamber without water supply, turn off humidity control in the "setpoints" menu (chap. 6.8) in order to avoid humidity alarms..



## 11.2 State of alarm

1. Visual indications in Normal display: alarm message, screen header flashing in red color
2. Audible alert, if the buzzer is enabled (chap. 11.4).
3. In some cases: Switching the zero-voltage relay alarm output (chap. 11.6) to transmit the alarm e.g., to a central monitoring system.




Normal display in state of alarm (example).

(a) Screen header flashing in red color and showing the alarm message

(b) **Alarm** icon on the bottom of the screen: change to the list of active alarms and alarm acknowledgement


(c) If applicable, information icon in the screen header. Indication of a certain condition

## 11.3 Resetting an alarm, list of active alarms

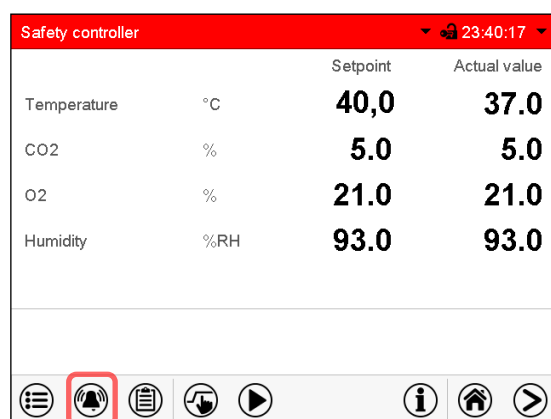

Normal display in state of alarm (example).

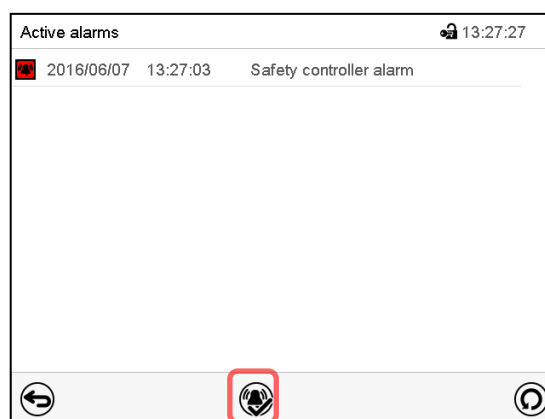Press the **Alarm** icon



List of active alarms.

Press the **Reset alarm** icon.

Pressing the **Reset alarm** icon mutes the buzzer for all active alarms. The icon then disappears.

- Acknowledging while the alarm condition persists: Only the buzzer turns off. The visual alarm indication remains on the controller display. The alarm remains in the list of active alarms.


When the alarm condition has ended, the visual alarm indication is automatically cleared. The alarm is then no longer in the list of active alarms.

- Acknowledging after the alarm condition has ended: The buzzer and the visual alarm indication are reset together. The alarm is then no longer in the list of active alarms.
- The zero-voltage relay alarm output resets together with the alarm.

For appropriate actions in the event of an alarm, please refer to chap. 24 "Troubleshooting".

## 11.4 Activating / deactivating the audible alarm (alarm buzzer)

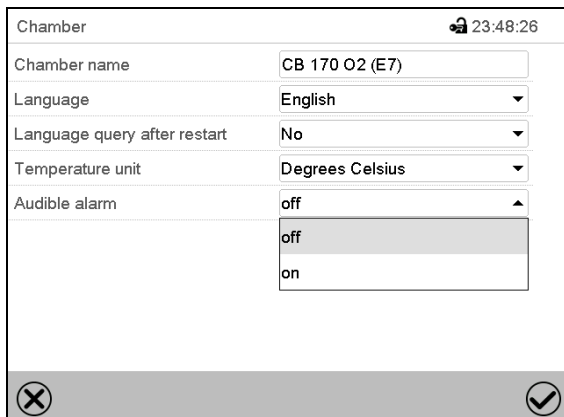Path: [Main menu](#) > [Settings](#) > [Chamber](#)




“Chamber” submenu (example).

In the field “Audible alarm” select the desired setting “off” or “on” and press the **Confirm** icon.

## 11.5 Tolerance range settings and alarm delay times

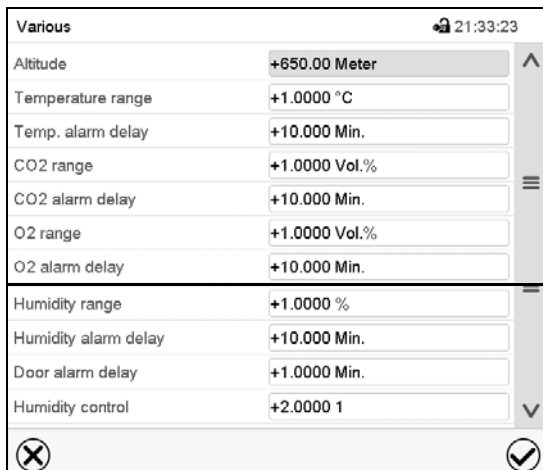Path: [Main menu](#) > [Settings](#) > [Various](#)



Submenu “Various”.

Select the desired function.

In this menu you can set the deviation between the actual value and setpoint for temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification) which that shall cause a tolerance range alarm, and you can define the delay time after which the tolerance alarm will be triggered.

Also you can set the delay time after opening the door will trigger an alarm.

The entered value defines the tolerance range around the setpoint value. Example: Temperature setpoint: 37 °C, tolerance range value: +/- 2 °C, i.e. this defines a tolerance range from 35 °C up to 39 °C.

### Alarms following deviations from the tolerance range or door opening

If the actual value, after having reached the setpoint, deviates by more than the set tolerance range value and longer than 10 min. from the setpoint **or** the temperature doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door, after the configured interval (e.g., “Temp range alarm delay”) an alarm is triggered. It is visually indicated in Normal display. If the alarm buzzer is activated (chap. 11.4) there is an audible alert. The zero-voltage relay alarm output (chap. 11.6) switches to transmit the alarm. The alarm is shown in the list of active alarms (chap. 11.3).

Alarm message	Start after condition occurred
"Temperature range"	after configurable time
"CO2 range"	after configurable time
"O2 range"	after configurable time
"Humidity range"	after configurable time
"Door open"	after configurable time

If the condition (deviation from the tolerance range or door opening) ends within the configured alarm delay time, no alarm will be triggered after the delay time.



When operating the chamber with deactivated CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control) or humidity (chamber with active humidification) control (chap. 6.8), the corresponding tolerance range alarms will not occur.

### Setting:

- Select the field "Temperature range" and enter the desired value for the temperature range. Setting range: 1 °C to 10 °C. Confirm entry with **Confirm** icon.
- Select the field "Temp alarm delay" and enter the time in minutes, after which the range alarm shall be triggered. Setting range: 1 min to 120 min. Confirm entry with **Confirm** icon.
- Select the field "CO2 range" and enter the desired value for the CO<sub>2</sub> range. Setting range: 1 vol.-% to 10 vol.-%. Confirm entry with **Confirm** icon.
- Select the field "CO2 alarm delay" and enter the time in minutes, after which the range alarm shall be triggered. Setting range: 1 min to 120 min. Confirm entry with **Confirm** icon.
- Select the field "O2 range" and enter the desired value for the O<sub>2</sub> range. Setting range: 1 vol.-% to 10 vol.-%. Confirm entry with **Confirm** icon.
- Select the field "O2 alarm delay" and enter the time in minutes, after which the range alarm shall be triggered. Setting range: 1 min to 120 min. Confirm entry with **Confirm** icon.
- Select the field "Humidity range" and enter the desired value for the humidity range. Setting range: 5% r.h. to 20% r.h. Confirm entry with **Confirm** icon.
- Select the field "Humid alarm delay" and enter the time in minutes, after which the range alarm shall be triggered. Setting range: 1 min to 120 min. Confirm entry with **Confirm** icon.
- Select the field "Door alarm delay" and enter the time in minutes, after which the range alarm shall be triggered. Setting range: 1 min to 120 min. Confirm entry with **Confirm** icon.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

## 11.6 Zero-voltage relay alarm output

The chamber is equipped with a zero-voltage relay output which permits the transmission of some alarm messages to a central monitoring system.

The connection is realized as a DIN socket (3) on the chamber rear. A suitable DIN plug is enclosed.


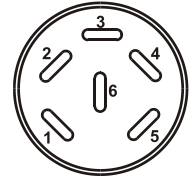


Figure 20: Pin configuration of the DIN socket (3) on the chamber rear

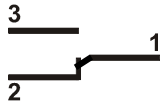



ALARM

Pin 1: Pole, Pin 2: Break relay, Pin 3: Make contact

In case there is no alarm, contact 1 closes with contact 3.

Closing contact 1 with contact 2 switches the zero-voltage relay alarm output.

**Maximum loading capacity of the switching contacts: 24V AC/DC – 2.5A**

	 <b>DANGER</b>
<p><b>Electrical hazard.</b>  <b>Danger of death.</b>  <b>Damage to switching contacts and connection socket.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT exceed the maximum switching load of 24V AC/DC – 2.5A.</li> <li>⊘ Do NOT connect any devices with a higher loading capacity.</li> </ul>	


The zero-voltage relay alarm output switches at the following events:

Alarm message	Condition	Switching the alarm contact
---	Power failure	immediately
"Temperature range"	Temperature tolerance range alarm (see chap. 11.1.2)	After configurable delay time (chap. 11.5). Factory setting: 10 minutes
"CO2 range"	CO <sub>2</sub> tolerance range alarm (see chap. 11.1.2)	After configurable delay time (chap. 11.5). Factory setting: 10 minutes
"O2 range"	O <sub>2</sub> tolerance range alarm (chamber with O <sub>2</sub> control) (see chap. 11.1.2)	After configurable delay time (chap. 11.5). Factory setting: 10 minutes
"Humidity range"	Humidity tolerance range alarm (chamber with active humidification) (see chap. 11.1.2)	After configurable delay time (chap. 11.5). Factory setting: 10 minutes
"Door open"	Chamber door open alarm (see chap. 11.1.2)	After configurable delay time (chap. 11.5). Factory setting: 1 minutes

In case of a tolerance range alarm or a door open alarm, the alarm message on the controller display remains on during the alarm transmission via the zero-voltage relay outputs.

As soon as the cause of the alarm is identified and resolved, you can reset the alarm transmission via the zero-voltage relay outputs together with the alarm message in the "Active alarms" menu by pressing the **Reset alarm** icon.

In case of a power failure, transmission of the alarm via zero-voltage relay outputs remains active for the duration of the power failure. After power returns, contact 1 closes automatically with contact 3.

	<p>When using the communication software APT-COM™ 3 DataControlSystem (option, chap. 18.1) for data acquisition, the alarm messages are only recorded in the protocol.</p> <ul style="list-style-type: none"> <li>➤ Set the tolerance limits for limit alarms by APT-COM™ 3 separately in the APT-COM™ 3 measuring window.</li> </ul>
---	---

## 12. Temperature safety devices

### 12.1 Over temperature protective device (class 1)

The chamber is equipped with an internal temperature safety device, class 1 acc. to DIN 12880:2007. It serves to protect the chamber and prevents dangerous conditions caused by major defects.




If the actual temperature exceeds the nominal temperature by approx. 10 °C, the over temperature protective device permanently turns off the chamber. The user cannot restart the device again. The protective cut-off device is located internally. Only a service specialist can replace it. Therefore, please contact an authorized service provider or BINDER Service.

### 12.2 Overtemperature safety controller class 3.1

The chambers are regularly equipped with an electronic overtemperature safety controller (temperature safety device class 3.1 according to DIN 12880:2007). The safety controller is functionally and electrically independent of the temperature control system. If an error occurs, it performs a regulatory function.

Please observe the DGUV guidelines 213-850 on safe working in laboratories (formerly BGI/GUV-I 850-0, BGR/GUV-R 120 or ZH 1/119, issued by the employers' liability insurance association) (for Germany).

The overtemperature safety controller serves to protect the chamber, its environment and the contents from exceeding the maximum permissible temperature. In the case of an error, it limits the temperature inside the chamber to the entered safety controller set-point. This condition (state of alarm) is indicated visually and additionally with an audible alert if the buzzer is enabled (chap. 11.4). The alarm persists until the chamber cools down below the configured safety controller setpoint.

	Check the setting regularly and adjust it following changes of the set-point or charge.
	The safety controller only activates after the set-point has been reached once.
	During a running sterilization (chap. 21.3) the safety controller is non-functional

## 12.2.1 Safety controller modes

You can select between “Limit (absolute)” and “Offset (relative)” safety controller mode

- **Limit:** Absolute maximum permitted temperature value

This setting offers high safety as a defined temperature limit will not be exceeded. It is important to adapt the safety controller set-point after each modification of the temperature set-point. Otherwise, the limit could be too high to ensure efficient protection, or, in the opposite case, it could prevent the controller from reaching an entered set-point outside the limit range.

- **Offset:** Maximum overtemperature above any active temperature set point. The maximum temperature changes internally and automatically with every set-point change.


This setting is recommended for program operation. It is important to check the safety controller set-point and safety controller mode occasionally, as it does not offer a fix, independent limit temperature value, which would never be exceeded.

**Example:** Desired temperature value: 37 °C, desired safety controller value: 39 °C.

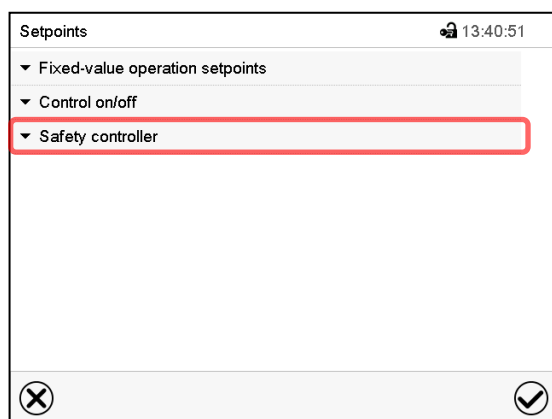
Possible settings for this example:

Temperature set point	Safety controller mode	Safety controller set-point
37 °C	Limit (absolute)	39 °C
	Offset (relative)	2 °C

## 12.2.2 Setting the safety controller




Press the **Setpoint setting** icon to access the “Setpoint” setting menu from Normal display.




“Setpoints” menu.

Select the field “Safety controller” to access the settings.

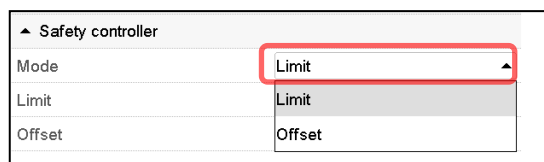- In the field “Mode” select the desired setting “Limit” or “Offset”.



- Select the corresponding field “Limit” or “Offset” according to the selected mode and enter the desired safety controller setpoint. Confirm entry with **Confirm** icon.

	<p>Regularly check the safety controller setting for set-point type "Limit" or "Offset"</p> <ul style="list-style-type: none"> <li>• in Fixed value operating mode according to the entered set-point temperature value</li> <li>• in program mode according to the highest temperature value of the selected temperature program</li> </ul> <p>Set the safety controller set-point by approx. 2 °C above the desired temperature set-point.</p>
---	--

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

### 12.2.3 Message and measures in the state of alarm


The state of alarm is indicated visually in Normal display by the alarm message "Safety controller alarm" and the screen header flashing in red color. If the buzzer is enabled (chap. 11.4) there is an additional audible alert (chap. 11.2). The alarm remains active until it is acknowledged on the controller and the inner temperature falls below the set safety controller setpoint. Then the heating is released again.

Safety controller				23:40:17	
		Setpoint	Actual value		
Temperature	°C	40,0	37.0		
CO2	%	5.0	5.0		
O2	%	21.0	21.0		
Humidity	%RH	93.0	93.0		

Normal display with safety controller alarm.

Press the **Alarm** icon



Active alarms			13:27:27
	2016/06/07 13:27:03	Safety controller alarm	

List of active alarms.

Press the **Reset alarm** icon.

### 12.2.4 Function check

Check the safety controller at appropriate intervals for its functionality. It is recommended that the authorized operating personnel should perform such a check, e.g., before starting a longer work procedure.

## 13. User management

### 13.1 Authorization levels and password protection

The available functions depend on the current authorization level “Master”, “Service”, “Admin” or “User”.

The authorization levels are hierarchical: Every authorization includes all functions of the next lower level.

#### **“Master” authorization level**

- Highest authorization level, only for developers
- Extensive authorization for controller operation and configuration, outputs/inputs, alarm settings, parameter sets and operating ring display
- All passwords can be changed in the “log out” submenu (chap. 13.3).

#### **“Service” authorization level**

- Authorization level only for BINDER service
- Extensive authorization for controller operation and configuration, access to service data
- The passwords for “Service”, “Admin” and “User” authorization levels can be changed in the “log out” submenu (chap. 13.3).

#### **“Admin” authorization level**

- Expert authorization level, for the administrator
- Authorization for controller configuration and network settings and for operating those controller functions required for operating the chamber. Restricted access to service data.
- Password (factory setting): “2”.
- The passwords for “Admin” and “User” authorization levels can be changed in the “log out” submenu (chap. 13.3).

#### **“User” authorization level**

- Standard authorization level for the chamber operator
- Authorization for operating the controller functions required for operating the chamber.
- No authorization for controller configuration and network settings. The “Settings” and “Service” submenus of the main menu are not available.
- Password (factory setting): “1”
- The password for the “User” authorization level can be changed in the “log out” submenu (chap. 13.3).

As soon as a password has been assigned for an authorization level, the access to this level and the related controller functions are only available after log-in with the appropriate password.

If for an authorization level no password is assigned, the related controller functions of this level are available for every user without login.

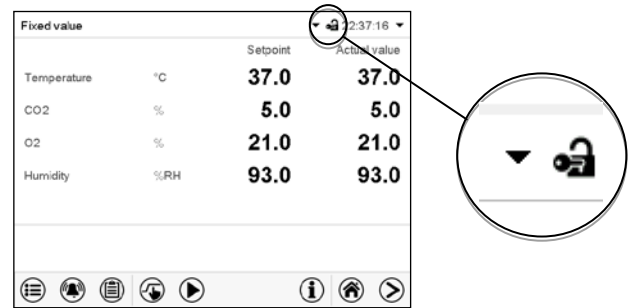
If passwords have been assigned for all authorization levels, access to the controller functions is locked without login.



## Operation after user login

At user login, the authorization level is selected and confirmed by entering the respective password.


Following user login, controller operation is available, recognizable by the open-lock icon in the header. The available controller functions correspond to the user's authorization level.



## Password protection activated for all levels: operation without user login is locked

If passwords have been assigned for all authorization levels, the controller is locked without registration of a user.

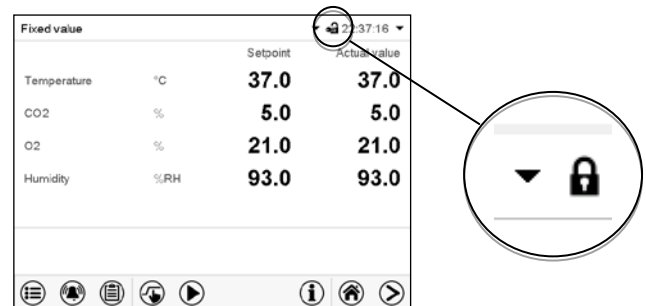As long as no user is registered, controller operation is locked, recognizable at the closed-lock icon in the header. This requires that the user management has been activated by the assignment of passwords for the individual authorization levels.



## Password protection for at least one level deactivated: operation without user login is possible

If passwords have not been assigned for all authorization levels, after turning on the chamber there are those controller functions available, which correspond to the highest authorization level without password protection.

No lock icon is shown in the display header.

User login is neither required nor possible.

To activate the password protection and user login, perform new password assignment (chap. 13.5.3).



## Information window

To check the authorization level of the user currently logged-in, select in Normal display the arrow far right in the display header.

Fixed value		14:19:52	
	Setpoint	Actual value	
Temperature	°C	40.0	40.0

The information window shows date and time, the controller's free memory space and under "Authorization" the authorization level of the current user.

If passwords have been assigned for all authorization levels, a user without login (password entry) has no authorization. There are only viewing functions available.

Fixed value		14:32:10	
Tuesday, 2018/05/24			
Authorization:		Free storage: 98%	

Display when all authorization levels are password protected and no user has logged in:

No authorization level is displayed.

If passwords have been assigned only for some of the authorization levels, a user without login (password entry) has access to the functions of the highest authorization level without password protection.

Fixed value		14:29:26	
Tuesday, 2018/05/24			
Authorization: Admin		Free storage: 98%	

Display when only some of the authorization levels are password protected (example: no protection for the "User" and "Admin" levels) and no user has logged in:

The user's effective authorization (due to lack of password protection) is shown.

Example: user with "Admin" authorization.

If passwords have been assigned for some or all of the authorization levels, user login (password entry) provides the authorization for the corresponding password-protected level.

Fixed value		14:29:26	
Tuesday, 2018/05/24			
Authorization: Admin		Free storage: 98%	

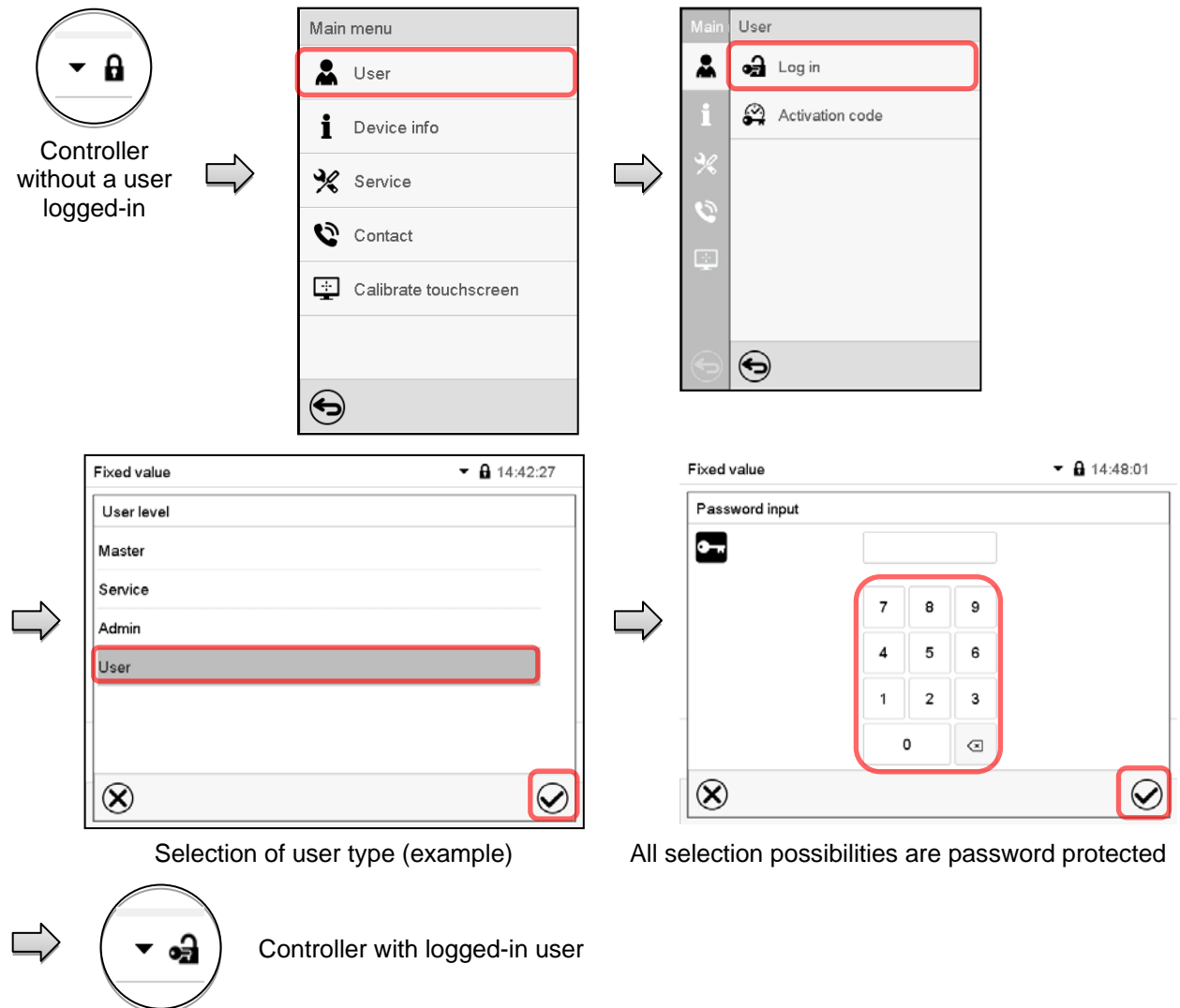
Display when at least some of the authorization levels are password protected and a user has logged in.

The user's authorization (by password entry) is shown.


Example: user with "Admin" authorization.

## 13.2 Log in

Path: **Main menu > User > Log in**



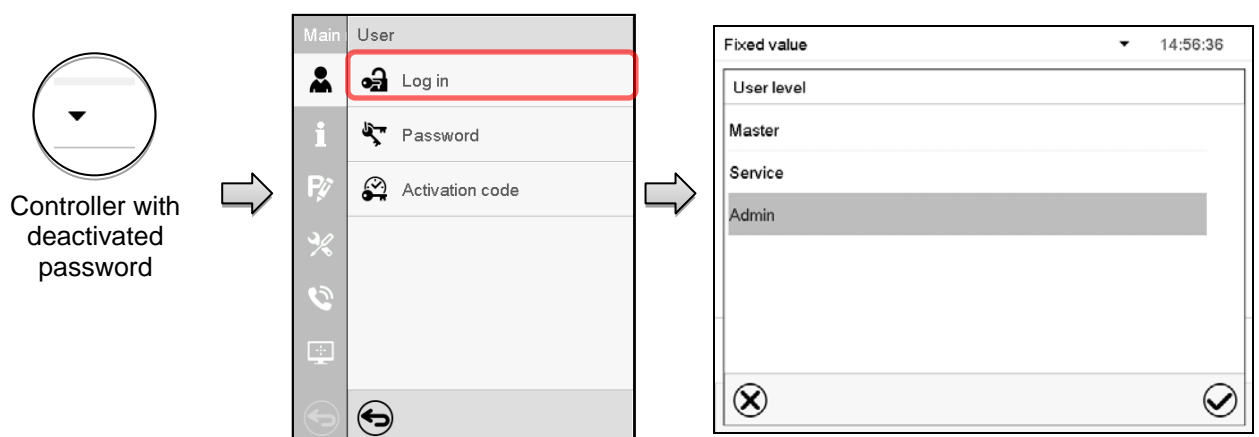
After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.




### 13.3 Log out

Path: [Main menu > User > Log out](#)

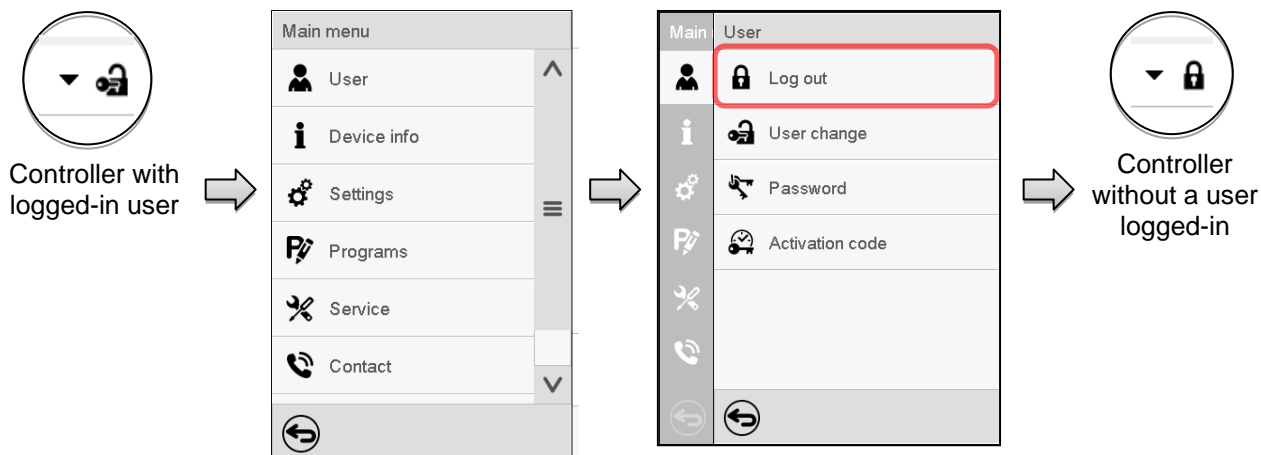User logoff with “Admin” authorization




User logoff with “Admin” authorization

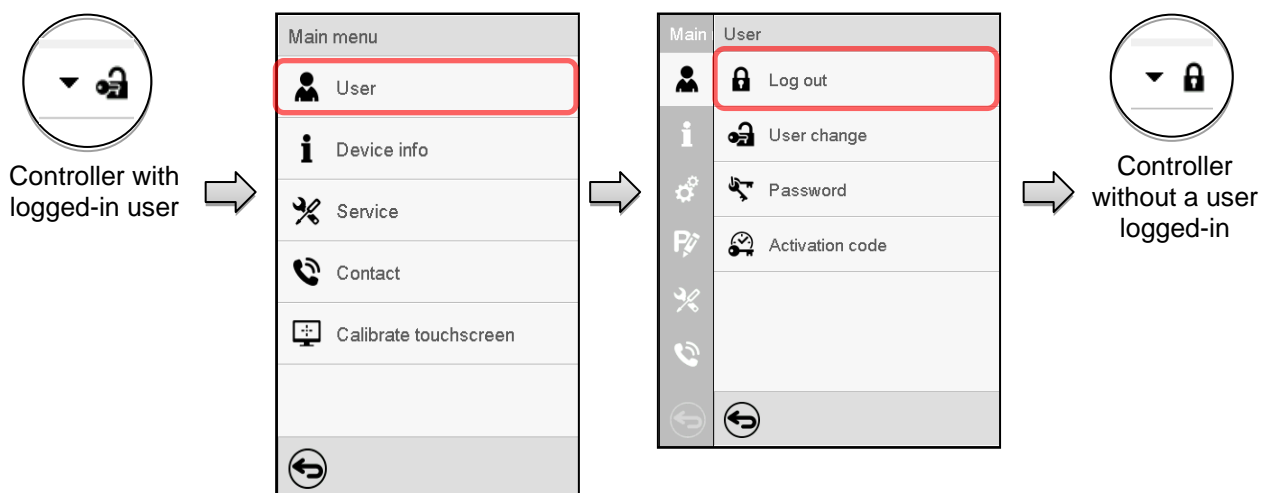


### 13.4 User change

If the password function has been deactivated (chap.13.5.2) this function is not available.

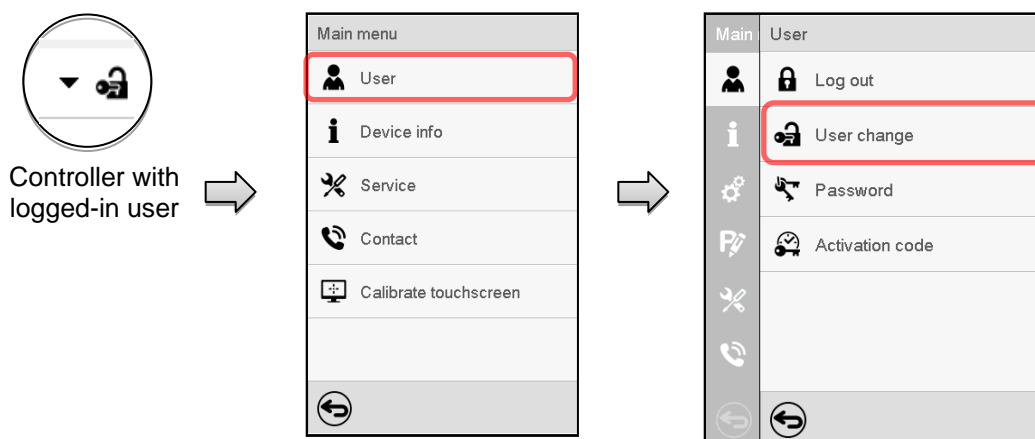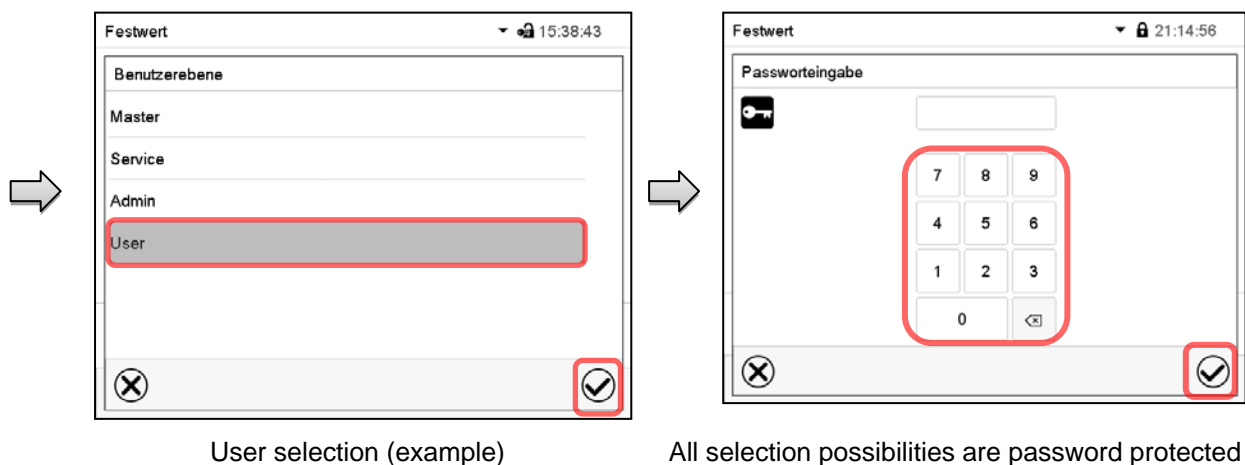Path: [Main menu > User > User change](#)





## 13.5 Password assignment and password change

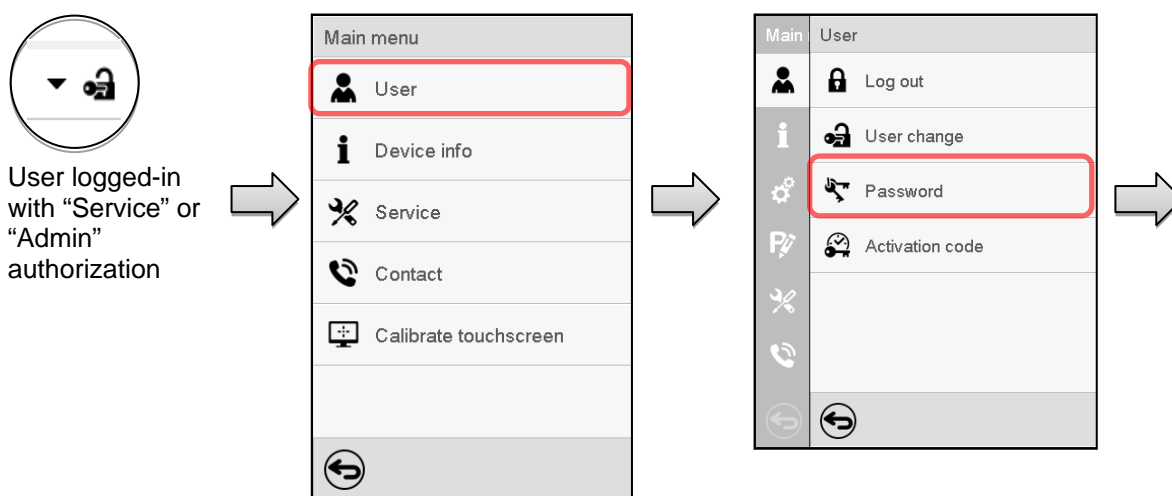
This function is not available for a user logged-in with "User" authorization.

### 13.5.1 Password change

A logged-in user can change the passwords of his current level and of the next lower level(s).


**Example:** A user with "Admin" authorization can change the passwords for the "Admin" and "User" authorization levels.

Path: [Main menu](#) > [User](#) > [Password](#)






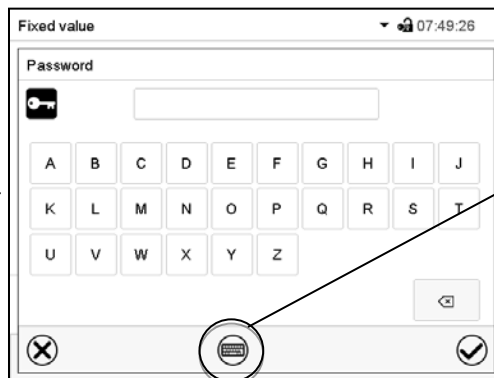
Selection of the authorization level  
(example: view with "Admin" authorization)



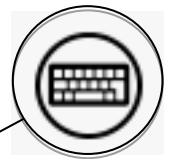
Enter desired password. If desired, press the **Change keyboard** icon to access other entry windows.



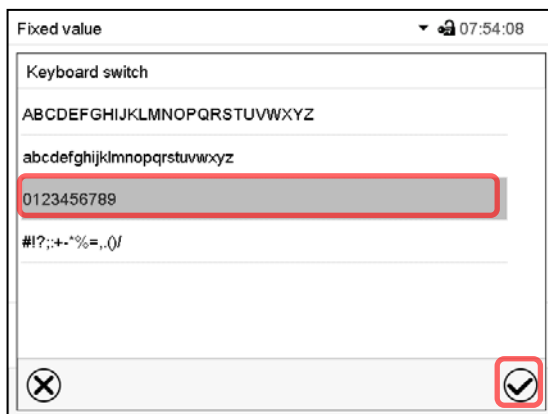In the "Keyboard switch" window you can select different keyboards to enter uppercase and lowercase letters, digits, and special characters. All types of characters can be combined within one single password.




Example: access the digit entry window

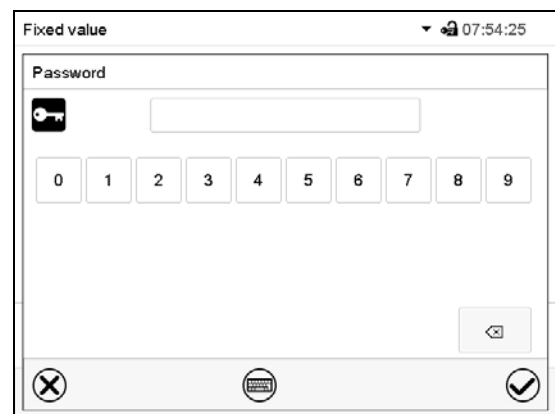

Entry of digits

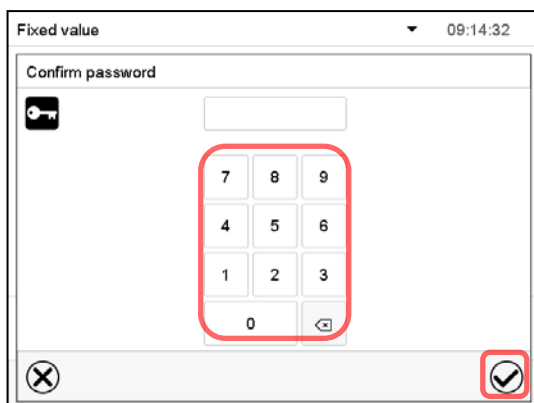To confirm the entry, press the **Confirm** icon.





Repeat the password entry for confirmation (sample picture). For each character of the password, the required keyboard appears automatically. Then press the **Confirm** icon.

### 13.5.2 Deleting the password for an individual authorization level

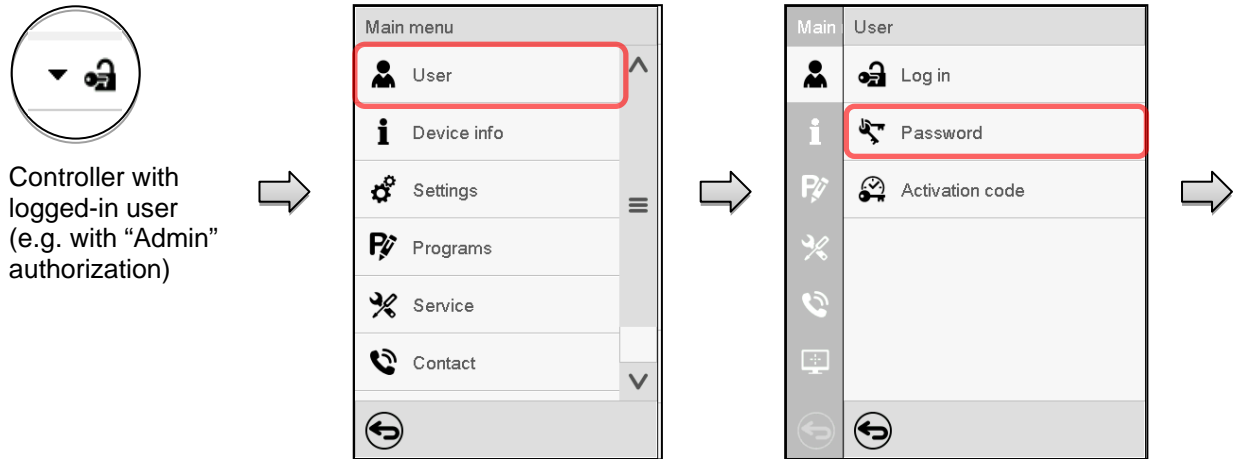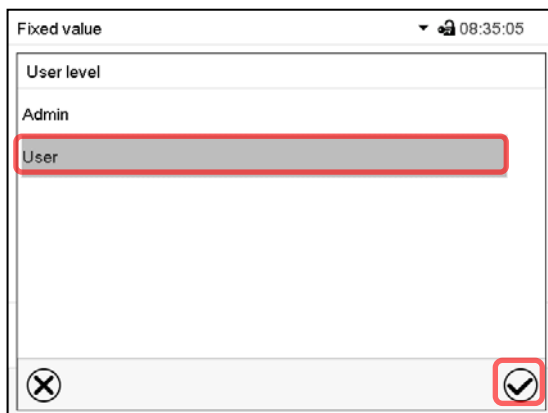A user logged-in with “Service” or “Admin” authorization can delete the passwords of his current level and of the next lower level(s). To do this no password is entered during a password change.

Path: [Main menu](#) > [User](#) > [Password](#)

Fixed value 08:35:05

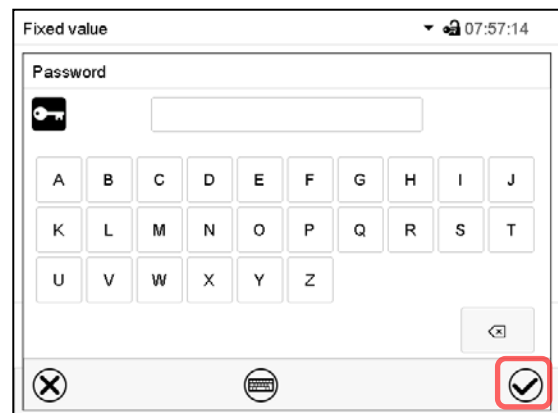
User level

Admin

User

Navigation icons: back, cancel (X), confirm (checkmark).

Select the authorization level for which the password shall be deleted.





Fixed value 07:57:14

Password

Navigation icons: back, cancel (X), confirm (checkmark).

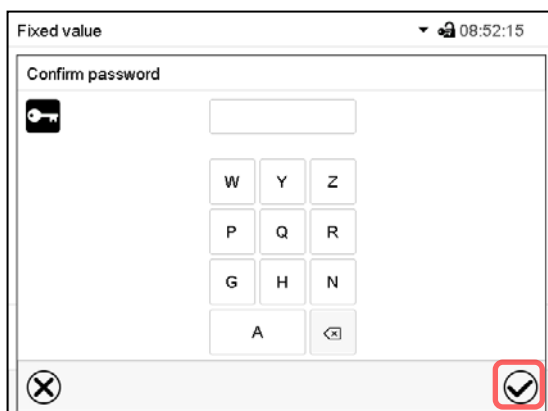Do NOT enter anything in the “Password” screen. Press the **Confirm** icon.

Fixed value 08:52:15

Confirm password

Navigation icons: back, cancel (X), confirm (checkmark).

Do NOT enter anything in the “Confirm password” screen. Press the **Confirm** icon.

The password is deleted.


### 13.5.3 New password assignment for “service” or “admin” authorization level when the password function was deactivated

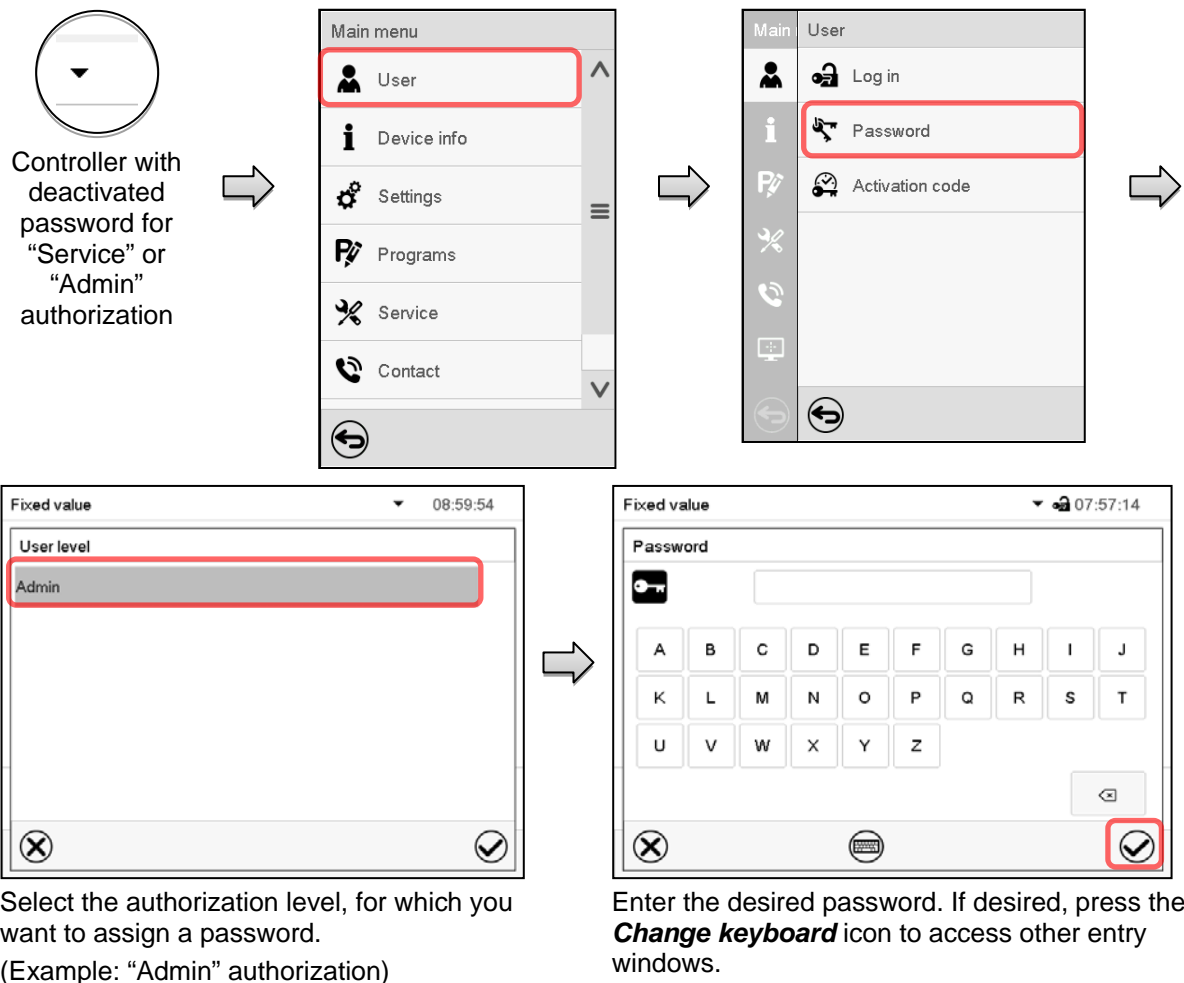If the password protection for an authorization level has been deactivated, i.e., no password is assigned, no login for this level is possible. Therefore access to this authorization level is available without login.

If the password for the “Service” or “Admin” authorization has been deleted (chap. 13.5.2), a new password can be assigned for the current level and the next lower level(s) without user login.

**Example:** The password for the “Admin” authorization level was deleted, therefore every user without login has full access to the functions of the “Admin” authorization level. If access to this level shall become password protected again, the user can assign a new password for the “Admin” authorization level with the “Password” function.

Path: [Main menu](#) > [User](#) > [Password](#)



To confirm the entry, press the **Confirm** icon.

Repeat the password entry for confirmation. For each character of the password, the required keyboard appears automatically. Then press the **Confirm** icon.




## 13.6 Activation code

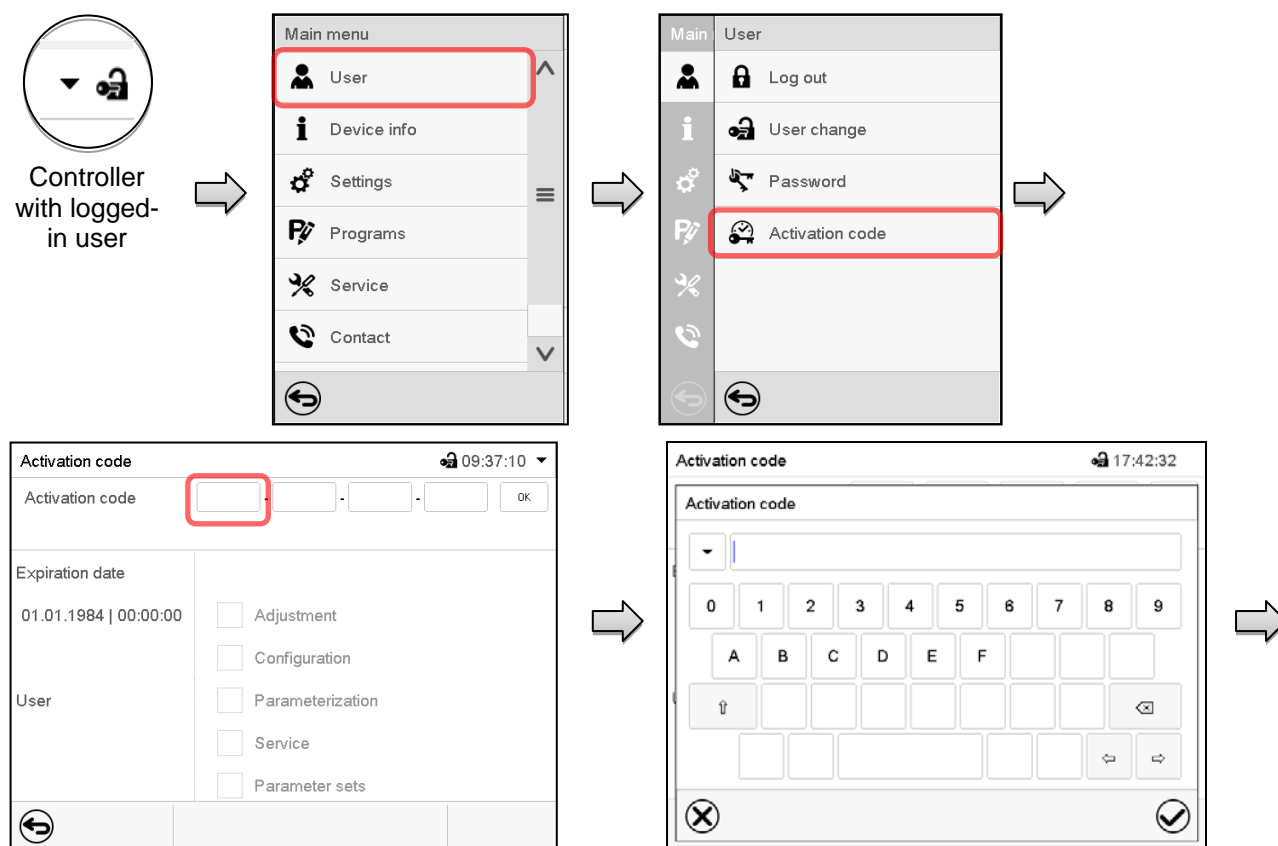Certain functions of the controller can be unlocked with a previously generated activation code.

The activation code enables access to functions available only in the “Service” authorization level by users without a “Service” authorization. Such functions include e.g., adjustment or extended configurations.

The activation code is available in authorization levels.

Path: **Main menu > User > Activation code**




Activation code menu.

Select the first of the four entry fields.

Activation code entry window.

Enter the first four characters of the activation code and press the **Confirm** icon.

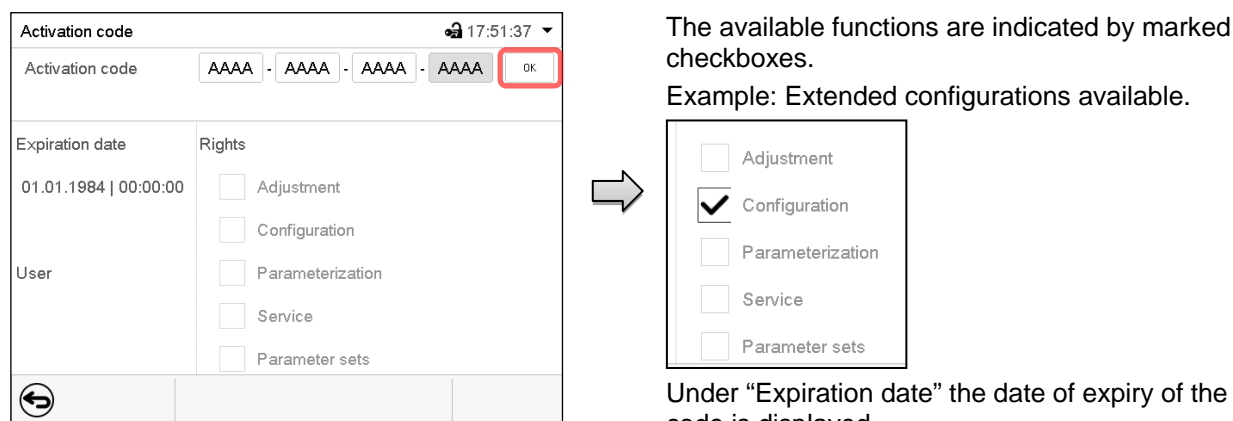Select the next of the four entry fields and proceed accordingly until the entire code has been entered.



“Activation code” menu with entered code (sample view).

Press **OK** to take over the entry


## 14. General controller settings

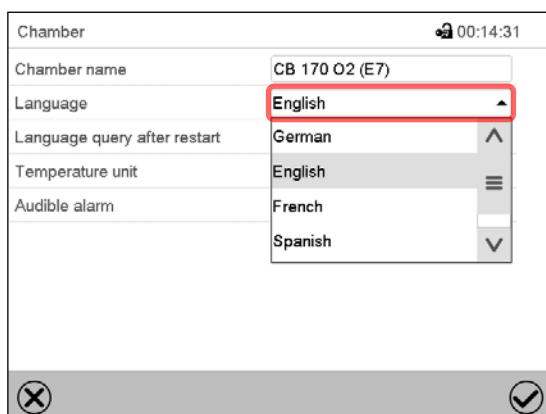Most of the general settings can be accessed in the “Settings” submenu, which is available for users with “Service” or “Admin” authorization level. It serves to enter date and time, select the language for the controller menus and the desired temperature unit and to configure the controller’s communication functions.

### 14.1 Selecting the controller’s menu language

The MB2 program controller communicates by a menu guide using real words in German, English, French, Spanish, and Italian.

Path: [Main menu](#) > [Settings](#) > [Chamber](#)




The screenshot shows the 'Chamber' submenu with the following settings:

Chamber name	CB 170 O2 (E7)
Language	English
Language query after restart	German
Temperature unit	English
Audible alarm	French
	Spanish

The 'Language' dropdown menu is highlighted with a red box, showing the selection of 'English'.

“Chamber” submenu.

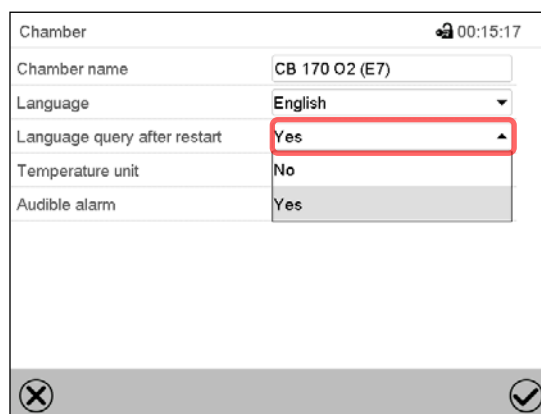Select the desired language.



The screenshot shows the 'Chamber' submenu with the following settings:

Chamber name	CB 170 O2 (E7)
Language	English
Language query after restart	Yes
Temperature unit	No
Audible alarm	Yes

The 'Language query after restart' dropdown menu is highlighted with a red box, showing the selection of 'Yes'.


“Chamber” submenu.

Select if there shall be a language query after restarting the chamber and press the **Confirm** icon.

Return to Normal display with the **Back** icon to take over the entries.

### 14.2 Setting date and time

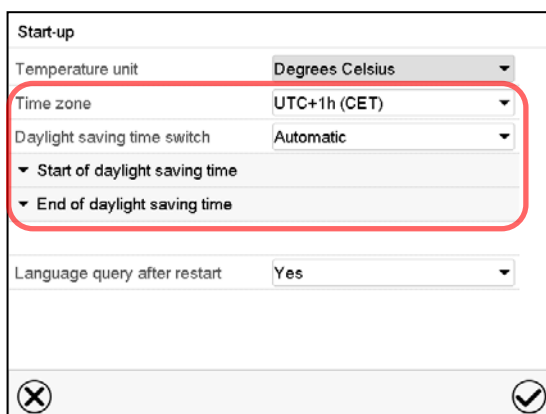Following start-up of the chamber after language selection:



The screenshot shows the 'Start-up' submenu with the following settings:


Temperature unit	Degrees Celsius
Time zone	UTC+1h (CET)
Daylight saving time switch	Automatic
Start of daylight saving time	
End of daylight saving time	
Language query after restart	Yes

The 'Time zone' and 'Daylight saving time switch' dropdown menus are highlighted with a red box, showing the selection of 'UTC+1h (CET)' and 'Automatic' respectively.


Select the time zone and configure the daylight saving time switch.

Or later:

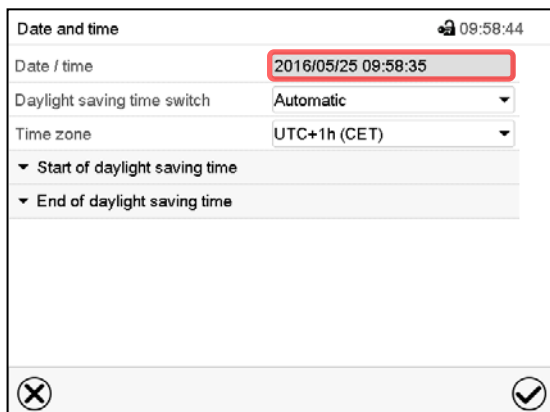
Path: [Main menu](#) > [Settings](#) > [Date and time](#)



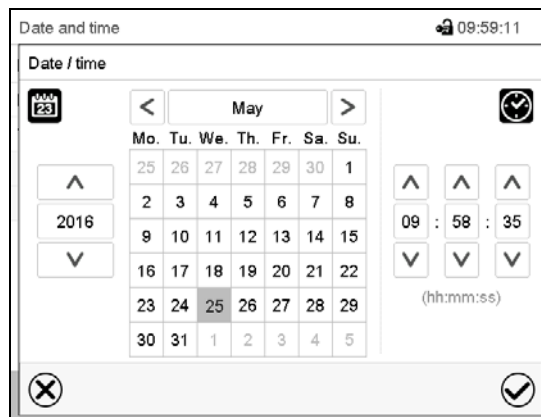
“Date and time” submenu.  
Select the field “Date / time”.

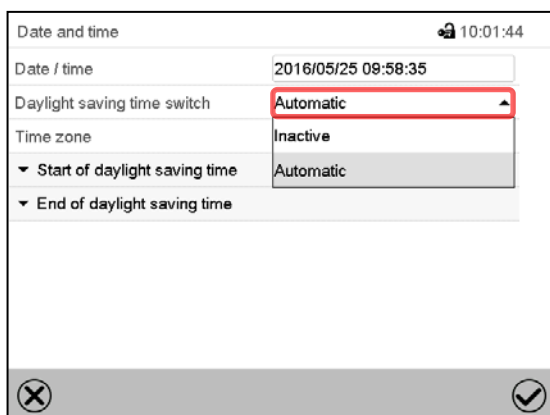
“Date / time” entry menu.  
Enter date and time and press the **Confirm** icon.




“Date and time” submenu.  
In the field “Daylight saving time switch” select the desired setting “Automatic” or “Inactive”.

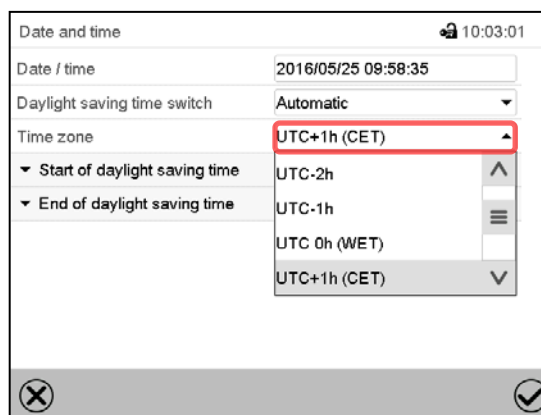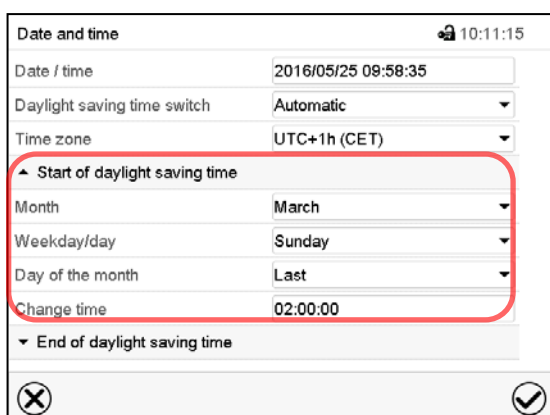“Date and time” submenu.  
Select the desired time zone and press the **Confirm** icon.



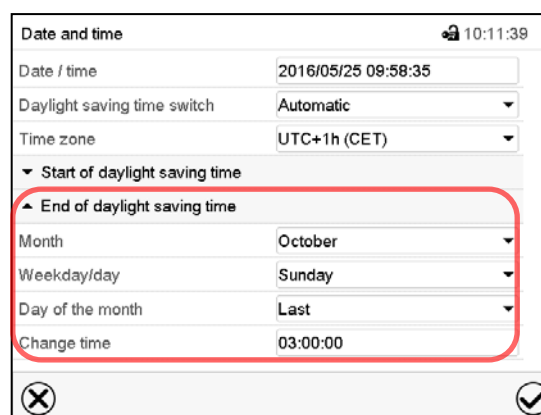“Date and time” submenu.  
Select the desired start of the daylight saving time.





“Date and time” submenu.  
Select the desired end of the daylight saving time and press the **Confirm** icon.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

## 14.3 Selecting the temperature unit

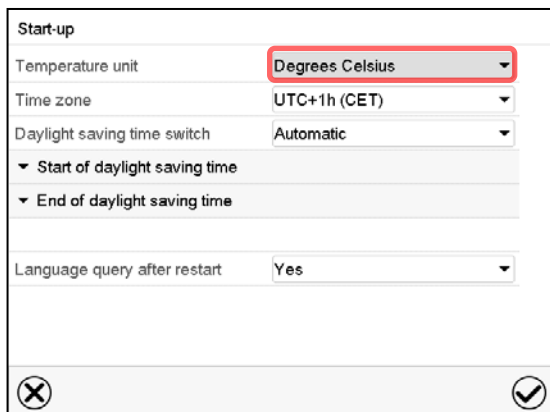Following start-up of the chamber:



Start-up

Temperature unit: Degrees Celsius

Time zone: UTC+1h (CET)

Daylight saving time switch: Automatic

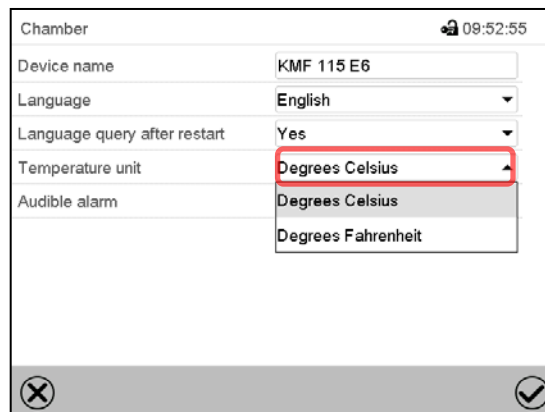
Start of daylight saving time

End of daylight saving time

Language query after restart: Yes

Or later:

Path: [Main menu](#) > [Settings](#) > [Chamber](#)



Chamber 09:52:55

Device name: KMF 115 E6

Language: English

Language query after restart: Yes


Temperature unit: Degrees Celsius

Audible alarm: Degrees Celsius

Select the desired temperature unit and press the **Confirm** icon.

Change of the temperature unit between °C and °F.

If the unit is changed, all values are converted accordingly


	C = degree Celsius	0 °C = 32°F	Conversion: [value in °F] = [value in °C] * 1,8 + 32
	F = degree Fahrenheit	100 °C = 212°F	

## 14.4 Display configuration
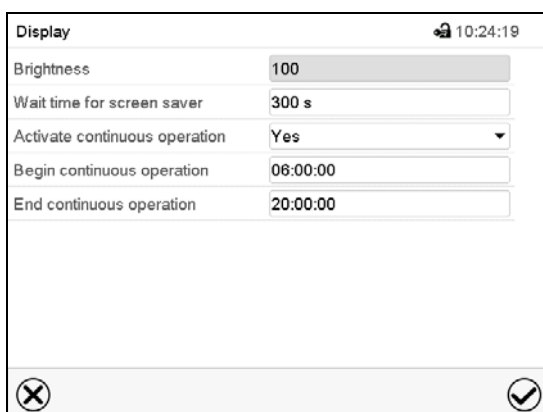
### 14.4.1 Adapting the display parameters

This function serves to configure parameters like display brightness and operating times.

Path: [Main menu](#) > [Settings](#) > [Display](#) > [Display](#)



Display 10:24:19

Brightness: 100

Wait time for screen saver: 300 s

Activate continuous operation: Yes

Begin continuous operation: 06:00:00

End continuous operation: 20:00:00

“Display” submenu.

- Select the field “Brightness”.

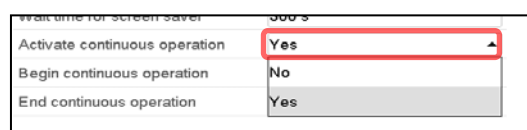
Move the grey slide to the left or right to define the brightness of the display

- left = darker (minimum value: 0)
- right = brighter (maximum value: 100)

Press the **Confirm** icon.



- Select the field “Wait time for screen saver” and enter the desired waiting time for the screen saver in seconds. Setting range: 10 sec up to 32767 sec. During the waiting time the display is off. Confirm entry with **Confirm** icon.
- In the field “Activate continuous operation” select the desired setting “Yes” or “No”.




- Select the field “Begin continuous operation” (possible only if continuous operation is activated) and enter the time with the arrow keys. Confirm entry with **Confirm** icon.
- Select the field “End continuous operation. (only possible if continuous operation is activated) and enter the time with the arrow keys. Confirm entry with **Confirm** icon.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.


#### 14.4.2 Touchscreen calibration

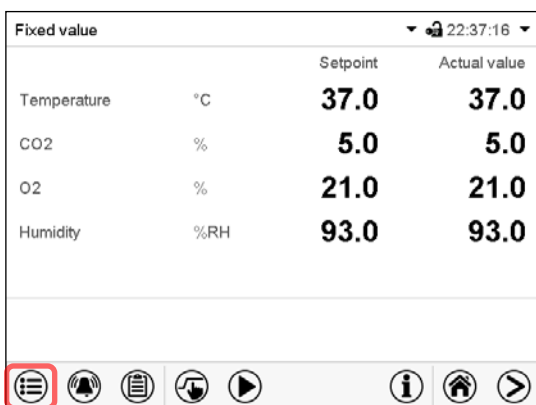This function serves to optimize the display for the user's individual angular perspective.

Path: [Main menu](#) > [Calibrate touchscreen](#)




Normal display.

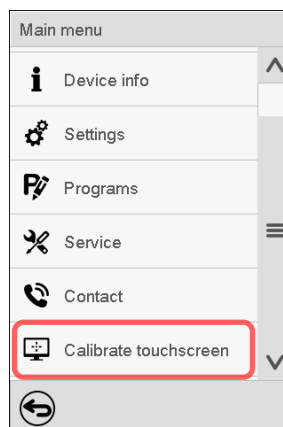

Select “Calibrate touchscreen” and follow the instructions on the display.

You need to touch all four corners of the touchscreen to calibrate it. Appropriate boxes appear successively in each corner.

	<p>The waiting icon shows how much time there is left to touch the currently activated box. If the box is not touched within this period, calibration is aborted and the display changes to Normal display.</p>
---	---

After completing the calibration, i.e., touching all four boxes, the display changes to Normal display.

## 14.5 Network and communication


For these settings at least the “Admin” authorization level is required.

### 14.5.1 Serial RS485 interface (available via BINDER INDIVIDUAL Customized Solutions)

This menu allows to configure the communication parameters of the RS485 interface.

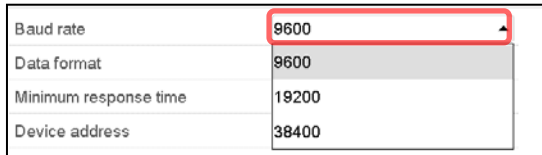
The device address is required to recognize chambers with this interface type in a network, e.g. when connecting it to the optional communication software BINDER APT-COM™ 3 DataControlSystem (chap. 18.1). In this case do not change the other parameters.

Path: [Main menu](#) > [Settings](#) > [Serial interfaces](#)




“Serial interfaces” submenu.

- Select the desired setting in the field “Baud rate”.



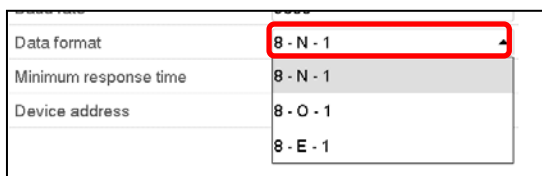- Select the desired setting in the field “Data format”.




- Select the field “Minimum response time” and enter the desired minimum response time. Confirm entry with **Confirm** icon.
- Select the field “Device address” and enter the device address. Factory setting is “1”. Confirm entry with **Confirm** icon.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

## 14.5.2 Ethernet
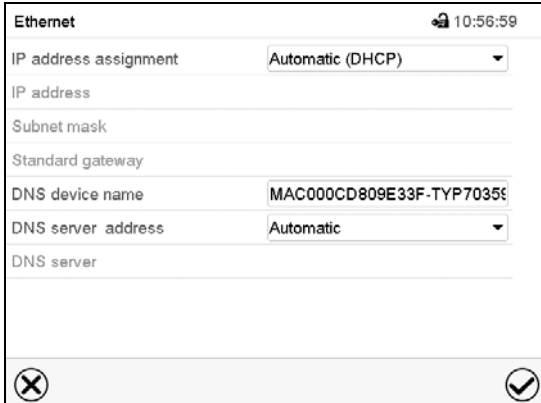
### 14.5.2.1 Configuration



Path: [Main menu](#) > [Settings](#) > [Ethernet](#)



“Ethernet” submenu.



- In the field “IP address assignment” select the desired setting “Automatic (DHCP)” or “Manual”.

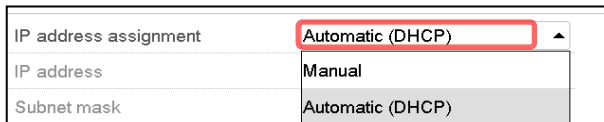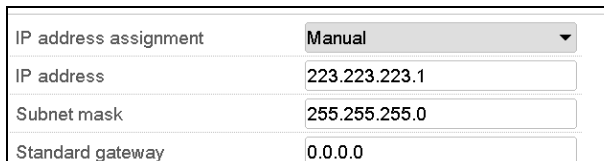With selection “Manual” you can enter the IP-address, the subnet mask and the standard gateway manually.

- Select “DNS device name” and enter the DNS device name. Confirm entry with **Confirm** icon.
- In the field “DNS server address” select the desired setting “Automatic” or “Manual”.


With selection „Manual“ you can enter the DNS server address manually.

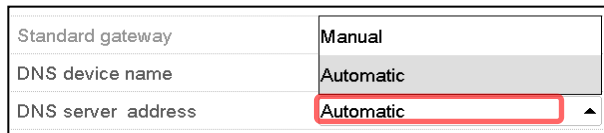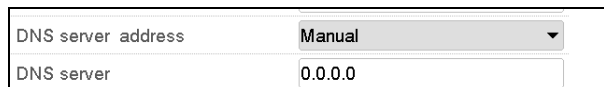After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

### 14.5.2.2 Display of MAC address

Path: [Main menu](#) > [Device info](#) > [Ethernet](#)




Ethernet	
Ethernet	Yes
MAC address	00-0C-D8-09-E3-3F
IP address	192.168.14.87
Subnet mask	255.255.255.0
Standard gateway	192.168.14.1
DNS server	192.168.10.5
DNS device name	MAC000CD809E33F-TYP703596

“Ethernet” submenu (example).

### 14.5.3 Web server

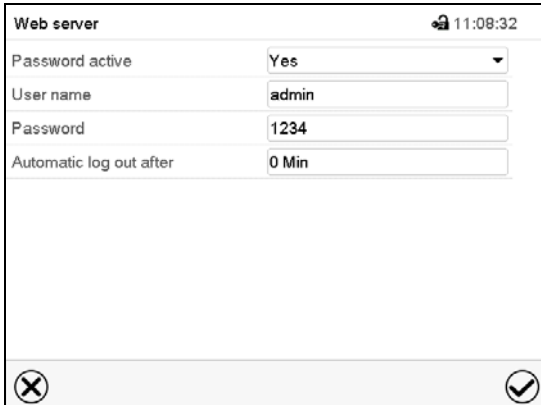This controller menu serves to configure the web server. Then you can enter the chamber's IP-address in the Internet. The IP address is available via [Chamber information](#) > [Ethernet](#). The BINDER web server opens. Enter the user name and password which have been assigned for the web server in the controller menu. This enables online access to the controller display, to see e.g., the event list or error messages. In this view no settings can be changed.


Path: [Main menu](#) > [Settings](#) > [Web server](#)



Web server	
Password active	Yes
User name	admin
Password	1234
Automatic log out after	0 Min

“Webserver” submenu.

- In the field “Password active” select the desired setting “Yes” or “No”.



Password active	Yes
User name	No
Password	Yes

- Select the field “User name” and enter the desired user name. Confirm entry with **Confirm** icon.
- Select the field “Password” and enter the desired password. Confirm entry with the **Confirm** icon.
- Select the field “Automatic log out after” and enter the time in minutes after which the webserver shall log out automatically. Setting range: 0 min to 65535 min. Confirm entry with **Confirm** icon.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.




## 14.5.4 E-Mail

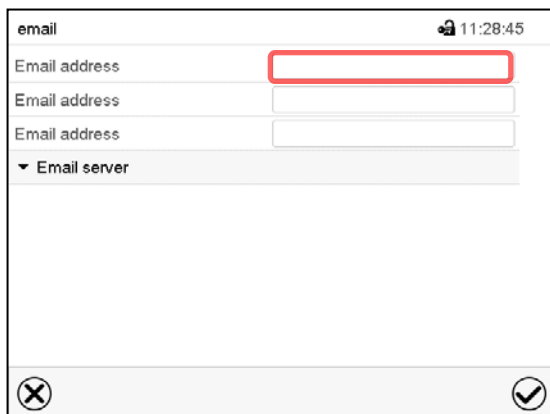As soon as an alarm was triggered, an e-mail is sent to the configured e-mail address.

Path: [Main menu](#) > [Settings](#) > [Email](#)


### E-mail address entry:



“Email” submenu.

Select the desired e-mail address field and enter the e-mail address. You can use the **Keyboard change** icon for entry. Confirm entry with **Confirm** icon.

### E-mail server settings:

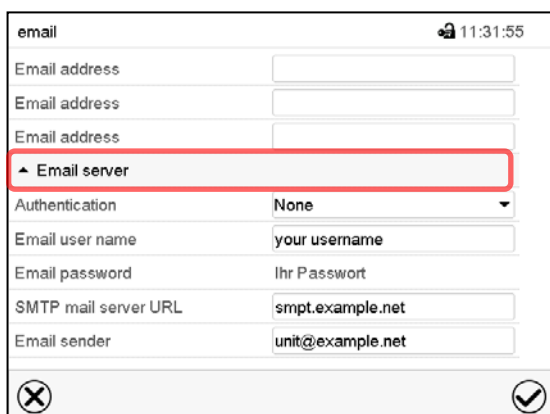

“Email” submenu.

Select the field “Email server” to access the settings

- In the field “Authentication” select the desired setting “None” or “SMTP” auth“.

With the setting “SMTP auth”, you can enter a password under “Email password”.

Authentication	None
Email user name	None
Email password	SMTP auth
SMTP mail server URL	192.168.10.45

- Select the field “Email user name” and enter the desired user name. Confirm entry with **Confirm** icon.
- Select the field “SMTP mail server URL” and enter the SMPT mail server URL. Confirm entry with **Confirm** icon.
- Select the field “Email sender” and enter the desired Email sender. Confirm entry with **Confirm** icon.


After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

## 14.6 USB menu: Data transfer via USB interface

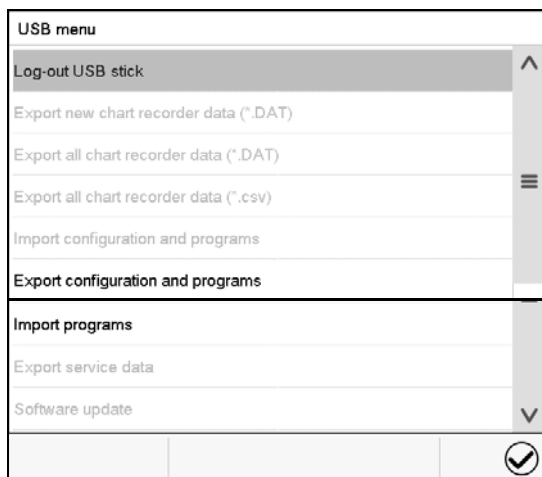
The USB port is located in the instrument box.

When you insert a USB-stick, the “USB” menu opens.

Depending on the user’s authorization level, different functions (highlighted in black) are available for the logged-in user.



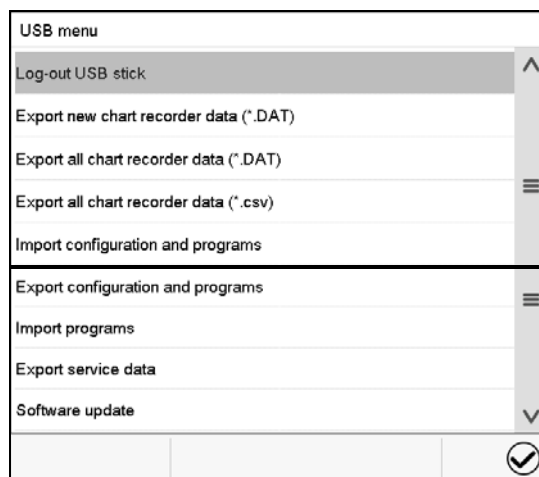Available functions  
with “User” authorization level




Available functions  
with “Admin” authorization level

Function	Explanation
Log-out USB stick	Log-out USB stick bevor pulling it
Export new chart recorder data (*.DAT)	Export chart recorder data, which have been added since last export, in .dat format
Export all chart recorder data (*.DAT)	Export all chart recorder data in .dat format
Export all chart recorder data (*.csv)	Export all chart recorder data in .csv format
Import configuration and programs	Import configuration and timer / time / week programs
Export configuration and programs	Export configuration and timer / time / week programs
Import programs	Import timer / time / week programs
Export service data	Export service data
Software update	Controller firmware update

## 15. General information
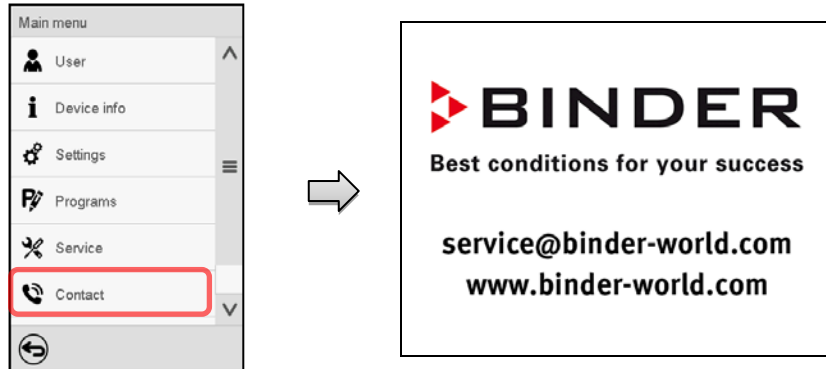
### 15.1 Service contact page


Path: [Main menu](#) > [Contact](#)



### 15.2 Current operating parameters



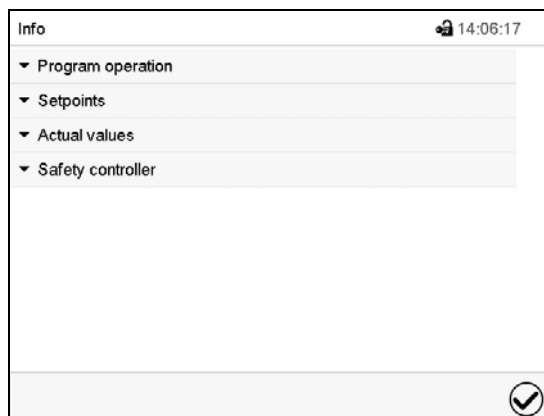Press the **Information** icon to access the “Info” menu from Normal display.




“Info” menu.  
Select the desired information.




- Select “Program operation” to see information on a currently running program.
- Select “Setpoints” to see information on the entered setpoints and operation lines.
- Select “Actual values” to see information on the current actual values.
- Select “Safety controller” to see information on the safety controller status.

## 15.3 Event list


The “Event list” displays status information and errors of the current day. It enables to view the last 100 events or defective conditions of the chamber.




Press the **Event list** icon to access the event list from Normal display.

Event list			13:18:52
2016/06/07	09:09:53	Login Service (Touch)	^
2016/06/07	09:09:53	Automatic log out Admin	
2016/06/07	07:47:25	Login Admin (Touch)	
2016/06/07	07:46:15	Automatic log out Admin	
	2016/06/07	07:46:15	Power on
	2016/06/06	16:08:09	Power off
	2016/06/06	10:50:25	Login Admin (Touch)
	2016/06/06	10:49:44	Automatic log out Admin
	2016/06/06	10:49:44	Power on
			v

Event list
















Press the **Update** icon to update the event list.



**Attention:** Following a modification of the language setting (chap. 14.1) or the storage interval of the chart recorder (chap. 16.2) the Event list is cleared.

## 15.4 Technical chamber information


Path: [Main menu](#) > [Device info](#)

Main	Device info		
	 General	Chamber name and setup	
	 v1.x Versions	Versions of CPU, I/O module and safety controller	for BINDER Service
	 In-/Outputs	Information on digital and analog inputs and outputs and phase angle outputs	for BINDER Service
	 Modbus inputs	Information on modbus analog and digital inputs	for BINDER Service
	 Ethernet	Information on Ethernet connection, MAC address display	chap. 14.5.2
			
		Back to main menu	



## 16. Chart recorder display

This view offers graphic representation of the measurement course. Data representation imitates a chart recorder and allows recalling any set of measured data at any point of time taken from the recorded period.

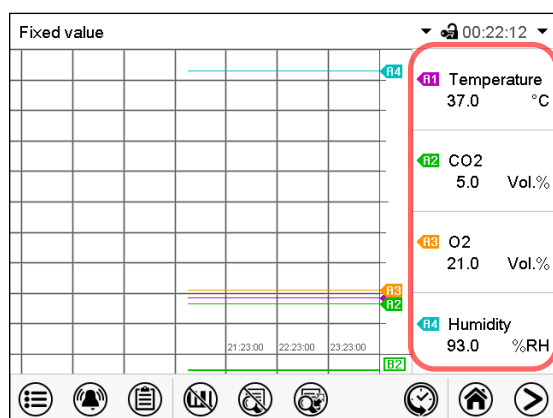
### 16.1 Views

	Press the <b>Change view</b> icon to access the pen recorder display.
---	---

#### 16.1.1 Show and hide legend


	<b>Show legend</b>		<b>Hide legend</b>
---	--------------------	---	--------------------

Press the **Show legend** icon to display the legend on the right side of the display




Legend shown on the right side of the display

#### 16.1.2 Switch between legend pages



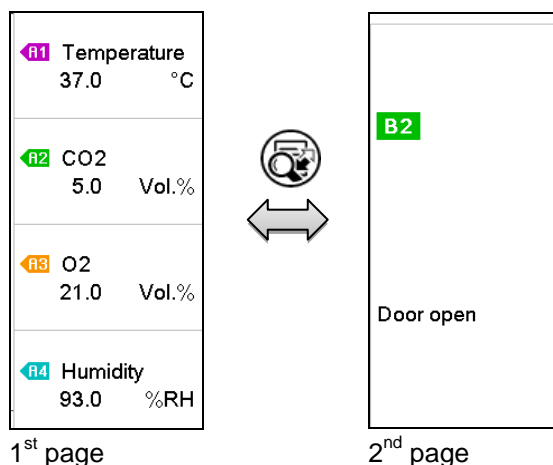	<b>Switch legend</b>
---	----------------------

Press the **Switch legend** icon to switch between the legend pages




Switching between the legend pages

### 16.1.3 Show and hide specific indications

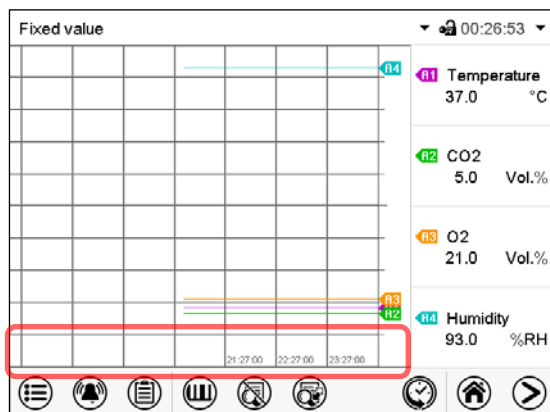
	<b>Show indications</b>		<b>Hide indications</b>
---	-------------------------	---	-------------------------

Press the **Show indications** icon to display the indication "Door open" (B2).




Indication "Door open" displayed.

### 16.1.4 History display

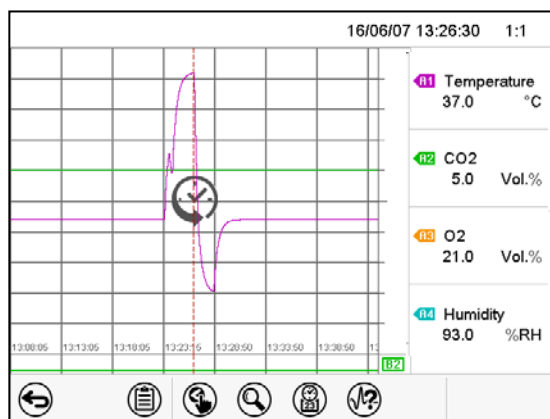	<b>History display</b>
---	------------------------

Press the **History display** icon to change to the history display.



History display.

The chart recorder is paused. Data recording continues in the background.

Move the central red line by tapping and holding to the desired position.


The legend at the right side shows the values of the current line position.

Then further icons appear:

## History display: Curve selection



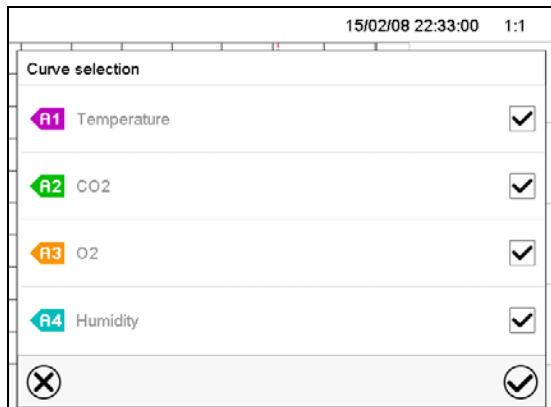
Press the **Curve selection** icon to access the “Curve selection” submenu.




“Curve selection” submenu.

Select the curves to be displayed by checking the checkbox of the corresponding parameter. Press the **Confirm** icon

## History display: Search the required instant

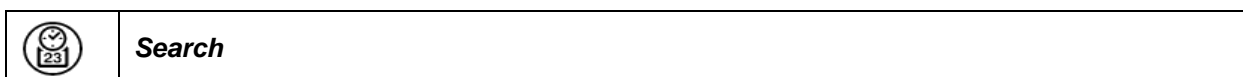

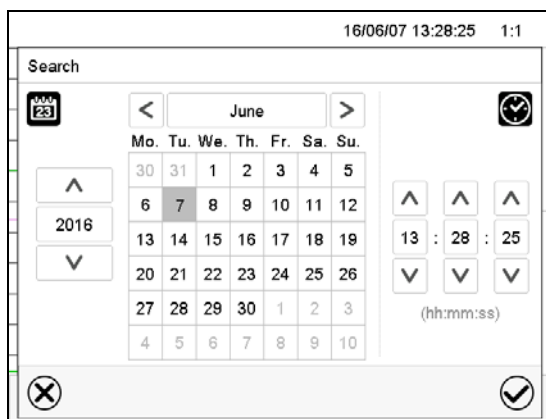Press the **Search** icon to access the “Search” submenu.




“Search” submenu.

Select the required instant by entering its date and time and press the **Confirm** icon

## History display: Zoom function



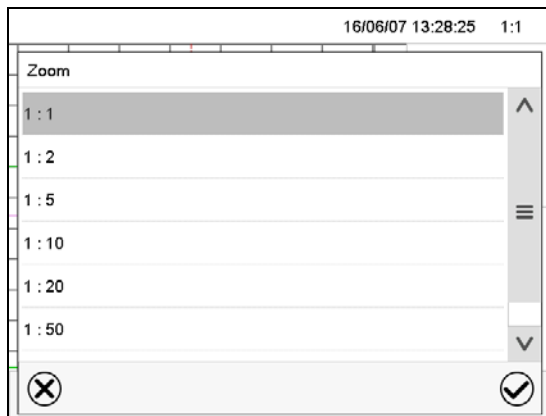Press the **Zoom** icon to access the “Zoom” submenu.




“Zoom” submenu.

Select the zoom factor and press the **Confirm** icon

## History display: Show and hide scroll buttons to scroll to an instant




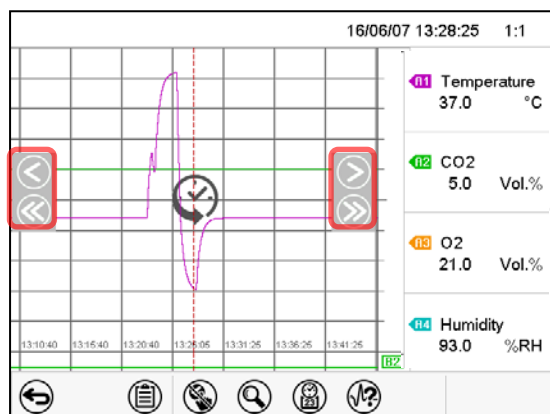Press the **Show scroll buttons** icon to access the “Page selection” submenu.



“Page selection” submenu.

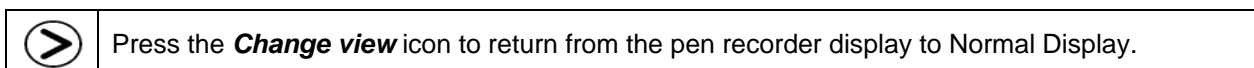Scroll buttons are shown on the left and on the right.  
Use them to move along the timeline.






## 16.2 Setting the parameters

This menu allows setting the storage interval, the type of values to be shown and the scaling of the temperature and humidity charts.

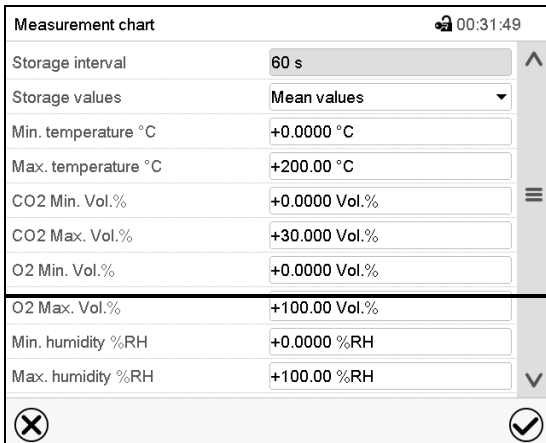Path: **Main menu > Settings > Measurement chart**




Measurement chart	
Storage interval	60 s
Storage values	Mean values
Min. temperature °C	+0.0000 °C
Max. temperature °C	+200.00 °C
CO2 Min. Vol.%	+0.0000 Vol.%
CO2 Max. Vol.%	+30.000 Vol.%
O2 Min. Vol.%	+0.0000 Vol.%
O2 Max. Vol.%	+100.00 Vol.%
Min. humidity %RH	+0.0000 %RH
Max. humidity %RH	+100.00 %RH

“Measurement chart” submenu.

- Select the field “Storage interval” and enter the desired storage interval. Confirm entry with **Confirm** icon.

The available presentation depends on the pre-selected storage rate. Factory setting: 60 seconds. This means the higher the storage rate, the more precisely but shorter the data representation will be.

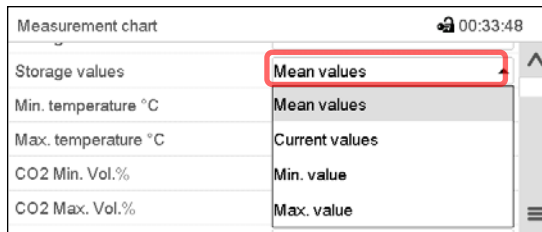- In the field “Storage values” select the desired value type to be displayed.



Measurement chart	
Storage values	Mean values
Min. temperature °C	Mean values
Max. temperature °C	Current values
CO2 Min. Vol.%	Min. value
CO2 Max. Vol.%	Max. value

- For scaling the representation select the desired minimum and maximum temperature or humidity value and enter the desired values. Confirm each entry with **Confirm** icon.

### Display ranges:

- Temperature: 0 °C up to 200 °C
- CO<sub>2</sub>: 0 Vol.-% up to 100% Vol.-%.
- O<sub>2</sub>: 0 Vol.-% up to 100% Vol.-%.
- Humidity: 0% r.h. up to 100% r.h.



### CAUTION

**Setting the storage rate or rescaling (minimum and/or maximum) will clear the measured-value memory and the event list.**

#### Danger of information loss.

- Change the storage rate or scaling **ONLY** if the previously registered data is no longer needed.

After completing the settings, press the **Confirm** icon to take over the entries and exit the menu, **or** press the **Close** icon to exit the menu without taking over the entries.

## 17. Humidification system (chamber with active humidification)

The chamber is equipped with a capacitive humidity sensor. This results in a control accuracy of up to  $\pm 3\%$  r.H. of the set point.

You can turn humidity control on or off in the controller (chap. 6.8).

With humidity control turned off, the humidification module cools down. After activation it will take up to 10 minutes until the humidification function is fully available again. This setting is required when operating the chamber with the Permadry water pan in order to avoid humidity alarms.



Heat emission of electrical devices connected inside the chamber may modify the temperature and humidity range.

The chambers are equipped with a door heating system to avoid condensation in the door area.

### 17.1 Function of the humidifying system

The humidifying system is located in the humidity generation module. In a cylindrical container with a volume of approx. 0,07 liters an electrical resistance heating evaporates water. The water content is kept close to the boiling point, and thus steam can be immediately generated in sufficient quantity for rapid humidity increases or for compensation of humidity losses, e.g. by door openings.

Freshwater is supplied by manually filling a freshwater bag.



**In order to ensure accurate humidifying, observe the following points with regard to the freshwater supply:**

- Water type: deionized (demineralized) water
- To ensure humidification during 24 hours even at high humidity set-points, we recommend filling the freshwater bag at the end of each day.
- Water intake temperature NOT below  $+5\text{ }^{\circ}\text{C}$  /  $41\text{ }^{\circ}\text{F}$  and not exceeding  $40\text{ }^{\circ}\text{C}$  /  $104\text{ }^{\circ}\text{F}$ .



BINDER GmbH is NOT responsible for the water quality provided by the customer. Any problems and malfunctions that might arise following use of water of deviating quality is excluded from liability by BINDER GmbH.

The humidity system is functional only if the freshwater bag is sufficiently filled. Check the filling level daily. The water reserve is sufficient for one up to two weeks, depending on the humidity demand (set humidity set-point, number of door openings).

For error indications concerning water supply and humidity system, see chap. 11 and 24.

## 18. Options

### 18.1 Communication software APT-COM™ 3 DataControlSystem (option)

The chamber is regularly equipped with an Ethernet interface (5) that can connect the BINDER communication software APT-COM™ 3 DataControlSystem. The actual values of temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control) and humidity (chamber with active humidification) are given at adjustable intervals. Programming can be performed graphically via PC. Up to 40 chambers can be cross-linked. The MAC Address is indicated in the "Device info" controller menu (chap. 14.5.2.2). For further information on networking please refer to the operating manual of the BINDER communication software APT-COM™ 3.

### 18.2 Silicone access ports 30 mm / 1.18 in, closable from both sides with silicon plugs (option)









Figure 21: Positions of the optional silicon access ports left (a), rear (b), and right (c)

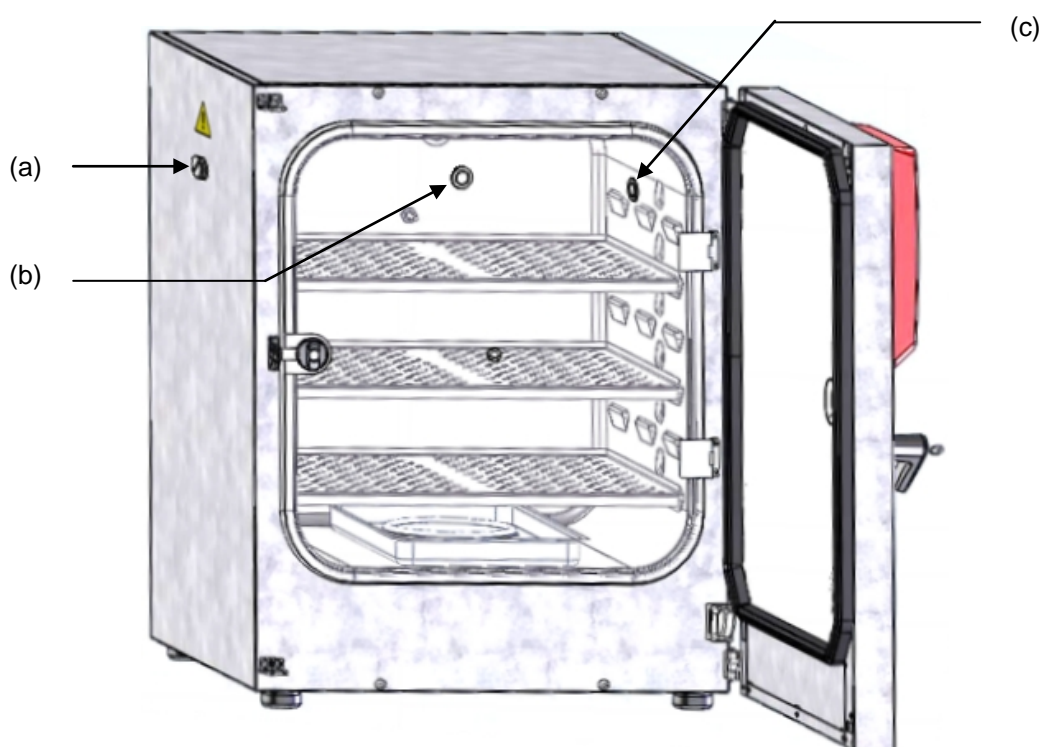A warning sticker is located above each access port on the outside of the incubator.



When operating a CB with silicon access ports, both silicon plugs must tightly close the access ports. If the plugs are inserted in a not-gastight manner, or if plugs are missing, CO<sub>2</sub> / O<sub>2</sub> / N<sub>2</sub> gas (chamber with O<sub>2</sub> control) may escape into the environment. The CO<sub>2</sub> control and the O<sub>2</sub> control (chamber with O<sub>2</sub> control) only turn off when the chamber door is opened.

	<div style="background-color: orange; text-align: center; padding: 5px;">  <b>WARNING</b> </div> <p><b>High concentration of CO<sub>2</sub> (&gt; 4 Vol.-%).</b>  <b>Risk of death by suffocation.</b>  <b>Danger of poisoning.</b></p> <p>➤ Tightly close each access port with two plugs during operation.</p>
---	---

**Note for chambers with O<sub>2</sub> control:**

 	<div style="background-color: orange; text-align: center; padding: 5px;">  <b>WARNING</b> </div> <p><b>High concentration of O<sub>2</sub> (&gt; 21 % O<sub>2</sub>).</b>  <b>Fire and explosion hazard through contact of combustible materials with O<sub>2</sub>.</b>  <b>Risk of burns and other injuries.</b></p> <p>➤ Tightly close each access port with two plugs during operation.</p>
--	--

### 18.3 Interior socket 230V (option, available via BINDER INDIVIDUAL Customized Solutions)

	<p>The interior socket is activated and deactivated in the controller via the switching state of an operation line (chap. 7.3, 9.7.3, 10.6.5). When the interior socket is activated, the notification "Interior socket" is shown highlighted in black next to the information icon in the screen header in Normal display (chap. 11.1.1)</p>	
---	---	---

The interior socket (G) is located at the upper left corner on the rear wall of the inner chamber. It is closed with a waterproof lid. The interior socket is suitable to supply electrical devices inside the chamber with 230V AC voltage.

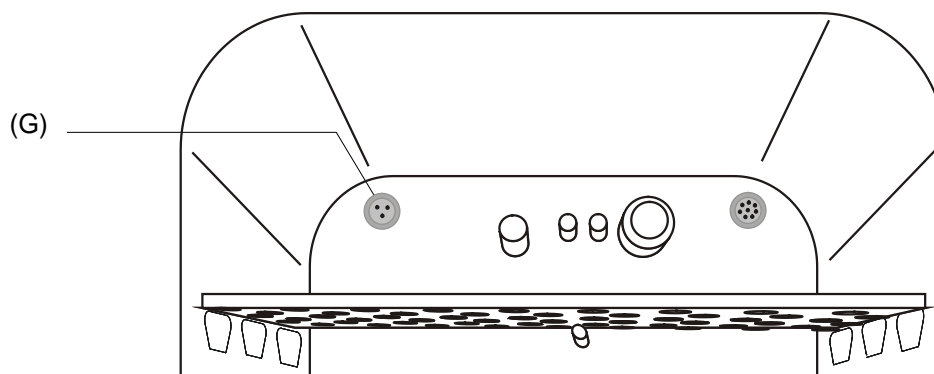




Figure 22: Position of the interior socket 230 V



(G) Interior socket

The maximum head load must not exceed a capacity of 20 W.

	<b>CAUTION</b>
	<p><b>Exceeding the setpoint temperature.</b>  <b>Damage to the cultures.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT exceed the maximum head load of 20 W.</li> <li>⊘ Do NOT connect equipment with a nominal capacity &gt; 20 W.</li> </ul>

	<p>Heat emission of electrical devices connected inside the CO<sub>2</sub> incubator may modify the temperature range.</p>
---	--

The maximum load must not exceed 3 Amp.

	 <b>WARNING</b>
	<p><b>Overload of contacts.</b>  <b>Damage to contacts and connection socket.</b>  <b>Electrical hazard.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT exceed the maximum load of 3 A.</li> <li>⊘ Do NOT connect any devices with a nominal current &gt; 3 A.</li> </ul>

The interior socket's voltage is turned off at the chamber controller via operation line (chap. 7.3, 9.7.3, 10.6.5).

Turning off the chamber at the main power switch also switches the interior socket voltage-free.

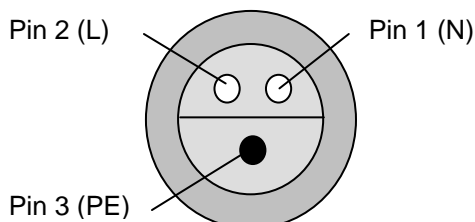


Figure 23: Interior socket (G) (front view)

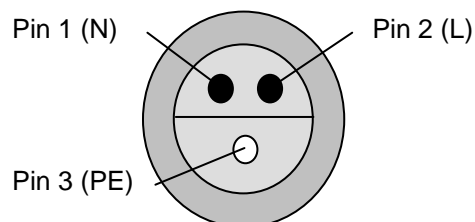




Figure 24: Supplied plug (front view)

Electrical data socket and plug: IP system of protection 65, 230 V 1N ~ 50-60 Hz

	<b>CAUTION</b>
	<p><b>Risk of short circuit.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>➤ Use delivered plug only (IP protection type 65).</li> <li>➤ If the socket is not used, close it with the waterproof lid.</li> </ul>


When inserting a heat load into the inner chamber by introducing electrical devices such as shakers or rollers, the ambient temperature of the incubator must not exceed 25 °C / 77°F. Otherwise, temperature control to 37 °C / 98.6°F cannot be assured. When operating electrical devices inside the incubator, always check the desired incubation temperature **before** introducing any cells.

	<b>CAUTION</b>
	<p><b>Excessive inner temperature &gt; 37 °C / 98.6°F when operating electrical devices inside the incubator.</b></p> <p><b>Damage to the cell cultures.</b></p> <ul style="list-style-type: none"> <li>➤ When operating electrical devices inside the incubator: Ambient temperature ≤ 25 °C / 77 °F.</li> <li>➤ Check the temperature observation before introducing any cells.</li> </ul>

## 18.4 Analog outputs for temperature and CO<sub>2</sub> (option)

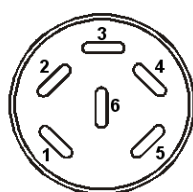With this option, the chamber is equipped with analog outputs 4-20 mA for temperature and CO<sub>2</sub>. These outputs allow transmitting data to external data registration systems or devices.

The connection is realized as a DIN socket (4) on the chamber rear. A suitable DIN plug is enclosed.



### ANALOG OUTPUT 4-20 mA DC

PIN 1: Temperature +

PIN 2: Temperature -

PIN 3: CO<sub>2</sub> +

PIN 4: CO<sub>2</sub> -

CO<sub>2</sub> range: 0 vol.-% up to 20 vol.-%

Temperature range: 0 °C / 32 °F up to +200 °C / 392 °F

Figure 25: Pin configuration of the DIN socket (4) on the chamber rear

## 18.5 Data Logger kits (option)


BINDER Data Logger Kits offer an independent long-term measuring system for temperature and humidity, available for different temperature ranges. According to the selected kit, the Data Logger can measure and record also the ambient temperature and humidity values via a second multi-function sensor.

BINDER Data Loggers are equipped with a keyboard and a large LCD display, alarm functions and a real-time function. Measurement data are recorded in the Data Logger and can be read out after the measurement via the RS232 interface of the Data Logger. It offers a programmable measuring interval and permits storing up to 64000 measuring values. Reading out is done with the Data Logger evaluation software. You can give out a combined alarm and status protocol directly to a serial printer.

**Data Logger Kit T 220:** Sensor for chamber temperature: Temperature range -90 °C / -130 °F up to 220 °C / 428 °F.

**Data Logger Kit TH 70:** Multi-function sensor for chamber temperature and humidity: Temperature range -40 °C / -40 °F up to 70 °C / 158 °F, humidity range 0% r.h. up to 100% r.h.

**Data Logger Kit TH 70/70:** Multi-function sensor for chamber temperature and humidity: Temperature range -40 °C / -40 °F up to 70 °C / 158 °F, humidity range 0% r.h. up to 100% r.h. Multi-function sensor for ambient temperature and humidity: Temperature range -40 °C / -40 °F up to 70 °C / 158 °F, humidity range 0% r.h. up to 100% r.h.

	<p>For detailed information on installation and operation of the BINDER Data Logger, please refer to the mounting instructions Art. No. 7001-0204 and to the original user manual of the manufacturer, supplied with the data logger.</p>
---	---

## 18.6 Access port for extra-low voltage (option)

The access port (B) (8-pin) for extra-low voltage (ELV) consists of a LEMO socket (which can be covered) and a LEMO connector. It is suitable to connect two electrical devices, one inside and one outside the incubator. You can use it e.g. for devices with a monitoring and control part remaining outside the incubator, whereas its mechanical component like roller or shaker systems is located inside the chamber.

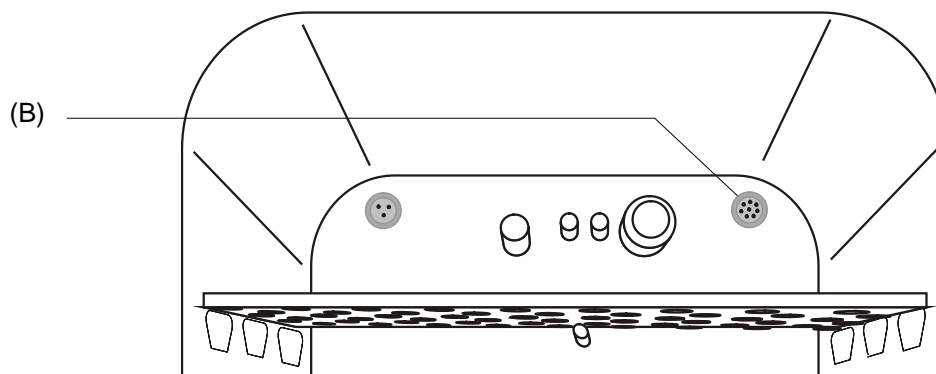




Figure 26: Position of the access port for extra-low voltage

**Maximum power rating 24V AC/DC 2Amp.**

	 <b>WARNING</b>
	<p><b>Overload of contacts.</b>  <b>Damage to contacts and connection socket.</b>  <b>Electrical hazard.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>⊗ Do NOT exceed the maximum load of 24V AC/DC 2Amp.</li> <li>⊗ Do NOT connect any devices with a nominal current &gt; 2 Amp.</li> <li>⊗ Do NOT connect any devices with a nominal voltage &gt; 24 V DC.</li> </ul>

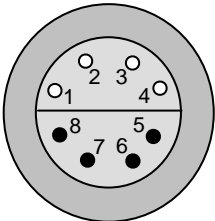
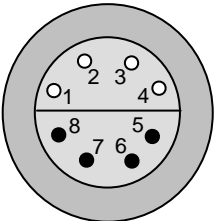
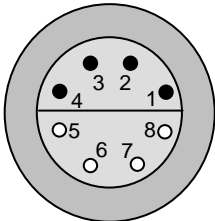

		
<p>On chamber rear: Exterior Lemo socket (2) on the rear of the chamber</p>	<p>Interior: LEMO socket (B) located at the upper right corner of the rear wall of the inner chamber</p>	<p>Supplied LEMO connector: on external and internal side</p>


Figure 27: Pin allocation (front view) of the Lemo sockets and plug

When introducing a heat load into the inner chamber by inserting electrical devices like shakers or rollers, the ambient temperature of the incubator must not exceed 25 °C / 77°F. Otherwise, temperature control to 37 °C / 98.6°F cannot be assured. When operating electrical devices inside the incubator, always check the desired incubation temperature **before** introducing in any cells.

	<b>CAUTION</b>
	<p><b>Excessive inner temperature &gt; 37 °C / 98.6°F when operating electrical devices inside the incubator.</b></p> <p><b>Damage to the cell cultures.</b></p> <ul style="list-style-type: none"> <li>➤ When operating electrical devices inside the incubator: Ambient temperature ≤ 25 °C / 77°F.</li> <li>➤ Before introducing any cells, check the correct ambient temperature.</li> </ul>

## 18.7 BINDER Gas Supply Service – External bottle changer for CO<sub>2</sub>, N<sub>2</sub> or O<sub>2</sub> (option)

The external bottle changer permits automatic switching to a second gas cylinder as soon as the first cylinder is empty. It can be used for a maximum of two CB chambers.

	<p>Instructions 7001-0196 delivered with the external bottle changer (Art. no. 8012-0408) describe its installation and operation.</p>
---	--


## 18.8 Stands

### 18.8.1 Stacking stand (option)

We recommend not stacking CB incubators directly on top of one another in order to avoid transmission of shocks and vibrations from one chamber to the other. This could happen e.g. while opening or closing the door, cleaning, charging and discharging the chamber. BINDER offers stable, vibration-free stands with castors (2 lockable by brakes) for the safe stacking of two chambers.

The stacking stand ensures that the set incubation parameters are precisely maintained also during sterilization of the other chamber in the same stacking stand (chap. 21.3) by thermal decoupling.


Using the stacking stand offers more advantages: You can pull the lower chamber forward separately (e.g., for access to the rear), and not have to place it directly on the floor (important for sanitary purpose).

	<p>The mounting instructions 7001-0194 delivered with the stacking stand describe its installation (Art. No. 9051-0020 for CB 170, Art. No. 9051-0023 for CB 260).</p>
---	--

### 18.8.2 Stacking adapter for direct thermal decoupled stacking (option)

We recommend not stacking CB incubators directly on top of one another in order to avoid transmission of shocks and vibrations from one chamber to the other. This could happen e.g. while opening or closing the door, cleaning, loading and unloading the chamber. BINDER offers a stacking adapter for direct thermal decoupled stacking of two chamber.

The stacking adapter ensures the exact maintenance of the set incubation parameters also during sterilization of the other chamber in the same stacking stand (chap. 21.3).

	<p>The mounting instructions 7001-0145 delivered with the stacking adapter describe its installation (Art. No. 9051-0032 for CB 56, Art. No. 9051-0026 for CB 170, Art. No. 9051-0030 for CB 260).</p>
---	--



### 18.8.3 Base on castors (option)

In order to obtain easy access to the chamber and to avoid contamination of the chamber caused by soil pollution, BINDER recommends using the base on castors.



The mounting instructions 7001-0147 delivered with the base on castors describe its installation (Art. No. 9051-0031 for CB 56, Art. No. 9051-0028 for CB 170, Art. No. 9051-0029 for CB 260).

## 19. Reference measurements

Reference measurements of the temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control), and humidity can be performed via the silicone measuring port (N) located on the inner glass door. Reference temperature measurements always take place under equilibrated conditions with both doors closed.

### 19.1 CO<sub>2</sub> reference measuring

There are three possibilities to perform CO<sub>2</sub> test measurements between the recommended annual maintenance procedures. To test the CO<sub>2</sub> concentration inside an incubator, see chapters 19.1.1 to 19.1.3.

#### 19.1.1 Measuring CO<sub>2</sub> concentration indirectly via the pH of the cell medium

By using the indirect determination of CO<sub>2</sub> concentration via the pH-value of the nutrient, it is possible to check the CO<sub>2</sub> concentration inside the chamber. This is a simple method to test the correct CO<sub>2</sub> concentration without any special CO<sub>2</sub> measuring equipment. You need only use an accurate pH indicator or a pH-measuring electrode, which are standard equipment in cell culture laboratories.

This method is based on the acid base equilibrium of the buffer system in the culture media. NaHCO<sub>3</sub> buffers the common media. From the pH value of the medium, it is possible to conclude its CO<sub>2</sub> concentration. Figure 28 shows the relationship between CO<sub>2</sub> concentration in vol.-% and the pH of different NaHCO<sub>3</sub> buffered media.



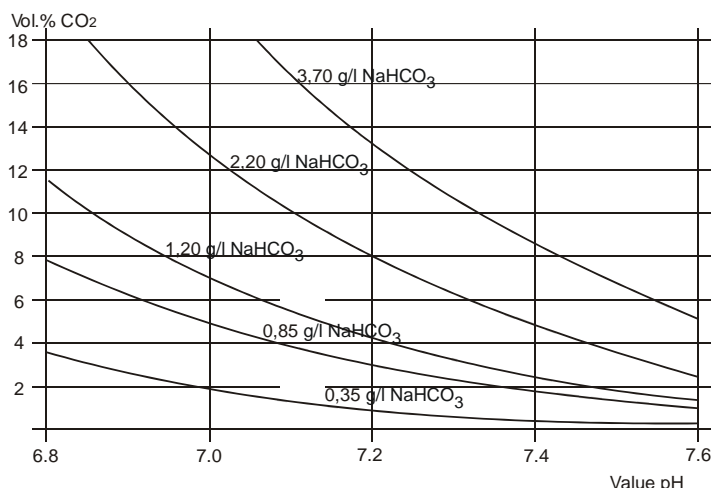
These test systems are not suitable for calibrating the BINDER sensor system.

#### Recommended procedure:

- Incubate an empty sample with medium for 1/2 day under the same conditions as the cells. You can perform the incubation in a cell culture cylinder or in a 50 ml Falcon tube with open lid.
- After gassing, remove the empty sample from the incubator and within 5 minutes measure the pH-value with a glass electrode.

During the measurement, the medium should have the least possible surface contact with the ambient air, so that the CO<sub>2</sub> can evaporate only slightly. A significant downward movement will happen only after 5 minutes, permitting sufficient time for measurement.

- In addition, you can of course use pH-test strips (pH range 6 to 8, not bleeding).



Trade names of common media:

	NaHCO <sub>3</sub> [g/l]
DMEM	3.70
BME	2.20
MEM	2.20
Medium 199	2.20
Mc Coy	2.20
F10	1.20
F12	1.20

Figure 28: Value pH of NaHCO<sub>3</sub> buffered media as a function of the CO<sub>2</sub> concentration:

Example: If a pH of 7.2 is measured in a medium buffered with 2.20 g NaHCO<sub>3</sub> per liter, there must be 8 vol.-% CO<sub>2</sub> surrounding this medium.

### 19.1.2 Measuring CO<sub>2</sub> directly via chemical indicator tubes

This is a common “do-it-yourself” test for many users. A chemical color reaction in a glass tube shows the CO<sub>2</sub> concentration. A standardized volume of air from inside the incubator has to be suctioned through this glass tube to get a quantitative test result. Therefore, use a special hand pump with a standardized suction volume.

#### Procedure (example):

All the necessary equipment must be supplied by one manufacturer only and in a defined test system.

Note: These test systems are not very accurate. A typical accuracy is around 10 % of the full-scale value.



These test systems are not suitable for calibrating the BINDER sensor system.

### 19.1.3 Measuring CO<sub>2</sub> directly with an electronic infrared measuring device

Another possible way of measuring the CO<sub>2</sub> concentration is by electronic sensor systems. Portable measuring devices are suitable for measuring temperature and CO<sub>2</sub> concentration and can be used both for reference measurements in certified laboratories, and for service purposes.

## 19.2 Temperature reference measurement

When performing a temperature reference measurement using an electronic measuring, and temperature display device, it is important to use a device traceable to an acknowledged standards/calibration institution (DKD, PTB for Germany) with a valid calibration certificate.

Note: The cable of the sensor must be thin enough to lay it over the door gasket of the incubator without causing any leakages.

## **20. Avoiding microbial contamination**

The main types of microbial contaminants in cell and tissue culture are bacteria, fungi, yeast, mycoplasma, and viruses. This chapter gives an overview of potential sources of contamination and precautions and measures to eliminate them.

### **20.1 Cells and media**

- Primary cultures from the original tissue
- Cells / cell lines from unknown sources or from cell banks: Use only cells of known and tested origin. Monitoring and routine screening of new cultures.
- Media and sera: Use only sera of known and tested origin (mycoplasma free, e.g., UV or  $\gamma$  radiated).
- Virus suspensions, antibody solutions etc. Use only reagents of known and tested origin.
- Laboratory instruments, media and reagents that were exposed to possible contaminated cultures must be sterilized / autoclaved / disposed.
- Antibiotics in the cell culture media may prevent bacteria detection: Use antibiotics selectively and economically.

### **20.2 Laboratory conditions / equipment around the incubator**

Possible sources of contamination in the cell culture lab are airborne germs, lab equipment, building features, and the lab personnel.

- Keep pipettes and instruments sterile after autoclaving.
- Bio safety cabinets (laminar air flow) should have a minimum of items apart from aspirator tube and burner. Items shall be positioned within easy reach and separate from each other. Disinfect surfaces with an alcohol-based disinfectant before and after use, clean the space underneath the bench, and carry out regular sterility tests of the filters.
- Regular cleaning / disinfection of laboratory equipment such as a centrifuge, microscope, water bath, refrigerator, and telephone.
- No equipment should be placed on the floor.
- Rough or humid walls are not suitable.
- Identify leaking doors and windows and make them airtight.
- Use air conditioning with special filters.
- Reduce the number of personnel and their movements in the lab by careful positioning all relevant equipment. For practical reasons, install the chamber close to the laminar air flow bench.
- Regular microbiological monitoring of the cell culture laboratory.

### **20.3 Working and behavior in the lab**

Sources of contamination are often the laboratory personnel themselves (surface germs, oral flora droplet transfer) and handling the equipment and cultures. We recommend staff training in aseptic techniques, laboratory safety and good laboratory practice (GLP).

#### **Examples of general rules to reduce the contamination risks**

- Reduce hand germ count (wash hands with antimicrobial soap, dry with paper tissues, and rub dry hands with alcohol-based solution).
- Wear appropriate clothing (work coat, shoes, face mask)
- Keep as few personnel as possible in the cell culture lab.

### **Examples of sterile working method**

- Work "clean-to-dirty", i.e., handle confirmed uncontaminated cells first, unknown or untested cells next, and lastly, if necessary, cells suspected to be contaminated.
- Perform daily microscopic observations of cultures and specific tests for the bacteria and fungi as part of a controlled routine. Test cultures for sterility before starting work.
- Keep working surfaces clean. Immediately wipe spilled liquids with alcohol solutions.
- No mouth contact on pipettes.
- Never work on top of open sterile containers.

## **20.4 Chamber design and equipment of the CO<sub>2</sub> incubator**

The design concepts behind the CO<sub>2</sub> incubator CB considerably reduce the risk of contamination. Among them are:

### **Even surfaces for easy manual cleaning**

- The inner surfaces are smooth and therefore easy to clean. The inner chamber is deep-drawn from a single piece, polished (suitable for pharmaceutical work) and has no welds or inaccessible corners. The hinges and the seal of the inner glass door are glued from the outside, which also aids cleaning of the inner chamber.

### **Removable parts for cleaning and autoclaving**

- The shelves are easily removed without screws. It is possible to autoclave the shelves. But this is generally not necessary because they can remain inside the chamber during sterilization.

### **Door gasket**

- The inner door gasket is removable and autoclavable.

### **Gas fine filter**

- The incoming gas used in the operation passes through a fine filter (aseptic filter, filtration efficiency 99.99 %, particle size 0.45 µm) with a high filtration efficiency, which can also filter the smallest particles.

### **CO<sub>2</sub> measuring system in the inner chamber**

- The CO<sub>2</sub> sensor can be hot-air sterilized inside the chamber.

### **O<sub>2</sub> measuring system in the inner chamber (chamber with O<sub>2</sub> control)**

- The O<sub>2</sub> sensor can be hot-air sterilized inside the chamber.

### **Humidity measuring system in the inner chamber (chamber with active humidification)**

- The humidity sensor can be hot-air sterilized inside the chamber.

### **Condensation prevention**

- Condensation in the inner chamber represents a particular risk of contamination. The Permadyr™ system developed by BINDER is an effective and easy way to ensure high humidity (≥95 % r.h.) inside the chamber without any condensation forming on the inner surfaces.

### **Hot air sterilization at 180 °C / 356 °F**

- The heating system of the CO<sub>2</sub> incubator permits hot-air auto-sterilization at a setpoint of 187.5 °C / 369.5 °F. Thus, a temperature of 180 °C / 356 °F is maintained for at least 30 minutes on all internal surfaces, resulting in sterilization of the entire inner chamber. Therefore, this procedure meets all international guidelines regarding hot air sterilization, e.g. AAMI ST63, DIN 58947, European Pharmacopoeia.

## 20.5 Handling the CO<sub>2</sub> incubator

Any manipulation of the CO<sub>2</sub> incubator involves some contamination risks, from installation to opening the doors and regular cleaning.

### Installation away from sources of contamination

- Do not place the chamber on the floor or close to windows and doors. Use the optional stand, if appropriate.

### Reduce the periods in which the door is open

- Do not open the door too frequently.
- Placing items in order inside the chamber results in shorter door opening times.

### Permadyr™ water pan

- Fill the outer pan with distilled, sterilized water (chap. 4.2). Never use ion exchange water; ion exchangers are propagation sites for bacteria.
- Clean and refill the pans 2 to 3 times a week. For evacuation, remove the Permadyr™ water pan. It is autoclavable.
- If desired, you can add microbiologically inhibiting substances as copper chips, copper sulfate or ethylene diamine tetra-vinegar acid (EDTA) in a concentration of 1 to 5 mmol/l.

### Avoiding condensation caused by ambient conditions

Ambient room conditions have an effect on condensation inside the chamber, which can be caused by insufficient wall clearances, preventing even dissipation of heat, air movement or direct sunlight. If the temperature distribution inside the chamber becomes uneven, condensation may form on the cooler surfaces.

- Maintain distances from the wall: rear 100 mm / 3.94 in, sides 50 mm / 1.97 in
- Do not place the chamber in front of a window. No direct sunlight. No air movement.
- Permissible ambient temperature range for operation: +18 °C / 64.4°F to +30 °C / 86°F.

Ideal ambient temperature: by at least 7 °C / 12.6 °F below the intended working temperature. E.g., working temperature 37 °C / 98.6 °F = ambient temperature 30 °C / 86 °F and less.

- The chamber should be calibrated / precisely adjusted.
- Let BINDER Service adjust the setting of the door heating in reaction to critical ambient conditions.






### Regular cleaning, decontamination and sterilization

- Clean the shelves, glass door, gaskets, and inner chamber weekly (for cleaning see chap. 21.1, for decontamination see chap. 21.2. You can clean the shelves in a laboratory dishwasher and, if needed, individually autoclave them.
- Regularly use the hot air sterilization function (chap. 21.3) following cleaning. The shelves and the emptied water pan can remain inside the chamber during this operation.
- Have replaced the CO<sub>2</sub> sterile filter (gas fine filter) once or twice a year.

### What to do in case of contamination?


- Throw away / autoclave contaminated cultures.
- Carefully inspect cultures that seem to be uncontaminated.
- Clean the chamber as described. Wipe the inner chamber and the doors with a disinfectant and allow drying. Autoclave the shelves. Empty the water pan and autoclave it.
- Perform a hot air sterilization.

## 21. Cleaning, decontamination / disinfection, and sterilization

   	<div style="background-color: red; color: white; text-align: center; padding: 5px;">  <b>DANGER</b> </div> <p><b>Electrical hazard.</b>  <b>Danger from cleaning agents on hot surfaces.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT spill water or cleaning agents over the inner and outer surfaces.</li> <li>➤ Before cleaning, turn off the chamber at the main power switch and disconnect the power plug. Let the chamber cool down to ambient temperature.</li> <li>➤ Completely dry the chamber before turning it on again.</li> </ul>
---	---

### 21.1 Cleaning


Disconnect the chamber from the power supply before cleaning. Disconnect the power plug.


	<p>The interior of the chamber must be kept clean. Thoroughly remove any residues of the charging material.</p>
---	---

Wipe the surfaces with a moistened towel. In addition, you can use the following cleaning agents:


Exterior surfaces, instrument panel	Standard commercial cleaning detergents free from acid or halides. Alcohol-based solutions. We recommend using the neutral cleaning agent Art. No. 1002-0016.
Inner chamber, shelves	Standard commercial cleaning detergents free from acid or halides. Copper sulfate solutions or alcohol-based solutions. We recommend using the neutral cleaning agent Art. No. 1002-0016.
Silicone door gasket	Alcohol-based solutions or the neutral cleaning agent Art. No. 1002-0016.
Zinc coated hinge parts rear chamber wall	Standard commercial cleaning detergents free from acid or halides. Do NOT use a neutral cleaning agent on zinc coated surfaces.

Do not use cleaning agents that may cause a hazard due to reaction with components of the device or the charging material. If there is doubt regarding the suitability of cleaning products, please contact BINDER service.


	<p>We recommend using the neutral cleaning agent Art. No. 1002-0016 for thorough cleaning.  Any corrosive damage that may arise following use of other cleaning agents is excluded from liability by BINDER GmbH.  Any corrosive damage caused by a lack of cleaning, is excluded from liability by BINDER GmbH.</p>
---	--


	<div style="background-color: yellow; text-align: center; padding: 5px;"> <b>CAUTION</b> </div> <p><b>Danger of corrosion.</b>  <b>Damage to the chamber.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT use acidic or chlorine cleaning detergents.</li> <li>Ø Do NOT use a neutral cleaning agent on other kind of surfaces e.g., the zinc coated hinge parts or the rear chamber wall.</li> </ul>
---	---


To maintain the proper function of the CO<sub>2</sub> sensor, never spray the sensor with any cleaning detergents or disinfectants. Generally perform any cleaning only when the chamber is turned off.

	<b>CAUTION</b>
	<p><b>Damage to the CO<sub>2</sub> sensor.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT spray cleaning detergents directly on the CO<sub>2</sub> sensor.</li> <li>➤ Wipe the sensor with a lint-free cloth soaked with cleaning agent.</li> </ul>


Use only the detergents or disinfectants recommended by BINDER. Other products are not approved, since they could in particular damage the CO<sub>2</sub> sensor.

	<p>For surface protection, perform cleaning as quickly as possible. After cleaning completely remove any cleaning agents from the surfaces by using a moistened towel. Let the chamber dry.</p>
---	---





	<p>Soapsuds may contain chlorides and must therefore NOT be used for cleaning.</p>
---	--


	<p>With every decontamination method, always use adequate personal safety controls.</p>
---	---

Following cleaning, leave the chamber door open or remove the access port plugs.

	<p>The neutral cleaning agent may cause health problems in contact with skin and if ingested. Follow the operating instructions and safety hints labeled on the bottle of the neutral cleaning agent.</p>
---	---

Recommended precautions: To protect the eyes use sealed protective goggles. Suitable protective gloves with full contact: butyl or nitrile rubber, penetration time >480 minutes.

	<table border="1"> <tr> <td data-bbox="541 1245 1481 1323" style="background-color: yellow; text-align: center;">  <b>CAUTION</b> </td> </tr> <tr> <td data-bbox="541 1323 1481 1635"> <p><b>Contact with skin, ingestion.</b> <b>Skin and eye damage due to chemical burns.</b></p> <ul style="list-style-type: none"> <li>⊘ Do not ingest. Keep away from food and beverages.</li> <li>⊘ Do NOT empty into drains.</li> <li>➤ Wear protective gloves and goggles.</li> <li>➤ Avoid skin contact.</li> </ul> </td> </tr> </table>	 <b>CAUTION</b>	<p><b>Contact with skin, ingestion.</b> <b>Skin and eye damage due to chemical burns.</b></p> <ul style="list-style-type: none"> <li>⊘ Do not ingest. Keep away from food and beverages.</li> <li>⊘ Do NOT empty into drains.</li> <li>➤ Wear protective gloves and goggles.</li> <li>➤ Avoid skin contact.</li> </ul>
 <b>CAUTION</b>			
<p><b>Contact with skin, ingestion.</b> <b>Skin and eye damage due to chemical burns.</b></p> <ul style="list-style-type: none"> <li>⊘ Do not ingest. Keep away from food and beverages.</li> <li>⊘ Do NOT empty into drains.</li> <li>➤ Wear protective gloves and goggles.</li> <li>➤ Avoid skin contact.</li> </ul>			

	<p>Following use of the neutral cleaning agent and prior to hot-air sterilization, remove any agent residues by using a moistened towel in order to avoid formation of permanent residues.</p>
---	--



## 21.2 Decontamination / chemical disinfection of the chamber


The operator must ensure that proper decontamination is performed in case a contamination of the chamber by hazardous substances has occurred.

Disconnect the chamber from the power supply prior to chemical decontamination. Disconnect the power plug.


Do not use decontamination agents that may cause a hazard due to reaction with components of the device or the charging material. If there is doubt regarding the suitability of cleaning products, please contact BINDER service.

You can use the following disinfectants:


Inner chamber	Standard commercial surface disinfectants free from acid or halides. Alcohol based solutions. We recommend using the disinfectant spray Art. No. 1002-0022.
---------------	---

	For chemical disinfection, we recommend using the disinfectant spray Art. No. 1002-0022. Any corrosive damage that may arise following use of other disinfectants is excluded from liability by BINDER GmbH.
---	---

Important: To maintain the proper function of the CO<sub>2</sub> sensor, never spray the sensor with any cleaning detergents or disinfectants. Generally perform any disinfection only when the chamber is turned off. The sensor may be only superficially disinfected with a damp cloth. The hot air sterilization routine is intended for a thorough sterilization of the entire chamber.

	<p style="text-align: center;"><b>CAUTION</b></p> <p><b>Damage to the CO<sub>2</sub> sensor.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT spray the disinfectant directly on the CO<sub>2</sub> sensor.</li> <li>➤ Wipe the sensor with a lint-free cloth soaked with the disinfectant.</li> </ul>
---	---

Use only the detergents or disinfectants recommended by BINDER. Other products are not approved, since they could in particular damage the CO<sub>2</sub> sensor.


	With every decontamination method, always use adequate personal safety controls.
---	--

In case of contamination of the interior by biologically or chemically hazardous material, there are two possible procedures depending on the type of contamination and charging material.

(1) Spray the inner chamber with an appropriate disinfectant.



Before start-up, the chamber must be absolutely dry and ventilated, as explosive gases may form during the decontamination process.


(2) If necessary, have strongly contaminated inner chamber parts removed by an engineer for cleaning, or have them exchanged. Sterilize the inner chamber parts in a sterilizer or autoclave.


	In case of eye contact, the disinfectant spray may cause eye damage due to chemical burns. Follow the operating instructions and safety hints labeled on the bottle of the disinfectant spray.
---	--




Recommended precautions: To protect the eyes use sealed protective goggles.

	<div style="background-color: yellow; text-align: center; padding: 5px;">  <b>CAUTION</b> </div> <p><b>Eye contact.</b>  <b>Eye damage due to chemical burns.</b></p> <p>Ø Do NOT empty into drains.          ➤ Wear protective goggles.</p>
---	---

	<p>Following frequent use of the disinfectant spray and prior to hot-air sterilization, remove any agent remainder by using the neutral cleaning agent and then a moistened towel to avoid formation of permanent residues.</p>
---	---

	<p>After using the disinfectant spray, allow the chamber to dry thoroughly, and aerate it sufficiently.</p>
---	---

## 21.3 Hot-air sterilization at 180 °C / 356 °F

	<p>The very first sterilization after operation may cause an odor. This is not a quality defect. We recommend ventilating well the room during sterilization.</p>
---	---

### 21.3.1 Overview

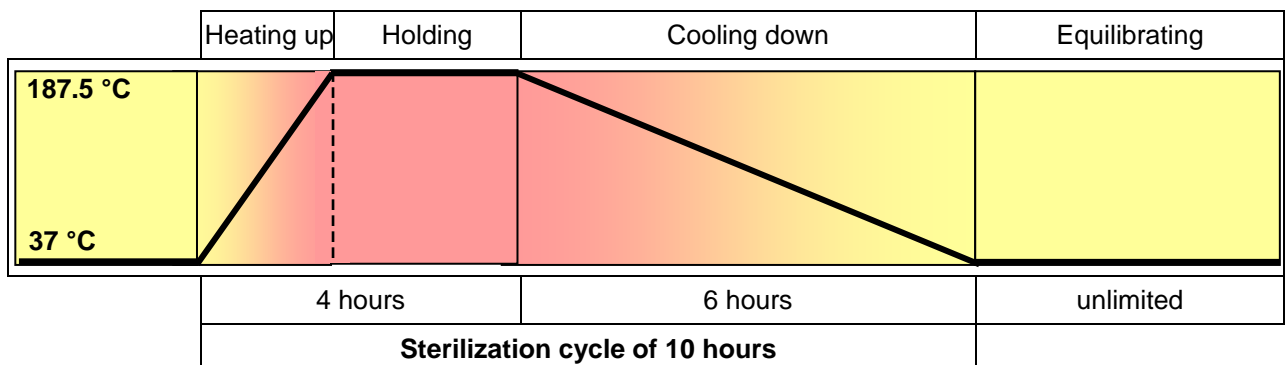


Figure 29: Setpoint profile during the sterilization cycle


The chamber can perform an automatically controlled hot-air sterilization cycle. This procedure will take approx. 10 hours and consists of the following steps:

- **Heating up phase:** The chamber heats up the inner chamber as fast as possible to the sterilization setpoint temperature
- **Holding phase:** Constant sterilization setpoint temperature  
 The sterilization setpoint temperature is pre-set in factory to 187.5 °C / 369.5 °F. The duration of the heating up and holding phase is in total 4 hours. This ensures that 180 °C / 356 °F is maintained on all internal surfaces for at least 30 minutes
- **Cooling down phase:** Programmed duration of 6 hours until 37 °C / 98.6 °F is reached
- Then follows an **equilibration phase:** After the 10-hour sterilization cycle the controller changes to fixed value operation mode and equilibrates to the set-point which has been entered there (e.g., 37 °C / 98.6 °F).


Note: in the case when two chambers are directly stacked on top of each other **without** using the original BINDER stacking stand or the BINDER stacking adapter, the exact maintenance of the incubation set parameters in one of the chambers while carrying out sterilization in the other one cannot be assured. Therefore, without using the stacking stand or the stacking adapter, safe incubation is not possible during sterilization of the other chamber. We recommend using a stacking stand (chap. 18.8.1) or a stacking adapter (chap. 18.8.2).


CO<sub>2</sub> / O<sub>2</sub> control is deactivated during the sterilization cycle. To activate gas control, the following conditions are required:


- Completing or prematurely terminating the sterilization cycle and confirming the corresponding information message
- Inner chamber temperature below 80 °C


	<p>If the temperature exceeds &lt; 80 °C, gas control will remain deactivated even after completing or prematurely terminating the sterilization cycle.</p> <p>Only after the temperature has decreased below 80 °C, gas control will turn on again.</p>
---	--

### 21.3.2 Preparation for a hot-air sterilization



	<p>Before carrying out the first hot-air sterilization, remove any protective lamination sheet from the inner metal surfaces.</p>
--	---

	<p>When starting a hot-air sterilization, all gas controls automatically deactivate.</p>
---	--


	<p>The safety controller settings are inactive during sterilization. They become functional again following completing or prematurely terminating of the sterilization cycle and/or restart of the chamber at the main power switch.</p>
---	--

	<p>The O<sub>2</sub> sensor must be plugged in during a hot-air sterilization.</p>
---	--


- Empty the Permadyr™ water pan.

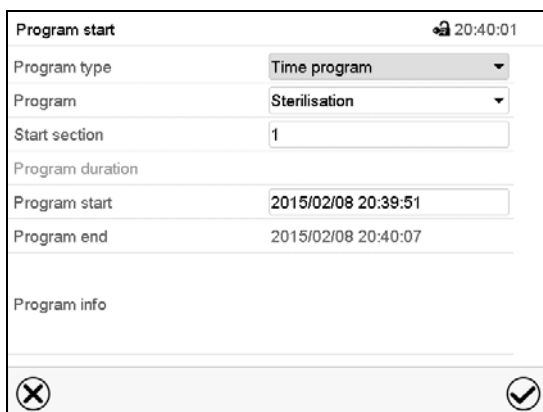
	 <b>WARNING</b>
	<p><b>Danger of implosion.</b>  <b>Destruction of the chamber.</b></p> <p>➤ Empty the Permadyr™ water pan before starting the hot-air sterilization.</p>

- Water pans and shelves must be inside the chamber.
- Close the inner glass door and the outer chamber door.

	<p>Before starting a hot-air sterilization, the entire interior must be clean and dry. No residue of e.g., water, medium or plastic must remain inside the chamber.</p>
---	---

### 21.3.3 Starting the sterilization cycle and running the hot-air sterilization


 In Normal display press the **Program start** icon to access the “Program start” menu.



“Program start” menu

- In the field “Program type” select the setting “Time program”.
- In the field “Program” select the program “Sterilization”.
- Select the field “Program start” and enter the desired program start time in the “Program start” entry menu. Press the **Confirm** icon.

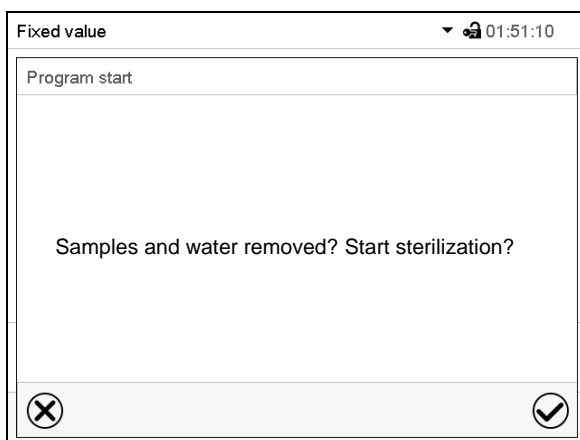A confirmation prompt is displayed.




Make sure that the water pan was emptied and all samples were removed from the chamber and press the **Confirm** icon to confirm the message.



The controller returns to the initial view and the program delay time until program start begins.

When the program delay time is over, the hot-air sterilization cycle begins.

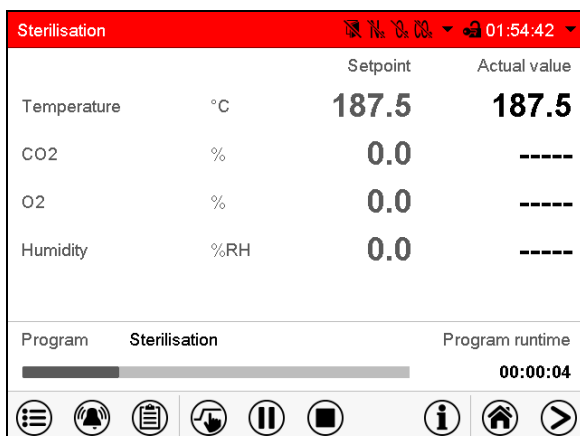Confirmation prompt.



	Setpoint	Actual value
Temperature °C	187.5	187.5
CO2 %	0.0	----
O2 %	0.0	----
Humidity %RH	0.0	----


Program: Sterilisation      Program runtime: 00:00:04



Normal display with alarm message “Sterilization”.

The sterilization is running.

The zero-voltage relay alarm output and the buzzer are not activated.


Opening the outer door leads to prematurely terminating the running sterilization cycle. If this happens during the heating up or holding phase (during 4 hours after starting the sterilization cycle), sterilization will be ineffective (chap. 21.3.4.1).

	<b>CAUTION</b>
	<p><b>Opening the outer door during the heating up or holding phase.</b>  <b>Interruption of the temperature reaction time.</b>  <b>Ineffective sterilization.</b></p> <p>⊘ Do NOT open the chamber doors during the heating up or holding phase of the sterilization cycle</p>

	 <b>CAUTION</b>
	<p><b>Glass door and inner chamber become hot during sterilization.</b>  <b>Danger of burning.</b></p> <p>⊘ Do NOT touch the glass door and inner surfaces during sterilization</p>

Note: Since the stored sterilization program cannot be changed, it is not accessible via the program path [Main menu > Programs > Time program](#)

The sterilization program must not be interrupted using the **Program pause** icon.

	<b>CAUTION</b>
	<p><b>Pausing the sterilization program during the heating up or holding phase.</b>  <b>Interruption of the temperature reaction time.</b>  <b>Ineffective sterilization.</b></p> <p>⊘ DO NOT interrupt a running sterilization program during the heating up or holding phase of the sterilization cycle.</p> <p>➤ If necessary, repeat the sterilization.</p>

#### 21.3.4 Prematurely terminating the sterilization cycle

Three events lead to terminating the sterilization cycle prematurely:

- Cancelling the program via the controller menu  
 To prematurely terminate the sterilization cycle via the controller menu, press the **Program cancelling** icon and confirm the confirmation prompt with the **Confirm** icon (chap. 9.2.2).
- Opening the outer door
- Turning off the chamber at the main power switch, or a power failure

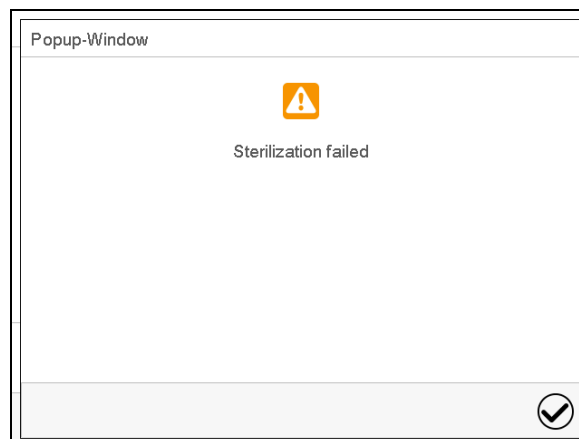
Prematurely terminating the sterilization cycle means terminating it before the entire 10-hour duration is completed. Whether effective sterilization has occurred depends on the time that has elapsed before.

#### 21.3.4.1 Prematurely terminating the sterilization cycle after less than 4 hours: Ineffective sterilization

After prematurely terminating the sterilization cycle after less than 4 hours, the information "Sterilization failed" is shown on the controller display.

The interior may still be hot. Do not touch the glass door and inner surfaces!

Confirm the message with the **Confirm** icon.



After confirming the message the controller changes to Fixed value operation mode. The setpoints of Fixed value operation mode are then equilibrated.



CO<sub>2</sub> / O<sub>2</sub> control is deactivated and will only be activated after confirming the information message "Sterilization failed" and when the temperature has cooled down to a value below < 80 °C.

When prematurely terminating the sterilization cycle, it may be that the cells/pathogens inside the chamber have not all been killed. If necessary, you should repeat the sterilization.



#### CAUTION

**Interruption of temperature reaction time.**

**Ineffective sterilization.**

➤ If necessary, repeat the sterilization.



#### CAUTION

**Glass door and inner chamber become hot during sterilization.**

**Danger of burning.**

Ø Do NOT touch the glass door and inner surfaces for approx. 7 hours after prematurely terminating the sterilization cycle.

#### 21.3.4.2 Prematurely terminating the sterilization cycle after more than 4 hours, i.e., during the cooling-down phase: successful sterilization

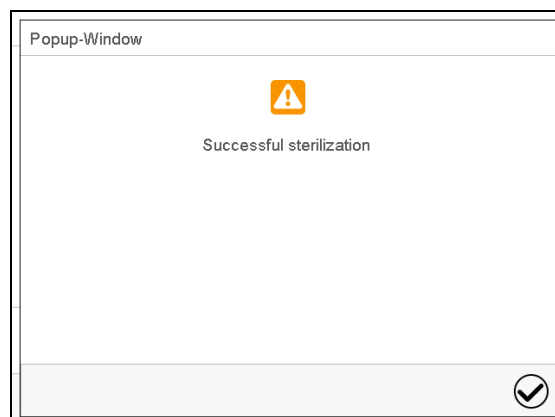
The duration of the entire sterilization is 10 hours. If you want to shorten the sterilization procedure in order to save time, you can prematurely terminate it during the cooling-down phase, i.e. no sooner than after 4 hours. At this point, the inner temperature is still approx. 140 °C / 284 °F.

The effective sterilization phase (heating up and holding phase) is automatically finished after 4 hours. The chamber is in the cooling-down phase, meaning that the necessary duration for the proper sterilization has occurred.

The information "Successful sterilization" on the controller display indicates a successful hot-air sterilization.

The interior may still be hot. Do not touch the glass door and inner surfaces!

Confirm the message with the **Confirm** icon.



After confirming the message the controller changes to Fixed value operation mode. The setpoints of Fixed value operation mode are then equilibrated.



CO<sub>2</sub> / O<sub>2</sub> control is deactivated and will only be activated after confirming the information message "Successful sterilization" and when the temperature has cooled down to a value below < 80 °C.

If you prefer accelerating cooling-down, proceed as follows:

- If required, open the outer door.
- Cooling-down time:
  - Front door open: at least 1 hour
  - Front door closed: at least 4 hours
- If required, open the glass door.



## CAUTION

**Glass door and inner chamber become hot during sterilization.**  
**The glass door handle reaches a temperature of approx. 150 °C / 302 °F.**  
**Danger of burning.**

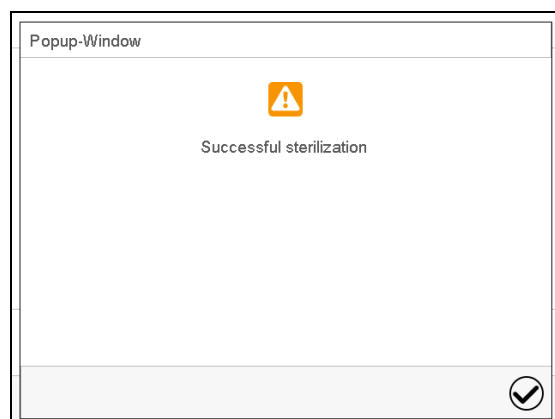
- ⊘ Do NOT touch the glass door and inner surfaces.
- Use gloves or a tool (e.g. pincers) to open the glass door.


### 21.3.5 Completing the entire sterilization cycle

The effective sterilization phase (heating up and holding phase) is automatically finished after 4 hours. It is followed by a defined 6-hours cooling down phase until 37 °C / 98.6°F is reached. After this sterilization cycle of 10 hours in total, the controller changes to fixed value operation mode and equilibrates to the set-point which has been entered there (e.g., 37 °C / 98.6 °F).

The information "Successful sterilization" on the controller display indicates a successful hot-air sterilization.




Confirm the message with the **Confirm** icon.




	CO <sub>2</sub> / O <sub>2</sub> control is deactivated and will only be activated after confirming the information message "Successful sterilization" and when the temperature has cooled down to a value below < 80 °C.
---	---

## 22. Maintenance and service


### 22.1 Maintenance intervals, service

 	 <b>DANGER</b>
	<p><b>Electrical hazard.</b>  <b>Danger of death.</b></p> <ul style="list-style-type: none"> <li>⊘ The chamber must NOT become wet during operation or maintenance works.</li> <li>⊘ Do NOT remove the rear panel of the chamber.</li> <li>➤ Before conducting maintenance work, turn off the chamber at the main power switch and disconnect the power plug.</li> <li>➤ Any maintenance work must be conducted by licensed electricians or experts authorized by BINDER.</li> </ul>

Ensure regular maintenance work is performed at least once a year and that the legal requirements are met regarding the qualifications of service personnel, scope of testing and documentation.

	The warranty becomes void if maintenance work is conducted by non-authorized personnel.
---	---


*Chamber with active humidification:* Have conducted regular maintenance work on the steam humidifier at least once a year. The operating behavior and the maintenance intervals of the humidifier essentially depend on the available water quality and the amount of steam produced in the meantime.

	Replace the door gasket only when cold. Otherwise, the door gasket may become damaged.
---	--

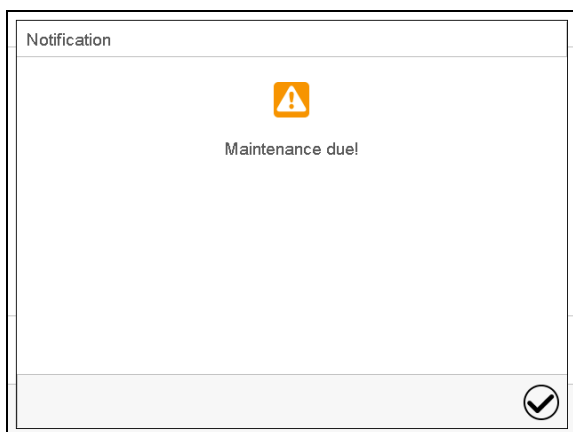
We recommend taking out a maintenance agreement. Please consult BINDER Service.

BINDER telephone hotline:	+49 (0) 7462 2005 555
BINDER fax hotline:	+49 (0) 7462 2005 93555
BINDER e-mail hotline:	service@binder-world.com
BINDER service hotline USA:	+1 866 885 9794 or +1 631 224 4340 x3 (toll-free in the USA)
BINDER service hotline Asia Pacific:	+852 390 705 04 or +852 390 705 03
BINDER service hotline Russia and CIS	+7 495 988 15 16
BINDER Internet website	<a href="http://www.binder-world.com">http://www.binder-world.com</a>
BINDER address	BINDER GmbH, post office box 102, 78502 Tuttlingen, Germany

International customers, please contact your local BINDER distributor.

	The gas sensors are especially adjusted for the specific incubator. When exchanging a sensor, you must repeat control adjustment for CO <sub>2</sub> and O <sub>2</sub> (chamber with O <sub>2</sub> control).
---	--

After 8760 operating hours or two years the following message appears:



After confirmation with the **Confirm** icon, the message window will pop up again every two weeks until it is reset by BINDER Service.

## 22.2 Checking the humidity system fan

The user should regularly perform the following checks:

Hold a sheet of paper in front of the ventilation silts at the bottom rear of the chamber. If the sheet of paper is suctioned inwards, the fan operates correctly. When opening the door, the fan automatically turns off.

## 22.3 Gas inlet fine filter

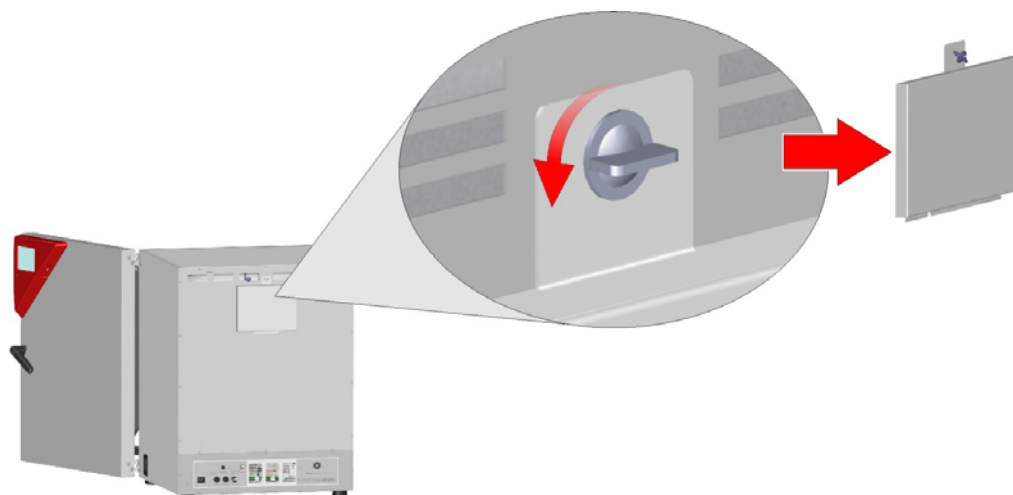
When the chamber is in operation, the incoming gas passes through a fine gas filter (aseptic filter, filtration efficiency 99.99 %, particle size 0.45 µm). This fine filter prevents dirt accumulating in the gas inlet valves and the tubes leading into the inner chamber, which could be in the gas cylinder or in the supply tubes.

When using gas with a technical grade of 99.5 %, we recommend changing the fine gas filter once a year. Please consult BINDER Service. When using gases with less pureness, the changing intervals should be shorter.

## 22.4 Replacing the CO<sub>2</sub> sensor


Pull the power plug.

Remove the cover on the chamber rear:






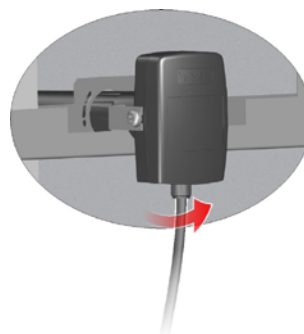
Unscrew the cable from the sensor.  
Make sure that the sensor cable does not fall into the inner chamber.



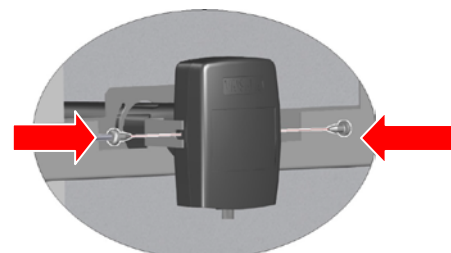Remove the fixing screws on both sides of the sensor.



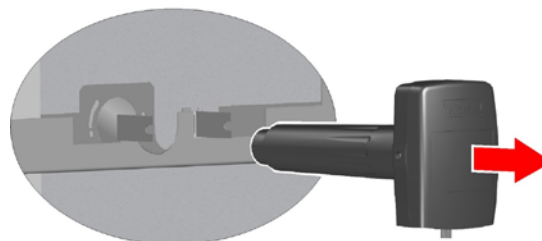Pull the sensor backwards out of the holder.



Then mount the new sensor proceeding in reverse order



After replacing the CO<sub>2</sub> sensor repeat the CO<sub>2</sub> calibration.

## 22.5 Sending the chamber back to BINDER GmbH

If you return a BINDER product to us for repair or any other reason, we will only accept the product upon presentation of an **authorization number** (RMA number) that has previously been issued to you. An authorization number will be issued after receiving your complaint either in writing or by telephone **prior** to your sending the BINDER product back to us. The authorization number will be issued following receipt of the information below:

- BINDER product type and serial number
- Date of purchase
- Name and address of the dealer from which you bought the BINDER product
- Exact description of the defect or fault
- Complete address, contact person and availability of that person
- Exact location of the BINDER product in your facility
- A contamination clearance certificate (chap. 28) must be faxed in advance

The authorization number must be applied to the packaging in such a way that it can be easily recognized or be recorded clearly in the delivery documents.



For security reasons we cannot accept a chamber delivery if it does not carry an authorization number.

**Return address:**


BINDER GmbH  
Abteilung Service

Gänsäcker 16  
78502 Tuttlingen, Germany

## 23. Disposal

### 23.1 Disposal of the transport packing
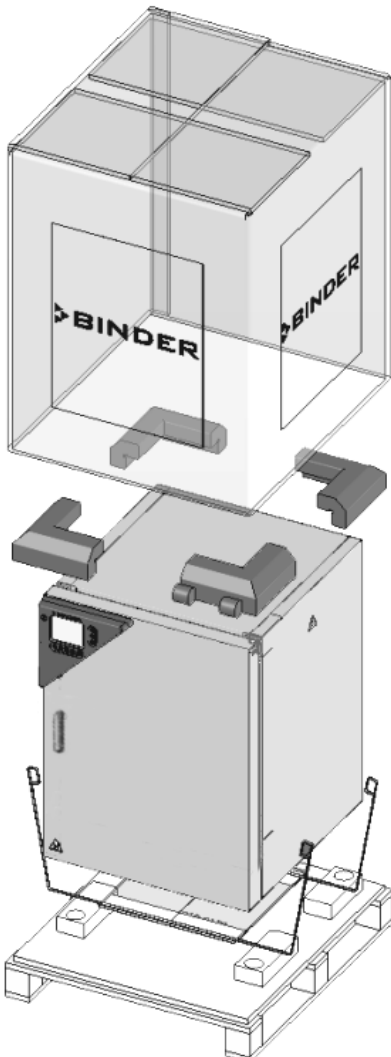
#### 23.1.1 Outer chamber packing

Packing element		Material	Disposal
	Straps to fix packing on pallet (no image)	Plastic	Plastic recycling
	Shipping box	Cardboard	Paper recycling
	Edge stuffing, top	PE foam	Plastic recycling
	Pallet with foamed plastic stuffing	PE foam	Plastic recycling
		Solid wood (IPPC standard)	Wood recycling

#### 23.1.2 Packing inside the chamber and equipment

Packing element	Material	Disposal
Door protection	PE foam	Plastic recycling
Packing box equipment	Cardboard	Paper recycling
Insulating air cushion foil	PE foil	Plastic recycling
Bag for operating manuals	PE foil	Plastic recycling



If recycling is not possible, all packing parts can also be disposed of with normal waste.

## 23.2 Decommissioning

- Turn off the chamber at the main power switch and disconnect it from the power supply (pull the power plug).



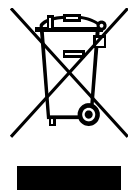
When switching off the main power switch, the stored parameters remain saved.

- Turn off the CO<sub>2</sub> supply and the O<sub>2</sub> / N<sub>2</sub> supplies (chamber with O<sub>2</sub> control). Remove the gas connections.
- *Chamber with active humidification*: Empty the freshwater bag.
- Let the inner chamber sufficiently cool down after a hot-air sterilization before removing any parts.
- The Permadyr™ water pan must not remain filled while the chamber is out of operation. Otherwise condensation on the inner surfaces may occur. In this case, clean and dry the chamber with doors open for at least one hour before restarting the chamber. BINDER recommends performing a hot air sterilization of the chamber before commissioning.
- Temporal decommissioning: See indications for appropriate storage, chap. 3.3.
- Final decommissioning: Dispose of the chamber as described in chap. 23.3 to 23.5.
- When restarting the chamber, please pay attention to the corresponding safety information.


## 23.3 Disposal of the chamber in the Federal Republic of Germany

According to Annex I of Directive 2012/19/EU of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE), BINDER devices are classified as “monitoring and control instruments” (category 9) only intended for professional use“. They must not be disposed of at public collecting points.


The chambers bear the symbol for the marking of electrical and electronic equipment manufactured / placed on the market in the EU after 13 August 2005 and be disposed of in separate collection according to Directive 2012/19/EU on waste electrical and electronic equipment (WEEE) and German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG). WEEE marking: crossed-out wheeled bin with solid bar under. A significant part of the materials must be recycled in order to protect the environment.





At the end of the device's service life, have the chamber disposed of according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG from 20 October 2015, BGBl. I p. 1739) or contact BINDER service who will organize taking back and disposal of the chamber according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG from 20 October 2015, BGBl. I p. 1739).

	<b>CAUTION</b>
	<p><b>Violation against existing law.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT dispose of BINDER devices at public collecting points.</li> <li>➤ Have the device disposed of professionally at a recycling company which is certified according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG from 20 October 2015, BGBl. I p. 1739).</li> <li style="text-align: center;">or</li> <li>➤ Instruct BINDER Service to dispose of the device. The general terms of payment and delivery of BINDER GmbH apply, which were valid at the time of purchasing the chamber.</li> </ul>

Certified companies disassemble waste (used) BINDER equipment in primary substances for recycling according to Directive 2012/19/EU. The devices must be free from toxic, infectious or radioactive substances in order to eliminate any health hazards to the employees of the recycling companies.


	<ul style="list-style-type: none"> <li>• Prior to handing the chamber over to a recycling company, it is the user's responsibility that it is free from toxic, infectious or radioactive substances.</li> <li>• Prior to disposal, clean all introduced or residual toxic substances from the chamber.</li> <li>• Prior to disposal, disinfect the chamber from all sources of infection. Be aware that sources of infection may also be located outside the inner chamber.</li> <li>• If you cannot safely remove all toxic substances and sources of infection from the chamber, dispose of it as special waste according to national law.</li> <li>• Fill out the contamination clearance certificate (chap. 28) and enclose it with the chamber.</li> </ul>
---	---

	<div style="background-color: orange; text-align: center; padding: 5px;">  <b>WARNING</b> </div> <p><b>Contamination of the chamber with toxic, infectious or radioactive substances.</b>  <b>Danger of intoxication.</b>  <b>Danger of infection.</b></p> <ul style="list-style-type: none"> <li>⊘ NEVER take a chamber contaminated with toxic substances or sources of infection for recycling according to Directive 2012/19/EU.</li> <li>➤ Prior to disposal, remove all toxic substances and sources of infection from the chamber.</li> <li>➤ A chamber from which all toxic substances or sources of infection cannot be safely removed must be considered as "special" waste according to national law. Dispose of it accordingly.</li> </ul>
---	---


## 23.4 Disposal of the chamber in the member states of the EU except for the Federal Republic of Germany

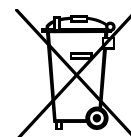According to Annex I of Directive 2012/19/EU of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE), BINDER devices are classified as "monitoring and control instruments" (category 9) only intended for professional use". They must not be disposed of at public collecting points.


The chambers bear the symbol for the marking of electrical and electronic equipment manufactured / placed on the market in the EC after 13 August 2005 and be disposed of in separate collection according to the Directive 2012/19/EU on waste electrical and electronic equipment (WEEE). WEEE marking: crossed-out wheeled bin with solid bar under.





At the end of the device's service life, notify the distributor who sold you the device, who will take back and dispose of the chamber according to the Directive 2012/19/EU on waste electrical and electronic equipment (WEEE).


	<div style="background-color: yellow; text-align: center; padding: 5px;"> <b>CAUTION</b> </div> <p><b>Violation against existing law.</b></p> <ul style="list-style-type: none"> <li>⊘ Do NOT dispose of BINDER devices at public collecting points.</li> <li>➤ Have the device disposed of professionally at a recycling company that is certified according to conversion of the Directive 2012/19/EU into national law. or</li> <li>➤ Instruct the distributor who sold you the device to dispose of it. The agreements apply that were agreed with the distributor when purchasing the chamber (e.g. his general terms of payment and delivery).</li> <li>➤ If your distributor is not able to take back and dispose of the chamber, please contact BINDER service.</li> </ul>
---	--

Certified companies disassemble waste (used) BINDER equipment in primary substances for recycling according to Directive 2012/19/EU. The devices must be free from toxic, infectious or radioactive substances in order to eliminate any health hazards to the employees of the recycling companies.

	<ul style="list-style-type: none"> <li>• Prior to handing the chamber over to a recycling company, it is the user's responsibility that it is free from toxic, infectious or radioactive substances.</li> <li>• Prior to disposal, clean all introduced or residual toxic substances from the chamber.</li> <li>• Prior to disposal, disinfect the chamber from all sources of infection. Be aware that sources of infection may also be located outside the inner chamber.</li> <li>• If you cannot safely remove all sources of infection and toxic substances from the chamber, dispose of it as special waste according to national law.</li> <li>• Fill out the contamination clearance certificate (chap. 28) and enclose it with the chamber.</li> </ul>
---	---

	<div style="background-color: orange; text-align: center; padding: 5px;">  <b>WARNING</b> </div> <p><b>Contamination of the chamber with toxic, infectious or radioactive substances.</b>  <b>Danger of intoxication.</b>  <b>Danger of infection.</b></p> <ul style="list-style-type: none"> <li>Ø NEVER take a chamber contaminated with toxic substances or sources of infection for recycling according to Directive 2012/19/EU.</li> <li>➤ Prior to disposal, remove all toxic substances and sources of infection from the chamber.</li> <li>➤ A chamber from which all toxic substances or sources of infection cannot be safely removed must be considered as "special" waste according to national law. Dispose of it accordingly.</li> </ul>
---	---

### 23.5 Disposal of the chamber in non-member states of the EU

	<div style="background-color: yellow; text-align: center; padding: 5px;"> <b>CAUTION</b> </div> <p><b>Alteration of the environment.</b></p> <ul style="list-style-type: none"> <li>➤ For final decommissioning and disposal of the chamber, please contact BINDER service.</li> <li>➤ Follow the statutory regulations for appropriate, environmentally friendly disposal.</li> </ul>
---	--

The main board of the chamber includes a lithium cell. Please dispose of it according to national regulations.

## 24. Troubleshooting



Only qualified service personnel authorized by BINDER must perform repair.  
Repaired chambers must comply with the BINDER quality standards.

### 24.1 General

Fault description	Possible cause	Required measures
Chamber without function.	No power supply.	Check connection to power supply.
	Wrong voltage.	Check power supply for voltage of 100-120V or 200-240V.
	Chamber fuse has responded.	Check chamber fuse and replace it if appropriate. If it responds again, contact BINDER service.
	Controller defective.	Contact BINDER service.
	Nominal temperature exceeded by 10 °C due to chamber failure. Over temperature protective device (class 1) responds.	


### 24.2 Heating

Fault description	Possible cause	Required measures
<b>Excess temperature.</b> Having reached the setpoint, the temperature exceeds the setpoint by more than the set tolerance range value (more than 10 min.) Alarm message "Temp. range"	Site of installation too warm. Difference between the set temperature and the ambient temperature too low.	Acknowledge the alarm on the controller. Difference between the set temperature and the ambient temperature must be at least 7 °C / 12.6 °F.
	Controller defective.	Acknowledge the alarm on the controller. Contact BINDER service.
	Semiconductor relay defective	
	Temperature controller not adjusted.	Acknowledge the alarm on the controller. Calibrate and adjust the temperature controller.
<b>Excess temperature.</b> Overtemperature safety controller class 3.1 responds. Alarm message "Safety controller"	Safety controller setpoint value exceeded.	Acknowledge the alarm on the controller. Check setting of temperature setpoint and of the safety controller class 3.1 setpoint. If appropriate, select suitable value.
	Too much external heat load.	Acknowledge the alarm on the controller. Reduce heat load.
	Controller defective.	Acknowledge the alarm on the controller. Contact BINDER service.
	Semiconductor relay defective	
	Safety controller defective.	

Fault description	Possible cause	Required measures
<b>Too low temperature.</b> Having reached the setpoint, the temperature falls below the setpoint by more than the set tolerance range value (more than 10 min.) <b>or</b> the temperature doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door. Alarm message "Temp. range"	Chamber door not properly closed.	Acknowledge the alarm on the controller. Completely close chamber door.
	Door gaskets defective.	Acknowledge the alarm on the controller. Replace door gasket.
	Controller defective.	Acknowledge the alarm on the controller. Check the function of the temperature controller.
	Door gasket defective.	Replace door gasket.
	Temperature controller not adjusted.	Calibrate and adjust the temperature controller.
Chamber doesn't heat up. Alarm message "Safety controller"	Inner chamber temperature has reached the safety controller setpoint. Safety controller set too low.	Acknowledge the alarm on the controller. Check temperature setpoint setting. If appropriate, select suitable safety controller setpoint (chap. 12.2).
	Safety controller defective.	Acknowledge the alarm on the controller. Contact BINDER service.
Chamber doesn't heat up.	Pt 100 sensor defective.	Contact BINDER service.
	Heating element defective.	
	Semiconductor relay defective.	
	Controller defective.	
Very long heating-up times.	Chamber fully loaded.	Load the chamber less or consider longer heating-up times.
The displayed actual temperature value deviates largely compared with a reference method..	Temperature control not adjusted.	Calibrate and adjust the temperature control.
	Temperature sensor defective.	Transfer the cultures to another incubator and contact BINDER Service.

### 24.3 Gas cylinder pressure too low

The alarm messages indicate that the pressure in the supply lines of CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control) and N<sub>2</sub> (chamber with O<sub>2</sub> control) has dropped below 0.3 bar. If no gas cylinder changer is installed, the gas cylinder must be replaced. Observe the safety instructions and the maximum outlet pressure (chap. 4.4).

	Prerequisite for pressure alarm messages for O <sub>2</sub> and N <sub>2</sub> (chamber with O <sub>2</sub> control) is that the O <sub>2</sub> / N <sub>2</sub> control with the respective alarms are activated (chap. 6.8, 7.3).
---	---

**Note:** The recovery times of the gas concentrations inside the chamber following door opening, which are indicated in the technical data (chap. 25.4), refer to a connection pressure of 2.0 bar / 29 *psi*. Decreasing supply pressure down to the shift point (alarm point) of 0.3 bar / 4.4 *psi* results in longer recovery times. Check the pressure displays of your gas supply. If very short recovery times are required or the door is opened frequently, replace the gas cylinders promptly when the pressure decreases below 2.0 bar / 29 *psi*.



Fault description	Possible cause	Required measures
Low CO <sub>2</sub> outlet pressure (< 0.3 bar) Alarm message "CO <sub>2</sub> pressure"	CO <sub>2</sub> cylinder is not connected correctly.	Acknowledge the alarm on the controller. Correctly connect the gas cylinder.
	Connected gas cylinder is closed.	Acknowledge the alarm on the controller. Open the gas cylinder.
	Connected gas cylinder is empty.	Acknowledge the alarm on the controller. Replace the gas cylinder. Observe the safety instructions and the maximum outlet pressure (chap. 4.4).
	Gas hose is dirty or obstructed.	Acknowledge the alarm on the controller. Turn off the gas supply and remove the gas connections. Check the tube system for dirt accumulation or obstruction, clean or replace it.
	Controller malfunction.	Acknowledge the alarm on the controller. If necessary, contact BINDER Service.
	Pressure sensor system defective.	Acknowledge the alarm on the controller. Contact BINDER Service.
<i>Chamber with O<sub>2</sub> control:</i> Low O <sub>2</sub> outlet pressure (< 0.3 bar). Alarm message "O <sub>2</sub> pressure"	O <sub>2</sub> cylinder is not connected correctly.	Acknowledge the alarm on the controller. Correctly connect the gas cylinder.
	Connected gas cylinder is closed.	Acknowledge the alarm on the controller. Open the gas cylinder.
	Connected gas cylinder is empty.	Acknowledge the alarm on the controller. Replace the gas cylinder. Observe the safety instructions and the maximum outlet pressure (chap. 4.4).
	Gas hose is dirty or obstructed.	Acknowledge the alarm on the controller. Turn off the gas supply and remove the gas connections. Check the tube system for dirt accumulation or obstruction, clean or replace it.
	Controller malfunction	Acknowledge the alarm on the controller. If necessary, contact BINDER Service.
	Pressure sensor system defective.	Acknowledge the alarm on the controller. Contact BINDER Service.



Fault description	Possible cause	Required measures
<b>Chamber with O<sub>2</sub> control:</b> Low N <sub>2</sub> outlet pressure (< 0.3 bar). Alarm message "N <sub>2</sub> pressure"	N <sub>2</sub> cylinder is not connected correctly.	Acknowledge the alarm on the controller. Correctly connect the gas cylinder.
	Connected gas cylinder is closed.	Acknowledge the alarm on the controller. Open the gas cylinder.
	Connected gas cylinder is empty.	Acknowledge the alarm on the controller. Replace the gas cylinder. Observe the safety instructions and the maximum outlet pressure (chap. 4.4).
	Gas hose is dirty or obstructed.	Acknowledge the alarm on the controller. Turn off the gas supply and remove the gas connections. Check the tube system for dirt accumulation or obstruction, clean or replace it.
	Controller malfunction	Acknowledge the alarm on the controller. If necessary, contact BINDER Service.
	Pressure sensor system defective.	Acknowledge the alarm on the controller. Contact BINDER Service.

## 24.4 Gas concentration

Fault description	Possible cause	Required measures
<b>Excess CO<sub>2</sub> concentration.</b> Having reached the setpoint, CO <sub>2</sub> exceeds the setpoint by more than the set tolerance range value (more than 10 min.) Alarm message "CO <sub>2</sub> range"	Temporary disturbance of the CO <sub>2</sub> control.	Acknowledge the alarm on the controller. Turn off the chamber. Open the chamber doors for approx. 5 minutes. Observe the general information for safe handling of CO <sub>2</sub> (chap. 1.6). Turn on the chamber again. Upon renewed alarm, contact BINDER service.
	CO <sub>2</sub> sensor defective.	Acknowledge the alarm on the controller. Contact BINDER Service.
<b>Chamber with O<sub>2</sub> control:</b> <b>Excess O<sub>2</sub> concentration.</b> Having reached the setpoint, O <sub>2</sub> exceeds the setpoint by more than the set tolerance range value (more than 10 min.) (O <sub>2</sub> control option) Alarm message "O <sub>2</sub> range"	Temporary disturbance of the O <sub>2</sub> control.	Acknowledge the alarm on the controller. Turn off the chamber. Open the chamber doors for approx. 5 minutes. Observe the general information for safe handling of oxygen (chap. 1.6). Prevent oxygen enrichment in the ambience of the chamber. Turn on the chamber again. Upon renewed alarm, contact BINDER service.
	O <sub>2</sub> sensor defective.	Acknowledge the alarm on the controller. Contact BINDER Service.
Recovery time (up to 5 vol.-% CO <sub>2</sub> ) after doors were open for 2 minutes is < 2 minutes	CO <sub>2</sub> sensor system defective.	Contact BINDER Service.

Fault description	Possible cause	Required measures
<b>Too low CO<sub>2</sub> concentration.</b> Having reached the setpoint, CO <sub>2</sub> falls below the setpoint by more than the set tolerance range value (more than 10 min.) <b>or</b> CO <sub>2</sub> doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door. Alarm message "CO <sub>2</sub> range"	Temporary disturbance of the CO <sub>2</sub> control.	Acknowledge the alarm on the controller. Turn off the chamber. After approx. 5 minutes turn on the chamber again. Upon renewed alarm, contact BINDER service.
	CO <sub>2</sub> sensor defective.	Acknowledge the alarm on the controller. Contact BINDER Service.
<i>Chamber with O<sub>2</sub> control:</i> <b>Too low O<sub>2</sub> concentration.</b> Having reached the setpoint, O <sub>2</sub> falls below the setpoint by more than the set tolerance range value (more than 10 min.) <b>or</b> O <sub>2</sub> doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door. Alarm message "O <sub>2</sub> range".	Temporary disturbance of the O <sub>2</sub> control.	Acknowledge the alarm on the controller. Turn off the chamber. After approx. 5 minutes turn on the chamber again. Upon renewed alarm, contact BINDER service.
	O <sub>2</sub> sensor defective.	Acknowledge the alarm on the controller. Contact BINDER Service.
<b>Too low gas concentration.</b> The concentration of CO <sub>2</sub> or O <sub>2</sub> (chamber with O <sub>2</sub> control) does not reach the adjusted set value.	Door gaskets defective.	Replace door gaskets
	Doors not closed properly.	Close door properly.
	Connected gas cylinder is empty or not opened.	Open or replace gas cylinder.
	Gas cylinder is not connected correctly.	Correctly connect the gas cylinder.
	Gas hose is dirty or obstructed.	Check the tube system for dirt accumulation or obstruction, clean or replace it.
Recovery time (up to 5 vol.-% CO <sub>2</sub> ) after doors were open for 2 minutes is > 10 minutes.	Obstructed gas supply.	Check gas supply (cylinder, connections, hose system).
	Insufficient CO <sub>2</sub> input pressure	
Unusually high gas consumption.	Door gaskets defective.	Replace door gaskets
	Gas sensor not adjusted.	Calibrate the sensor.
	Gas sensor defective.	Contact BINDER Service
	Gas fine filter not connected correctly.	
The displayed actual value of CO <sub>2</sub> deviates largely compared with a reference method.. The pH indicator of the cell culture medium changes its normal color.	CO <sub>2</sub> control not adjusted.	Calibrate and adjust CO <sub>2</sub> control.
	CO <sub>2</sub> sensor defective.	Transfer the cultures to another incubator and contact BINDER Service.
<i>Chamber with O<sub>2</sub> control:</i> The displayed actual value of O <sub>2</sub> (option) deviates largely compared with a reference method.	O <sub>2</sub> control not adjusted.	Calibrate and adjust O <sub>2</sub> control.
	O <sub>2</sub> sensor defective.	Transfer the cultures to another incubator and contact BINDER Service.
No gas control after hot-air sterilization.	Sterilization cycle still running.	Terminate hot-air sterilization.
	Information message not confirmed.	Confirm the information message with "Ok".
	Inner chamber temperature above 80 °C	Let the inner chamber temperature decrease below 80 °C.

## 24.5 Sterilization

Indication	Possible cause	Required measures
The sterilization is running. Alarm message "Sterilization"	Sterilization cycle is running	Acknowledge the alarm on the controller. Do not open the door.

## 24.6 Humidity (chamber without active humidification)

Fault description	Possible cause	Required measures
Condensations inside the chamber.	Permadyr <sup>TM</sup> pan filled with water when the chamber is not operating.	Empty Permadyr <sup>TM</sup> pan when the chamber is not operating.
	Door gaskets defective.	Replace door gaskets
	Doors not closed properly.	Close door properly.
	Humidity set too high.	Reduce humidity (chap. 6.9)
Condensation on the doors or on the optional divided inner door, gas-proof.	Chamber placed on very cold floor.	Place the chamber on a BINDER stand to increase the distance to the floor.
	Door gaskets defective.	Replace door gaskets
	Doors not closed properly.	Close door properly.
	Humidity set too high.	Reduce humidity (chap. 6.9)
	Door heating set too low.	Contact BINDER Service to adjust the door heating
Too low humidity inside.	Humidity set too low.	Increase humidity (chap. 6.9)
	Chamber door not properly closed.	Completely close chamber door.
	Door gasket defective.	Replace door gasket.
	Permadyr <sup>TM</sup> pan empty.	Fill the outer Permadyr <sup>TM</sup> pan with water up to the marking on the inner pan with distilled, sterile water. The pan must have thorough contact to the bottom of the inner chamber, see chap. 4.2.

## 24.7 Humidity (chamber with active humidification)

Fault description	Possible cause	Required measures
Humidity fluctuation: Control accuracy of +/- 3 % r.h. is not reached.	Door gasket defective.	Replace door gasket.
	Door opened very frequently.	Open doors less frequently.
Humidity fluctuation, together with temperature fluctuation > 1 °C with a set-point approx. 3 °C above ambient temperature.	Place of installation too hot.	Select cooler place of installation or contact BINDER service.

## 24.8 Controller

Fault description	Possible cause	Required measures
No chamber function (dark display).	Display in standby mode.	Press on touchscreen.
	Main power switch turned off.	Turn on the main power switch.
Menu functions not available.	Menu functions not available with current authorization level.	Log in with the required higher authorization. or contact BINDER service to obtain an activation code (chap. 13.6).
No access to controller	Password incorrect.	Contact BINDER service.
Chart recorder function: measured-value memory cleared; information lost.	New setting of storage rate or scaling (minimum and/or maximum) (chap. 16.2).	Change the storage rate or scaling ONLY if the previously registered data are no longer required.
Controller does not equilibrate to setpoints entered in Fixed value operation mode	Controller is not in Fixed value operation mode.	Change to Fixed value operation mode.
	CO <sub>2</sub> control is deactivated.	Activate CO <sub>2</sub> control (chap. 6.8).
	<i>Chamber with O<sub>2</sub> control:</i> O <sub>2</sub> control is deactivated.	Activate O <sub>2</sub> control (chap. 6.8).
	<i>Chamber with active humidification:</i> Humidity control is deactivated.	Activate humidity control (chap. 6.8).
Controller does not equilibrate to program set-points.	Controller is not in program operation mode, or program delay time is running.	Start the program again. If appropriate, wait for the program delay time.
Program duration longer than programmed.	Tolerances have been programmed.	For rapid transition phases, do NOT program tolerance limits in order to permit maximum heating, refrigerating, or humidification speed.
Program keeps the last program setpoint constant while in setting "ramp".	Program line with setting "ramp" is incomplete.	When programming with setting "ramp", define the end value of the desired cycle by adding an additional section with a section time of at least one second.
Ramp temperature transitions are only realized as steps.	Setting "step" has been selected.	Select setting "ramp".
<i>Chamber with active humidification:</i> Humidity alarm when operating without water supply.	Humidity control is activated.	Deactivate humidity control (chap. 6.8).
Acknowledging the alarm does not cancel the alarm state.	Cause of alarm persists.	Remove cause of alarm. If the alarm state continues, contact BINDER service.
Alarm message: - - - - or <-<-< or >->->	Sensor rupture between sensor and controller or Pt 100 sensor defective.	Contact BINDER service.

## 24.9 Open door

Indication	Possible cause	Required measures
Outer door open (longer than the selected alarm delay time) Alarm message "Door open".	Outer door open or not properly closed.	Acknowledge the alarm on the controller. Close the door. The triggered zero-voltage relay alarm contact switches back.

## 25. Technical description

### 25.1 Factory calibration and adjustment

This chamber was calibrated and adjusted in the factory. Calibration and adjustment were performed using standardized test instructions, according to the QM DIN EN ISO 9001 system applied by BINDER (certified since December 1996 by TÜV CERT). All test equipment used is subject to the administration of measurement and test equipment that is also a constituent part of the BINDER QM DIN EN ISO 9001 systems. They are controlled and calibrated to a DKD-Standard at regular intervals.

A record of this calibration and adjustment is part of the BINDER test certificate of the chamber.

#### Adjustment in factory:

- **Temperature:** 37 °C / 98.6°F measured in the center of the usable volume
- **CO<sub>2</sub>:** 0 vol.-% CO<sub>2</sub> (100 vol.-% N<sub>2</sub>) and 5 vol.-% CO<sub>2</sub> (analyzed test gas directly exposed to the sensor head)
- **O<sub>2</sub>** (chamber with O<sub>2</sub> control): 0 vol.-% O<sub>2</sub> (100 vol.-% N<sub>2</sub>, analyzed test gas directly exposed to the sensor head) and 20.9 vol.-% O<sub>2</sub> (ambient air).
- **Humidity** (chamber with active humidification): 93 % r.h.



Repeated calibrations are recommended in periods of 12 months.

BINDER service uses an electronic measuring and display device for temperature traceable to an acknowledged standards/calibration institution (DKD or PTB for Germany) with valid calibration certificate.

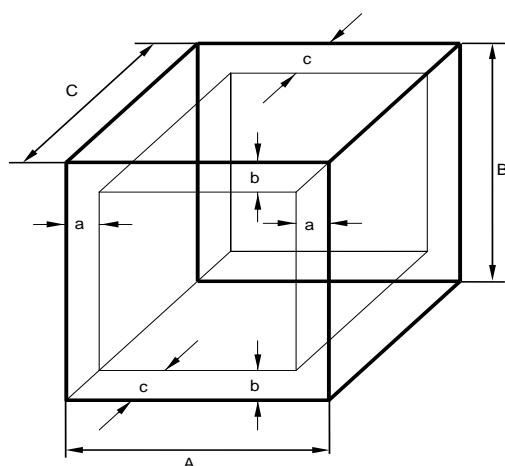
Test gases with an analyzed concentration serve to calibrate the sensor systems for CO<sub>2</sub> and O<sub>2</sub> (chamber with O<sub>2</sub> control). The sensor heads are exposed directly to the test gas.

### 25.2 Over current protection

The CB is protected by a chamber-protection against over current, accessible from the outside. It is located at the rear of the chamber below the strain relief of the power cord. The fuse holder is equipped with a fuse clip 5mm x 20 mm (cUL-Version 6.3 x32 mm). Replace this fuse only with a substitute of the same ratings. Refer to the technical data of the respective device type. If this fuse is blown, please inform an electronic engineer or BINDER service.

## 25.3 Definition of usable volume

The usable volume illustrated below is calculated as follows:



A, B, C = internal dimensions (W, H, D)

a, b, c = wall separation

$$a = 0.1 \cdot A$$

$$b = 0.1 \cdot B$$

$$c = 0.1 \cdot C$$

$$V_{USE} = (A - 2 \cdot a) \cdot (B - 2 \cdot b) \cdot (C - 2 \cdot c)$$

Figure 30: Determination of the usable volume

The technical data refers to the defined usable volume.



Do NOT place samples outside this usable volume.

Do NOT load this volume by more than half to enable sufficient airflow inside the chamber.

Do NOT divide the usable volume into separate parts with large area samples.

Do NOT place samples too close to each other in order to permit circulation between them and thus obtain a homogenous distribution of temperature, CO<sub>2</sub>, O<sub>2</sub> (chamber with O<sub>2</sub> control), and humidity.

## 25.4 CB technical data

Chamber size		56	170	260
<b>Exterior dimensions</b>				
Width, net	mm / inch	520 / 20.47	680 / 26.77	740 / 29.13
Height including feet	mm / inch	675 / 26.57	870 / 34.25	1025 / 40.35
Depth, net	mm / inch	550 / 21.65	715 / 28.15	775 / 30.51
Depth plus door handle, I-triangle	mm / inch	55 / 2.16	55 / 2.16	55 / 2.16
Depth plus power connection and gas connection	mm / inch	10 / 0.39	10 / 0.39	10 / 0.39
Wall clearance rear (minimum)	mm / inch	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (minimum)	mm / inch	50 / 1.97	50 / 1.97	50 / 1.97
<b>Doors</b>				
Quantity of outer doors		1	1	1
Quantity of inner doors (inner glass door or divided inner door)		1	1	1
Quantity of individual inner glass doors (with divided inner door)		4	6	8
<b>Interior dimensions</b>				
Width	mm / inch	400 / 15.75	560 / 22.05	620 / 24.41
Height	mm / inch	400 / 15.75	600 / 23.62	750 / 29.53
Depth	mm / inch	330 / 12.99	500 / 19.69	560 / 22.05
Interior volume	l / c.ft.	53 / 1.9	170 / 6.0	260 / 9.2

Chamber size		56	170	260	
Shelves					
Number of shelves, regular		2	3	4	
Number of shelves, max.		3	6	8	
Size of shelf (external) Width x Depth		mm x mm inch x inch	396 x 289 15.59 x 11.38	551 x 442 21.69 x 17.40	614 x 502 24.17 x 19.76
Maximum load per shelf		Kg / lbs.	10 / 22	10 / 22	10 / 22
Permitted total load		Kg / lbs	30 / 66	30 / 66	30 / 66
Weight					
Weight (empty)		Kg / lbs	***	***	***
Temperature data					
Temperature range	from ... degree above ambient	°C / °F	***	4 / 7.2	***
	up to	°C / °F	60 / 140	60 / 140	60 / 140
Temperature range, chamber with O <sub>2</sub> control	from ... degree above ambient	°C / °F	***	6 / 10.8	***
	up to	°C / °F	60 / 140	60 / 140	60 / 140
Temperature fluctuation at 37 °C / 98.6 °F		± K	0.1	0.1	0.1
Temperature uniformity (variation) at 37 °C/ 98.6 °F		± K	***	0.3	***
Recovery time after door was opened for 30 sec at 37 °C / 98.6 °F		minutes	***	6	***
CO <sub>2</sub> data					
CO <sub>2</sub> range		vol.-% CO <sub>2</sub>	0 to 20	0 to 20	0 to 20
Setting accuracy		vol.-% CO <sub>2</sub>	0.1	0.1	0.1
CO <sub>2</sub> fluctuation		vol.-% CO <sub>2</sub>	***	0.1	***
Recovery time after door was opened for 30 sec at 5 vol.-% CO <sub>2</sub>		minutes	***	***	***
CO <sub>2</sub> measurement			IR	IR	IR
Connection hose nozzle DN6 for CO <sub>2</sub> connection to the chamber for hose with internal diameter		mm / inch	6 / 0.24	6 / 0.24	6 / 0.24
Humidity data					
Humidity (without active humidification)		% r.h.	90 to 95	90 to 95	90 to 95
Humidity (chamber with active humidification)		% r.h.	60 to 93	60 to 93	60 to 93
Recovery time after door was opened for 30 sec at 93 % r.h. (chamber with active humidification)		minutes	***	***	***
O <sub>2</sub> Data (chamber with O <sub>2</sub> control)					
O <sub>2</sub> range (regular hypoxic control range)		vol.-% O <sub>2</sub>	0.2 to 20	0.2 to 20	0.2 to 20
O <sub>2</sub> range (optional alternative control range)		vol.-% O <sub>2</sub>	10 to 95	10 to 95	10 to 95
Setting accuracy with inlet pressure 2 bar		vol.-% O <sub>2</sub>	0.1	0.1	0.1
Recovery time after door was opened for 30 sec at 5 vol.-% O <sub>2</sub>		minutes	***	10	***
O <sub>2</sub> measurement			ZrO <sub>2</sub>	ZrO <sub>2</sub>	ZrO <sub>2</sub>
Connection hose nozzle DN6 for O <sub>2</sub> /N <sub>2</sub> connection to the chamber for hose with internal diameter		mm / inch	6 / 0.24	6 / 0.24	6 / 0.24




Chamber size			56	170	260
Electrical data					
IP system of protection acc. to EN 60529		IP	20	20	20
Nominal voltage (+/-10 %)	at 50 Hz power frequency	V	200-230	200-230	200-230
	at 60 Hz power frequency	V	200-230	200-230	200-230
Current type			1N~	1N~	1N~
Nominal power		kW	***	1.40	***
Power plug (IEC connector plug)			Grounded plug IEC 7/7		
Installation category acc. to IEC 61010-1			II	II	II
Pollution degree acc. to IEC 61010-1			2	2	2
Chamber fuse 4x20mm / 250V / time-lag (T), external		Amp	10	10	10
Additional temperature fuse class 1 (DIN 12880), internal		°C	240	240	240
Different electrical data for CB-UL constructed for the USA and Canada					
Nominal voltage (±10 %)	at 50 Hz power frequency	V	100-120	100-120	100-120
	at 60 Hz power frequency	V	100-120	100-120	100-120
Current type			1N~	1N~	1N~
Power plug		NEMA	5-15P	5-20P	5-20P
Chamber fuse 6.3x32 mm / 250V / time-lag (T), external		Amp	12.5	12.5	12.5
Environment-specific data					
Energy consumption at 37 °C/ 98.6 °F		Wh/h	***	60	***

\*\*\* Data not yet determined


**Note:** The recovery times of the gas concentrations inside the chamber after the door is opened coincide with a connection pressure of 2.0 bar / 29 psi. Decreasing supply pressure results in longer recovery times.

All technical data is specified for unloaded chambers with standard equipment at an ambient temperature of 22 +/- 3 °C / 71.6 +/- 5.4 °F and a power supply voltage fluctuation of +/-10%. Technical data is determined in accordance to BINDER Factory Standard Part 2:2015 and DIN 12880:2007.

**All indications are average values, typical for chambers produced in series. We reserve the right to change technical specifications at any time.**

	<i>Chamber with active humidification:</i> Bringing a source of humidity into the inner chamber will affect the minimum humidity specification
---	--

## 25.5 Equipment and Options (extract)

	To operate the chamber, use only original BINDER accessories or accessories / components from third-party suppliers authorized by BINDER. The user is responsible for any risk arising from using unauthorized accessories.
---	---


Regular equipment
Multifunction touchscreen controller MB2 for temperature, CO <sub>2</sub> , O <sub>2</sub> (chamber with O <sub>2</sub> control) and humidity (chamber with active humidification)
CO <sub>2</sub> infra-red absorption measuring system
Silicone jacket heating
Hot-air auto sterilization at 187.5 °C / 369.5 °F
Gas mixing head



<b>Regular equipment (continued)</b>
Permady™ humidity system
Weldless deep-drawn polished inner chamber made of stainless steel 1.4301/V2A, polished
FailSafe – Electronic error auto-diagnosis system with zero-voltage relay output
Zero-voltage relay alarm output with DIN socket (6 poles), DIN plug included
Lockable door
Tightly closing inner glass door
3 perforated shelves, stainless steel 1.4301/V2A
Overtemperature safety controller class 3.1 acc. to DIN 12880:2007
<i>Chamber with active humidification:</i> Microprocessor controlled humidifying system
<i>Chamber with O<sub>2</sub> control:</i> O <sub>2</sub> control in the hypoxic control range 0.2 to 20 vol. % O <sub>2</sub> .
Ethernet interface for computer communication

<b>Options / accessories</b>
Perforated shelf ,stainless steel
<i>Chamber with O<sub>2</sub> control:</i> Alternative control range 10 to 95 vol. % O <sub>2</sub> (available via BINDER INDIVIDUAL Customized Solutions)
Freshwater bag (0.9 liters / 0.32 cu.ft.)
Divided inner door, gas-proof, stainless steel, 4 times (CB 56), 6 times (CB 170), 8 times (CB 260)
Shelves for divided inner door, gas-proof, stainless steel
Stacking stand with castors lockable by breaks (CB 170 / CB 210)
Stacking adapter for direct, thermally decoupled stacking
Base with castors
Stacking adapter for combination C 150 on top of CB 170
Silicone access ports closable with 2 silicone plugs
BINDER Gas Supply Service: External CO <sub>2</sub> bottle changer CO <sub>2</sub> , O <sub>2</sub> or N <sub>2</sub>
Connection kit for CO <sub>2</sub> , O <sub>2</sub> or N <sub>2</sub> cylinder
Water tight internal socket 230 V (max. 3 A), with turn-off switch (available via BINDER INDIVIDUAL Customized Solutions)
Analog output temperature and CO <sub>2</sub> 4-20 mA 4-20mA with DIN socket (6 poles), DIN plug included
BINDER Data Logger kit for temperature: T 220 (chamber values), and for temperature / humidity: TH 70 (chamber values) or TH 70/70 (chamber and ambient values)
BINDER Pure Aqua Service
Cartridge for BINDER Pure Aqua Service
Cleaning kit (neutral cleaning agent, disinfection spray and lint-free disposable wipes, gloves and goggles)
Intelligent Fail-Safe CO <sub>2</sub> monitoring function
CELLROLL roller system with set of connection cables for extra-low voltage access port
Access port (8-pin) for extra-low voltage (ELV) with LEMO socket (can be covered) and LEMO connector, maximum power rating 24VAC/DC – 2.5 Amp
Data Logger Kit T 260
Qualification folder
Calibration of temperature including certificate
Calibration of CO <sub>2</sub> including certificate
Spatial temperature measurement including certificate
Spatial temperature measurement acc. to DIN 12880 including certificate
<i>Chamber with O<sub>2</sub> control:</i> Calibration of O <sub>2</sub> including certificate
<i>Chamber with active humidification:</i> Calibration of humidity including certificate
<i>Chamber with active humidification:</i> Spatial temperature and humidity measurement including certificate

## 25.6 Accessories and spare parts (extract)

	<p>BINDER GmbH is responsible for the safety features of the chamber only, provided skilled electricians or qualified personnel authorized by BINDER perform all maintenance and repair, and if components relating to chamber safety are replaced in the event of failure with original spare parts. The user is responsible for any risks arising from using unauthorized accessories/components.</p>
---	---

Description	Art. no.
Perforated shelf, stainless steel	6004-0137
Divided shelves (1 level) for divided inner door, stainless steel	6004-0226
Door gasket for outer chamber door	6005-0275
Door gasket for inner glass door or for divided inner door	6005-0508
Door gasket for individual glass door of divided inner door	6005-0293
Individual glass door of divided inner door, complete	8010-0111
Stacking stand with castors lockable by breaks	9051-0020
Stacking adapter for direct, thermally decoupled stacking	9051-0026
Base with castors	9051-0028
Stacking adapter for combination C 150/160/170 on top of CB 170	9051-0027
Flat stacking adapter	9051-0035
CELLROLL roller system with set of connection cables and extra-low voltage access port	8012-0571
Permadyr™ water pan	4022-0081
Plug for optional silicon access port d30	6016-0035
Gas cylinder connection kit for CO <sub>2</sub>	8012-0014
Gas cylinder connection kit for O <sub>2</sub>	8012-0015
Gas cylinder connection kit for N <sub>2</sub>	8012-0016
Chamber fuse 5x20mm 250 V 10 A time-lag (T)	5006-0088
Chamber fuse 6.3x32mm 250V 12.5 A time-lag (T) for CUL version only	5006-0096
CO <sub>2</sub> sensor	5002-0078
O <sub>2</sub> sensor for chamber with regular equipment: hypoxic control range 0.2 up to 20 vol.-% O <sub>2</sub>	5002-0084
O <sub>2</sub> sensor for chamber with O <sub>2</sub> control with optional alternative control range 10 up to 95 vol.-% O <sub>2</sub>	5002-0085
Humidity sensor for chamber with active humidification	5002-0077
Gas fine filter	8009-0369
Fresh water bag (0.9 liters / 0.32 cu.ft.)	6002-0626
BINDER Pure Aqua Service	8012-0759
Cartridge for BINDER Pure Aqua Service	6011-0165
Cleaning kit (neutral cleaning agent, disinfection spray and lint-free disposable wipes, protective gloves and goggles)	8012-0503
Neutral cleaning agent, 1 kg	1002-0016

Validation service	Art. no.
Qualification folder IQ-OQ	8012-0876
Qualification folder IQ-OQ-PQ	8012-0963
Execution of IQ-OQ	DL410200
Execution of IQ-OQ-PQ	DL440500

Calibration service	Art. no.
Calibration of temperature including certificate	8012-1132
Calibration of temperature and CO <sub>2</sub> including certificate	8012-1235
Calibration of O <sub>2</sub> including certificate (chamber with O <sub>2</sub> control)	8012-0229
Spatial temperature measurement including certificate (15-18 measuring points)	8012-1571
Spatial temperature measurement including certificate (9 measuring points)	8012-1550
Spatial temperature measurement including certificate (27 measuring points)	8012-1592

For information on components not listed here, please contact BINDER Service.

## 25.7 Important conversion data for non-SI units

1 ft = 0.305 m

1 m = 100 cm = 3.28 ft = 39.37 inch

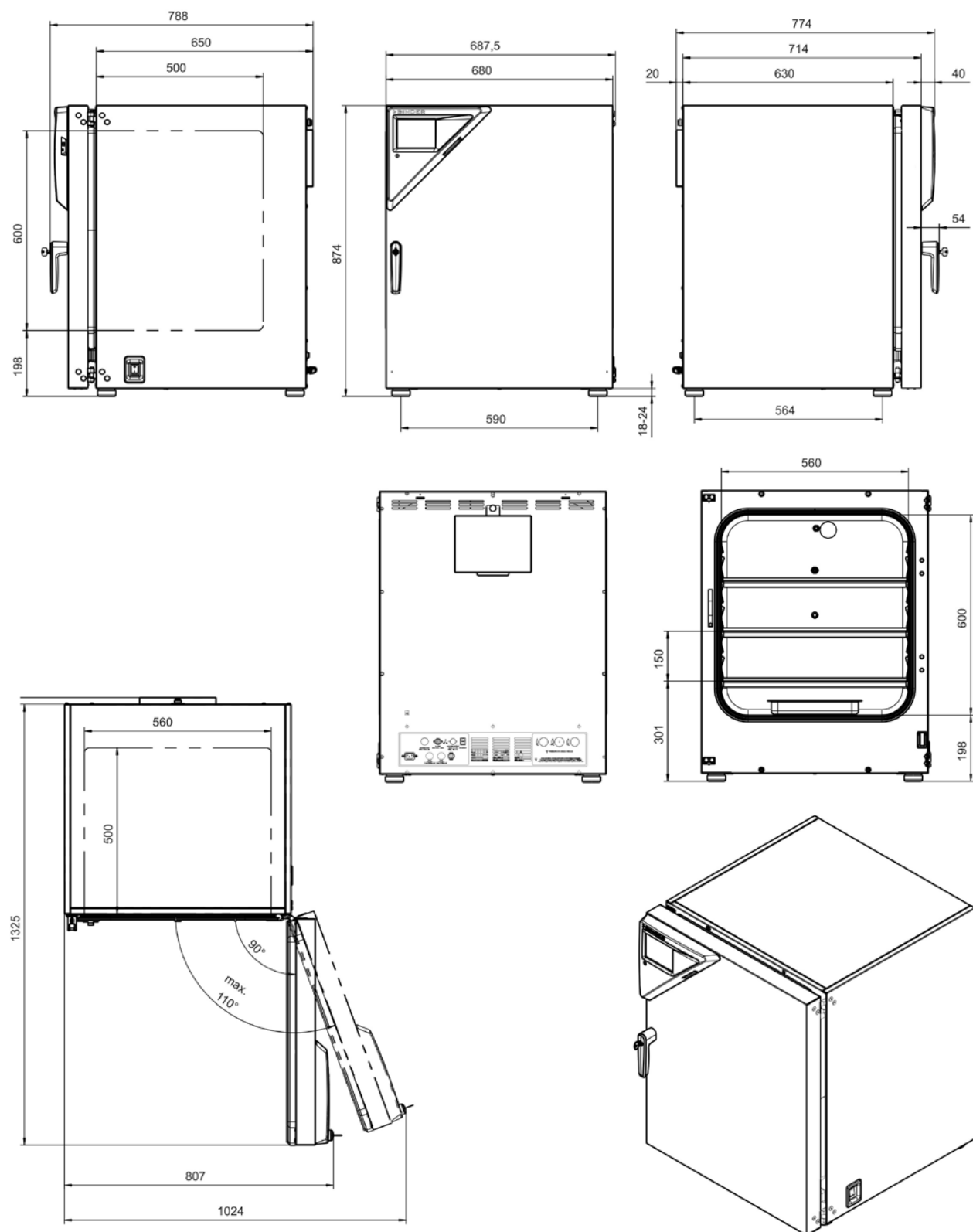
1 mbar = 0.0145 psi

## 25.8 Conversion table for gas inlet pressures, bar – psi

bar	psi	bar	psi	bar	psi
1	14.5	3	43.5	5	72.5
1.5	21.7	3.5	50.7	5.5	79.7
2	29.0	4	58.0	6	87.0
2.5	36.3	4.5	65.2		

## 25.9 Dimensions


### CB 170




(Dimensions in mm)

## 26. Certificates and declarations of conformity

### 26.1 EU Declaration of conformity

  
 Best conditions for your success

	<b>EU-Konformitätserklärung / EU Declaration of Conformity / Déclaration de conformité          UE / Declaración de conformidad UE / Dichiarazione di conformità UE / Декларация          соответствия EU</b>
---	---

Hersteller / Manufacturer / Fabricant / Fabricante / Fabbicante / Производитель	BINDER GmbH
Anschrift / Address / Adresse / Dirección / Indirizzo / Адрес	Im Mittleren Ösch 5, 78532 Tuttlingen, Germany
Produkt / Product / Produit / Producto / Prodotto / Продукт	CO <sub>2</sub> -Inkubatoren CO <sub>2</sub> Incubators Incubateurs à CO <sub>2</sub> Incubadoras de CO <sub>2</sub> Incubatori a CO <sub>2</sub> CO <sub>2</sub> инкубаторы
Typenbezeichnung / Type / Type / Tipo / Tipo / Тип	CB 170

Das oben beschriebene Produkt ist konform mit folgenden EU-Richtlinien:  
 The product described above is in conformity with the following EU Directives:  
 Le produit décrit ci-dessus est conforme aux directives UE suivantes:  
 El producto descrito arriba cumple con las siguientes directivas de la UE:  
 Il prodotto sopra descritto è conforme alle seguenti direttive UE:  
 Продукты, указанные выше, полностью соответствуют следующим EU руководствам:

- **2014/35/EU**  
 Niederspannungsrichtlinie 2014/35/EU / Low voltage directive 2014/35/EU / Directive basse tension  
 2014/35/UE / Directiva sobre baja tensión 2014/35/UE / Direttiva Bassa tensione 2014/35/UE /  
 Директива по низкому напряжению 2014/35/EU
- **2014/30/EU**  
 EMV-Richtlinie 2014/30/EU / EMC Directive 2014/30/EU / Directive CEM 2014/30/UE / Directiva CEM  
 2014/30/UE / Direttiva EMC 2014/30/UE / Директива ЭМС 2014/30/EU
- **2011/65/EU**  
 RoHS-Richtlinie 2011/65/EU / RoHS Directive 2011/65/EU / Directive RoHS 2011/65/UE / Directiva  
 RoHS 2011/65/UE / Direttiva RoHS 2011/65/UE / Директива RoHS 2011/65/EU

Die oben beschriebenen Produkte tragen entsprechend die Kennzeichnung CE.  
 The products described above, corresponding to this, bear the CE-mark.  
 Les produits décrits ci-dessus, en correspondance, portent l'indication CE.  
 Los productos descritos arriba, en conformidad, llevan la indicación CE.  
 I prodotti sopra descritti, conformi a quanto sopra, portano il marchio CE.  
 Данные продукты в соответствии с изложенным выше маркированы знаком CE.

1 / 2

BINDER GmbH Postfach 102 D-78502 Tuttlingen    Anschrift: BINDER GmbH Im Mittleren Ösch 5 D-78532 Tuttlingen  
 Kontakt: Telefon: +49 (0) 74 62 / 20 05 – 0 | Telefax: +49 (0) 74 62 / 20 05 – 100 | info@binder-world.com | www.binder-world.com  
 Geschäftsführung: Dipl.-Ing. Peter M. Binder | Amtsgericht Stuttgart, HRB 727150 | Sitz der Gesellschaft: Tuttlingen  
 Bankverbindung: Kreissparkasse Tuttlingen Konto-Nr.: 2266 BLZ: 643 500 70 | IBAN-Code: DE05643 500700 000002266 | SWIFT-Code: SOLA DE S1TUT  
 S-Account 2202 611 55 | IBAN-Code: DE7464350070 0220 261155 | SWIFT-Code: SOLA DE S1TUT  
 Deutsche Bank Tuttlingen Konto-Nr.: 2 138 709 BLZ: 653 700 75 | IBAN-Code: DE56653 70075 0213870900 | SWIFT-Code: DEUT DE SS603  
 Altgeräte-Entsorgung gemäß WEEE-Reg.-Nr. DE 37004983

Die oben beschriebenen Produkte sind konform mit folgenden harmonisierten Normen:  
 The products described above are in conformity with the following harmonized standards:  
 Les produits décrits ci-dessus sont conformes aux normes harmonisées suivantes:  
 Los productos descritos arriba cumplen con las siguientes normas:  
 I prodotti sopra descritti sono conformi alle seguenti normative armonizzate:  
 Продукты, указанные выше, полностью соответствуют следующим стандартам:

Sicherheit / Safety / Sécurité / Seguridad / Sicurezza / Нормативы по безопасности
<ul style="list-style-type: none"> <li>• EN 61010-1:2010</li> <li>• EN 61010-2-010:2014</li> </ul>
EMV / EMC / CEM / CEM / EMC / ЭМС
<ul style="list-style-type: none"> <li>• EN 61326-1:2013</li> </ul>
RoHS
<ul style="list-style-type: none"> <li>• EN 50581:2012</li> </ul>

78532 Tuttlingen, 28.05.2018  
 BINDER GmbH



P. M. Binder  
 Geschäftsführender Gesellschafter  
 Managing Director  
 Directeur général  
 Director general  
 Direttore Generale  
 Генеральный Директор



J. Bollaender  
 Leiter F & E  
 Director R & D  
 Chef de service R&D  
 Responsable I & D  
 Direttore R & D  
 Глава департамента R&D

2 / 2

**BINDER GmbH** Postfach 102 D-78502 Tuttlingen **Anschrift:** BINDER GmbH Im Mittleren Ösch 5 D-78532 Tuttlingen  
**Kontakt:** Telefon: +49 (0) 74 62 / 20 05 - 0 | Telefax: +49 (0) 74 62 / 20 05 - 100 | info@binder-world.com | www.binder-world.com  
**Geschäftsführung:** Dipl.-Ing. Peter M. Binder | Amtsgericht Stuttgart, HRB 727150 | Sitz der Gesellschaft: Tuttlingen  
**Bankverbindung:** Kreissparkasse Tuttlingen Konto-Nr.: 2266 BLZ: 643 500 70 | IBAN-Code: DE05643 500700 000002266 | SWIFT-Code: SOLA DE 31TUT  
 S-Account 2202 611 55 | IBAN-Code: DE7464350070 0220 261155 | SWIFT-Code: SOLA DE 31TUT  
 Deutsche Bank Tuttlingen Konto-Nr.: 2 138 709 BLZ: 653 700 75 | IBAN-Code: DE56653 70075 0213870900 | SWIFT-Code: DEUT DE 3303  
 Altgeräte-Entsorgung gemäß WEEE-Reg.-Nr. DE 37004983



## 27. Product registration

# Online Product Registration

Register your BINDER now!

[www.binder-world.com/register](http://www.binder-world.com/register)

The registration is free and takes just a few seconds

Advantages:

- ▶ Short response times if service is needed
- ▶ Fair prices when relocating or installing equipment
- ▶ Calibration as required at no charge in case of recalls
- ▶ Free information on news, product upgrades and accessories

Easy registered in 3 steps:



1. List serial number here:   -

2. Go online: [www.binder-world.com/register](http://www.binder-world.com/register)

3. Register serial number

## 28. Contamination clearance certificate

### 28.1 For chambers located outside the USA and Canada

#### Declaration regarding safety and health

Erklärung zur Sicherheit und gesundheitlichen Unbedenklichkeit

The German Ordinance on Hazardous Substances (GefStofV), and the regulations regarding safety at the workplace, require that this form be filled out for all products that are returned to us, so that the safety and the health of our employees can be guaranteed.

Die Sicherheit und Gesundheit unserer Mitarbeiter, die Gefahrstoffverordnung GefStofV und die Vorschriften zur Sicherheit am Arbeitsplatz machen es erforderlich, dass dieses Formblatt für alle Produkte, die an uns zurückgeschickt wird.



Note: A repair is not possible without a completely filled out form.  
Ohne Vorliegen des vollständig ausgefüllten Formblattes ist eine Reparatur nicht möglich.

- A completely filled out form must be transmitted via Fax (+49 (0) 7462 2005 93555) or by letter in advance, so that this information is available before the equipment/component part arrives. A second copy of this form must accompany the equipment/component part. In addition, the carrier should be informed.

Eine vollständig ausgefüllte Kopie dieses Formblattes soll per Telefax (Nr. +49 (0) 7462 2005 93555) oder Brief vorab an uns gesandt werden, so dass die Information vorliegt, bevor das Gerät/Bauteil eintrifft. Eine weitere Kopie soll dem Gerät/Bauteil beigelegt sein. Ggf. ist auch die Spedition zu informieren.

- Incomplete information or non-conformity with this procedure will inevitably lead to substantial delays in processing. Please understand the reason for this measure, which lies outside our area of influence and will help us to speed up this procedure.

Unvollständige Angaben oder Nichteinhalten dieses Ablaufs führen zwangsläufig zu beträchtlichen Verzögerungen in der Abwicklung. Bitte haben Sie Verständnis für Maßnahmen, die außerhalb unserer Einflussmöglichkeiten liegen und helfen Sie mit, den Ablauf beschleunigen.

- **Please print and fill out this form completely.**

Bitte unbedingt vollständig ausfüllen!

<b>1.</b>	<b>Chamber/ component part / type: / Gerät / Bauteil / Typ:</b>
<b>2.</b>	<b>Serial No./ Serien-Nr.:</b>
<b>3.</b>	<b>Details about utilized substances / biological substances / Einzelheiten über die eingesetzten Substanzen/biologische Materialien:</b>
<b>3.1</b>	<b>Designations / Bezeichnungen:</b>
a)	_____
b)	_____
c)	_____
<b>3.2</b>	<b>Safety measures required for handling these substances / Vorsichtsmaßnahmen beim Umgang mit diesen Stoffen:</b>
a)	_____
b)	_____
c)	_____



<b>3.3</b>	<b>Measures to be taken in case of skin contact or release into the atmosphere / Maßnahmen bei Personenkontakt oder Freisetzung:</b>
a)	_____
b)	_____
c)	_____
d)	_____
<b>3.4</b>	<b>Other important information that must be taken into account / Weitere zu beachtende und wichtige Informationen:</b>
a)	_____
b)	_____
c)	_____
<b>4.</b>	<b>Declaration on the risk of these substances (please checkmark the applicable items) / Erklärung zur Gefährlichkeit der Stoffe (bitte Zutreffendes ankreuzen) :</b>
<input type="checkbox"/> <b>4.1 For non-toxic, non-radioactive, biologically harmless materials / für nicht giftige, nicht radioaktive, biologisch ungefährliche Stoffe:</b>	
<b>We hereby guarantee that the above-mentioned chamber / component part... / Wir versichern, dass o.g. Gerät/Bauteil...</b>	
<input type="checkbox"/> Has not been exposed to or contains any toxic or otherwise hazardous substances / weder giftige noch sonstige gefährliche Stoffe enthält oder solche anhaften.	
<input type="checkbox"/> That eventually generated reaction products are non-toxic and also do not represent a hazard / auch evtl. entstandene Reaktionsprodukte weder giftig sind noch sonst eine Gefährdung darstellen.	
<input type="checkbox"/> Eventual residues of hazardous substances have been removed / evtl. Rückstände von Gefahrstoffen entfernt wurden.	
<input type="checkbox"/> <b>4.2 For toxic, radioactive, biologically harmful or hazardous substances, or any other hazardous materials / für giftige, radioaktive, biologisch bedenkliche bzw. gefährliche Stoffe oder anderweitig gefährliche Stoffe.</b>	
<b>We hereby guarantee that ... / Wir versichern, dass ...</b>	
<input type="checkbox"/> The hazardous substances, which have come into contact with the above-mentioned equipment/component part, have been completely listed under item 3.1 and that all information in this regard is complete / die gefährlichen Stoffe, die mit dem o.g. Gerät/Bauteil in Kontakt kamen, in 3.1 aufgelistet sind und alle Angaben vollständig sind.	
<input type="checkbox"/> That the chamber /component part has not been in contact with radioactivity / das Gerät/Bauteil nicht mit Radioaktivität in Berührung kam	
<b>5.</b>	<b>Kind of transport / transporter / Transportweg/Spediteur:</b>
Transport by (means and name of transport company, etc.) Versendung durch (Name Spediteur o.ä.)	
_____	
Date of dispatch to BINDER GmbH / Tag der Absendung an BINDER GmbH:	
_____	

**We hereby declare that the following measures have been taken / Wir erklären, dass folgende Maßnahmen getroffen wurden:**

- ☐ Hazardous substances were removed from the chamber including component parts, so that no hazard exists for any person in the handling or repair of these items / das Gerät/Bauteil wurde von Gefahrstoffen befreit, so dass bei Handhabung/Reparaturen für die betreffende Person keinerlei Gefährdung besteht
- ☐ The chamber was securely packaged and properly identified / das Gerät wurde sicher verpackt und vollständig gekennzeichnet.
- ☐ Information about the hazardousness of the shipment (if required) has been provided to the transporter / der Spediteur wurde (falls vorgeschrieben) über die Gefährlichkeit der Sendung informiert.

We hereby commit ourselves and guarantee that we will indemnify BINDER GmbH for all damages that are a consequence of incomplete or incorrect information provided by us, and that we will exempt BINDER GmbH from eventual damage claims by third parties./ Wir versichern, dass wir gegenüber BINDER für jeden Schaden, der durch unvollständige und unrichtige Angaben entsteht, haften und BINDER gegen eventuell entstehende Schadenansprüche Dritter freistellen.

We are aware that, in accordance with Article 823 of the German Civil Code (BGB), we are directly liable with regard to third parties, in this instance especially the employees of BINDER GmbH, who have been entrusted with the handling / repair of the chamber / component. / Es ist uns bekannt, dass wir gegenüber Dritten – hier insbesondere mit der Handhabung/Reparatur des Geräts/des Bauteils betraute Mitarbeiter der Firma BINDER - gemäß §823 BGB direkt haften

Name: \_\_\_\_\_

Position/Title: \_\_\_\_\_

Date / Datum: \_\_\_\_\_

Signature / Unterschrift: \_\_\_\_\_

Company stamp / Firmenstempel:



Equipment that is returned to the factory for repair must be accompanied by a completely filled out contamination clearance certificate. For service and maintenance on site, such a contamination clearance certificate must be submitted to the service technician before the start of any work. No repair or maintenance of the equipment is possible, without a properly filled out contamination clearance certificate.

## 28.2 For chambers in the USA and Canada

### Product Return Authorization Request

Please complete this form and the Customer Decontamination Declaration (next 2 pages) and attach the required pictures. E-mail to: IDL\_SalesOrderProcessing\_USA@binder-world.com

After we have received and reviewed the complete information we will decide on the issue of a RMA number. Please be aware that size specifications, voltage specifications as well as performance specifications are available on the internet at [www.binder-world.us](http://www.binder-world.us) at any time.

Take notice of shipping laws and regulations.

	Please fill:	
Reason for return request	<input type="radio"/> Duplicate order	
	<input type="radio"/> Duplicate shipment	
	<input type="radio"/> Demo	<i>Page one completed by sales</i>
	<input type="radio"/> Power Plug / Voltage	115V / 230 V / 208 V / 240V
	<input type="radio"/> Size does not fit space	
	<input type="radio"/> Transport Damage	Shock watch tripped? ( <i>pictures</i> )
	<input type="radio"/> Other (specify below)	
	_____	
Is there a replacement PO?	<input type="radio"/> Yes <input type="radio"/> No	
<i>If yes -&gt; PO #</i>		
<i>If yes -&gt; Date PO placed</i>		
Purchase order number		
BINDER model number		
BINDER serial number		
Date chamber was received		
Was the chamber unboxed?	<input type="radio"/> Yes <input type="radio"/> No	
Was the chamber plugged in?	<input type="radio"/> Yes <input type="radio"/> No	
Was the chamber in operation?	<input type="radio"/> Yes <input type="radio"/> No	
<i>Pictures of chamber attached?</i>	<input type="radio"/> Yes <input type="radio"/> No	Pictures have to be attached!
<i>Pictures of Packaging attached?</i>	<input type="radio"/> Yes <input type="radio"/> No	

	Customer Contact Information	Distributor Contact Information
Name		
Company		
Address		
Phone		
E-mail		

**Customer (End User) Decontamination Declaration****Health and Hazard Safety declaration**

To protect the health of our employees and the safety at the workplace, we require that this form is completed by the user for all products and parts that are returned to us. (Distributors or Service Organizations cannot sign this form)



NO RMA number will be issued without a completed form. Products or parts returned to our NY warehouse without a RMA number will be refused at the dock.

A second copy of the completed form must be attached to the outside of the shipping box.

<b>1. Chamber/ component part / type:</b>
<b>2. Serial No.</b>
<b>3. List any exposure to hazardous liquids, gasses or substances and radioactive material</b>
<b>3.1 List with MSDS sheets attached where available or needed (if there is not enough space available below, please attach a page):</b>
a) _____
b) _____
c) _____
<b>3.2 Safety measures required for handling the list under 3.1</b>
a) _____
b) _____
c) _____
<b>3.3 Measures to be taken in case of skin contact or release into the atmosphere:</b>
a) _____
b) _____
c) _____
d) _____
<b>3.4 Other important information that must be considered:</b>
a) _____
b) _____
c) _____

#### 4. Declaration of Decontamination

**For toxic, radioactive, biologically and chemically harmful or hazardous substances, or any other hazardous materials.**

**We hereby guarantee that**

- 4.1 Any hazardous substances, which have come into contact with the above-mentioned equipment / component part, have been completely listed under item 3.1 and that all information in this regard is complete.
- 4.2 That the chamber / component part has not been in contact with radioactivity
- 4.3 Any Hazardous substances were removed from the chamber / component part, so that no hazard exists for a persons in the shipping, handling or repair of these returned chamber
- 4.4 The chamber was securely packaged in the original undamaged packaging and properly identified on the outside of the packaging material with the chamber designation, the RMA number and a copy of this declaration.
- 4.5 Shipping laws and regulations have not been violated.

**I hereby commit and guarantee that we will indemnify BINDER Inc. for all damages that are a consequence of incomplete or incorrect information provided by us, and that we will indemnify and hold harmless BINDER Inc. from eventual damage claims by third parties.**

Name: \_\_\_\_\_

Position: \_\_\_\_\_

Company: \_\_\_\_\_

Address: \_\_\_\_\_

Phone #: \_\_\_\_\_

Email: \_\_\_\_\_

Date: \_\_\_\_\_

Signature: \_\_\_\_\_



Equipment returned to the NY warehouse for repair must be accompanied by a completed customer decontamination declaration. For service and maintenance works on site, such a customer decontamination declaration must be submitted to the service technician before the start of work. No repair or maintenance of the equipment is possible without a completed form.

