

# High Capacity Series for Storage or Transport (Round Canisters)



## Scope of Application

High Capacity Series for Storage or Transport (Round Canisters) provides two cryopreservation solutions for long-term static storage and transportation of biological samples.

## Innovative Design

- Numbered Index Canister Locations
- Round Canisters for Straw Storage
- Lockable
- LN<sub>2</sub> Dispenser
- High Thermal Efficiency
- Roller Bases Available
- High Capacity
- 5 Year Vacuum Warranty

# High Capacity Series for Storage or Transport (Round Canisters)



## Key Features

- Strong and lightweight aluminium construction
- Heavy duty lockable enclosure offers excellent security
- Numbered index location points for canisters
- Straw storage
- LN<sub>2</sub> dispense pump available
- Roller bases available
- High thermal efficiency
- 5 year vacuum warranty
- Ultra-low evaporation loss

## Product Advantages

Narrow neck design and excellent thermal conductivity reduces evaporation loss of liquid nitrogen to extend storage times.

Large-capacity models have wider necks for easy access to biological samples.

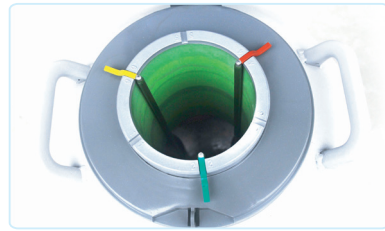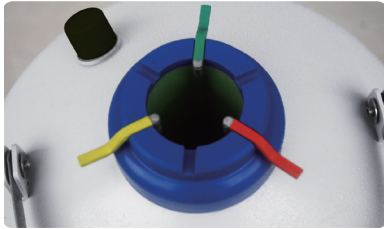

Long-term cryopreservation models have narrower necks with less heat intrusion.

## Technical Parameters

Model		YDS-2-30	YDS-2-35	YDS-3	YDS-6	YDS-10	YDS-10-80
Maximum Capacity	No. of Canister	3(3)	3(3)	6(0)	6(0)	6(0)	6(0)
	No. of Straws (Single Layer) (0.5ml each canister)	90	165	792	792	792	2244
	No. of Straws (Single Layer) (0.25ml each canister)	204	330	1788	1788	1788	5022
	No. of Straws (Double Layer) (0.5ml each canister)	-	-	-	-	-	-
	No. of Straws (Double Layer) (0.25ml each canister)	-	-	-	-	-	-
Performance	Volume of LN <sub>2</sub> (L)	2	2	3	6	10	10
	Static Evaporation* (L/Day)	0.07	0.08	0.12	0.12	0.12	0.21
	Static Holding Time** (Day)	28	24	26	52	86	48
Dimensions	Opening Diameter (mm)	30	35	50	50	50	80
	Height(mm)	415	435	445	495	565	567
	Outside Diameter (mm)	223	204	223	300	300	300
	Canister Outside Diameter (mm)	19	25	38	38	38	63
	Outside Height (mm)	120	120	120	120	120	120
	Empty Weight (kg)	2.8	2.6	3.1	4.8	6.1	6.2
	Full Weight (kg)	4.5	4.4	6	9.9	14.5	14.7
	Shipping Weight (kg)	5.1	4.9	5.8	8	9	9.3

# High Capacity Series for Storage or Transport (Round Canisters)

Model		YDS-10-125	YDS-13	YDS-15	YDS-15	YDS-30
Maximum Capacity	No. of Canister	1(0)	6(0)	6(0)	6(0)	6(0)
	No. of Straws (Single Layer) (0.5ml each canister)	854	-	792	792	792
	No. of Straws (Single Layer) (0.25ml each canister)	1940	-	1788	1788	1788
	No. of Straws (Double Layer) (0.5ml each canister)	-	1284	-	1284	1284
	No. of Straws (Double Layer) (0.25ml each canister)	-	2832	-	2832	2832
Performance	Volume of LN <sub>2</sub> (L)	10	13	15	20	30
	Static Evaporation* (L/Day)	0.42	0.12	0.11	0.12	0.12
	Static Holding Time** (Day)	24	109	134	168	254
Dimensions	Opening Diameter (mm)	125	50	50	50	50
	Height(mm)	625	635	610	700	725
	Outside Diameter (mm)	300	310	394	394	462
	Canister Outside Diameter (mm)	97	38	38	38	38
	Outside Height (mm)	120	276	120	120/276	120/276
	Empty Weight (kg)	6.3	6.3	8.5	9.5	12.9
	Full Weight (kg)	15.1	17.3	20.5	28.1	36.6
	Shipping Weight (kg)	8.6	11	13	15	18

Model		YDS-30-80	YDS-30-125	YDS-35	YDS-35-80	YDS-35-125	YDS-47-127
Maximum Capacity	No. of Canister	6(0)	6(0)	6(0)	6(0)	6(0)	10(0)
	No. of Straws (Single Layer) (0.5ml each canister)	2244	5124	792	2244	5124	-
	No. of Straws (Single Layer) (0.25ml each canister)	5022	11640	1788	5022	11640	-
	No. of Straws (Double Layer) (0.5ml each canister)	3624	-	1284	3624	9048	8200
	No. of Straws (Double Layer) (0.25ml each canister)	8460	-	2832	8460	20760	16400
Performance	Volume of LN <sub>2</sub> (L)	30	30	35	35	35	47
	Static Evaporation* (L/Day)	0.21	0.35	0.12	0.22	0.37	0.36
	Static Holding Time** (Day)	147	90	286	159	97	130
Dimensions	Opening Diameter (mm)	80	125	50	80	125	127
	Height(mm)	729	725	765	768	765	740
	Outside Diameter (mm)	462	462	462	462	462	508
	Canister Outside Diameter (mm)	63	97	38	63	97	72
	Outside Height (mm)	120/276	120	120/276	120/276	120/276	276
	Empty Weight (kg)	13	13	14.2	14.5	14.6	18.2
	Full Weight (kg)	37	38.9	42.7	42.9	43.2	56.2
	Shipping Weight (kg)	18.3	19.3	20	20.7	22	28

# High Capacity Series for Storage or Transport (Round Canisters)

Model		YDS-20B	YDS-30B	YDS-35B
Maximum Capacity	No. of Canister	6(0)	6(0)	6(0)
	No. of Straws (Single Layer) (0.5ml each canister)	792	792	792
	No. of Straws (Single Layer) (0.25ml each canister)	1788	1788	1788
	No. of Straws (Double Layer) (0.5ml each canister)	1284	1284	1284
	No. of Straws (Double Layer) (0.25ml each canister)	2832	2832	2832
Performance	Volume of LN <sub>2</sub> (L)	20	30	35
	Static Evaporation* (L/Day)	0.2	0.2	0.2
	Static Holding Time** (Day)	101	159	179
Dimensions	Opening Diameter (mm)	50	50	50
	Height(mm)	700	725	765
	Outside Diameter (mm)	394	462	462
	Canister Outside Diameter (mm)	38	38	38
	Outside Height (mm)	120/276	120/276	120/276
	Empty Weight (kg)	9.5	12.9	14.2
	Full Weight (kg)	27.9	37.1	42.8
	Shipping Weight (kg)	15.1	18.3	20.1

Model		YDS-35B-80	YDS-35B-125	YDS-50B	YDS-50B-125
Maximum Capacity	No. of Canister	6(0)	6(0)	6(0)	6(0)
	No. of Straws (Single Layer) (0.5ml each canister)	2244	5124	792	5124
	No. of Straws (Single Layer) (0.25ml each canister)	5022	11640	1788	11640
	No. of Straws (Double Layer) (0.5ml each canister)	3624	9048	1284	9048
	No. of Straws (Double Layer) (0.25ml each canister)	8460	20760	2832	20760
Performance	Volume of LN <sub>2</sub> (L)	35	35	50	50
	Static Evaporation* (L/Day)	0.3	0.41	0.24	0.45
	Duration of Normal Operation** (Day)	119	86	213	110
Dimensions	Opening Diameter (mm)	80	125	50	125
	Height(mm)	768	765	825	830
	Outside Diameter (mm)	462	462	462	462
	Canister Outside Diameter (mm)	63	97	38	97
	Outside Height (mm)	120/276	120/276	120/276	120/276
	Empty Weight (kg)	14.5	14.6	17.2	17.3
	Full Weight (kg)	43	43.5	60.3	62
	Shipping Weight (kg)	20.7	22	24	25.4

A(B): A for maximum number of canisters, B for number of canisters in standard configuration

\*Static Evaporation: The amount of liquid nitrogen lost by natural evaporation within 24 hours after the container is allowed to reach thermal equilibrium under the specified filling rate, which is converted into the evaporation value under the standard state (atmospheric pressure 101.325kPa, ambient temperature 20°C).

\*\*Static Holding Time: under the standard state (atmospheric pressure 101.325kPa, ambient temperature 20°C), the container is filled with liquid nitrogen pre-cooled and reaches thermal equilibrium, and the liquid nitrogen filling rate reaches 100%, the number of days when the liquid nitrogen is completely evaporated on standing.

# High Capacity Series for Storage or Transport (Round Canisters)

## Vials Cane Specification

Canister Height 120mm, Diameter 38mm (50mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	4	3	12
1.5ml	4	3	12
2ml	4	3	12
3ml	4	3	12
5ml	4	1	4

Canister Height 276mm, Diameter 38mm (50mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	4	6	24
1.5ml	4	6	24
2ml	4	6	24
3ml	4	6	24
5ml	4	3	12

Canister Height 120mm, Diameter 63mm (80mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	16	3	48
1.5ml	16	3	48
2ml	16	3	48
3ml	16	3	48
5ml	16	1	16

Canister Height 276mm, Diameter 63mm (80mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	16	6	96
1.5ml	16	6	96
2ml	16	6	96
3ml	16	6	96
5ml	16	3	48

Canister Height 120mm, Diameter 72mm (127mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	21	3	63
1.5ml	21	3	63
2ml	21	3	63
3ml	21	3	63
5ml	21	1	21

Canister Height 276mm, Diameter 72mm (127mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	21	6	126
1.5ml	21	6	126
2ml	21	6	126
3ml	21	6	126
5ml	21	3	63

Canister Height 120mm, Diameter 97mm (125mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	-	-	-
1.5ml	40	3	120
2ml	40	3	120
3ml	40	3	120
5ml	40	1	40

Canister Height 276mm, Diameter 97mm (125mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	-	-	-
1.5ml	40	6	240
2ml	40	6	240
3ml	40	6	240
5ml	40	3	120


Canister Height 120mm, Diameter 104mm (127mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	-	-	-
1.5ml	46	3	138
2ml	46	3	138
3ml	46	3	138
5ml	46	1	46

Canister Height 276mm, Diameter 104mm (127mm neck)			
Vials	Canes/Canister	Vials/Cane	Vials/Canister
0.5ml	-	-	-
1.5ml	46	6	276
2ml	46	6	276
3ml	46	6	276
5ml	46	3	138

## Accessories



Vial

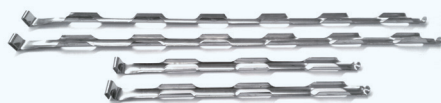

Vial cane



Canister

# High Capacity Series for Storage or Transport (Round Canisters)

## Product List

Model	Product	Product Model	Description and Accessories
YDS-2-30	Liquid Nitrogen container	YDS-2-30	Includes 3 canisters, lockable cap and PU bag
	Canister	TT2-30	φ19, 120mm single tier canister, No goblet, suitable for YDS-2-30
	Cork (30)	GS-30-125	Suitable for YDS-2-30
	Lockable cap	SG-30	Suitable for container with 30mm neck opening, suitable for YDS-2-30
YDS-2-35	Liquid Nitrogen container	YDS-2-35	Includes 3 canisters, lockable cap and Braces for carrying, no PU bag
	Canister	TT2-35	φ25, 120mm single tier canister, No goblet, suitable for YDS-2-35
	Cork (30)	GS-35-125 (L)	Suitable for YDS-2-35
	Lockable cap	SG-35	Suitable for container with 35mm neck opening, suitable for YDS-2-35
YDS-3	Liquid Nitrogen container	YDS-3	Includes 6 canisters (Height 120mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-3	Includes PU bag only
	Canister	TT3-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-3
	Cork (50)	GS-50-180	Suitable for YDS-3/YDH-3/YDS-6/YDS-10/YDS-13/YDS-25/YDS-50B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening
YDS-6	Liquid Nitrogen container	YDS-6	Includes 6 canisters (Height 120mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-6	Includes PU bag only
	Canister	TT6-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-6
	Cork (50)	GS-50-180	Suitable for YDS-3/YDH-3/YDS-6/YDS-10/YDS-13/YDS-25/YDS-50B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening
YDS-10	Liquid Nitrogen container	YDS-10	Includes 6 canisters (Height 120mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-10	Includes PU bag only
	Canister	TT10-50	φ38, 120mm single tier canister, No goblet, suitable for YDS-10
	Cork (50)	GS-50-180	Suitable for YDS-3/YDH-3/YDS-6/YDS-10/YDS-13/YDS-25/YDS-50B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening
YDS-10-80	Liquid Nitrogen container	YDS-10-80	Includes 6 canisters (Height 120mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-10-80	Includes PU bag only
	Canister	TT10-80	φ63, 120mm single tier canister, No goblet, suitable for YDS-10-80
	Cork (80)	GS-80-180	Suitable for YDH-6-80/YDS-10-80
	Lockable cap	SG-80	Suitable for container with 80mm neck opening
YDS-13	Liquid Nitrogen container	YDS-13	Includes 6 canisters (Height 276mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-13	Includes PU bag only
	Canister	TTS13-50	φ38, 276mm double-tier canister, 2 goblets included. Suitable for YDS-13
	Cork (50)	GS-50-180	Suitable for YDS-3/YDH-3/YDS-6/YDS-10/YDS-13/YDS-25/YDS-50B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening
YDS-15	Liquid Nitrogen container	YDS-15	Includes 6 canisters (Height 120mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-15	Includes PU bag only
	Canister	TT15-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-15
	Cork (50)	GS-50-210	Suitable for YDS-15/YDS-20/YDS-30/YDS-35/YDS-20B/YDS-30B/YDS-35B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening

# High Capacity Series for Storage or Transport (Round Canisters)

## Product List

Model	Product	Product Model	Description and Accessories
YDS-20	Liquid Nitrogen container	YDS-20	Includes 6 canisters (Height 276mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-20	Includes 6 canisters (Height 120mm), lockable cap and PU bag
	Liquid Nitrogen container	YDS-20	Includes PU bag only
	Canister	TT30-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-20/YDS-30
	Canister	TTS30-50	φ38,276mm double-tier canister, 2 goblets included. Suitable for YDS-20/YDS-30
	Cork (50)	GS-50-210	Suitable for YDS-15/YDS-20/YDS-30/YDS-35/YDS-20B/YDS-30B/YDS-35B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening
YDS-30	Liquid Nitrogen container	YDS-30	Includes 6 canisters (Height 276mm). lockable cap and PU bag. roller base is optional
	Liquid Nitrogen container	YDS-30	Includes 6 canisters(Height 120mm), lockable cap and PU bag, roller base is optional
	Liquid Nitrogen container	YDS-30	Includes PU bag only, roller base is optional
	Canister	TT30-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-20/YDS-30
	Canister	TTS30-50	φ38, 276mm double-tier canister, 2 goblets included. Suitable for YDS-20/YDS-30
	Roller Base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank. equipped with 5 universal wheels.
	Smart Cap	LT-50	Used for monitoring real time level and temperature for containers with 50mm neck opening, 6 grooves and 540mm height. suitable for YDS-20/YDS-30
	Cork (50)	GS-50-210	Suitable for YDS-15/YDS-20/YDS-30/YDS-35/YDS-20B/YDS-30B/YDS-35B
YDS-30-80	Liquid Nitrogen container	YDS-30-80	Includes 6 canisters (Height 276mm), lockable cap and PU bag. roller base is optional
	Liquid Nitrogen container	YDS-30-80	Includes 6 canisters(Height 120mm), lockable cap and PU bag, roller base is optional
	Liquid Nitrogen container	YDS-30-80	Includes PU bag only, roller base is optional
	Canister	TT30-80	φ63, 120mm single tier canister, No goblet. suitable for YDS-30-80
	Canister	TTS30-80	φ63,276mm double-tier canister, 2 goblets included. Suitable for YDS-30-80
	Roller Base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank. equipped with 5 universal wheels.
	Smart Cap	LT-80	Used for monitoring real time level and temperature for containers with 80mm neck opening, 6 grooves and 530mm height. suitable for YDS-20-80/YDS-30-80
	Cork (80)	GS-80-210	Suitable for YDS-30-80/YDS-35-80/YDS-35B-80
YDS-35	Liquid Nitrogen container	YDS-35	Includes 6 canisters (Height 276mm), lockable cap and PU bag, roller base is optional
	Liquid Nitrogen container	YDS-35	Includes 6 canisters(Height 120mm), lockable cap and PU bag, roller base is optional
	Liquid Nitrogen container	YDS-35	Includes PU bag only, roller base is optional
	Canister	TT35-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-25/YDS-35
	Canister	TTS35-50	φ38, 276mm double-tier canister, 2 goblets included. Suitable for YDS-35
	Roller Base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank. equipped with 5 universal wheels.
	Roller Base	LT-50	Used for monitoring real time level and temperature for containers with 50mm neck opening, 6 grooves and 570mm height. suitable for YDS-25/YDS-35
	Cork (50)	GS-50-210	Suitable for YDS-15/YDS-20/YDS-30/YDS-35/YDS-20B/YDS-30B/YDS-35B
YDS-35-80	Liquid Nitrogen container	YDS-35-80	Includes 6 canisters (Height 276mm). lockable cap and PU bag. roller base is optional
	Liquid Nitrogen container	YDS-35-80	Includes PU bag only, roller base is optional
	Canister	TT35-80	φ63, 120mm single tier canister, No goblet. suitable for YDS-35-80
	Canister	TTS35-80	φ63, 276mm double-tier canister, 2 goblets included. Suitable for YDS-35-80
	Roller Base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank. equipped with 5 universal wheels.
	Roller Base	LT-80	Used for monitoring real time level and temperature for containers with 80mm neck opening, 6 grooves and 580mm height. suitable for YDS-25-80/YDS-35-80
	Cork (80)	GS-80-210	Suitable for YDS-30-80/YDS-35-80/YDS-35B-80
	Lockable cap	SG-80	Suitable for container with 80mm neck opening

# High Capacity Series for Storage or Transport (Round Canisters)

## Product List

Model	Product	Product Model	Description and Accessories
YDS-20B	Liquid nitrogen container	YDS-20B	Includes PU bag only.
	Canister	TT30-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-20/YDS-30
	Canister	TTS30-50	φ38, 276mm double-tier canister, 2 goblets included. Suitable for YDS-20/YDS-30
	Cork (50)	GS-50-210	Suitable for YDS-15/YDS-20/YDS-30/YDS-35/YDS-20B/YDS-30B/YDS-35B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening
YDS-30B	Liquid nitrogen container	YDS-30B	Includes PU bag only, roller base is optional
	Canister	TT30-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-20/YDS-30
	Canister	TTS30-50	φ38, 276mm double-tier canister, 2 goblets included. Suitable for YDS-20/YDS-30
	Roller base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank, equipped with 5 universal wheels.
	Cork (50)	GS-50-210	Suitable for YDS-15/YDS-20/YDS-30/YDS-35/YDS-20B/YDS-30B/YDS-35B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening
YDS-35B	Liquid nitrogen container	YDS-35B	Includes PU bag only, roller base is optional
	Canister	TT35-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-25/YDS-35
	Canister	TTS35-50	φ38,276mm double-tier canister, 2 goblets included. Suitable for YDS-35
	Roller base	YSC-465-5W	Suitable for tanks with diameter less than φ465 Can be used for 30L, 35L and 50L tank, equipped with 5 universal wheels.
	Cork (50)	GS-50-210	Suitable for YDS-15/YDS-20/YDS-30/YDS-35/YDS-20B/YDS-30B/YDS-35B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening

# High Capacity Series for Storage or Transport (Round Canisters)

## Product List

Model	Product	Product Model	Description and Accessories
YDS-35B-80	Liquid Nitrogen container	YDS-35B-80	Includes PU bag only, roller base is optional
	Canister	TT35-80	φ63, 120mm single tier canister. No goblet, suitable for YDS-35-80
	Canister	TTS35-80	φ63, 276mm double-tier canister, 2 goblets included. Suitable for YDS-35-80
	Roller Base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank, equipped with 5 universal wheels.
	Cork (80)	GS-80-210	Suitable for YDS-30-80/YDS-35-80/YDS-35B-80
	Lockable cap	SG-80	Suitable for container with 80mm neck opening
YDS-35B-125	Liquid Nitrogen container	YDS-35B-125	Includes PU bag only
	Canister	TT35-125	φ97, 120mm single tier canister, No goblet, suitable for YDS-35-125
	Canister	TTS35-125	φ97,276mm double-tier canister, 2 goblets included. Suitable for YDS-35-125
	Rack	FTT-25-5H	Used for storing 5*5 cryo boxes(2ml), 5 layers in total, with 5*5 cryo boxes included, suitable for YDS-35-125
	Roller Base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank, equipped with 5 universal wheels.
	Cork (125)	GS-125-210	Suitable for YDS-30-125/YDS-35-125/YDS-47-127 /YDS-35B-125
YDS-50B	Liquid Nitrogen container	YDS-50B	Includes PU bag only, roller base is optional
	Canister	TT50-50	φ38, 120mm single tier canister, No goblet. suitable for YDS-50
	Canister	TTS50-50	φ38, 276mm double-tier canister, 2 goblets included. Suitable for YDS-50
	Roller Base	YSC-465-5W	Suitable for tanks with diameter less than φ465. Can be used for 30L, 35L and 50L tank, equipped with 5 universal wheels.
	Smart cap	LT-50-500-6H640	Used for monitoring real time level and temperature for containers with 50mm neck opening, 6 grooves and 640mm height. suitable for YDS-50
	Cork (50)	GS-50-180	Suitable for YDS-3/YDH-3/YDS-6/YDS-10/YDS-13/YDS-25/YDS-50B
	Lockable cap	SG-50	Suitable for container with 50mm neck opening

\*Haier Biomedical reserves the right to change products and specifications without prior notice.



# WolfLabs

**Pricing on any accessories shown can be found by keying the part number into the search box on our website.**

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

**[www.wolflabs.co.uk](http://www.wolflabs.co.uk)**

**Tel : 01759 301142**

**Fax : 01759 301143**

**[sales@wolflabs.co.uk](mailto:sales@wolflabs.co.uk)**

Please contact us if this literature doesn't answer all your questions.